Multimedia cameras - EDUCAM USB 64 BIT VERSION

INSTRUCTION MANUAL

Model
4083.4

Version: 1
Issued: 18, 08, 2014

Description

The **EDUCAM MULTIMEDIA USB** video camera is especially designed to meet the various requirements in the educational field.

It can be connected to a professional monitor, a TV set, or a PC through USB link. It can be used in the following ways:


- as an episcopo, for the reproduction of texts, documents, photographs, etc.
- as an enlarger, to enlarge small objects, insects, minerals, etc.
- for video-microscopy, connectable both to biological or stereo microscopes using the adapters included.
- as a overhead projector, for the projection of tracings
- as a camera for teleconferences, that can be connected to a computer and to the Internet
- as a camera for video-conferences, used in conferences, assemblies, meetings, etc.
- as a camera for filming, with the help of a video recorder.

This video camera enables you to record clearly even when there is little light available, thanks to its extremely sensitive electronic circuits. The special lens enables you to focus from 0,76 mm, up to an infinite distance.

At the camera base, there is an extremely sensitive microphone that, via the TV itself, or via a separate amplifying system, enables you to reproduce for example the teacher's voice during the lesson, or to record sounds and noises from the surrounding area.

The special flexible arm and the heavy weight of the base, gives the EDUCAM its versatile, sturdy and stabile characteristics. The micro-video camera placed on the end of the flexible arm can be placed in any position, even projecting from the base, without jeopardising the system's stability. A special joint allows the head to rotate avoiding connector breakages inside the flexible arm.

Overview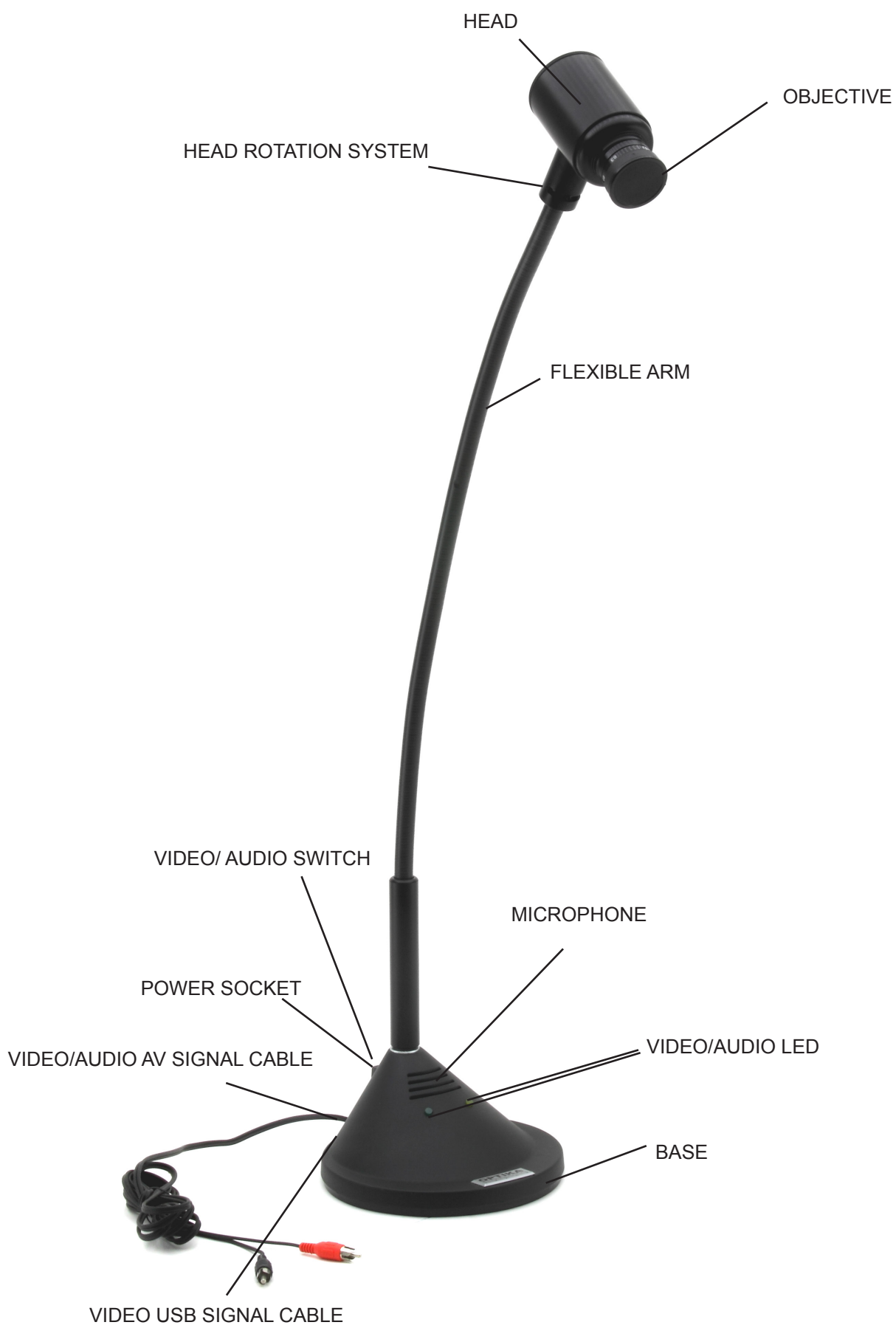

Product specifications

Sensor	1/3" interline transfer colour CCD element
Horizontal resolution	420 TV lines
Digital resolution	640x480
Total resolution	307k pixel
Objective	8 mm, F:1.2
Sensitivity	min. 0,08 lux at F:1.2
Picture controls	electronic shutter
Gain control	automatic
White balance	automatic
Video signal	PAL compatible - VHS
Working distance	from 0,76 mm to infinite
Magnifications	90x at 0,76 mm (with 21" monitor)
Microphone	built-in high sensitivity microphone
Audio signal	analogic
Power supply	15 V DC, 400 mA, with external power adapter
Switching on	via 3-position switch
PC link	standard USB 2.0
Length of flexible arm	65cm
Base diameter	17 cm
Weight	3.4 Kg

Assembly

Power supply

The Educam Multimedia video camera requires 15 V DC working voltage and comes with an external power pack. This power pack needs to be plugged directly into a 220 V mains sockets and its output cable should then be plugged directly into the video camera power supply socket.

Video and audio signals

The analog video signal from the Educam video camera is transmitted to a TV monitor and the audio signal to an amplifier or an amplified monitor. The small video/audio output signal cable, that can be found on the back of the camera base, should be connected as follows:

Red plug (video signal):

into the video input of your monitor/colour TV directly or via the SCART adapter provided (in the yellow video socket).

Black plug (audio signal):

into the audio input of your amplifier system or into the audio input of your monitor/colour TV or via the SCART adapter provided (in the white audio mono socket).

Using the SCART adapter, it has to be inserted into the SCART socket at the back of the TV set. See the TV manufacturer's instruction manual for details of how to switch on the TV and use its SCART circuit.

Connect the **USB** cable that can be found on the back of the camera base, to the PC in order to view the digital **USB** video signal (follow the instructions on next section to install the camera on your PC).

Switching on the video camera

The Educam video camera has a 3-position switch at the rear of its base. Each position corresponds to a specific function:

Central position "0" -> TV camera off.

Position "I" -> the yellow LED is lit and the Educam is operating with video signal only.

Position "II" -> both the yellow and green LEDs are lit and the Educam is operating with both video and audio signals.

Operation

Using the video camera

Position the camera head with the lens and the electronic sensor directly in front of the object that you wish to observe using its flexible arm. The lens provided with the camera has a manual focusing and must therefore be adjusted every time the distance between the object and the camera is changed. This can be done by simply turning the lens on the camera head. The object must be at least 0.76 mm away from the lens. There are no limits to the maximum distance. To use the microphone to capture audio, put the camera base with the microscope close to the audio source. It could happen that you hear a whistling noise while using the microphone. This whistle is not a defect of the Educam camera, but a normal audio effect ("Larsen effect") that occurs when the microscope is too close to the loudspeakers and the microphone is receiving the audio from the loudspeakers. The intensity also depends on the dimension and size of the room.


The best solution to avoid this effect is to place the Educam video camera far from the loudspeakers, to use a separated amplifier system and place it in a room without reverb. The "Larsen effect" will decrease in a room with many people and many pieces of furniture.

Capturing video from a microscope


The Educam video camera has been designed for effective use with microscopes.

The following special adapters (provided) are required for this application:

Adapter for biological microscopes: To be used directly in the optical tube of the microscope after having extracted the eyepiece. This adapter contains a WF10x eyepiece that needs to be replaced with the one belonging to the microscope.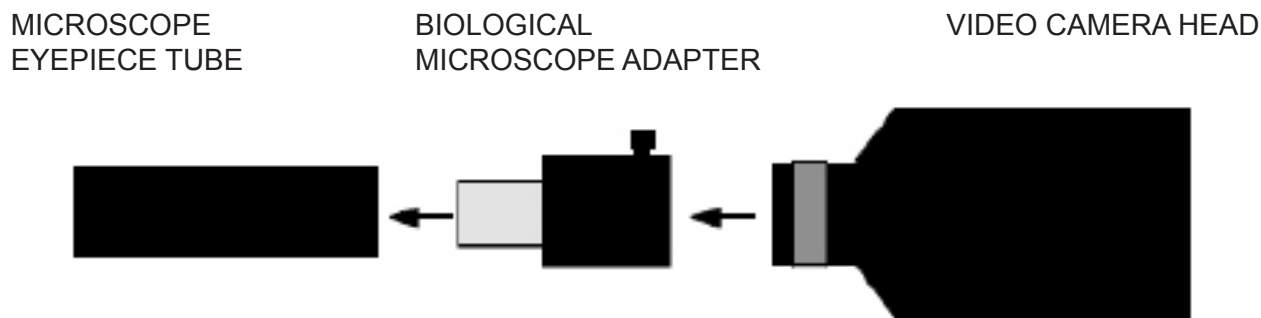

Adapter for stereo microscopes: To be used directly on one of the optical cylinders on a stereo microscope after having extracted the eyepiece. The biological microscope adapter needs to be inserted inside this adapter.

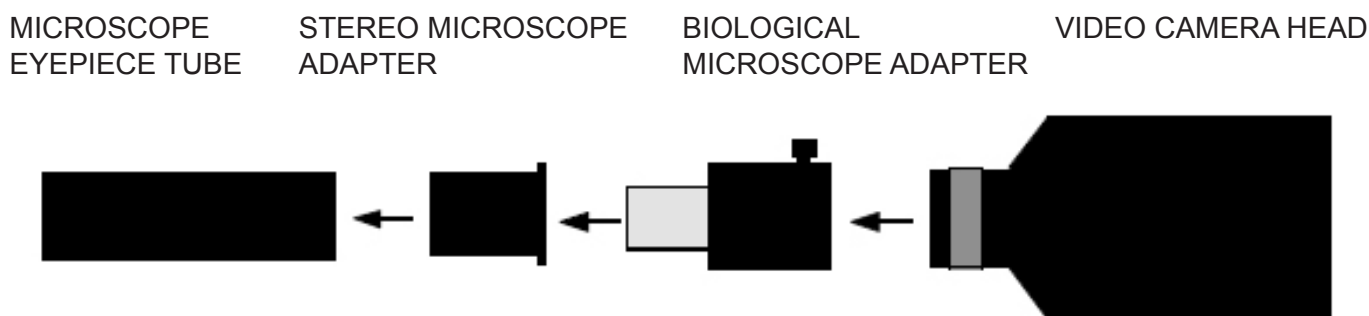Installation of the camera's drivers and pc

Drivers and Software Installation

Before connecting the camera to the PC, insert the CD for installation, choose the folder \ driver \ and run setup.exe.

When the Driver installation is completed choose the folder \ powerdirector8 \ and run setup.exe, wait for the software installation. The software's serial number can be found in the file CD \ SN.txt.

Connect the power supply and turn on the camera. Connect the USB cable to the PC. The camera will be recognized and the drivers installed.


Image and Video Recording:

Start the software by clicking on the icon:


Select **"Power director"**.


Select **"Capture"**.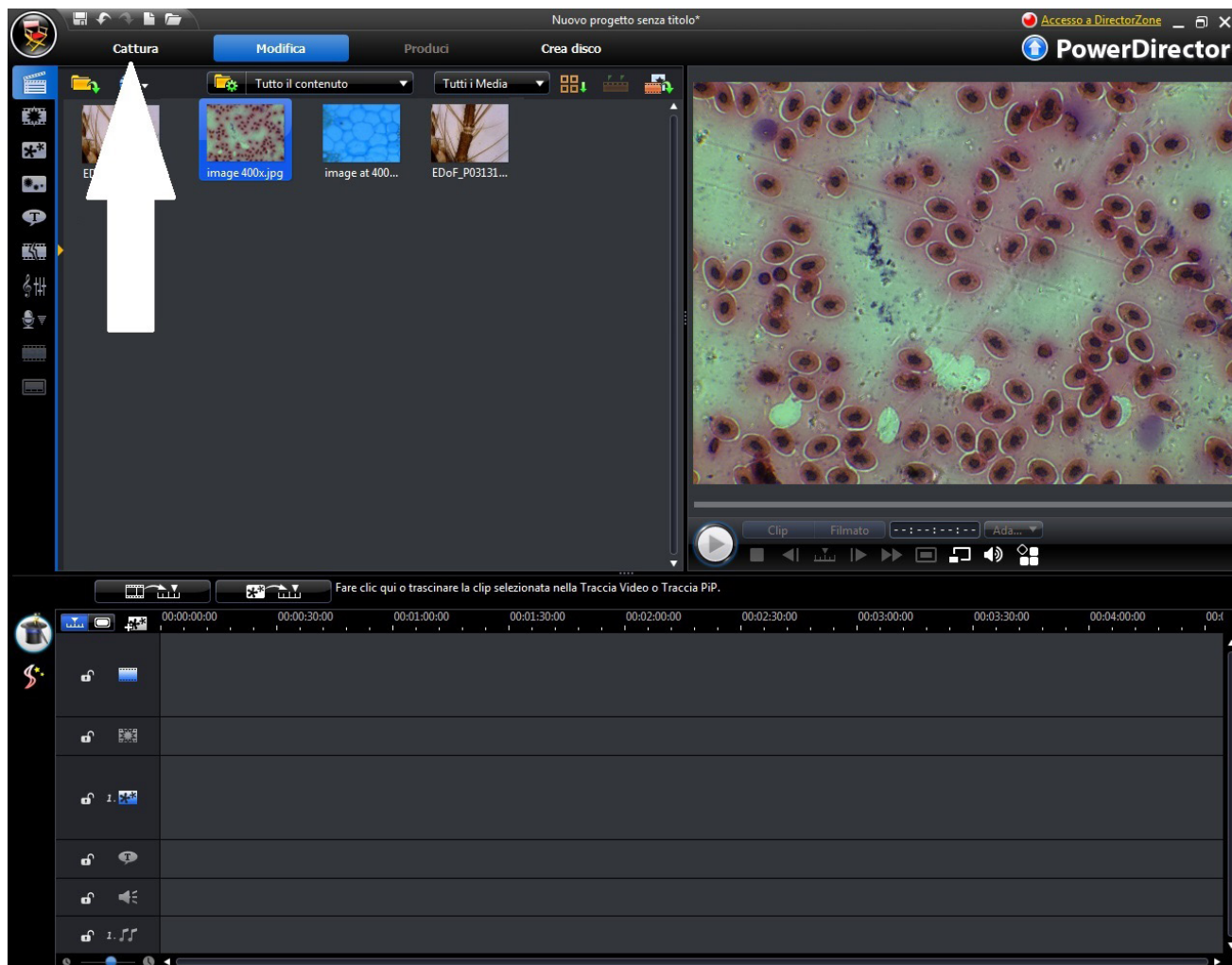

In order to adjust the image properties (colour, resolution) click on “Settings”.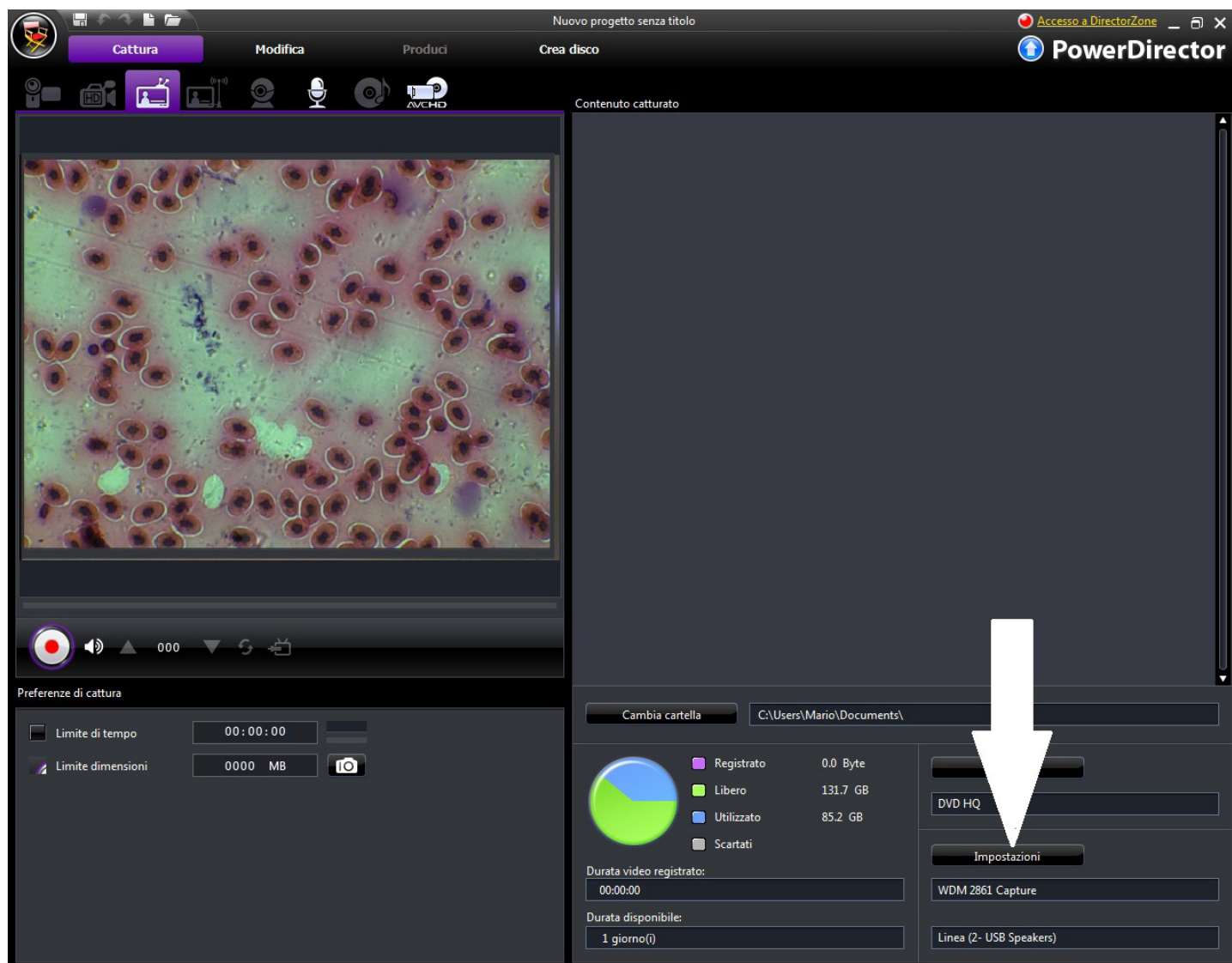

Modify the colour and the properties according to your requirements, then click on OK.

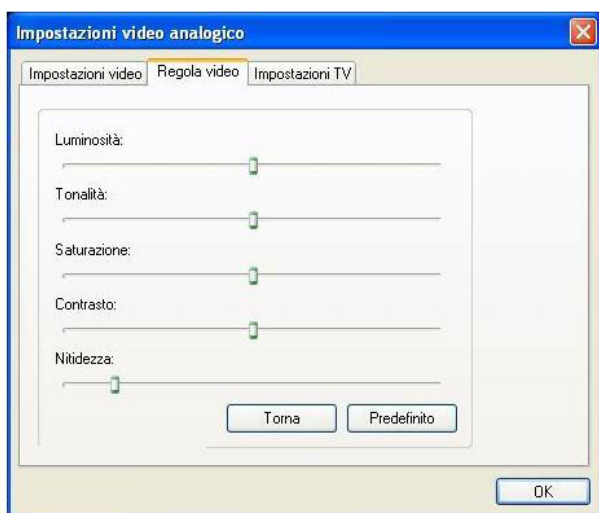Use the camera with Optika Vision Lite

Optika Vision Lite allows you to have a full screen view, to save images step by step, to save video, and to make line measurements on the captured image.

1. Please insert your CD and run \Optika Vision Lite 2.11\Software\setup.exe
2. Now Optika Vision Lite icon appears on your desktop.
3. Please install the additional Driver from the folder \Optika Vision Lite 2.11\Additional Driver\
4. Now you can run Optika Vision Lite and select the camera Driver USB2861Device from the list that appears when you select the microscope-shaped icon.

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection.

The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment.

Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.