B-510 Series

INSTRUCTION MANUAL

Model
B-510-2F
B-510-2
B-510-3
B-510-5

Version: 1

Issued: 14, 11, 2018

Table of Contents

1. **Warning**
 2. **Symbols and conventions**
 3. **Safety Information**
 4. **Intended use**
 5. **Overview**
 6. **Unpacking**
 7. **Assembling**
 8. **Use of the microscope**
 9. **Maintenance**
 10. **Troubleshooting**
- Equipment disposal**

1. Warning

This instruction manual is for the assembling, troubleshooting and maintenance of the multi-head discussion systems B-510-2F / B-510-2 / B-510-3 / B-510-5. The assembly sequence is identical for all systems.

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.

2. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution

ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock

3. Safety Information

Avoiding Electrical Shock


Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

4. Intended use


For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

5. Overview

5.1 B-510-2F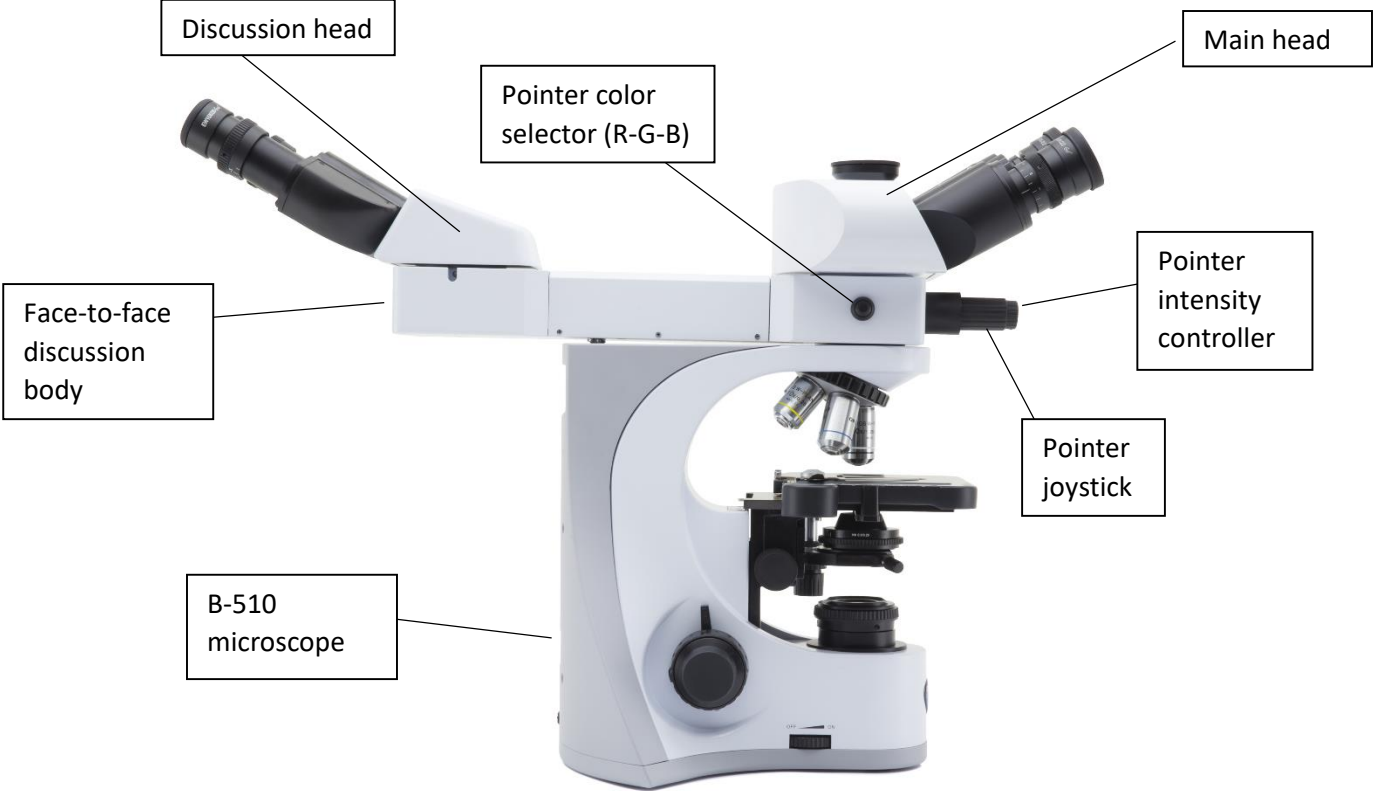

5.2 B-510-2 / B-510-3 / B-510-5

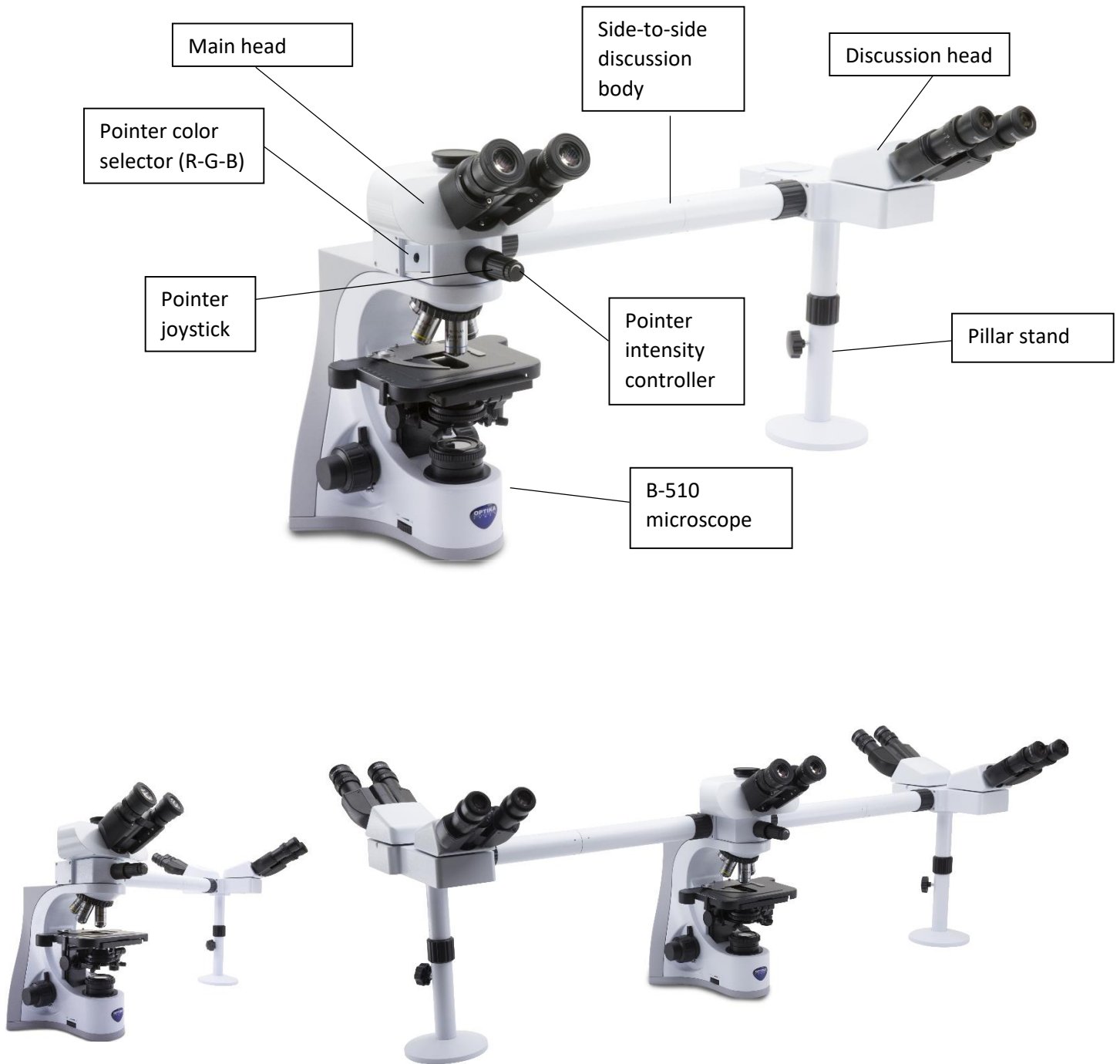6. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

7. Assembling

1. Install the microscope body as indicated in the specific instruction manual.
 - **Do not install the observation head.**
2. Place the splitter attachment of the multidiscussion system and tighten the lock screw on the right side. (Fig. 1).


3. In order to use the arrow which will be used as a pointer, connect the 5Vdc power supply to the rear socket of the splitter attachment. (Fig. 2). Now use the pointer joystick ① or the pointer intensity controller ② to adjust the position and intensity of the pointer. (Fig. 1)


4. Begin to place the connection tubes on both sides of the splitter attachment. Firmly screw the black ring nut. (Fig. 3-4).
 - **Every connection is identified by a letter printed on both sides of the connection. Make sure to match the letters in order to correctly assemble the microscope.**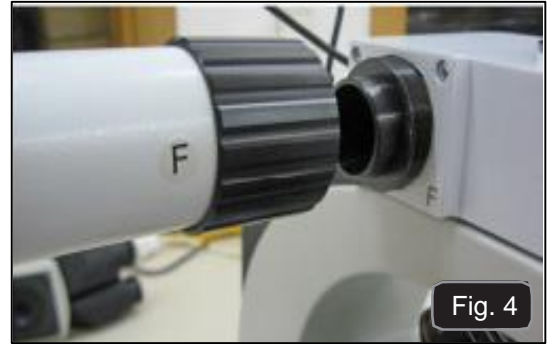

5. Insert the second part of the extension tube (Fig. 5).
6. Fully insert the second extension tube in the right position. Using the provided Allen wrench (small one) lock the fixing screws ① to block the extension tube

➤ **At the end of the first extension tube there is a lens (Fig. 6). Make sure it is free from dirt, dust or other contaminants before to proceed with the assembling of the second extension tube.**

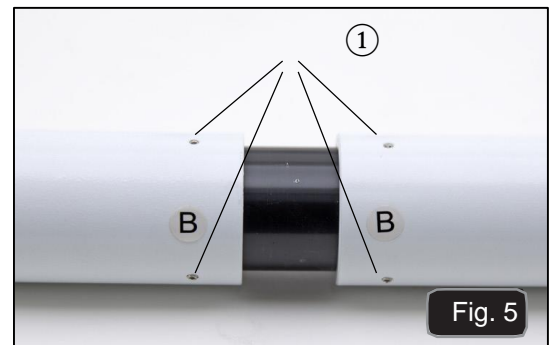7. Adjust the height of the multi-head holder. Loosen the base fixing knob ②, unscrew the base ③ in order to reach the desired height, then lock the knob. (Fig. 7). Make sure that each extension tube is perfectly horizontal.

8. Insert the binocular heads, matching the reference letters (Fig. 8).

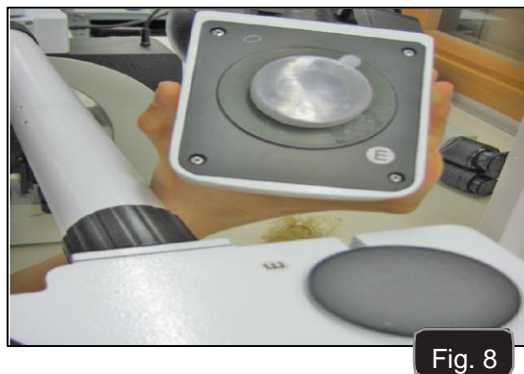
Fig. 8

9. Insert the provided eyepieces (WF10X/20) into binocular heads. (Fig. 9)

10. Repeat all the above operations for each observation point.

Fig. 9

11. Install the trinocular head over the splitter. (Fig. 10)

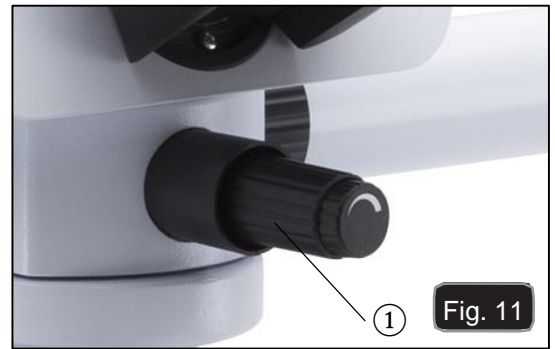
Fig. 10

8. Use of the microscope

Moving the pointer (Fig. 11)

1. By moving the pointer joystick ① it is possible to change the position of the light arrow visible inside the eyepieces.
2. This pointer is used by the teacher to indicate the issue under discussion.

Changing the pointer color (Fig. 12)

3. Press the color selection ② button placed on the left side of the splitter in order to change the color of the pointer. Several pressing allow the cyclical changing of the color in this sequence: RED → GREEN → BLUE → OFF. (Fig. 12).

Changing the intensity of the pointer (Fig. 13)

4. Rotate the pointer intensity controller ③ to change the intensity of the pointer. Adapt the intensity according to the specimen under observation.

9. Maintenance

Microscopy environment

This microscope is recommended to be used in a clean, dry and shock free environment with a temperature of 5°-40°C and a maximum relative humidity of 75 % (non condensing). Use a dehumidifier if needed.

To think about when and after using the microscope

- The microscope should always be kept vertically when moving it and be careful so that no moving parts, such as the eyepieces, fall out.
- Never mishandle or impose unnecessary force on the microscope.
- Never attempt to service the microscope yourself.
- After use, turn off the light immediately, cover the microscope with the provided dust-cover, and keep it in a dry and clean place.

Electrical safety precautions

- Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off-position.
- Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

Cleaning the optics

- If the optical parts need to be cleaned try first to: use compressed air.
- If that is not sufficient: use a soft lint-free piece of cloth with water and a mild detergent.
- And as a final option: use the piece of cloth moistened with a 3:7 mixture of ethanol and ether. Note: ethanol and ether are highly flammable liquids. Do not use them near a heat source, near sparks or near electric equipment. Use these chemicals in a well ventilated room.
- Remember to never wipe the surface of any optical items with your hands. Fingerprints can damage the optics.
- Do not disassemble objectives or eyepieces in attempt to clean them.

For the best results, use the OPTIKA cleaning kit (see catalogue).

If you need to send the microscope to Optika for maintenance, please use the original packaging.

10. Troubleshooting

Refer to the information listed in the microscope instruction manual to troubleshoot operating problems.

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA S.r.l.