HDMI Cameras – H Series

INSTRUCTION MANUAL

Model
C-H4K

Ver. 1.2 2020

Table of Contents

1. Warning	3
2. Symbols and conventions	3
3. Safety Information	3
4. Intended use	3
5. C-H4K Basic features	4
5.2 Quick Instructions for C-H4K camera	5
5.2.1 HDMI Mode	5
5.2.2 USB Mode	5
5.2.3 WLAN Mode	6
5.2.4 GE Ethernet Mode	7
5.2.5 Connecting multi-cameras to a router via GE port/WLAN for network application	9
6. C-H4K User Interface (UI) and its Functions	12
6.1 The Camera Control Panel	13
6.2 The Measurement Toolbar	14
6.3 Icons and Functions of the Camera Control Toolbar	15
6.3.1 Setting > Network > General	16
6.3.2 Setting > Network > LAN	16
6.3.3 Setting > Network > WLAN	18
6.3.4 Setting > Measurement	18
6.3.5 Setting > Magnification	19
6.3.6 Setting > Image Format	19
6.3.7 Setting > Storage	20
6.3.8 Setting > Time	20
6.3.9 Setting > Language	21
6.3.10 Setting > Miscellaneous	22
7. Micrometric Slide M-005	23
Equipment disposal	24

1. Warning

This device is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.

2. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

3. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

4. Intended use

Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

5. C-H4K Basic features

C-H4K is a multiple interfaces (HDMI + USB3.0 + GE + WiFi + SD card) CMOS camera. HDMI, USB3.0 and GE are used as the data transfer interface to HDMI display or computer.

For HDMI output, the *Camera Control Panel + Measurement Toolbar* and *Camera Control Toolbar* are overlaid on the HDMI screen when the mouse move to the related region. In this case, the USB mouse can be used to set the camera, browse and compare the captured image, play the video and perform the measurement.

For USB output, there are three possibilities:

1. Connect USB flash drive to save pictures and videos.
2. Connect 5G WLAN modules to transfer video wirelessly in real time (AP/STA)
3. Connect computer with USB connection to transfer video in real time.

For GE output, the camera can be connected with an Ethernet cable directly to a PC or to a router. From here the ProView software can control each camera function.

5.1 Package content

- | | |
|-------------------|---------------------|
| ① Camera | ⑤ Micrometric slide |
| ② HDMI cable | ⑥ SD card 16GB |
| ③ WLAN 5G adapter | ⑦ Power supply |
| ④ Wireless mouse | ⑧ USB3.0 cable |

5.2 Quick Instructions for C-H4K camera

Before turning on the camera, please connect the standard C-mount camera to the camera adapter and insert it into the microscope's trinocular port.

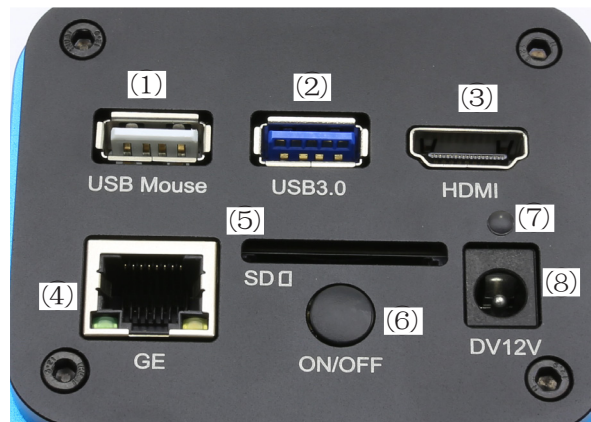
Figure 1 - C-H4K Rear Panel

- | | |
|-----------------------|---------------------|
| ① USB mouse connector | ⑤ SD card slot |
| ② USB3.0 connector | ⑥ ON-OFF button |
| ③ HDMI connector | ⑦ LED indicator |
| ④ GE connector | ⑧ Power supply plug |

You can use the C-H4K camera in 5 different ways. Each application requires different hardware environment.

5.2.1 HDMI Mode

1. Plug the HDMI cable into the HDMI port ③ to connect the C-H4K camera to HDMI display.
2. Plug the USB mouse into USB Mouse port ① to get control of the camera by using built-in software ProView.
3. Plug 12V/1A power adapter into DC 12V slot ⑧ to supply power for the camera. The LED Indicator ⑦ will turn into red.
4. Insert SD card into SD Card slot ⑤ or a USB flash disk into USB3.0 port ② for saving captured images and recorded videos.
5. Press ON/OFF Button ⑥ to turn on the camera. The LED indicator will turn into blue.
6. Move mouse cursor to the left side of the video window, a *Camera Control Panel* will appear. It includes Manual/ Automatic Exposure, White Balance, Sharpness and other functions, please refer to 6.1 for details.
7. Move mouse cursor to the bottom of the video window and a *Camera Control Toolbar* will appear. Operations like Zoom In, Zoom Out, Flip, Freeze, Cross Line, Comparison and so on can be achieved. Please refer to 6.3 for details.
8. Move mouse cursor to the upper side of the video window, a *Measurement Toolbar* with calibration and other measurement tools will appear, please refer to 6.2 for details. The measurement data can be output with *.CSV format.

5.2.2 USB Mode

1. Plug the USB cable provided with the camera into USB3.0 ② to connect the C-H4K camera to the computer.
2. Plug 12V/1A power cable into DC 12V slot ⑧ to supply power for the camera. The LED Indicator ⑦ will turn into red.
3. Press ON/OFF Button ⑥ to turn on the camera. The LED indicator will turn into blue.
4. Start the ProView software.
5. Clicking the camera name in the camera list starts the live video.
 - **USB3.0 video output function is implemented by UVC protocol based on ISO transfer. But USB3.0 UVC output is not supported by Windows 7 or below. It could work on Windows 8 and Windows 10. To use the USB video output function on Windows 7 or below, users must connect the USB3.0 cable to a USB2.0 port on the computer.**
 - **The video output is in MJPEG format and the frame rate of the video output for USB3.0 and USB2.0 will be the same.**

5.2.3 WLAN Mode

The PC should be a WLAN enabled one.

- For Windows user (Windows 7/8/10 [32/64 bit]), use ProView.
- For Mac OS and Linux user (Mac OS 10.10 or above or Linux distributions with kernel 2.6.27 or higher), use LiteView.
- When connecting the camera with a mobile device, the free App is required. Just make sure that the mobile device uses iOS 11 or higher/Android 5.1 or higher operating systems.

1. Plug the HDMI cable into the HDMI port ③ to connect the C-H4K camera to HDMI display.
2. Plug the USB mouse into USB Mouse port ① to get control of the camera by using built-in software ProView.
3. Plug 12V/1A power adapter into DC 12V slot ⑧ to supply power for the camera. The LED Indicator ⑦ will turn into red.
4. Press ON/OFF Button ⑥ to turn on the camera. The LED indicator will turn into blue.
5. Move the mouse to the bottom of the UI and click the  button on the *Camera Control Toolbar* at the bottom of the video window. A small window called *Settings* will pop up as shown below.



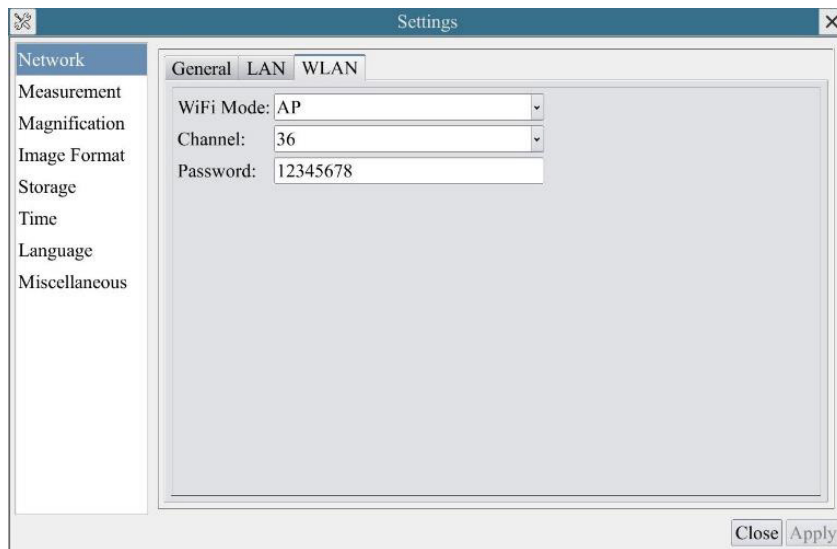
Figure 2 - WLAN configuration

6. Click *Network>WLAN* property page and choose the *AP* option in the *WiFi Mode* edit box (the factory default configuration is *AP* mode).
7. Install ProView or LiteView on your PC or install the app on the mobile device.
8. Plug the USB WLAN adapter into the camera's USB3.0 port.
9. Connect the PC or mobile device with the WLAN AP point that the camera provides.
 - The network name (*SSID*) and the WLAN password (default one is *12345678*) can be found on the camera's *Setting>Network>WLAN* page in *AP* mode.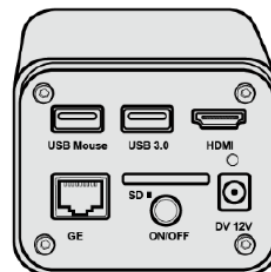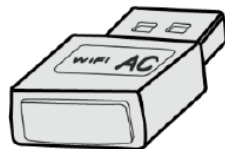

10. Start the ProView / LiteView software or the mobile App and check the configuration. Normally, active cameras are automatically recognized. The live image of each camera is displayed.
11. For the display, the *Camera List* tool window is used in the ProView/LiteView software, and the *Camera Thumbnail* is used in the mobile App.

5.2.4 GE Ethernet Mode

This application uses the camera as the network camera. User must configure the IP of the camera and PC manually and ensure their IP addresses are in the same net. The subnet mask and gateway of the camera and PC must be the same.

1. Plug the HDMI cable into the HDMI port ③ to connect the C-H4K camera to HDMI display.
2. Plug the USB mouse into USB Mouse port ① to get control of the camera by using built-in software ProView.
3. Plug 12V/1A power adapter into DC 12V slot ⑧ to supply power for the camera. The LED Indicator ⑦ will turn into red.
4. Press ON/OFF Button ⑥ to turn on the camera. The LED indicator will turn into blue.
5. Move the mouse to the bottom of the UI and clicking the  button on the *Camera Control Toolbar* at the bottom of the video window. A small window called *Settings* will pop up as shown below.
6. Click the *LAN* property page, uncheck the *DHCP* item. Input the *IP Address*, *Subnet Mask* and *Default Gateway* for the camera.

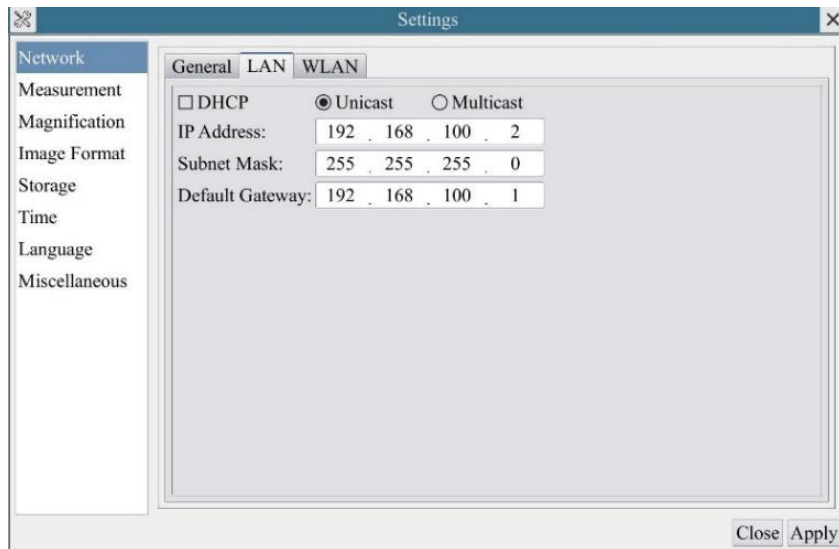
Figure 3 - IP Camera configuration

7. Designate the Internet Protocol Version 4 (TCP/IPv4) Settings page's IP address on the PC with similar configuration as shown below but with different IP address.



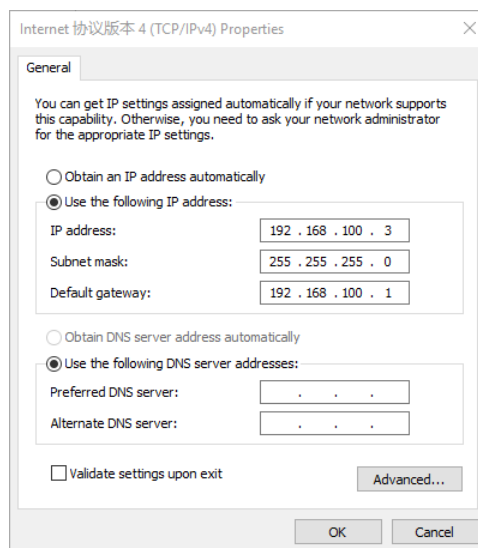
Figure 4 - PC's IP configuration


8. After the above configurations are finished, user can connect the camera to the computer through the USB to Ethernet adapter as shown below:
 - Connect the GE port with the Ethernet cable to the PC's network port.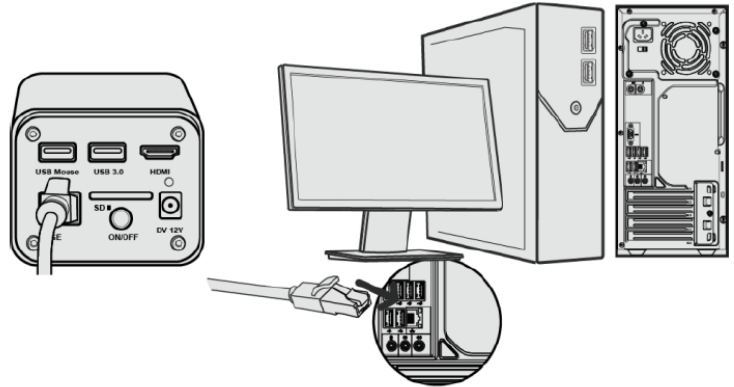

- Insert the SD card (supplied) or USB flash disk (not supplied) into the camera SD card slot / USB3.0 port.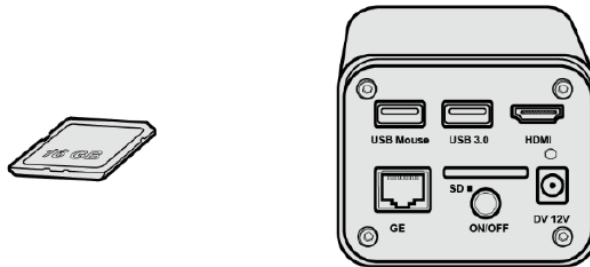

9. Start the ProView / LiteView software or the mobile App and check the configuration.
10. Clicking the camera name in the camera list starts the live video.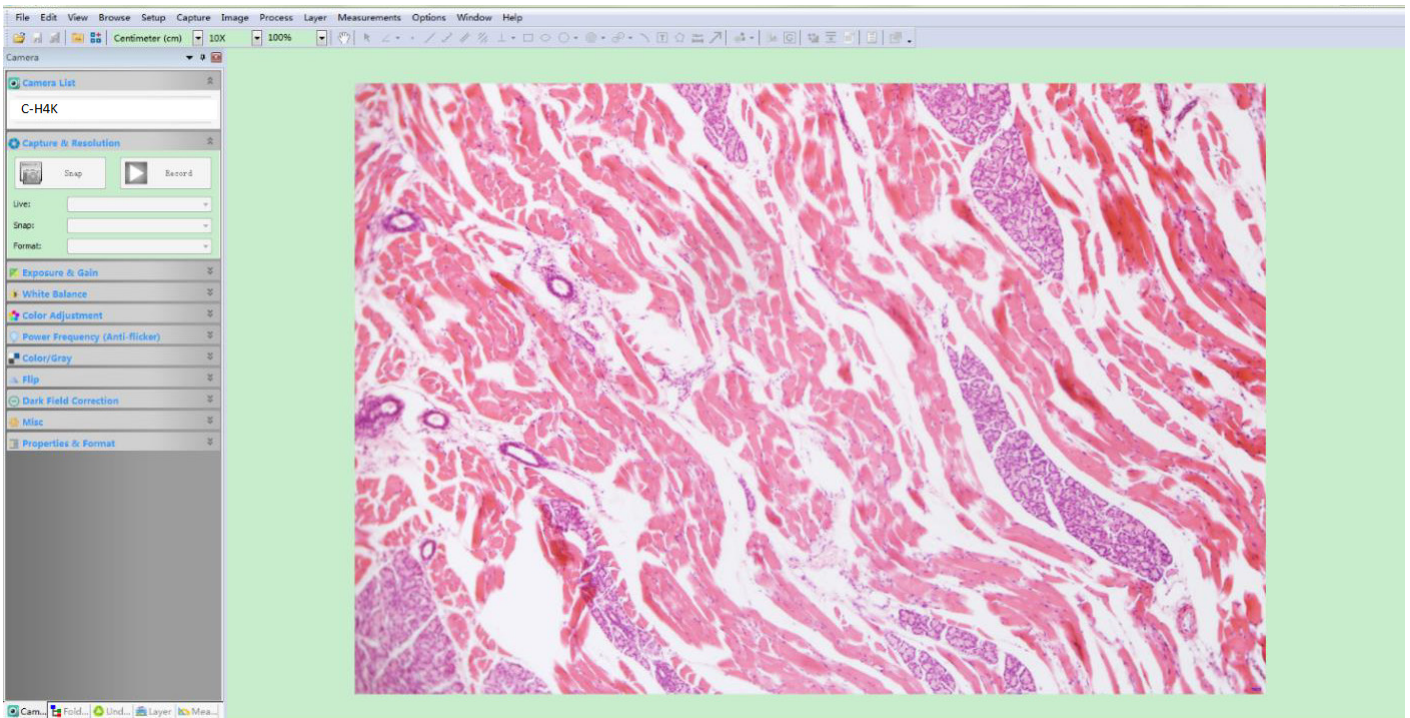

- When connecting via WLAN the connection and configuration are just the same as in Sec. 5.2.2 but here, users need to check *STA*.


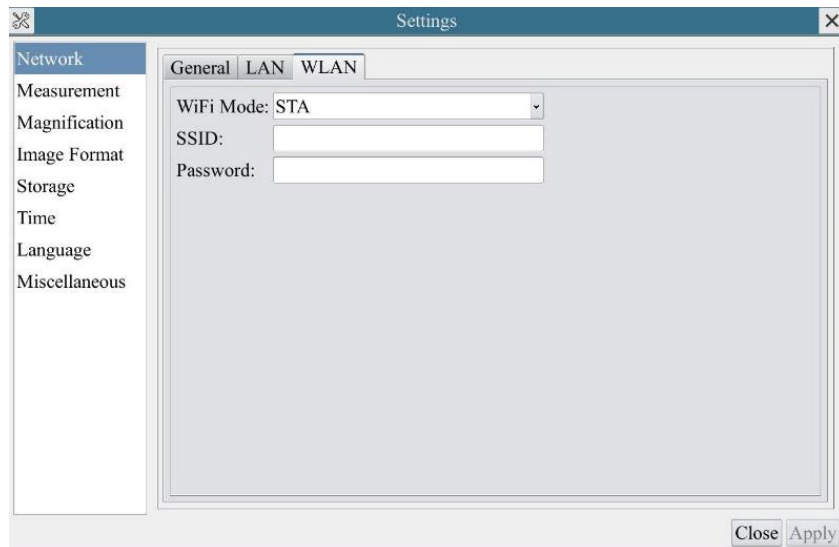
Figure 6 - WLAN configuration


- Install ProView or LiteView on your PC or install the app on the mobile device.
 - Connect the GE port with the Ethernet cable to the PC's network port or plug the USB WLAN adapter into the camera's USB3.0 port (for those connected to router with WLAN STA mode).

- Finally as shown below, 2 HDMI cameras are connected to the router with GE cable and 2 HDMI cameras are connected to the same router with WLAN STA mode (the number of the cameras, the connection mode (GE or WLAN STA)) connected to the router are determined by the router performance).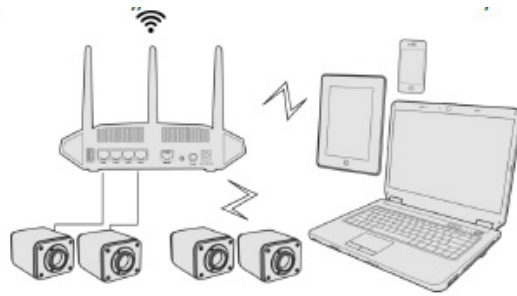

- Make sure that your PC or your mobile device is connected to the LAN or WLAN of the router.
- Start the ProView / LiteView software or the mobile App and check the configuration. Normally, active cameras are automatically recognized. The live image of each camera is displayed.
- For the display, the *Camera List* tool window is used in the ProView/LiteView software, and the *Camera Thumbnail* is used in the mobile App.

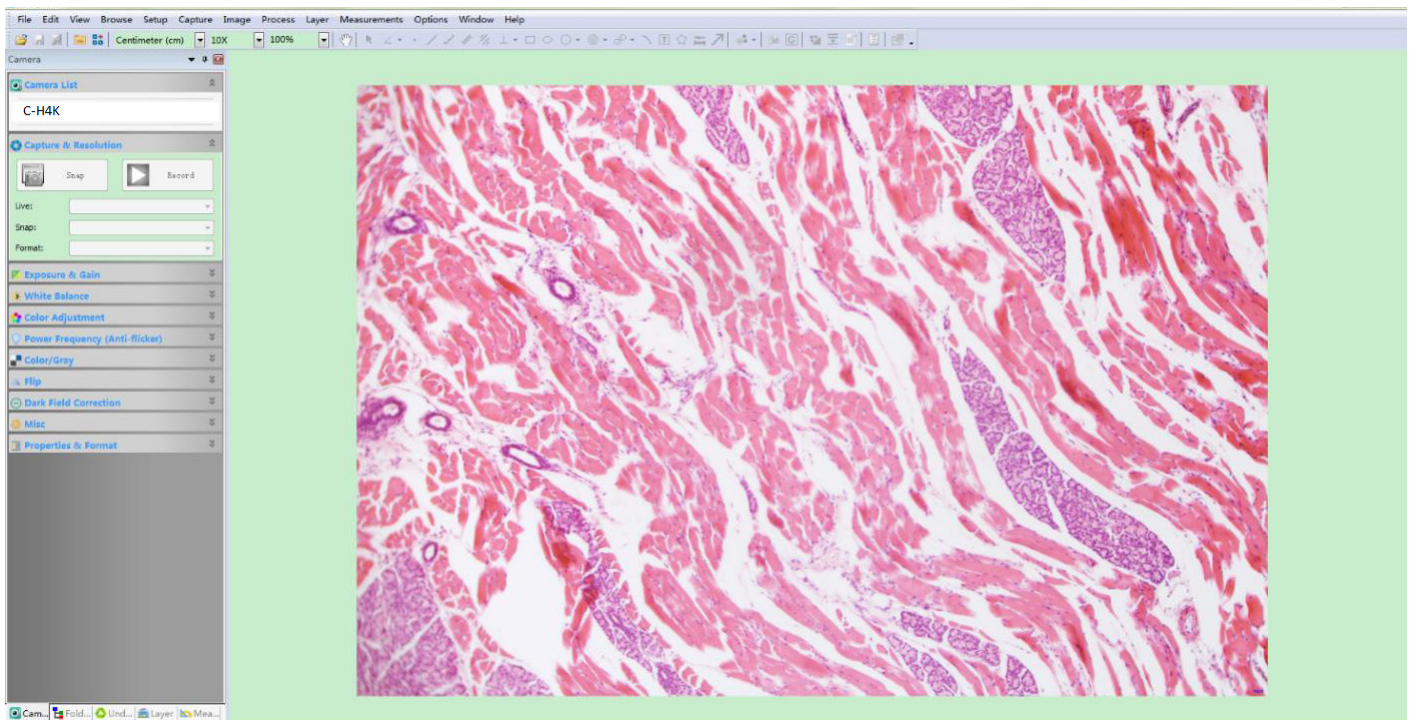- Note on data security**

The data transfer of the HDMI 4K camera in LAN or WLAN is not encrypted. Anyone who is connected to the network and has installed the Optika software or mobile App, can see the live image of all active cameras. Operate the camera with the ProView software, if you want to make sure that nobody in the network can see the camera's live image.

- About the routers/switches**

It is suggested that routers/switches supporting 802.11ac 5G segment should be selected to achieve better wireless connection experience.

6. C-H4K User Interface (UI) and its Functions

The C-H4K UI shown in Fig. 6 includes a *Camera Control Panel* on the left side of the video window, a *Measurement Toolbar* on the upper side of the video window, a *Camera Control Toolbar* on the bottom of the video window.

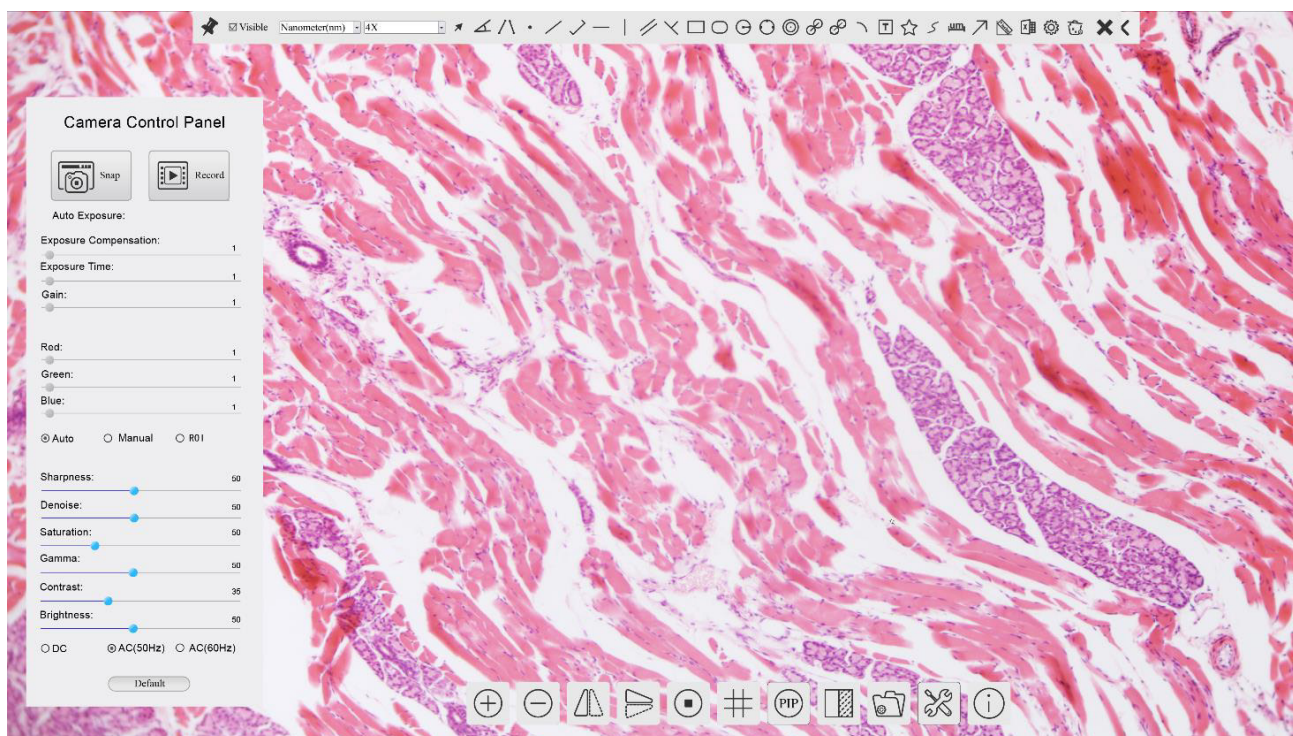

Figure 6 - Camera Control User Interface

1. Move the mouse cursor to the left side of the video window: the *Camera Control Panel* will pop up automatically.
2. Move the mouse cursor to the upper side of the video window: a *Measurement Toolbar* will pop up for the calibration and measurement operations.
 - When user left-clicks the *Float/Fixed* button  on the Measurement Toolbar, the Measurement Toolbar will be fixed. In this case the Camera Control Panel will not pop up automatically even if users move mouse cursor to left side of the video windows.
 - Only when user left-clicks the  button on the *Measurement Toolbar* to exit from measuring procedure it will be possible to perform other operations on the *Camera Control Panel*, or *Camera Control Toolbar*.
 - During the measuring process, when a specific measuring object is selected an *Object Location & Attributes Control Bar*  will appear for changing location and properties of the selected objects.
3. Move the mouse cursor to the bottom of the video window, the *Camera Control Toolbar* will pop up automatically.

6.1 The Camera Control Panel

The *Camera Control Panel* controls the camera to achieve the best image quality according to the specific applications. It will pop up automatically when mouse cursor is moved to the left side of the video window (in measurement status, the Camera Control Panel will not pop up. Only when measurement process is terminated will the Camera Control Panel pop up by moving mouse cursor to the left side of the video window). Left-clicking button  to achieve *Display/ Auto Hide* switch of the Camera Control Panel.

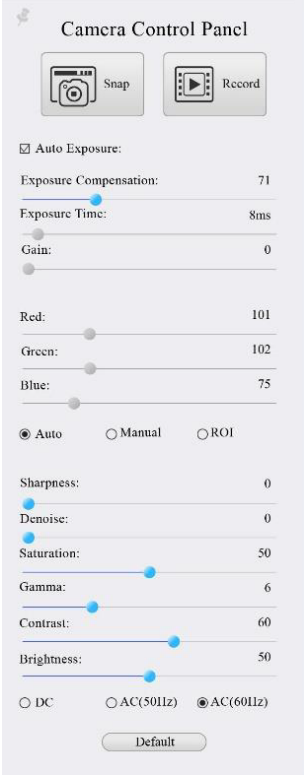Control Panel	Function	Description
 <p>The screenshot shows the Camera Control Panel with the following settings: Snap and Record buttons; Auto Exposure checked; Exposure Compensation: 71; Exposure Time: 8ms; Gain: 0; Red: 101; Green: 102; Blue: 75; Auto White Balance selected; Sharpness: 0; Denoise: 0; Saturation: 50; Gamma: 6; Contrast: 60; Brightness: 50; and AC(60Hz) selected for illumination.</p>	Snap	Capture image from the current video window
	Record	Record video from the current video window
	Auto Exposure	When <i>Auto Exposure</i> is checked, the system will automatically adjust exposure time according to the value of exposure compensation
	Exposure Compensation	Available when <i>Auto Exposure</i> is checked. Slide to left or right to adjust <i>Exposure Compensation</i> according to the current video brightness to achieve proper brightness value
	Exposure Time	Available when <i>Auto Exposure</i> is unchecked. Slide to left or right to reduce or increase exposure time, adjusting brightness of the video
	Gain	Adjust <i>Gain</i> to reduce or increase brightness of video. The Noise will be reduced or increased accordingly
	Red	Slide to left or right to decrease or increase the proportion of <i>Red</i> in RGB on video
	Green	Slide to left or right to decrease or increase the proportion of <i>Green</i> in RGB on video
	Blue	Slide to left or right to decrease or increase the proportion of <i>Blue</i> in RGB on the video
	Auto White Balance	<i>White Balance</i> adjustment according to the window video every time the button is clicked
	Manual White Balance	Slide the <i>Red</i> or <i>Blue</i> to manually set the video <i>White Balance</i>
	Sharpness	Adjust <i>Sharpness</i> level of the video window
	Denoise	Slide left or right to <i>denoise</i> the image
	Saturation	Adjust <i>Saturation</i> level of the video window
	Gamma	Adjust <i>Gamma</i> level of the video. Slide to the right side to increase gamma and to the left to decrease gamma
	Contrast	Adjust <i>Contrast</i> level of the video. Slide to the right side to increase contrast and to the left to decrease contrast
	DC	For <i>DC</i> illumination, there will be no fluctuation in light source so no need for compensating light flickering
	AC(50HZ)	Check <i>AC(50HZ)</i> to eliminate flickering “strap” caused by 50Hz illumination
	AC(60HZ)	Check <i>AC(60HZ)</i> to eliminate flickering “strap” caused by 60Hz illumination
	Default	Restore all the settings in the <i>Camera Control Panel</i> to default values

6.2 The Measurement Toolbar

The *Measurement Toolbar* will pop up when moving mouse cursor to any place near the upper side of the video window.

Figure 7 - The Measurement Toolbar

Icon	Function
	<i>Float/ Fix</i> switch of the <i>Measurement Toolbar</i>
<input checked="" type="checkbox"/> Visible	Define measuring object in <i>Show up/ Hide</i> mode
Nanometer(nm)	Select the desired <i>Measurement Unit</i>
4X	Select <i>Magnification for Measurement after Calibration</i>
	Object <i>Select</i>
	<i>Angle</i>
	<i>4 Points Angle</i>
	<i>Point</i>
	<i>Arbitrary Line</i>
	<i>3 Points Line</i>
	<i>Horizontal Line</i>
	<i>Vertical Line</i>
	<i>3 Points Vertical line</i>
	<i>Parallel</i>
	<i>Rectangle</i>
	<i>Ellipse</i>
	<i>Circle</i>
	<i>3 Points Circle</i>
	<i>Annulus</i>
	<i>Two Circles and Center Distance</i>
	<i>3 Points Two Circles and Center Distance</i>
	<i>Arc</i>
	<i>Text</i>
	<i>Polygon</i>

	Curve
	Scale Bar
	Arrow
	Execute <i>Calibration</i> to determine the corresponding relation between magnification and resolution, which will establish the corresponding relationship between measurement unit and the sensor pixel size. <i>Calibration</i> needs to be done with the help of a micrometer. For detailed steps of carrying out <i>Calibration</i> please refer to ProView help manual.
	Export the <i>Measurement</i> to CSV file(*.csv)
	<i>Measurement Setup</i>
	Delete All the Measurement Objects
	Exit from <i>Measurement</i> mode
	When the measurement ends, left-click on a single measuring object and the <i>Object Location & Properties Control Bar</i> will show up. User could move the object by dragging the object with the mouse. But more accurate movement could be done with the control bar. The icons on the control bar mean <i>Move Left, Move Right, Move Up, Move Down, Color Adjustment</i> and <i>Delete</i> .

- Note:
1. When user left-clicks *Display/Hide* button  on the *Measurement Toolbar*, it will be fixed. In this case the *Camera Control Panel* will not pop up automatically even if moving the mouse cursor to the left edge of the video window. Only when user left-click the  button on the *Measurement Toolbar* to exit from the measurement mode it will be possible to perform other operations on the *Camera Control Panel* or on the *Camera Control Toolbar*.
 2. When a specific item is selected during the measurement process, the *Object Location & Attributes Control Bar*  will appear for changing the object location and properties of the selected objects.

6.3 Icons and Functions of the Camera Control Toolbar

Icon	Function	Icon	Function
	Zoom In the Video Window		Zoom Out the Video Window
	Horizontal Flip		Vertical Flip
	Video Freeze		Display Cross Line
	Picture in Picture		Compare Image with the current video
	Browse Images and Videos in the SD Card		Settings
	Check the ProView Version		

The  *Setting* function is relatively more complicated than the other functions. Here are more info about it:

6.3.1 Setting > Network > General

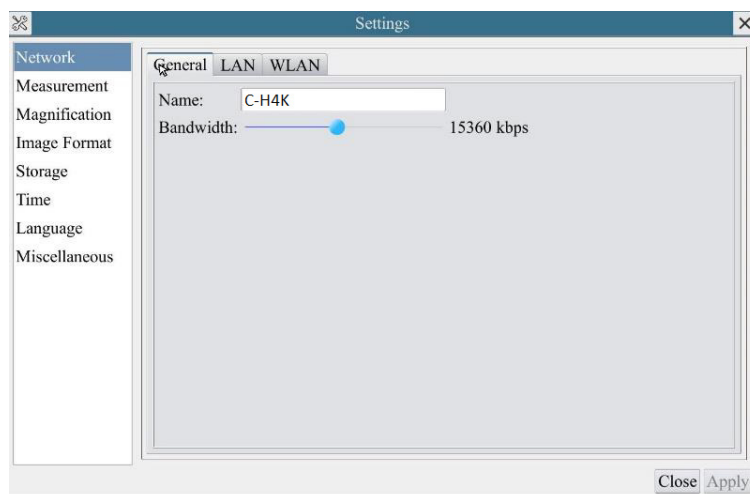
Figure 7 - General Network Setting Page

Name The current camera name recognized as the network name.

Bandwidth The encoding bandwidth for the video transmission stream. The larger the bandwidth, the higher quality the video.

6.3.2 Setting > Network > LAN

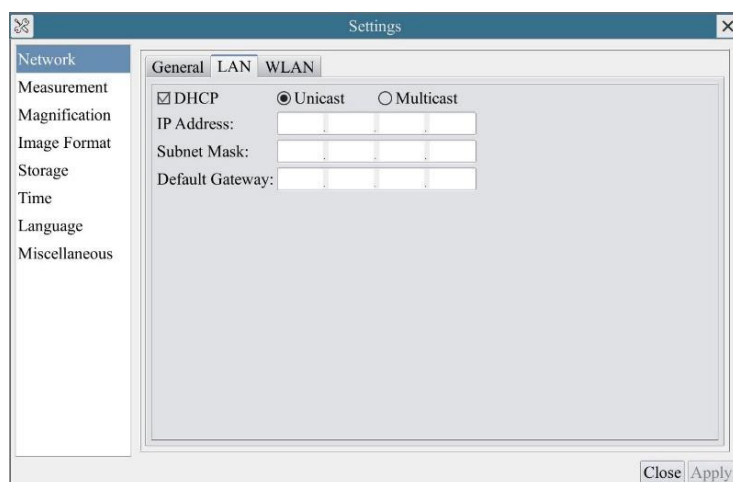
Figure 8 - LAN Network Setting Page

DHCP Dynamic host control protocol allows DHCP server to automatically assign IP information to the camera. Only in Sec. 5.2.2 LAN networking this item should be checked, so that cameras can automatically get IP information from routers/switches to facilitate networking operation.

Unicast/Multicast By default, *Unicast* function is used. Only in Sec. 5.2.3 networking environment, when the router/switch has *Multicast* function, camera can switch to multicast mode, which can save the network bandwidth consumed by the camera and facilitate the connection of more cameras in the same network.

Every machine on a network has a unique identifier. Just as you would address a letter to send in the mail, computers use the unique identifier to send data to specific computers on a network. Most networks today, including all computers on the Internet, use the TCP/IP protocol as the standard to communicate on the network. In the TCP/IP protocol, the unique identifier for a computer is called *IP address*.

There are two standards for *IP address*: IP Version 4 (IPv4) and IP Version 6 (IPv6). All computers with IP addresses have an IPv4 address, and many are starting to use the new IPv6 address system as well.

IP Address

Users must manually configure their *IP addresses* on the camera and computer side. The camera side and computer side should be in the same network segment. The specific settings are shown.

This is usually a private address. Private address is a non-registered address used exclusively within an organization. The internal private addresses retained are listed below: Class A 10.0.0-10.255.255; Class B 172.16.0-172.31.255.255; Class C 192.168.0-192.168.255.255. The suggested *IP address* is Class C.

Subnet Mask

Subnet Mask is used to distinguish network domain from host domain in 32-bit IP address.

Default Gateway

A default gateway allows computers on a network to communicate with computers on another network. Without it, the network is isolated from the outside. Basically, computers send data that is bound for other networks (one that does not belong to its local IP range) through the default gateway.

Network administrators configure the computer's routing capability with an IP range's starting address as the default gateway and point all clients to that IP address.

Uncheck the *DHCP* and select the *Unicast* item, user still need to set the *IP address*, *Subnet mask* and *Default Gateway* as shown below:

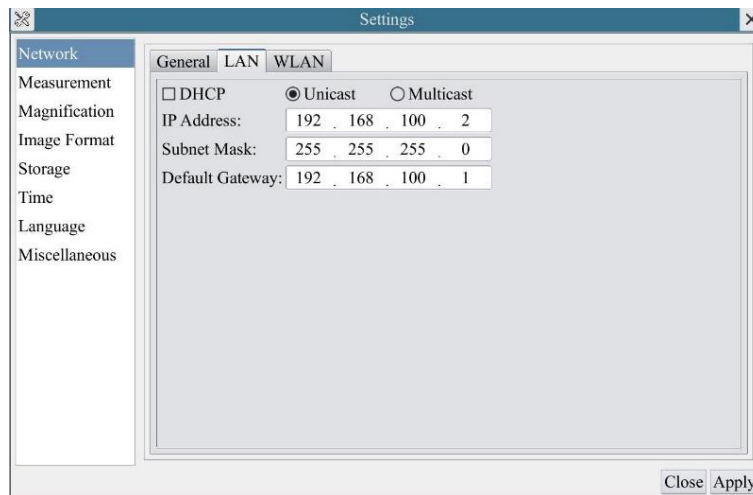
Figure 9 - Manual DHCP in Unicast

Uncheck the *DHCP* and select the *Multicast* item, user still need to set the *IP address*, *Subnet Mask* and *Default Gateway* as shown below:

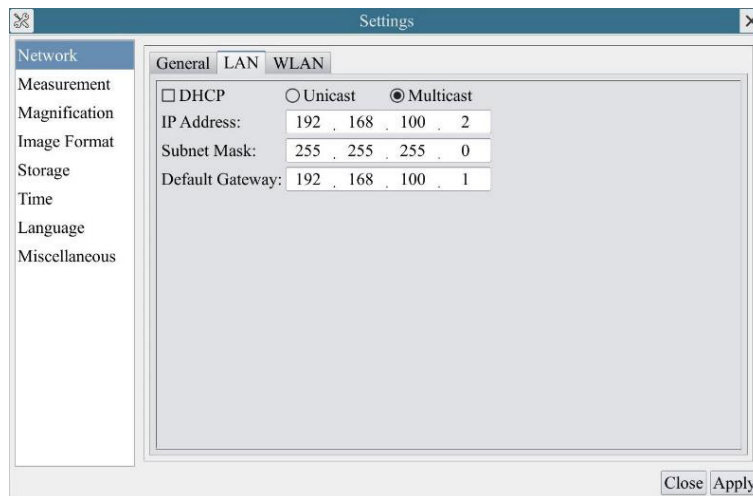
Figure 10 - Manual DHCP in Multicast

6.3.3 Setting > Network > WLAN

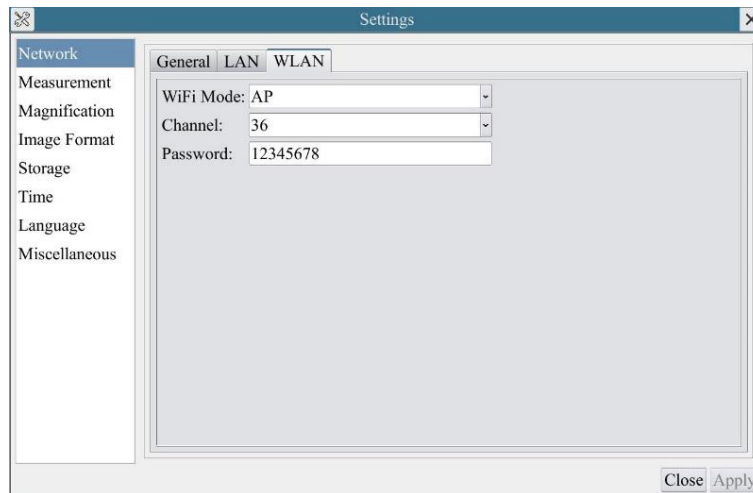
Figure 11 - WLAN Network Setting Page

WiFi Mode AP/STA mode to select.

Channel/SSID Channel for the AP mode and SSID for the STA mode. Here, the SSID is the router's SSID.

Password Camera Password for the AP mode. Router Password for the STA mode.

6.3.4 Setting > Measurement

This page is used for the definition of the *Measurement Object* properties

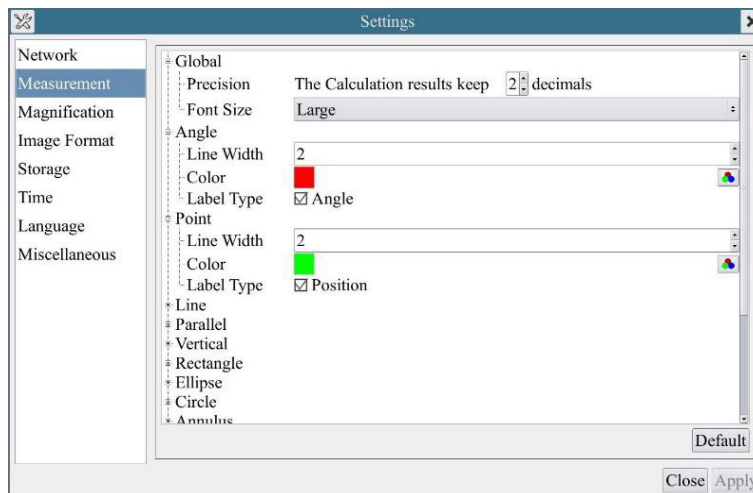
Figure 12 - Measurement Page

Global: Used for setting digits behind the decimal point for measurement results.

Calibration *Line Width* Used for defining width of the lines for calibration.

Color Used for defining color of the lines for calibration.

EndPoint *Type:* Used for defining shape of the endpoints of lines for calibration.

Null: Means no endpoints.

Rectangle: Means rectangle type of endpoints. It makes alignment more easily.

Point, Angle, Line, Horizontal Line, Vertical Line, Rectangle, Circle, Ellipse, Annulus, Two Circles, Polygon, Curve

Left-click the  beside the measuring patterns mentioned above will unfold the corresponding attribute settings to set the individual property of the measuring object.

6.3.5 Setting > Magnification

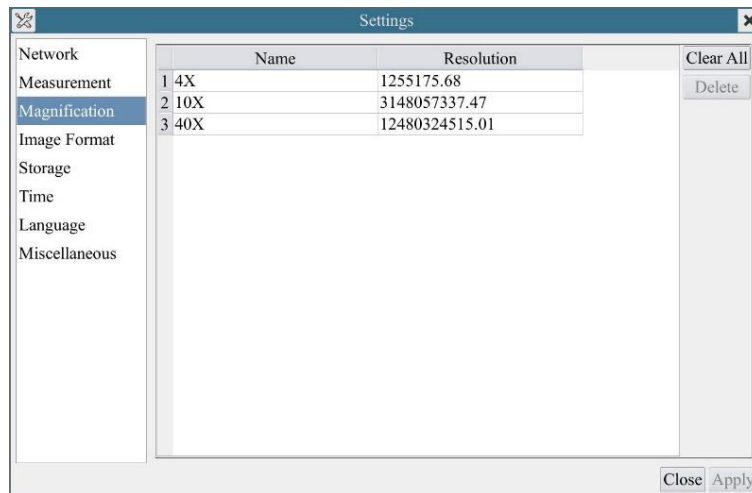
Figure 13 - Magnification Settings Page

- Name** Names such as 10X, 40X, 100X are based on magnification of the microscopes. For continuous zoom microscopes, ensure that the selected magnification coincides with the scale alignment line on the microscope zoom knob.
- Resolution** Pixels per meter. Devices like microscopes have high resolution value.
- Clear All** Click the *Clear All* button will clear the calibrated magnifications and resolutions.
- Delete** Click *Delete* to delete the selected item for specific resolution.

6.3.6 Setting > Image Format

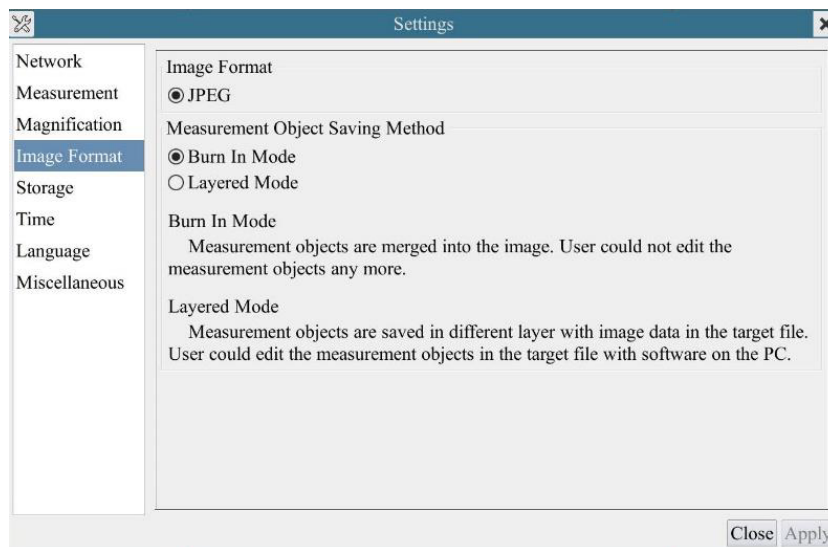
Figure 14 - Image Format Settings Page

- Image Format** *JPEG*: The extension of JPEG file can get very high compression rate and display very rich and vivid images by removing redundant images and color data. In other words, it can get better image quality with the least disk space. If measurement objects are available, the measurement objects will be burned into the image and the measurement cannot be edited.
- Measurement Object** *Burn in Mode*: The measurement objects are merged into the current image. User could not edit the measurement objects any more. This mode is not reversible.
- Save Method** *Layered Mode*: The measurement objects are saved in different layer with current image data in the target file. User could edit the measurement objects in the target file with some software on the PC. This mode is reversible.

6.3.7 Setting > Storage

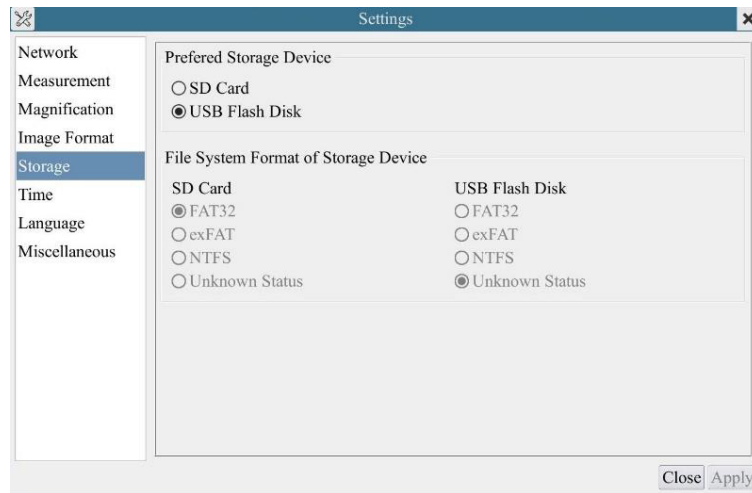
Figure 15 - SD Card Settings Page

Preferred Storage Device

SD Card: Select it to save the video and image to the *SD card*
USB Flash Disk: Select it to save the video and image to the *Flash Disk*

File System

List the file system format of the current storage device

Format of the

FAT32: The file system of SD card is FAT32. The maximum video file size of single file in FAT32 file system is 4G Bytes

Storage Device

NTFS: The file system of SD card is NTFS. The maximum video file size of single file is 2T Bytes

Use PC to format the SD cards and switch between FAT32 and NTFS.

Unknown Status:

SD card not detected or the file system is not identified

6.3.8 Setting > Time

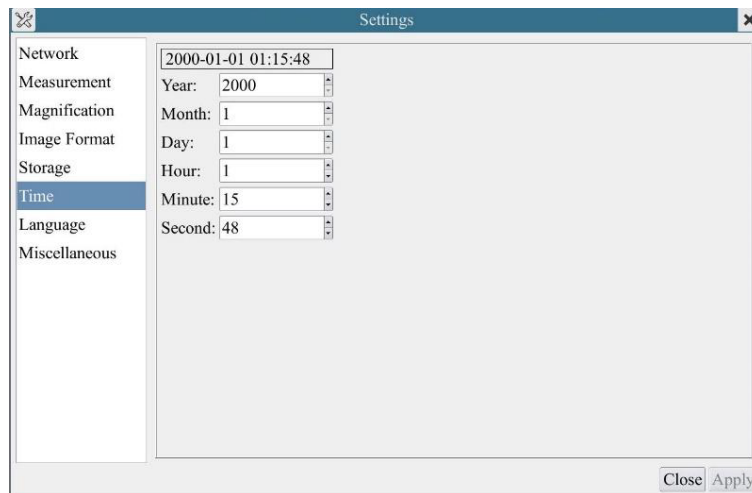
Figure 16 - Time Settings Page

Time

User can set *Year, Month, Day, Hour, Minute* and *Second* in this page

6.3.9 Setting > Language

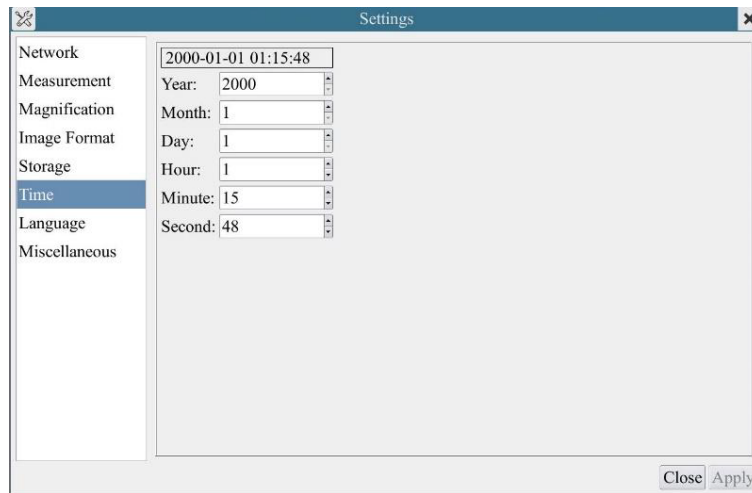
Figure 17 - Language Selection Settings Page

<i>English</i>	Set language of the whole software into English
<i>Simplified Chinese</i>	Set language of the whole software into Simplified Chinese
<i>Traditional Chinese</i>	Set language of the whole software into Traditional Chinese
<i>Korean</i>	Set language of the whole software into Korean
<i>Thailand</i>	Set language of the whole software into Thai
<i>French</i>	Set language of the whole software into French
<i>German</i>	Set language of the whole software into German
<i>Japanese</i>	Set language of the whole software into Japanese
<i>Italian</i>	Set language of the whole software into Italian

6.3.10 Setting > Miscellaneous


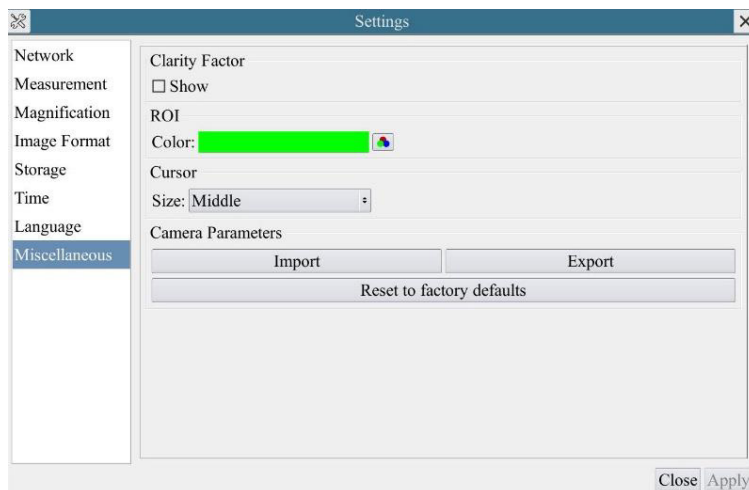
Figure 18 - Miscellaneous Settings Page

<i>Clarity Factor Show</i>	Check this to show the <i>Clarity Factor</i> on the video window screen and inform if the camera is correctly focused or not.
<i>ROI Color</i>	Choosing the <i>ROI</i> rectangle line color.
<i>Cursor</i>	Choosing the <i>Cursor</i> size according to the screen resolution or personal preference.
<i>Camera Parameters Import</i>	Import the <i>Camera Parameters</i> from the SD card or USB flash disk to use the previously exported <i>Camera Parameters</i> .
<i>Camera Parameters Export</i>	Export the <i>Camera Parameters</i> to the SD card or USB flash disk to use the previously imported <i>Camera Parameters</i> .
<i>Reset to factory defaults</i>	Restore camera parameters to its factory status.


7. Micrometric Slide M-005

Micrometric slide, 26x76mm, with 2 scales
(1mm/100div. for biological microscopes / 10mm/100div. for stereomicroscopes)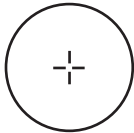

1 DIV=0.01mm

For biological microscopes calibration

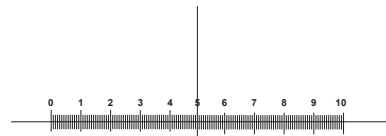1 DIV=0.1mm

For stereomicroscopes calibration

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.