HDMI Autofocus camera, 2 MP CMOS

INSTRUCTION MANUAL

Model
C-HA

Ver. 1.0 2019

Summary

1. Warning	3
2. Symbols and conventions	3
3. Safety Information	3
4. Intended use	3
6. Unpacking	5
7. Assembling	5
7.1 Camera installation	6
7.1.1 Use of “C” mount adapters	6
7.1.2 Use of projection lens	7
7.2 Camera connection	7
7.3 Parfocality adjustment	8
8. Use of the camera	9
8.1 “Control” Menu	9
8.1.1 “Focus” Section	10
8.1.2 “AF ROI” Section	10
8.1.3 “Exposure” Section	12
8.1.4 “Color” Section	13
8.1.5 “Function” Section	13
8.2 “Crosshair” Menu	16
8.3 “System” Menu	17
9. Maintenance	18
Equipment disposal	19

1. Warning

This camera is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.

2. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

3. Safety Information

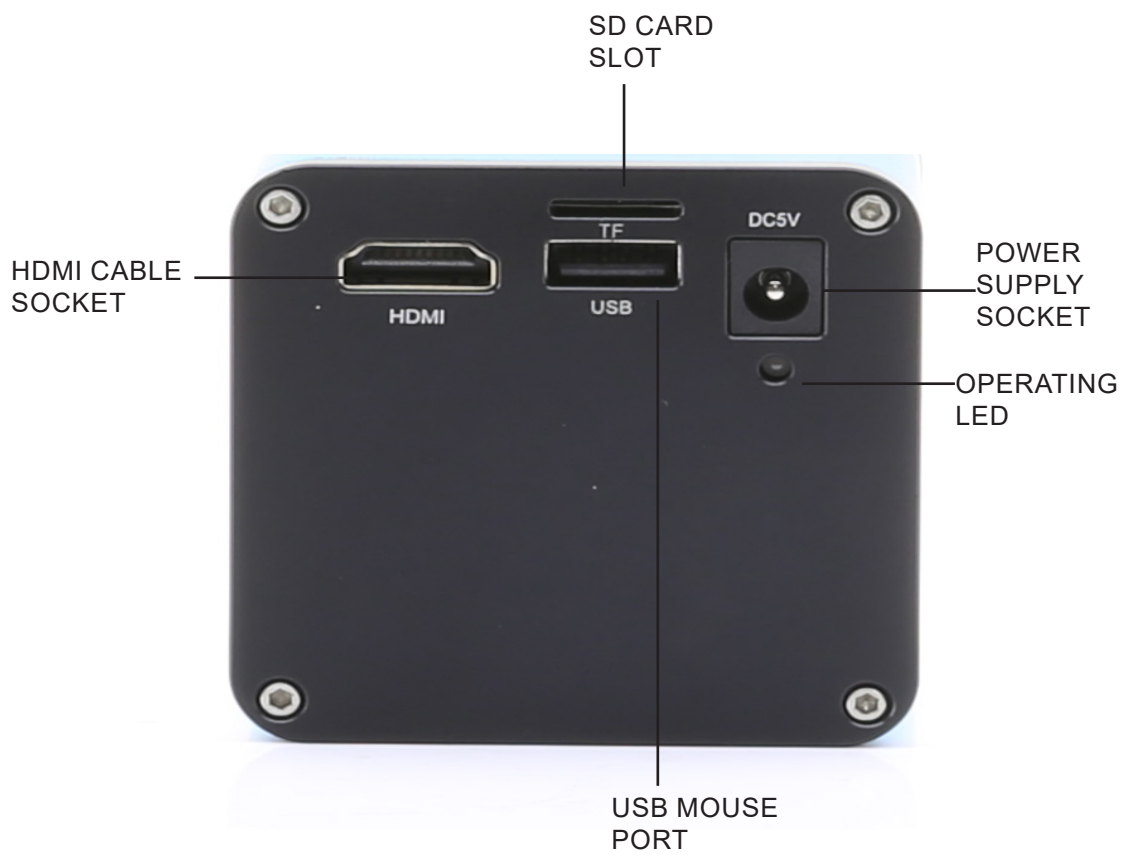
Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

4. Intended use

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

5. Overview

6. Unpacking

The device is housed in a moulded Styrofoam container. Take some care to avoid that the optical items fall out and get damaged. Carefully remove the camera from the container and put it on a stable desk.

 Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

7. Assembling

Once opened the box, the camera parts are the following:

- ① Camera
- ② 16GB micro SD card
- ③ HDMI cable

- ④ Wireless mouse + battery
- ⑤ Power supply

7.1 Camera installation

7.1.1 Use of “C” mount adapters

1. Remove dust caps from camera and “C” mount.
2. Screw the “C” mount to camera thread. (Fig. 1)

3. Loosen the clamping screw from the trinocular head. (Fig. 2)

4. Install the previously assembled C-mount + camera on the trinocular head. (Fig. 3)

5. Tighten the clamping screw firmly. (Fig. 4)

7.1.2 Use of projection lens

1. Remove dust caps from camera and projection lens.
2. Screw the projection lens to camera thread. (Fig. 1)
3. Insert the projection lens on the trinocular port. (Fig. 2)

7.2 Camera connection

1. Connect one side of the HDMI cable to the HDMI socket ①. (Fig. 7)
2. Open the mouse battery cover ②. (Fig. 8)

3. Remove the wireless mouse receiver ③. (Fig. 9)

4. Plug the mouse receiver into the USB port ④ of the camera. (Fig. 10)
5. Insert the battery (respect polarity) and put back the battery cover.

6. Plug the power supply into the DC5V socket ⑤ to power the camera. (Fig. 11)
 - **Once the power supply is connected, the camera takes about 10 seconds to switch on and the LED ⑥ lights up blue. This is not a defect.**

7.3 Parfocality adjustment

In order to have the same focus when observing the specimen through the eyepieces and on the screen/monitor, please verify the microscope is properly installed and set and follow the instructions below.

In case of a biological microscope:

1. Use a low power objective and focus the specimen
2. Switch to the highest dry objective available on the microscope (40x or 60x) and focus the specimen again
3. Turn on the live-view on the camera, without changing the focus on the microscope
4. While observing the image on the screen/monitor, adjust the focus by turning the knurled knob on the C-mount adapter

In case of a stereomicroscope:

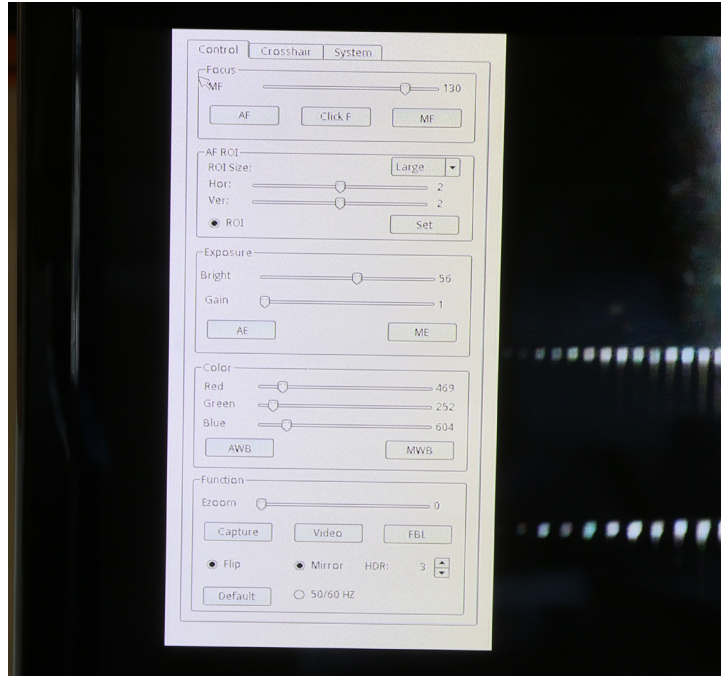
5. Using a low power magnification and focus the specimen
6. Reach the highest magnification available using the zoom knob and then focus the specimen again
7. Turn on the live-view on the camera, without changing the focus on the microscope
8. While observing the image on the screen/monitor, adjust the focus by turning the knurled knob on the C-mount adapter

The proper parfocality adjustment is obtained when the same focus is reached when looking into the eyepieces and on the screen/monitor.


8. Use of the camera

Camera is supplied with an “on screen” menu, which is activated when the mouse cursor is positioned in the upper left corner of the monitor.


Moving the cursor to a different area causes the menu to disappear.

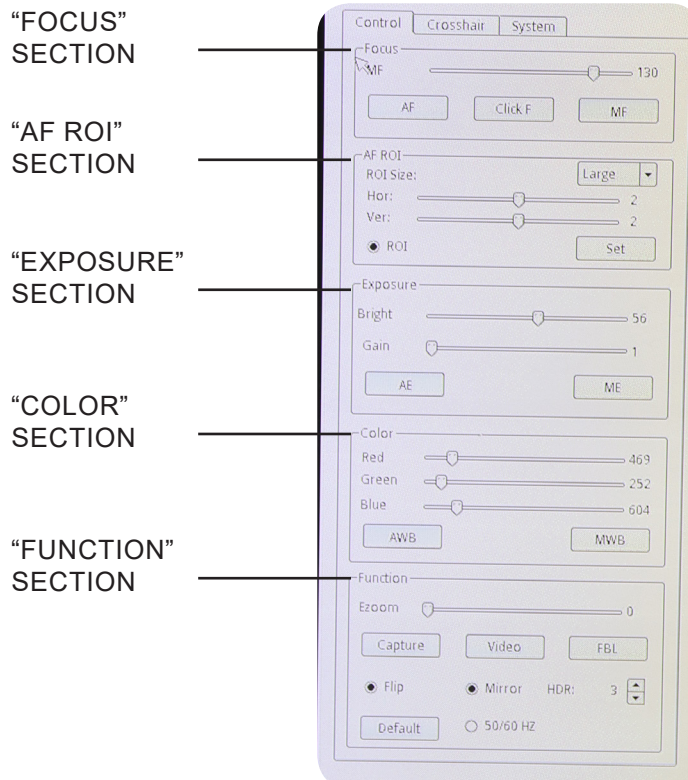8.1 “Control” Menu

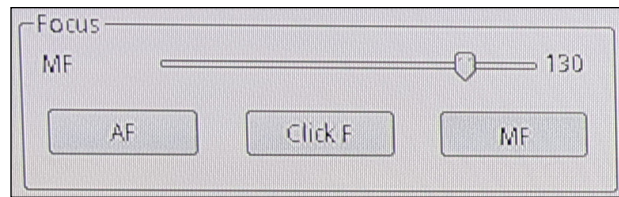8.1.1 “Focus” Section


It is possible to work either in autofocus (AF) or manual focus (MF).

Clicking on “**AF**” button will allow the camera to work in autofocus mode. It will be sufficient to position the target to be observed in front of the camera lens and the focus will be obtained. The autofocus will be independent of the optical or digital zoom value of the camera: the camera must have enough contrast in the sample to be observed.

In the case of objects with low contrast, it is possible that the camera can not get the focus. In this case we recommend using manual focus.

- **It is important to remind that the camera doesn't have an infinity autofocus capability. The user must approximately focus the specimen with the microscope focusing knobs. From this moment the camera will be able to automatically focus the specimen when the stage is moved to another position or the slide is changed on the stage.**

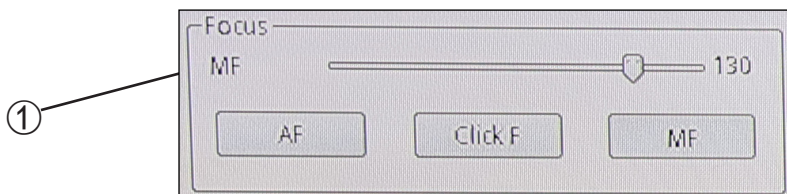
Clicking on “**MF**” button, the camera will work in manual focus mode. Activation of the MF button automatically activate the “**MF**” slider ①, with which you can operate to focus the camera without having to act on the microscope focusing knobs.

Clicking on “**Click F**” the camera is taught which is the optimal focus for the sample that is being observed. From now on, any movement that will be made on the sample or any change in zoom will have to be compensated again with a new focus.


If you work in “AF” mode, this operation will not be necessary.

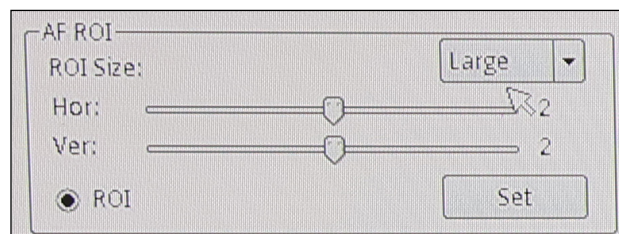8.1.2 “AF ROI” Section


In the “AF ROI” section it is possible to set the size and the position of the ROI (Region Of Interest) where the autofocus will be detected.

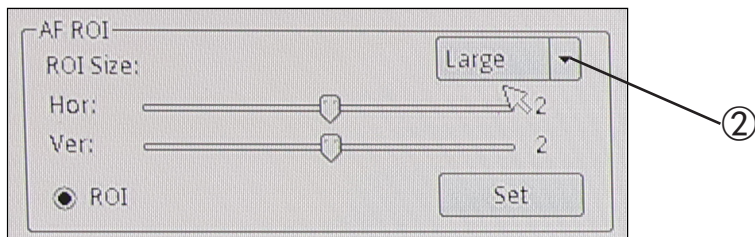
ROI size can be either “Large” or Small”. This can be selected by clicking on the arrow ②.

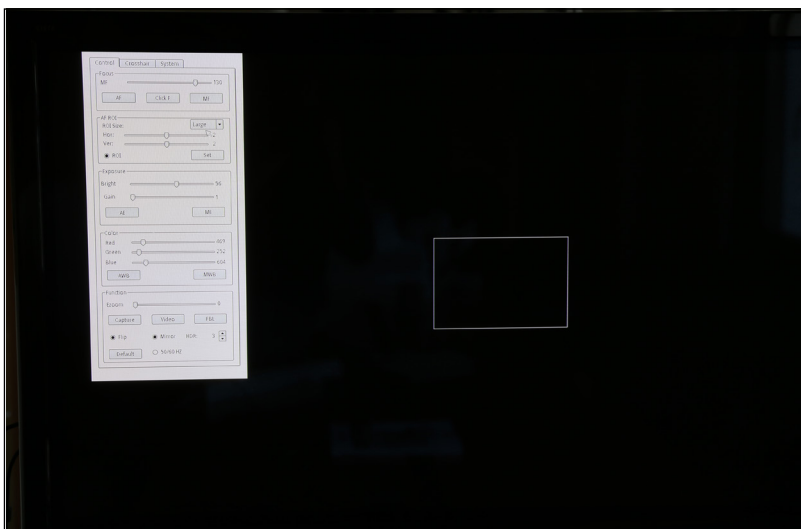A drop down list will appear. Here it is possible to select the size of the ROI area.


LARGE

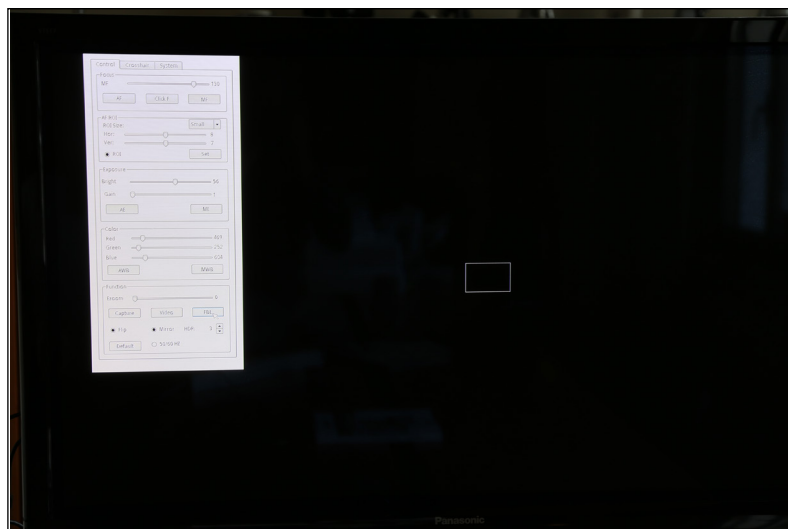
SMALL

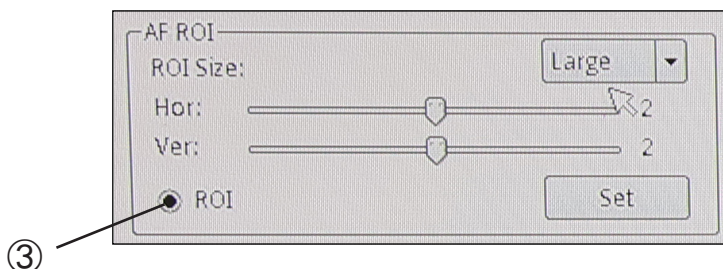Clicking on the radio button “ROI” ③ will activate or deactivate the ROI.

ROI can also be moved horizontally or vertically by using the two sliders “**Hor**” and “**Ver**” ④, in order to move the ROI in the position where the autofocus will be detected.

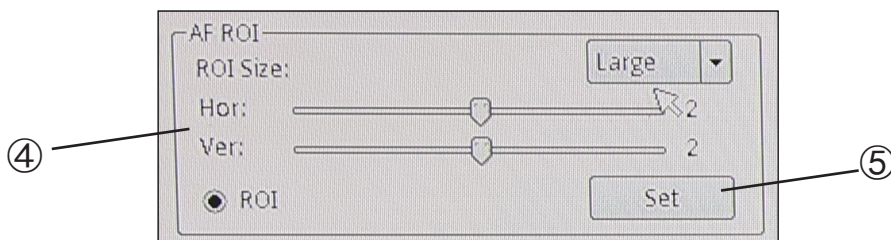
Horizontal movement ranges from 0 to 16, while vertical movement ranges from 0 to 14.

Clicking on “**Set**” ⑤ will apply all the settings.

8.1.3 “Exposure” Section

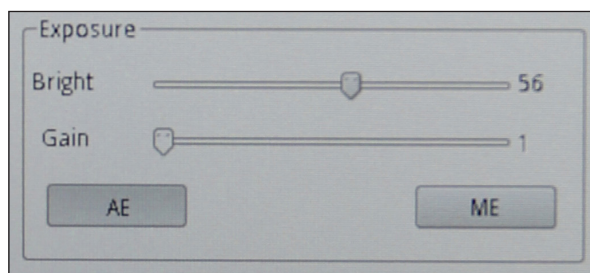“**Exposure**” section controls the exposure time and the brightness of the camera.

The camera can work in automatic exposure.

By activating the “**AE**” function, the camera works completely automatically. In this case it is not possible to act on the parameters of the camera, therefore in situations of insufficient lighting it is possible that the camera can not provide optimal image quality.

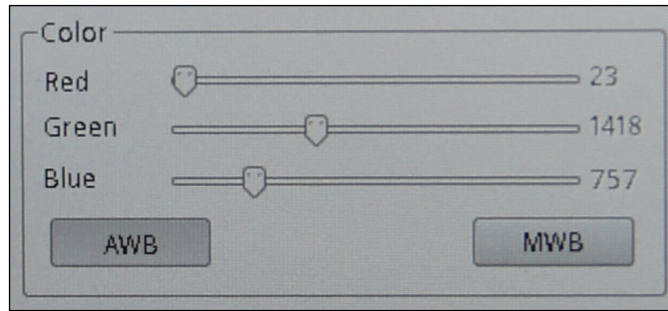
By activating the “**ME**” the camera works manually. It is therefore necessary to act on the camera parameters to adapt the exposure time to the lighting condition of the sample.

Moving the slider relative to “**Bright**” increases or decreases the exposure time of the camera. Depending on the amount of light that is used to illuminate the sample under examination, you will have to move to the right (increase) or to the left (decrease) the cursor to get a perfectly lit image avoiding under or overexposure.

The cursor relative to the “**Gain**” commands the so-called “gain” of the camera. The gain is a function that allows you to increase the sensitivity of the camera for shooting in low light. This obviously increases the background noise.

8.1.4 “Color” Section

The “**Color**” section controls the color rendering of the camera.


Activating the “**AWB**” function, the camera uses automatic white balance. Place an empty slide on the microscope front of the camera sensor and click “**AWB**”. The camera adjusts the colors output from the camera to the color temperature of the scene being shot.

Activating the “**MWB**” function, the camera uses manual white balance. In this case it is necessary to move the sliders related to Red, Green and Blue to obtain a faithful reproduction of what is being observed.

8.1.5 “Function” Section

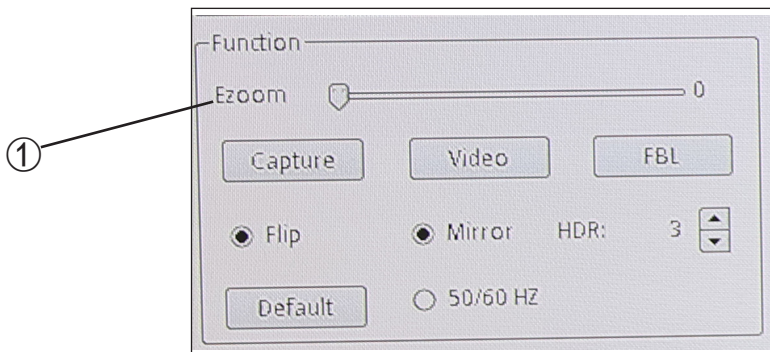
“**Function**” section controls the image display on the monitor.

The “**Ezoom**” cursor ① allows to vary the magnification of the projection to the monitor of the framed object.

It is possible to change the zoom factor (optical zoom) from 0 to 50 operating on the on the cursor in the on-screen panel.

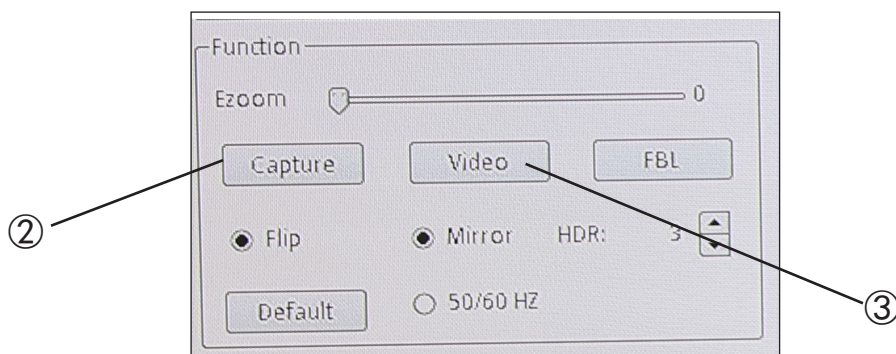Camera allows to capture single images or videos simply by clicking on the “**Capture**” ② button for single images or on “**Video**” ③ for film sequences of desired length.


Still images or video will be stored on the SD card.

To view images or video the user must remove the SD card, simply pushing down the card. The internal locking mechanism will release the SD card that can be freely removed from its slot.


To re-insert the SD card: fully insert the card and push it downward until the locking mechanism is engaged.

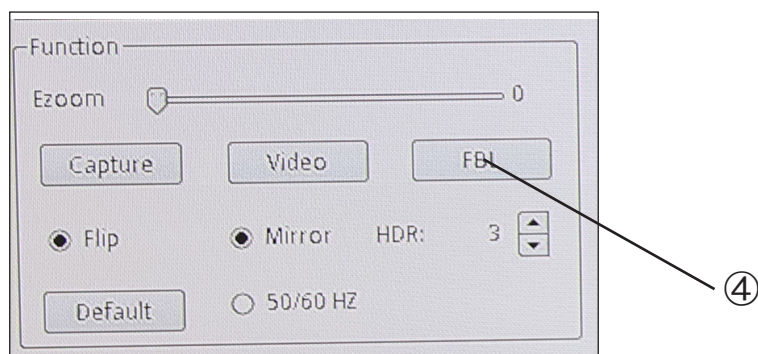“**FBL**” button ④ fully disable the autofocus function of the camera, putting back the sensor motor to the so-called “zero position”.


By clicking “**AF**” button or “**MF**” button again, the camera will allow to obtain an automatic or manual focus.

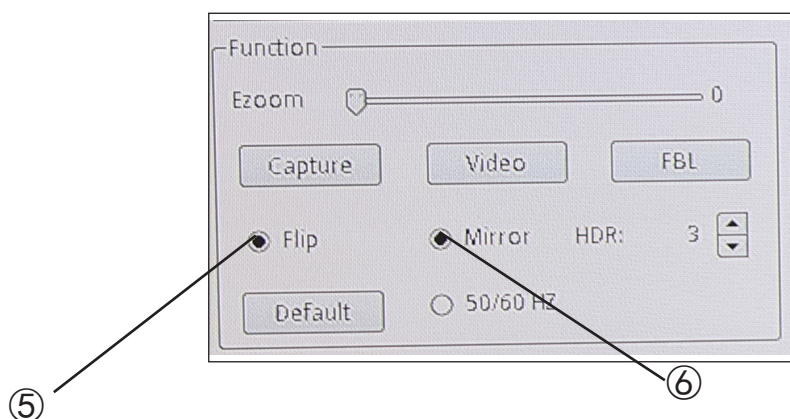The “**Flip**” ⑤ function flips the framed image vertically, while the “**Mirror**” ⑥ function horizontally flips it. This is useful for making what is observed in the camera consistent with the real object.

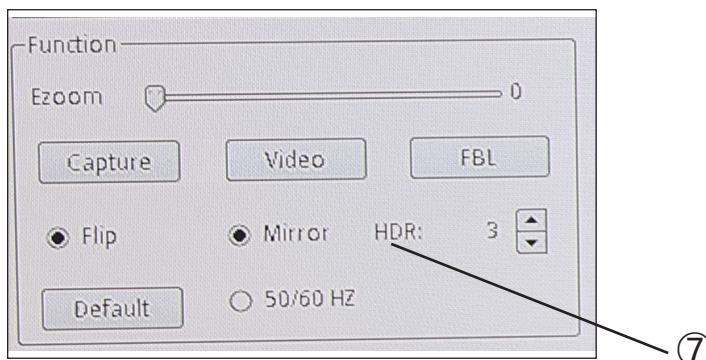
The “**HDR**” function ⑦ (High Dynamic Range) optimizes the image quality in the case of subjects with very clear framed areas together with very dark areas. Often in these conditions the camera favors or clear areas (omitting the dark areas making them black) or dark areas (with a consequent overexposure of the light areas, making them almost white).


An “**HDR**” setting of 0 indicates that the camera will acquire a single image and store it. Increasing the number (max 3) the camera acquires 3 images at different exposures (usually an underexposure, one with optimal exposure and one overexposed), combines them to optimize the different zones with different lighting and creates a single image at exposure “virtually correct”.

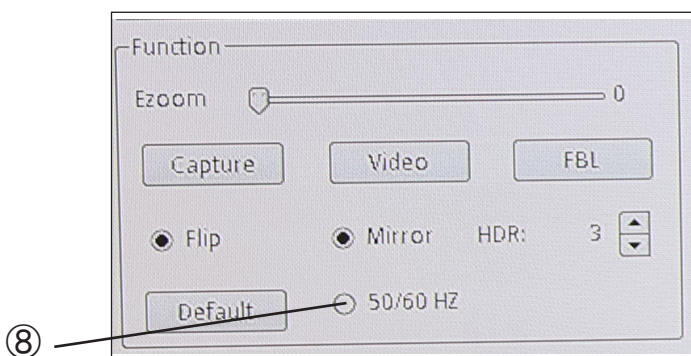
The “**50 / 60Hz**” function ⑧ is used to eliminate flickering of the image due to the power supply frequency.

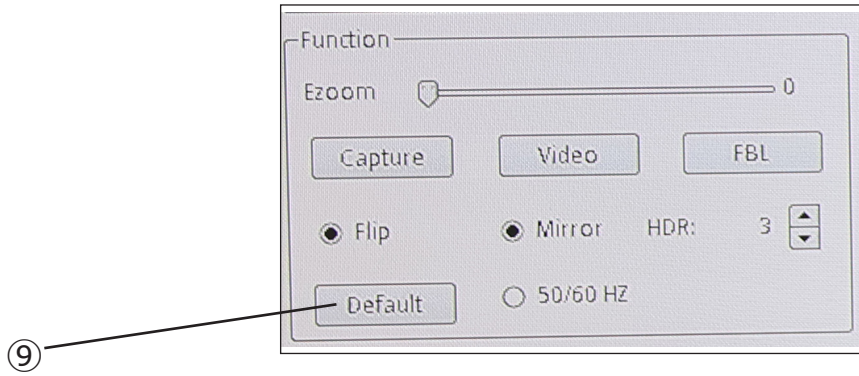The **“Default”** key ⑨ resets all the settings made and restores the factory settings.

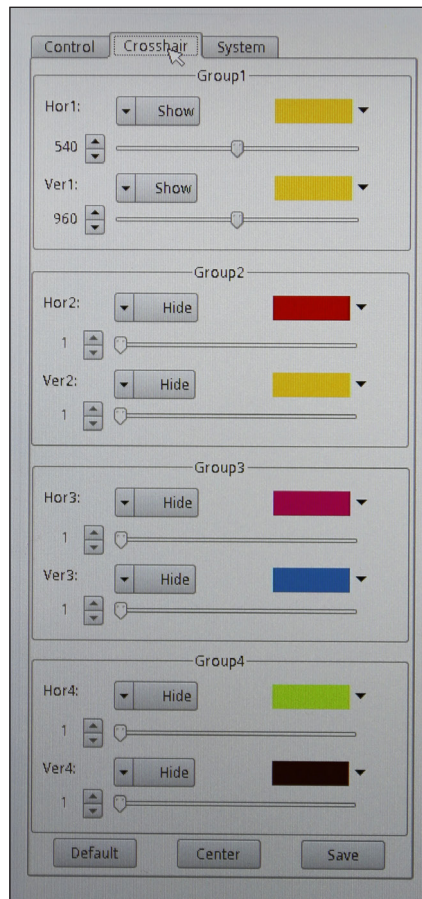8.2 “Crosshair” Menu


Up to 4 sets of horizontal or vertical rows can be superimposed on the monitor.


These lines can be placed anywhere on the monitor by simply dragging the cursor relative to the desired line.

The lines can be projected with different colors in order to be easily visible on the monitor, depending on the color of the sample in question.

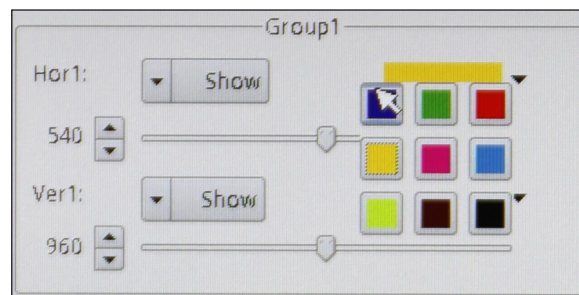Clicking with the mouse on the color palette of the desired row opens the context menu, with which you select the color of choice.

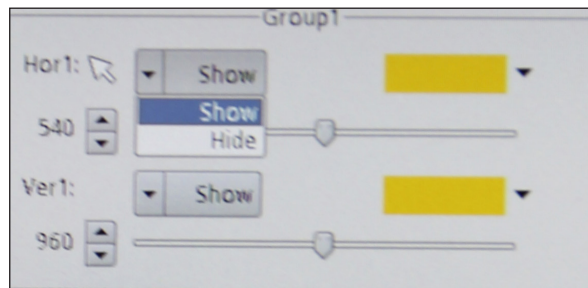The lines, once traced, can be visible or hidden. Clicking on the “**Show / Hide**” button opens the context menu with which you can view a “**Show**” row or hide it “**Hide**”. The lines, if hidden, will not be deleted from memory but simply will not be visible.


The “**Default**” key resets all the settings made and restores the factory settings.

The “**Center**” button places the lines exactly in the center of the monitor (possibly overlapping the different lines drawn).

The “**Save**” button saves the settings of the placed rows.

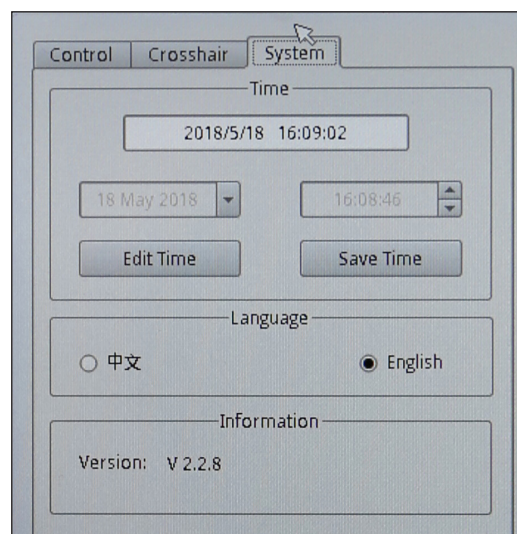8.3 “System” Menu

With this menu, the system information of the camera is changed.

The “**Edit Time**” button changes the current date and time, while the “**Save Time**” button saves the settings made just recently.

It is possible to change the menu language from English to Chinese.

The “**Information**” field displays the current version of the software.

9. Maintenance

To think about when and after using the system

- The system should always be kept vertically when moving it and be careful so that no moving parts, such as the eyepieces, fall out.
- Never mishandle or impose unnecessary force on the device.
- Never attempt to service the device yourself.
- After use, cover the microscope with the provided dust-cover, and keep it in a dry and clean place.

Electrical safety precautions

- Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off-position.
- Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

Cleaning the optics

- If the optical parts need to be cleaned try first to: use compressed air.
- If that is not sufficient: use a soft lint-free piece of cloth with water and a mild detergent.
- And as a final option: use the piece of cloth moistened with a 3:7 mixture of ethanol and ether.
- Note: ethanol and ether are highly flammable liquids. Do not use them near a heat source, near sparks or near electric equipment. Use these chemicals in a well ventilated room.
- Remember to never wipe the surface of any optical items with your hands. Fingerprints can damage the optics.
- Do not disassemble objectives or eyepieces in attempt to clean them.

For the best results, use the OPTIKA cleaning kit (see catalogue).

If you need to send the microscope to Optika for maintenance, please use the original packaging.

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.