HDMI Cameras – H Series

INSTRUCTION MANUAL

Model
C-HE
C-HESC
C-WH5
C-WH5SC

Ver. 4.1 2021

Table of Contents

1.	Warning	3
2.	Safety Information	3
3.	Package content	4
3.1	C-HE	4
3.2	C-HESC	4
3.3	C-WH5	5
3.4	C-WH5SC	5
4.	Unpacking	6
5.	Intended use	6
6.	Symbols and conventions	6
7.	C-HE Basic features	7
7.1	Quick Instructions for C-HE camera	7
7.1.1	HDMI Mode	7
8.	C-HE User Interface (UI) and Its Functions	8
8.1	The Camera Control Panel	9
8.2	Icons and Functions of the Camera Control Toolbar	9
8.3	The Measurement Toolbar	10
8.4	Camera calibration	11
9.	C-WH5 Basic features	13
9.1	Quick Instructions for C-WH5 camera	13
9.1.1	WiFi mode ProView controlled	14
9.1.2	WiFi mode App controlled	14
9.1.3	HDMI mode	14
10.	C-WH5 User Interface (UI) and Its Functions	15
10.1	The Camera Control Panel	16
10.2	Icons and Functions of the Camera Control Toolbar	17
10.3	The Measurement Toolbar	22
10.4	Camera calibration	23
11.	C-HESC / C-WH5SC	25
11.1	Instrument description	25
11.2	Assembling	26
12.	C-HESC / C-WH5SC Use of the system	27
13.	Micrometric Slide M-005	27
	Equipment disposal	28

1. Warning

This device is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.

2. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

3. Package content

3.1 C-HE

- ① Camera
- ② HDMI cable
- ③ Wireless mouse

- ④ SD card 16GB
- ⑤ Micrometric slide
- ⑥ Power supply

3.2 C-HESC

- ① C-HE camera
- ② HDMI Monitor
- ③ Mounting bracket
- ④ HDMI cable

- ⑤ Wireless mouse
- ⑥ SD card 16GB
- ⑦ Micrometric slide
- ⑧ Power supply (2 pcs)

3.3 C-WH5

- ① Camera
- ② WIFI adapter
- ③ HDMI cable
- ④ Wireless mouse

- ⑤ Micrometric slide
- ⑥ SD card 16GB
- ⑦ Power supply

3.4 C-WH5SC

- ① C-WH5 camera
- ② WiFi adapter
- ③ HDMI Monitor
- ④ Mounting bracket
- ⑤ HDMI cable

- ⑥ Wireless mouse
- ⑦ SD card 16GB
- ⑧ Micrometric slide
- ⑨ Power supply (2 pcs)

4. Unpacking

The system is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items fall out and get damaged.

5. Intended use

Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

6. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

7. C-HE Basic features

C-HE is a HDMI + SD card CMOS camera.
HDMI is used as the data transfer interface to HDMI display.

For HDMI output, the Camera Control Panel+ Measurement Toolbar and Camera Control Toolbar are overlaid on the HDMI screen when the mouse move to the related region.

The USB mouse is used to set the camera.

The C-HE's basic features are as follows:

For HDMI output:

- All in 1(HDMI + SD card) C-mount camera with Aptina high sensitivity CMOS sensor.
- Resolution 2MP (720p):
- Built-in mouse control.
- Built-in image capture to SD card.
- Built-in camera control panel, including exposure (manual/auto) / gain, color adjustment, sharpness control.
- Built-in toolbar including zoom, mirror, freeze, crosshair:
- Built-in toolbar for measurement.

7.1 Quick Instructions for C-HE camera

Before turning on the camera, please connect the standard C-mount camera to the camera adapter and insert it into the microscope's trinocular port.

Figure 1 - C-HE Rear Cover

- | | |
|---------------------|-----------------|
| ① HDMI plug | ④ SD card slot |
| ② USB mouse plug | ⑤ LED indicator |
| ③ Power supply plug | |

7.1.1 HDMI Mode

1. Plug the HDMI cable into the HDMI port ① to connect the C-HE camera to HDMI display.
2. Plug a USB mouse into USB Mouse port ② to get control of the camera by using built-in software.
 - **The camera cannot be connected to a PC in any way. The available USB port is used to connect the mouse.**
3. Insert SD card into SD Card Slot ④ for saving captured images.
4. Plug 12V/1A power adapter into DC 12V slot ③ to supply power for the camera. The LED Indicator ⑤ will turn into blue.
5. Move mouse cursor to the left side of the video window, a *Camera Control Panel* will appear. It includes Manual/ Automatic Exposure, White Balance, Sharpness and other functions, please refer to 8.1 for details.
6. Move mouse cursor to the bottom of the video window and a *Camera Control Toolbar* will appear. Operations like Zoom In, Zoom Out, Flip, Freeze, Cross Line, and etc. can be achieved. Please refer to 8.2 for details.
7. Move mouse cursor to the upper side of the video window, a *Measurement Toolbar* with calibration and other measurement tools will appear, please refer to 6.3 for details. The measurement data can be output with *.CSV format.

8. C-HE User Interface (UI) and Its Functions

The C-HE UI shown in Fig. 2 includes a *Camera Control Panel* on the left side of the video window, a *Camera Control Toolbar* on the bottom of the video window and a *Measurement Toolbar* on the top of the video.

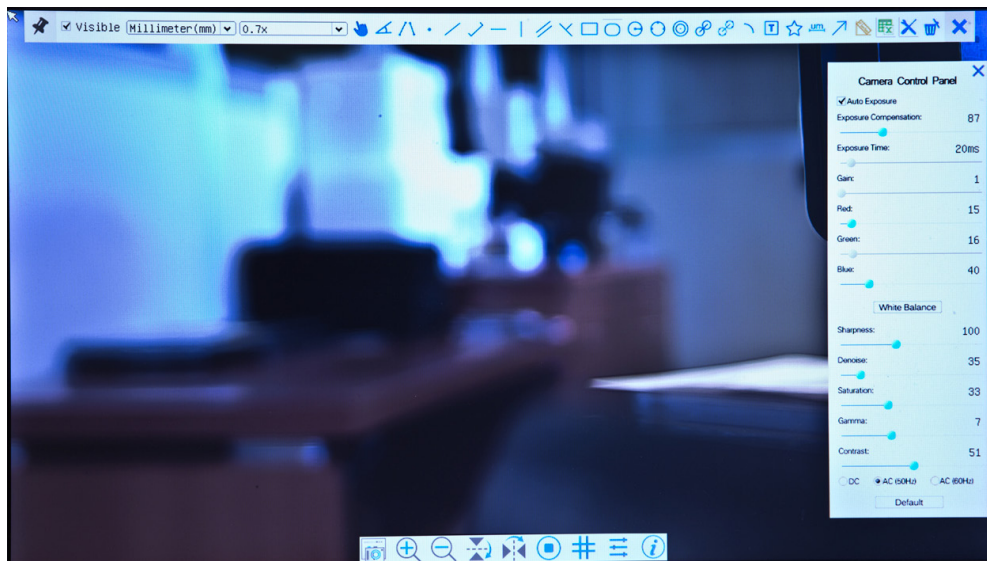
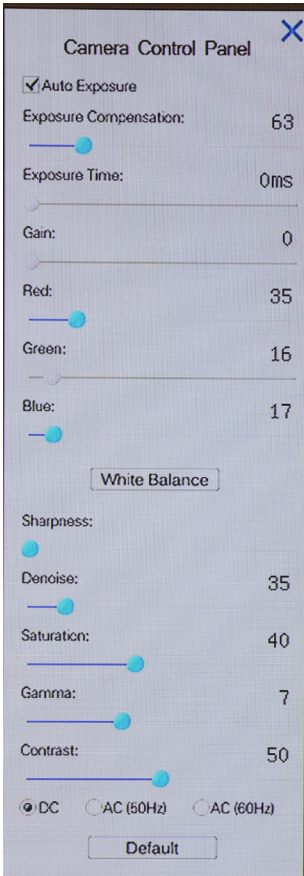
Figure 2 - Camera Control User Interface

Notes:

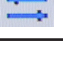
1. When users click on the  button, the *Camera Control Panel* will pop up automatically.
2. When users move mouse cursor to the bottom of the video window, the *Camera Control Toolbar* will pop up automatically.
3. Move the mouse cursor to the upper side of the video window: a *Measurement Toolbar* will pop up for the calibration and measurement operations. When user left-clicks the *Float/Fixed* button  on the Measurement Toolbar, the Measurement Toolbar will be fixed. In this case the Camera Control Panel will not pop up automatically even if users move mouse cursor to left side of the video windows. Only when user left-clicks the  button on the Measurement Toolbar to exit from measuring procedure it will be possible to perform other operations on the *Camera Control Panel*.

8.1 The Camera Control Panel

The Camera Control Panel controls the camera to achieve the best image quality according to the specific applications. It will pop up automatically when mouse cursor is moved to the left side of the video window.

Camera Control Panel	Function	Description
	Auto Exposure	When Auto Exposure is checked, the system will automatically adjust exposure time according to the value of exposure compensation
	Exposure Compensation	Available when Auto Exposure is selected. Slide left or right to adjust the exposure compensation according to the current video brightness to get the correct brightness value.
	Exposure Time	Available when Auto Exposure is unchecked. Slide to left or right to reduce or increase exposure time, adjusting brightness of the video
	Gain	Adjust Gain to reduce or increase brightness of video. The Noise will be reduced or increased accordingly
	Red	Slide to left or right to decrease or increase the proportion of Red in RGB on video
	Green	Green is base for reference and cannot be adjusted
	Blue	Slide to left or right to decrease or increase the proportion of Blue in RGB on the video
	White Balance	Adjusting the White Balance according to the monitor display
	Sharpness	Adjust Sharpness level of the video window
	Denoise	Adjust Noise level of the video window
	Saturation	Adjust Saturation level of the video window
	Gamma	Adjust Gamma level of the video. Slide to the right side to increase gamma and to the left to decrease gamma
	Contrast	Adjust Contrast level of the video. Slide to the right side to increase contrast and to the left to decrease contrast
	DC	For DC illumination, there will be no fluctuation in light source so no need for compensating light flickering
	AC(50HZ)	Check AC(50HZ) to eliminate flickering "strap" caused by 50Hz illumination
AC(60HZ)	Check AC(60HZ) to eliminate flickering "strap" caused by 60Hz illumination	
Default	Set all the settings in the Camera Control Panel to default values	

8.2 Icons and Functions of the Camera Control Toolbar

Icon	Function	Icon	Function
	Snap the image and save it to SD card		Horizontal Flip
	Zoom In the Video Window		Video Freeze
	Zoom Out the Video Window		Display Cross Line
	Vertical Flip		Toggle ON/OFF Camera Control Panel

8.3 The Measurement Toolbar

The *Measurement Toolbar* will pop up when moving mouse cursor to any place near the upper side of the video window.

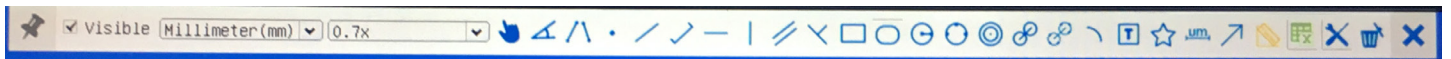
Figure 3 - The Measurement Toolbar

Icon	Function
	Float/ Fix switch of the Measurement Toolbar
<input checked="" type="checkbox"/> Visible	Define measuring object in Show up/ Hide mode
Nanometer (nm)	Select the desired Measurement Unit
4x	Choose the same Magnification as the microscope to ensure accuracy of measurement result when measurement unit is not in Pixel units
	Object Select
	Point
	Angle
	4 point angle
	Arbitrary Line
	3 points Line
	Vertical 3 points line
	Parallel
	Horizontal Line
	Vertical Line
	Rectangle
	Circle
	3 points Circle
	Ellipse
	Annulus
	3 points Circle and Center Distance
	Two Circles and Center Distance
	Arc
	Polygon

	Curve
	Arrow
	Ruler
	Make Calibration to determine the corresponding relation between magnification and resolution, this will establish the corresponding relationship between measurement unit and the sensor pixel size. Calibration needs to be done with the help of a micrometer. For detailed steps of carrying out calibration please refer to chapter 8.4
	Export the measurement information to CSV file(*.csv)
	Delete All the Measurement Objects
	Settings
	Exit from Current Measurement Mode

Notes:

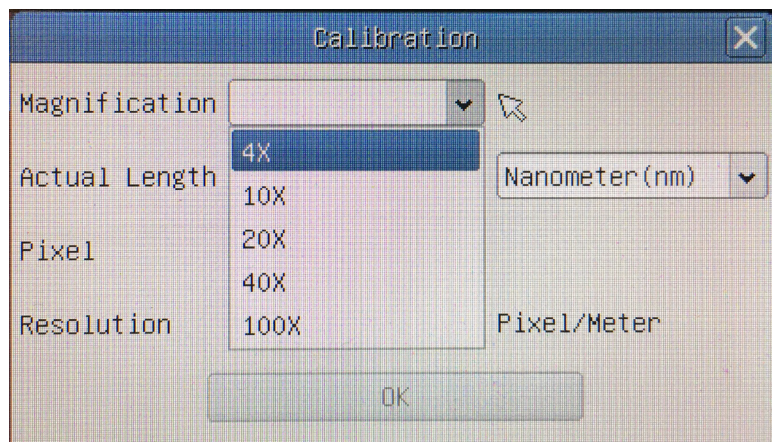
1. When user left-clicks *Display/Hide* button  on the *Measurement Toolbar*, the *Measurement Toolbar* will be fixed. In this case the *Camera Control Panel* will not pop up automatically even if moving mouse cursor to the left side of the video window. Only when users left click the  button on the *Measurement Toolbar* to exit from the measurement mode they will be able to do other operations in the *Camera Control Panel*, the *Auto Focus Control Panel* or the *Synthesis Camera Control Toolbar*.

8.4 Camera calibration


The camera must be calibrated before any measurements can be taken. To do this it is necessary to use the micrometer slide supplied with the camera.

Detailed procedure is the following:


1. Insert the objective with the lowest magnification available and put the stage micrometer on the stage.
2. Click on the “Calibrate”  button in the *Measurement toolbar*.
3. The following dialog box appears:

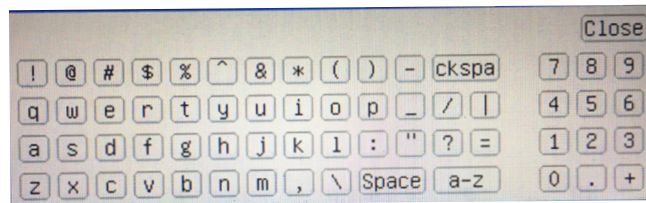4. In the **Magnification** field enter the current magnification.
 - Some magnification values (default ones 4x, 10x, 20x, 40x and 100x) are preset in the drop-down list.
 - If user must insert a different magnification value, double click in the **Magnification** field. The following dialog box appears:

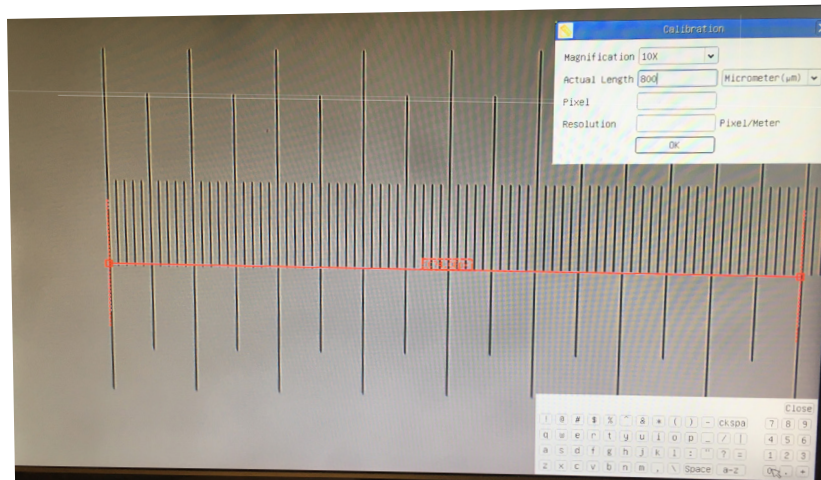- From here user can enter any desired magnification value.
5. Focus the ruler under the microscope.
 - Along with the previous dialog box a red line appears on the image. This line is the calibration line.

6. Align the two ends of the red line with the ruler scale (try to use as many notches as possible to keep the calibration accuracy).
7. Set the desired unit (micrometer, millimeter or nanometer) in the **Calibration** dialog box.
8. Insert the length of the drawn line in the **Actual Length** field (using the same procedure described above).
 - Micrometric slide has two rulers (see chapter 11). One is 1 mm/100 (for biological microscopes) and the other is 10mm/100 (for stereomicroscopes).
 - When using the ruler 1mm/100 the interval between two long lines is 100 μm , while using the ruler 10mm/100 the interval between two long lines is 1000 μm .
9. If everything is fine, click **OK** to end the calibration. The new magnification (the number you entered in the Magnification edit box) will be available in the Magnification drop-down list.
10. Repeat the same procedure for each magnification available in your microscope.

9. C-WH5 Basic features

C-WH5 is a multiple interfaces (HDMI + WiFi + SD card) CMOS camera. HDMI and WIFI are used as the data transfer interface to HDMI display or computer.

For HDMI output, the *Camera Control Panel + Measurement Toolbar* and *Camera Control Toolbar* are overlaid on the HDMI screen when the mouse move to the related region. in this case, the USB mouse can be used to set the camera, browse and compare the captured image, play the video and perform the measurement.

In the WiFi interface, connect the WiFi adapter to the camera's USB port. The camera generates a WiFi signal that is recognized by a computer with a WLAN connection or by a mobile device (IOS or Android).

The camera can be managed via a PC using the ProView software. With ProView, you can control the camera, process video and images in the same way as with all other cameras in the OPTIKA series.

Managing the camera via a mobile device requires an App to be downloaded on your device, allowing multiple users to view and capture images from the camera.

The C-WH5's basic features are as follows:

For HDMI output:

- C-mount camera with Sony high sensitivity CMOS sensor.
- Resolution 2MP (1080p).
- HDMI output.
- Built-in mouse control.
- Built-in image capture & video record to SD card.
- Built-in camera control panel, including exposure (manual/auto) / gain, white balance (lockable), color adjustment, sharpness control.
- Built-in toolbar for measurement.
- Built-in toolbar including zoom, mirror, comparison, freeze, cross, browser function.
- Built-in image & video browsing, display & play.
- Real time clock (RTC).

For WiFi output:

- Ultra-Fine color engine with perfect color reproduction capability.
- Resolution 5MP.
- With advanced video & image processing application ProView, which including professional image processing such as 2D measurement, HDR, image stitching, EDF (Extended Depth of Focus), image segmentation & count, image stacking, color composite and denoising.
- Supports standard UVC protocol for Windows.

9.1 Quick Instructions for C-WH5 camera

Before turning on the camera, please connect the standard C-mount camera to the camera adapter and insert it into the microscope's trinocular port.

Figure 4 - C-WH5 Rear Cover

- | | |
|------------------------------------|-----------------|
| ① HDMI connector | ④ SD card slot |
| ② USB mouse/WiFi adapter connector | ⑤ ON-OFF button |
| ③ Power supply plug | ⑥ LED indicator |

9.1.1 WiFi mode ProView controlled

1. Connect the 12V/1A power cable to the DC 12V connector ③ to power the camera. The ⑥ LED indicator will turn red.
2. Connect the provided WiFi adapter to the WIFI/USB Mouse port ②.
3. Press the ON/OFF button ⑤ to turn on the camera. the LED indicator ⑥ will turn blue.
4. Start the computer (equipped with WiFi connection) and in the “WiFi” settings of the computer select **C-WH5-xxxxx**.
 - The password for connecting the camera is “12345678”.
5. Start the ProView software, activate C-WH5 by clicking the camera name in the Camera List. For more details, refer to the ProView instruction manual.

9.1.2 WiFi mode App controlled

1. Connect the 12V/1A power cable to the DC 12V connector ③ to power the camera. The ⑥ LED indicator will turn red.
2. Connect the provided WiFi adapter to the WIFI/USB Mouse port ②.
3. Press the ON/OFF button ⑤ to turn on the camera. the LED indicator ⑥ will turn blue.
4. In the Settings of your mobile device select the WiFi device **C-WH5-xxxxx**.
 - The password for connecting the camera is “12345678”.
5. Download and launch the free App (available on AppStore for IOS and PlayStore for Android).
6. Manage the camera using the App to control all functions (Exposure Control, White Balance, Image or Movies Acquisition).

NOTE: The maximum number of users who can connect simultaneously is limited to 3. This is to ensure that the camera does not lose speed when many users are connected simultaneously. When 3 users are connected, the camera will not allow a fourth user to connect.

9.1.3 HDMI mode

1. Plug the HDMI cable into the HDMI port ① to connect the C-WH5 camera to HDMI display.
2. Plug a USB mouse into USB Mouse port ② to get control of the camera by using built-in software.
3. Plug 12V/1A power adapter into DC 12V slot ③ to supply power for the camera. The LED Indicator ⑥ will turn into red.
4. Insert SD card into SD Card slot ④ for saving captured images and recorded video.
5. Press ON/OFF Button ⑤ to turn on the camera. The LED indicator ⑥ will turn into blue.
6. Move mouse cursor to the left side of the video window, a *Camera Control Panel* will appear. It includes Manual/ Automatic Exposure, White Balance, Sharpness and other functions, please refer to 10.1 for details.
7. Move mouse cursor to the bottom of the video window and a *Camera Control Toolbar* will appear. Operations like Zoom In, Zoom Out, Flip, Freeze, Cross Line, Comparison and etc. can be achieved. Please refer to 10.2 for details.
8. Move mouse cursor to the upper side of the video window, a *Measurement Toolbar* with calibration and other measurement tools will appear, please refer to 10.3 for details. The measurement data can be output with *.CSV format.

10. C-WH5 User Interface (UI) and Its Functions

The C-WH5 UI shown in Fig. 5 includes a *Camera Control Panel* on the left side of the video window, a *Measurement Toolbar* on the upper side of the video window, a *Camera Control Toolbar* on the bottom of the video window.


Figure 5 - Camera Control User Interface


Notes:

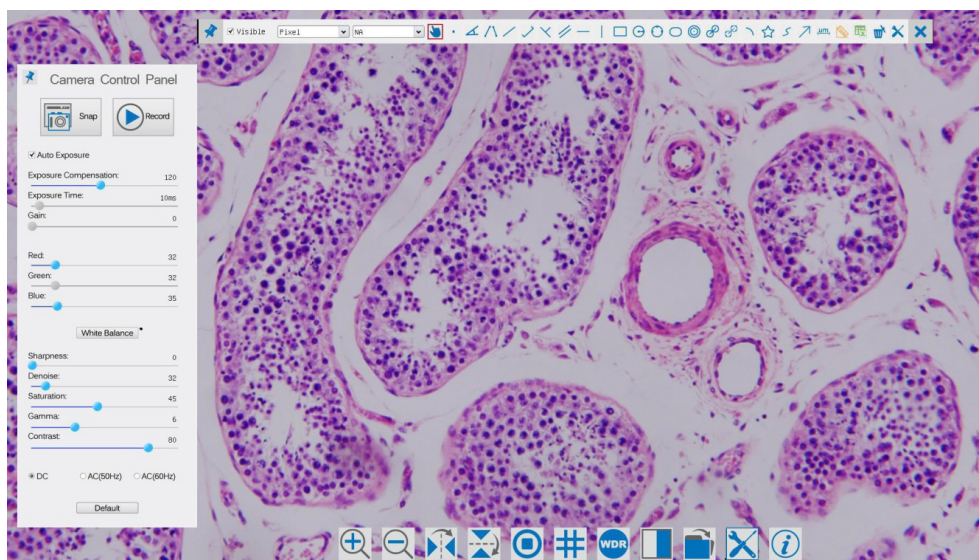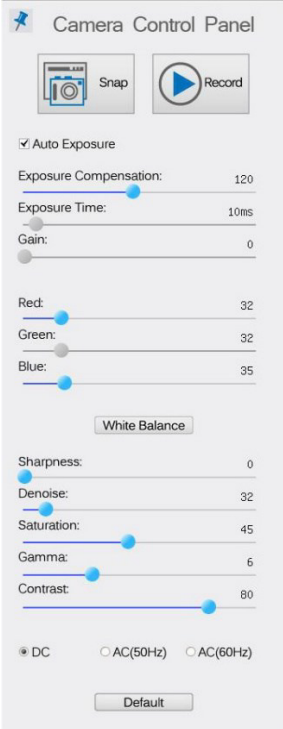1. When users move mouse cursor to the left side of the video window, the *Camera Control Panel* will pop up automatically.
2. When user moves mouse cursor to the bottom of the video window, the *Camera Control Toolbar* will pop up automatically.
3. Move the mouse cursor to the upper side of the video window: a *Measurement Toolbar* will pop up for the calibration and measurement operations. When user left-clicks the *Float/Fixed* button  on the Measurement Toolbar, the Measurement Toolbar will be fixed. In this case the Camera Control Panel will not pop up automatically even if users move mouse cursor to left side of the video windows. Only when user left-clicks the  button on the Measurement Toolbar to exit from measuring procedure it will be possible to perform other operations on the *Camera Control Panel*, or *Synthesis Camera Control Toolbar*. During the measuring process, when a specific measuring object is selected an *Object Location & Attributes Control Bar*  will appear for changing location and properties of the selected objects.

10.1 The Camera Control Panel

The *Camera Control Panel* controls the camera to achieve the best image quality according to the specific applications. It will pop up automatically when mouse cursor is moved to the left side of the video window (in measurement status, the Camera Control Panel will not pop up. Only when measurement process is terminated will the Camera Control Panel pop up by moving mouse cursor to the left side of the video window).

Left-clicking button  to enable *Display / Auto Hide* switch of the Camera Control Panel.

Control Panel	Function	Description
 <p>The screenshot shows the Camera Control Panel interface. It includes a 'Snap' button, a 'Record' button, and a 'White Balance' button. There are several sliders for 'Exposure Compensation' (120), 'Exposure Time' (10ms), 'Gain' (0), 'Red' (32), 'Green' (32), 'Blue' (35), 'Sharpness' (0), 'Denoise' (32), 'Saturation' (45), 'Gamma' (6), and 'Contrast' (80). At the bottom, there are radio buttons for 'DC', 'AC(50Hz)', and 'AC(60Hz)', and a 'Default' button.</p>	Snap	Capture image from the current video window
	Record	Record video from the current video window
	Auto Exposure	When Auto Exposure is checked, the system will automatically adjust exposure time according to the value of exposure compensation
	Exposure Compensation	Available when Auto Exposure is selected. Slide left or right to adjust the exposure compensation according to the current video brightness to get the correct brightness value.
	Exposure Time	Available when Auto Exposure is unchecked. Slide to left or right to reduce or increase exposure time, adjusting brightness of the video
	Gain	Adjust Gain to reduce or increase brightness of video. The Noise will be reduced or increased accordingly
	Red	Slide to left or right to decrease or increase the proportion of Red in RGB on video
	Green	Green is base for reference and cannot be adjusted
	Blue	Slide to left or right to decrease or increase the proportion of Blue in RGB on the video
	White Balance	Adjusting the White Balance according to the monitor display
	Sharpness	Adjust Sharpness level of the video window
	Denoise	Adjust Noise level of the video window
	Saturation	Adjust Saturation level of the video window
	Gamma	Adjust Gamma level of the video. Slide to the right side to increase gamma and to the left to decrease gamma
	Contrast	Adjust Contrast level of the video. Slide to the right side to increase contrast and to the left to decrease contrast
	DC	For DC illumination, there will be no fluctuation in light source so no need for compensating light flickering
	AC(50HZ)	Check AC(50HZ) to eliminate flickering "strap" caused by 50Hz illumination
	AC(60HZ)	Check AC(60HZ) to eliminate flickering "strap" caused by 60Hz illumination
	Default	Set all the settings in the Camera Control Panel to default values

10.2 Icons and Functions of the Camera Control Toolbar

Icon	Function	Icon	Function
	Zoom In the Video Window		Zoom Out the Video Window
	Vertical Flip		Horizontal Flip
	Video Freeze		Display Cross Line
	Enable Wide Dynamic Range mode		Compare Image with the current video
	Browse Images and Videos in the SD Card		Settings
	Informations on SW version		

The  *Setting* function is relatively more complicated than the other functions. Here are more info about it:

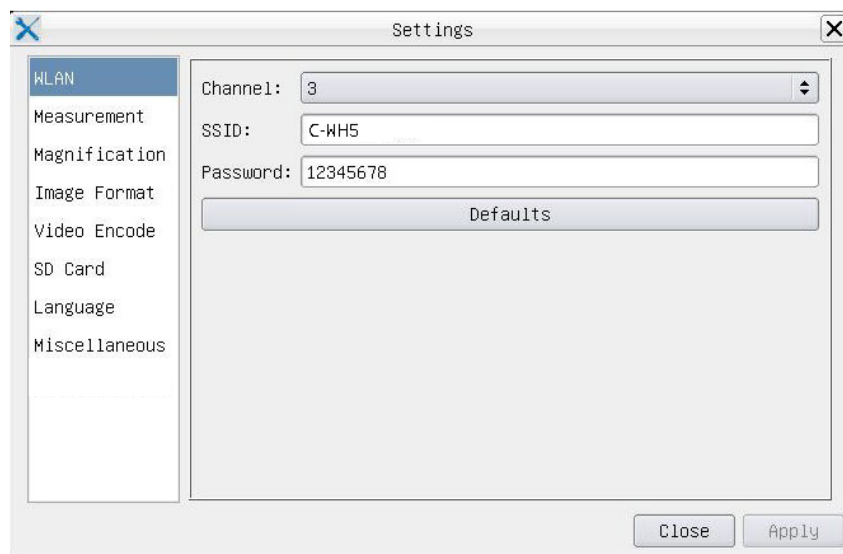
Figure 6 - WIFI Settings Page

Channel: Wireless communication channel, is a data transmission channel with wireless signals as transmission media.

SSID: Camera name sent by WIFI adapter.

Password: WIFI password. default is 12345678.

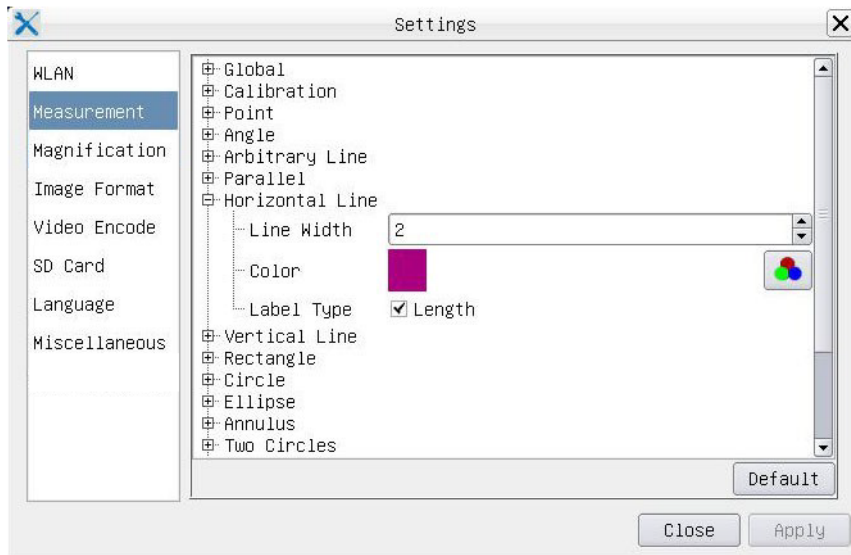
Figure 7 - Measurement Settings Page

Global: Used for setting digits behind the decimal point for measurement results.

Calibration Line Width Used for defining width of the lines for calibration.

Color Used for defining color of the lines for calibration.

EndPoint Type: Used for defining shape of the endpoints of lines for calibration.

Null: Means no endpoints

Rectangle: Means rectangle type of endpoints. It makes alignment more easily

Point, Angle, Line, Horizontal Line, Vertical Line, Rectangle, Circle, Ellipse, Annulus, Two Circles, Polygon, Curve: Left-click the  beside the measuring patterns mentioned above will unfold the corresponding attribute settings to set the individual property of the measuring objects.

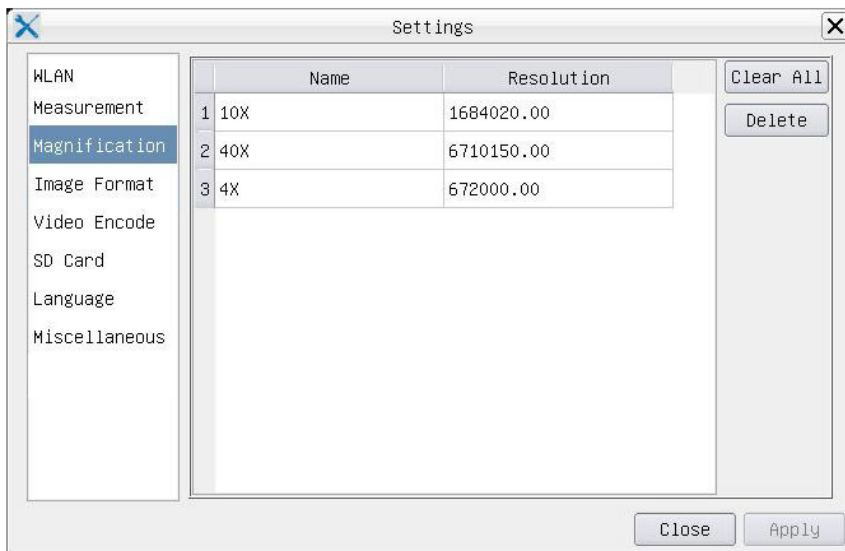
Figure 8 - Magnification Calibration Management Settings Page

Name: Names such as 10X, 40X, 100X are based on magnification of the microscopes. For continuous zoom microscopes, ensure that the selected magnification coincides with the scale alignment line on the microscope zoom knob.

Resolution: Pixels per meter. Devices like microscopes have high resolution value.

Clear All: Click the Clear All button will clear the calibrated magnifications and resolutions.

Delete: Click Delete to delete the selected item for specific resolution.

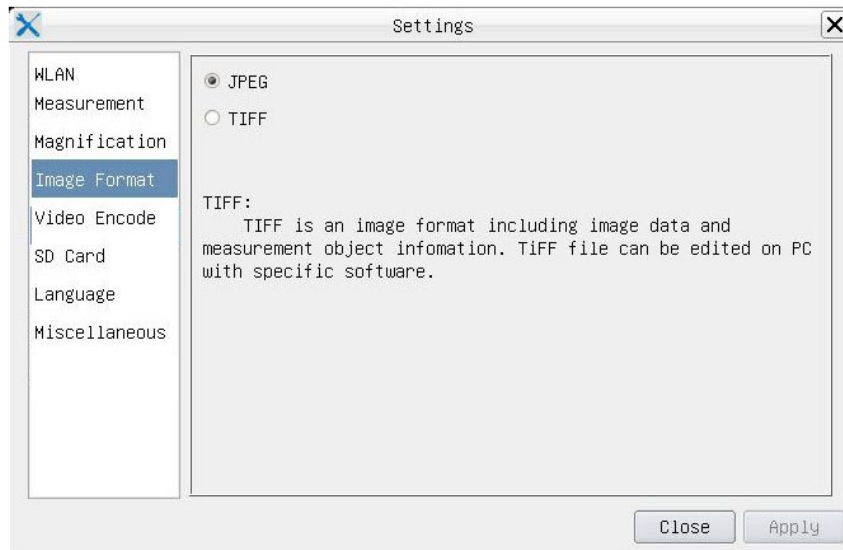
Figure 9 - Image Format Settings Page

JPEG The JPEG file can achieve a very high compression rate and display very rich and vivid images by removing redundant images and color data. In other words, you can get better image quality with minimal disk space. If measuring objects are available, they are burned in the image and the measurement cannot be changed.

TIFF TIFF is an image format that can include both the image data and the measurement object. The measurement objects are saved separately from the image data and can be edited later with specific software.

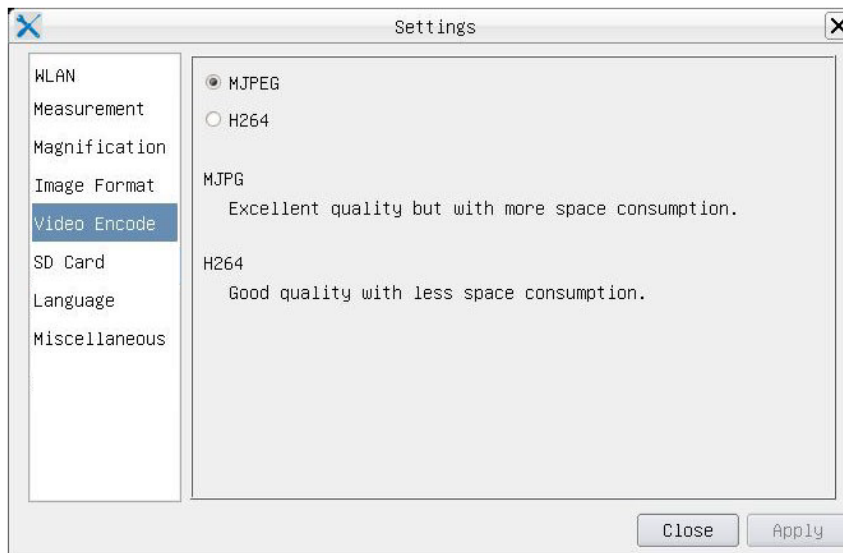
Figure 10 - Video Encode Settings Page

MJPEG Codec format for video is MJPEG. Excellent quality but with more space consumption than H264.

H264 Codec format for video is H264. Good quality with less space consumption.

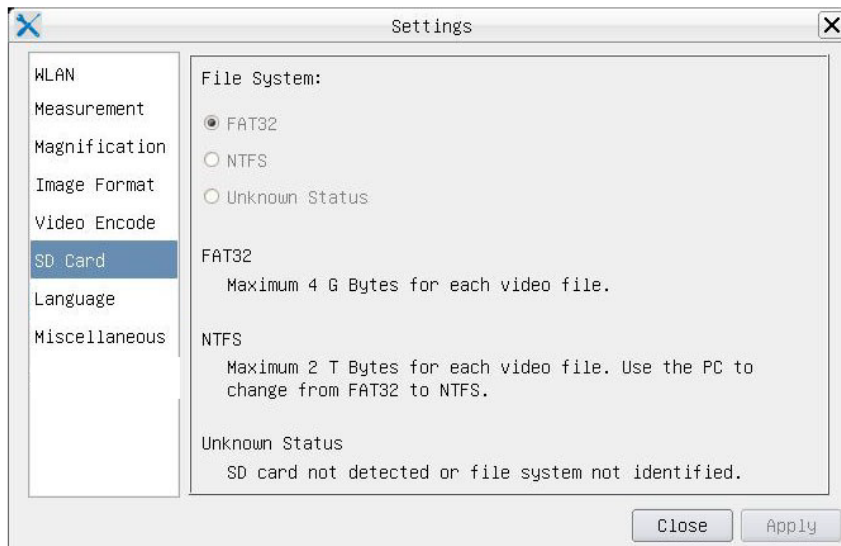
Figure 11 - SD Card Settings Page

- File System* **FAT32:** The file system of SD card is FAT32. The maximum video file size of single file in FAT32 file system is 4GB.
- NTFS:** The file system of SD card is NTFS. The maximum video file size of single file is 2TB. Use PC to format the SD cards and switch between FAT32 and NTFS.
- Unknown Status:** SD card not detected or the file system is not identified.

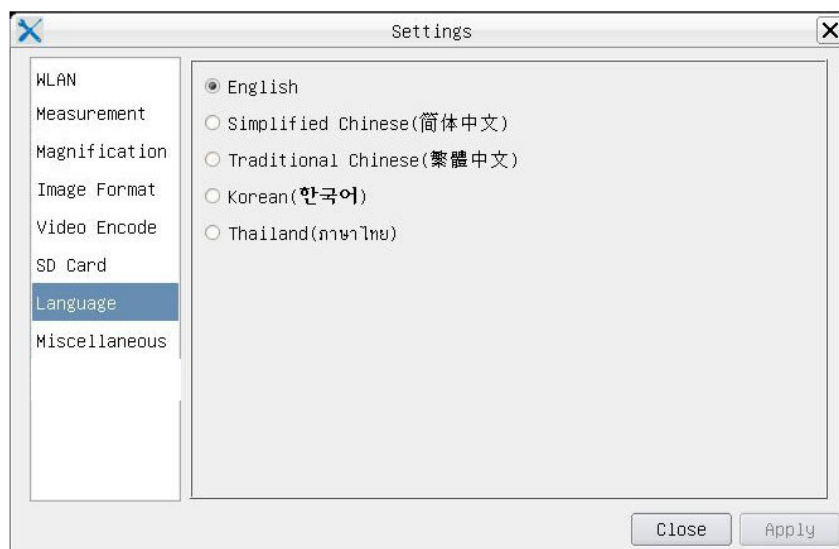
Figure 12 - Language Selection Settings Page

- English:** Set language of the whole software into English.
- Simplified Chinese:** Set language of the whole software into Simplified Chinese.
- Traditional Chinese:** Set language of the whole software into Traditional Chinese.
- Korean:** Set language of the whole software into Korean.
- Thailand:** Set language of the whole software into Thailand.

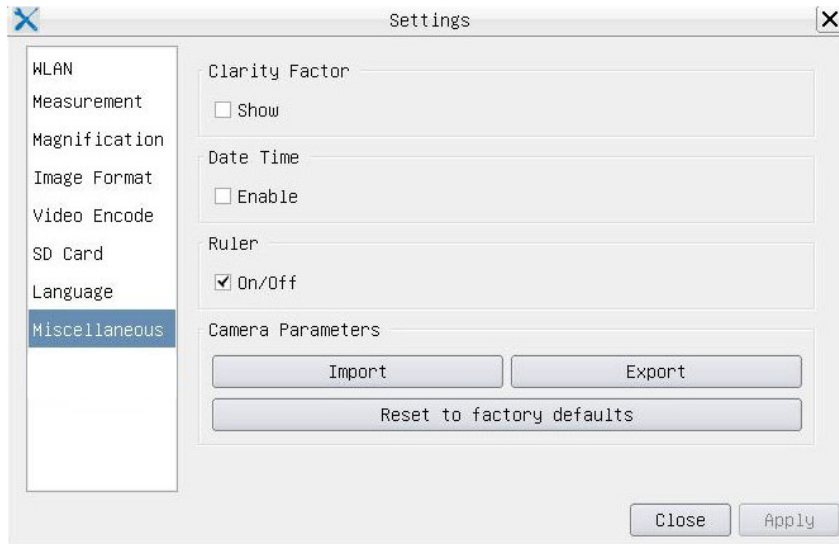
Figure 13 - Miscellaneous Settings Page

- Clarity Factor** The Clarity Factor informs if the sample is in good focus state or not. The larger the Clarity Factor, the better the focus of the sample. Enable *Show* to see if the video is fully focused or not.
- Date Time** For cameras without RTC (Real Time Clock) configuration, users can configure the system date after installation of the camera, which is not saved because there is no built-in RTC.
- Ruler (On/Off)** Enable to show the Ruler on the video window.
- Import** Import parameter settings saved from SD card.
- Export** Export current camera parameters to SD card for later use.
- Reset to factory defaults** Reset camera parameters to factory conditions.

10.3 The Measurement Toolbar

The *Measurement Toolbar* will pop up when moving mouse cursor to any place near the upper side of the video window.

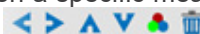
Figure 14 - The Measurement Toolbar

Icon	Function
	Float/ Fix switch of the Measurement Toolbar
<input checked="" type="checkbox"/> Visible	Define measuring object in Show up/ Hide mode
Nanometer (nm) dropdown	Select the desired Measurement Unit
4x dropdown	Choose the same Magnification as the microscope to ensure accuracy of measurement result when measurement unit is not in Pixel units
	Object Select
	Point
	Angle
	4 points angle
	Arbitrary Line
	3 points line
	Vertical 3 points line
	Parallel
	Horizontal Line
	Vertical Line
	Rectangle
	Circle
	3 points circle
	Ellipse
	Annulus
	Two Circles and Center Distance
	3 points circle and e Center Distance
	Arc
	Polygon

	Curve
	Arrow
	Ruler
	Make Calibration to determine the corresponding relation between magnification and resolution, this will establish the corresponding relationship between measurement unit and the sensor pixel size. Calibration needs to be done with the help of a micrometer. For detailed steps of carrying out calibration please refer to chapter 10.4
	Export the measurement information to CSV file(*.csv)
	Delete All the Measurement Objects
	Settings
	Exit from Current Measurement Mode
	When the measurement ends, left-click on a single measuring object and the Object Location & Properties Control Bar will show up. The icons on the control bar mean Move Left, Move Right, Move Up, Move Down, Color Adjustment and Delete

Note:

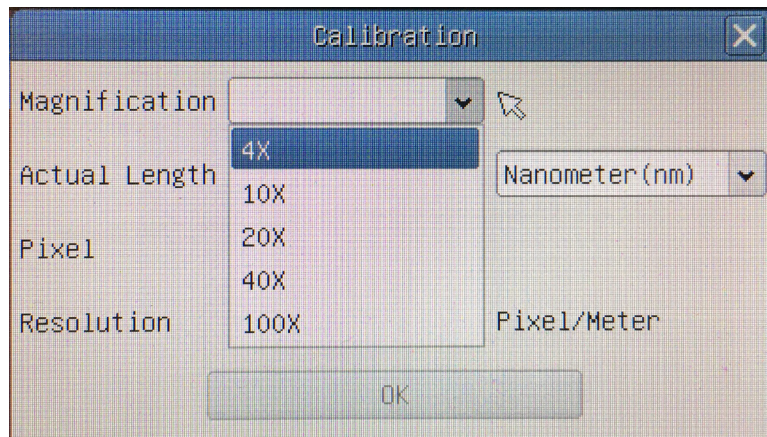
1. When user left-clicks *Display/Hide* button  on the *Measurement Toolbar*, the *Measurement Toolbar* will be fixed. In this case the *Camera Control Panel* will not pop up automatically even if moving mouse cursor to the left side of the video window. Only when users left click the  button on the *Measurement Toolbar* to exit from the measurement mode they will be able to do other operations in the *Camera Control Panel*, the *Auto Focus Control Panel* or the *Synthesis Camera Control Toolbar*.
2. When a specific measuring object is selected during the measuring process, the *Object Location & Attributes Control Bar*  will appear for changing the object location and properties of the selected objects.

10.4 Camera calibration


The camera must be calibrated before any measurements can be taken. To do this it is necessary to use the micrometer slide supplied with the camera.

Detailed procedure is the following:


1. Insert the objective with the lowest magnification available and put the stage micrometer on the stage.
2. Click on the “Calibrate”  button in the *Measurement toolbar*.
3. The following dialog box appears:

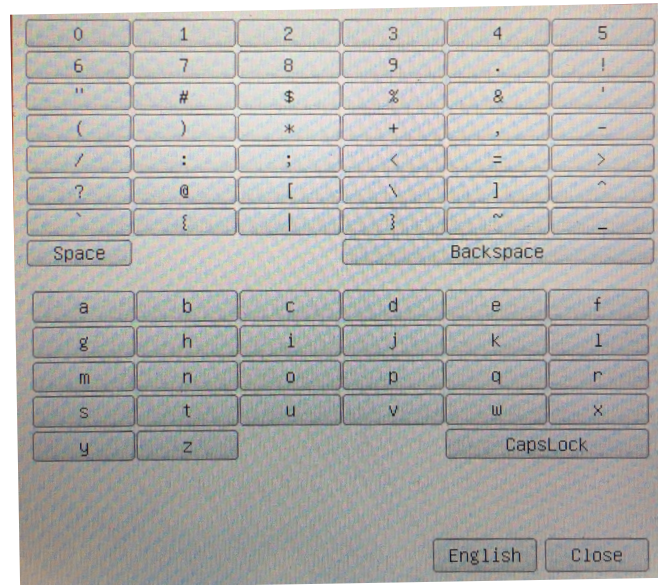4. In the **Magnification** field enter the current magnification
 - Some magnification values (default ones 4x, 10x, 20x, 40x and 100x) are preset in the drop-down list.
 - If user must insert a different magnification value, double click in the **Magnification** field. The following dialog box appears:

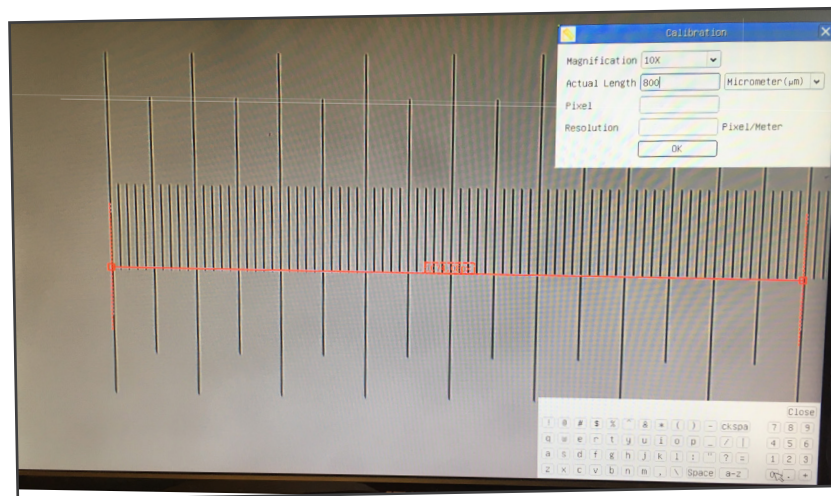- From here user can enter any desired magnification value.
5. Focus the ruler under the microscope.
 - Along with the previous dialog box a red line appears on the image. This line is the calibration line.

6. Align the two ends of the red line with the ruler scale (try to use as many notches as possible to keep the calibration accuracy).
7. Set the desired unit (micrometer, millimeter or nanometer) in the **Calibration** dialog box.
8. Insert the length of the drawn line in the **Actual Length** field (using the same procedure described above).
 - Micrometric slide has two rulers (see chapter 11). One is 1 mm/100 (for biological microscopes) and the other is 10mm/100 (for stereomicroscopes).
 - When using the ruler 1mm/100 the interval between two long lines is 100 μm , while using the ruler 10mm/100 the interval between two long lines is 1000 μm .
9. If everything is fine, click **OK** to end the calibration. The new magnification (the number you entered in the Magnification edit box) will be available in the Magnification drop-down list.
10. Repeat the same procedure for each magnification available in your microscope.

11. C-HESC / C-WH5SC

11.1 Instrument description

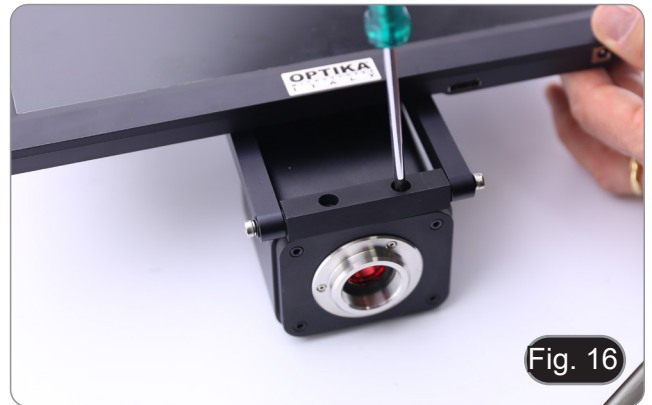
11.2 Assembling

1. Screw the four provided Allen screws on the back side of the monitor. (Fig. 15)

2. Screw the two provided cross screws on the back side of the camera. (Fig. 16)
3. Install the "C" mount and connect HDMI and power cables to monitor and camera.
4. System is ready.


12. C-HESC / C-WH5SC Use of the system

For the operating instructions of the C-HE camera, refer to chapter 8) and for the operating instructions of the C-WH5 camera, refer to chapter 10).


13. Micrometric Slide M-005

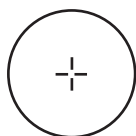Micrometric slide, 26x76mm, with 2 scales
(1mm/100div. for biological microscopes / 10mm/100div. for stereomicroscopes)

1 DIV=0.01mm

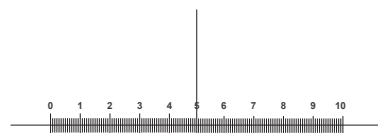For biological microscopes calibration

1 DIV=0.1mm

For stereo microscopes calibration

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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3.1 C-HE

- ① Telecamera
- ② Cavo HDMI
- ③ Mouse wireless

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3.3 C-WH5

- ① Telecamera
- ② Adattatore WIFI
- ③ Cavo HDMI
- ④ Mouse wireless

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OPTIKA® S.r.l.