HDMI Cameras – H Series

INSTRUCTION MANUAL

Model
C-HP
C-HPSC
C-HE
C-HESC
C-WH5

Ver. 3.1 2019

Table of Contents

1. Warning	3
2. Symbols and conventions	3
3. Safety Information	3
4. Intended use	3
5. C-HP Basic features	4
5.1 Package content	5
5.2 Quick Instructions for C-HP camera	5
5.2.1 USB Video Mode	5
5.2.2 HDMI Mode	6
6. C-HP User Interface (UI) and its Functions	7
6.1 The Camera Control Panel	8
6.2 Icons and Functions of the Camera Control Toolbar	9
6.3 The Measurement Toolbar	12
7. C-HE Basic features	14
7.1 Package content	14
7.2 Quick Instructions for C-HE camera	15
7.2.1 HDMI Mode	15
8. C-HE User Interface (UI) and Its Functions	16
8.1 The Camera Control Panel	17
8.2 Icons and Functions of the Camera Control Toolbar	17
9. C-HPSC / C-HESC	18
9.1 Overview	18
9.2 Package content	18
9.2.1 C-HPSC	18
9.2.2 C-HESC	19
9.3 Assembling	19
10. C-HPSC / C-HESC Use of the system	20
11. C-WH5 Basic features	20
11.1 Package content	21
11.2 Quick Instructions for C-WH5 camera	21
11.2.1 WiFi mode ProView controled	21
11.2.2 WiFi mode App controled	22
11.2.3 HDMI mode	22
12. C-WH5 User Interface (UI) and Its Functions	23
12.1 The Camera Control Panel	24
12.2 Icons and Functions of the Camera Control Toolbar	25
12.3 The Measurement Toolbar	30
13. Micrometric Slide M-005	32
Equipment disposal	33

1. Warning

This device is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.

2. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

3. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

4. Intended use

Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

5. C-HP Basic features

C-HP is a multiple interfaces (HDMI + USB2.0 + SD card) CMOS camera. HDMI + USB2.0 are used as the data transfer interface to HDMI display or computer.

For HDMI output, the *Camera Control Panel + Measurement Toolbar* and *Camera Control Toolbar* are overlaid on the HDMI screen when the mouse move to the related region. In this case, the USB mouse can be used to set the camera, browse and compare the captured image, play the video and perform the measurement.

For USB Video camera mode, plug in the micro USB host cable to the camera USB video port and computer USB port, then the video stream can be transfer to computer with the advanced software ProView. With ProView, you can control the camera, process the video and image as OPTIKA's other USB series camera.

The C-HP's basic features are as follows:

For HDMI output:

- All in 1(HDMI + USB + SD card) C-mount camera with Sony high sensitivity CMOS sensor;
- Resolution 2MP (1080p);
- Simultaneous HDMI & USB output;
- Built-in mouse control;
- Built-in image capture & video record to SD card;
- Built-in camera control panel, including exposure (manual/auto) / gain, white balance (lockable), color adjustment, sharpness control;
- Built-in video and image measurement;
- Built-in toolbar including zoom, mirror, comparison, freeze, cross, browser functions;
- Built-in image & video browsing, display & play;
- Real time clock (RTC).

For USB Video output:

- Ultra-Fine color engine with perfect color reproduction capability (USB);
- Resolution 2MP (USB);
- With advanced video & image processing application ProView, which including professional image processing such as 2D measurement, HDR, image stitching, EDF (Extended Depth of Focus), image segmentation & count, image stacking, color composite and denoising (USB);
- Supports standard UVC protocol for Windows (USB).

5.1 Package content

- ① Camera
- ② HDMI cable
- ③ USB-Micro USB cable
- ④ Wireless mouse
- ⑤ Micrometric slide
- ⑥ SD card 16GB
- ⑦ Power supply

5.2 Quick Instructions for C-HP camera

Before turning on the camera, please connect the standard C-mount camera to the camera adapter and insert it into the microscope's trinocular port.

Figure 1 - C-HP Rear Cover

- ① HDMI connector
- ② USB mouse connector
- ③ USB PC connector
- ④ SD card slot
- ⑤ Power supply plug
- ⑥ ON-OFF button
- ⑦ LED indicator

5.2.1 USB Mode

1. Plug the USB cable provided with the camera into USB Video port ③ to connect the video to the computer
2. Plug 12V/1A power cable into DC 12V slot ⑤ to supply power for the camera. The LED Indicator ⑦ will turn into red;
3. Press ON/OFF Button ⑥ to turn on the camera. The LED indicator will turn into blue;
4. Start ProView software, select "C-HP" by clicking the camera model name listed in Camera List. For more details please refer to the ProView help manual.

5.2.2 HDMI Mode

1. Plug the HDMI cable into the HDMI port ① to connect the C-HP camera to HDMI display;
2. Plug a USB mouse into USB Mouse port ② to get control of the camera by using built-in software ProView;
3. Plug 12V/1A power adapter into DC 12V slot ⑤ to supply power for the camera. The LED Indicator ⑦ will turn into red;
4. Insert SD card into SD Card slot ④ for saving captured images and recorded videos;
5. Press ON/OFF Button ⑥ to turn on the camera. The LED indicator will turn into blue;
6. Move mouse cursor to the left side of the video window, a *Camera Control Panel* will appear. It includes Manual/ Automatic Exposure, White Balance, Sharpness and other functions, please refer to 6.1 for details;
7. Move mouse cursor to the bottom of the video window and a *Camera Control Toolbar* will appear. Operations like Zoom In, Zoom Out, Flip, Freeze, Cross Line, Comparison and etc. can be achieved. Please refer to 6.2 for details;
8. Move mouse cursor to the upper side of the video window, a *Measurement Toolbar* with calibration and other measurement tools will appear, please refer to 6.3 for details. The measurement data can be output with *.CSV format.

6. C-HP User Interface (UI) and its Functions

The C-HP UI shown in Fig. 2 includes a *Camera Control Panel* on the left side of the video window, a *Measurement Toolbar* on the upper side of the video window, a *Camera Control Toolbar* on the bottom of the video window.


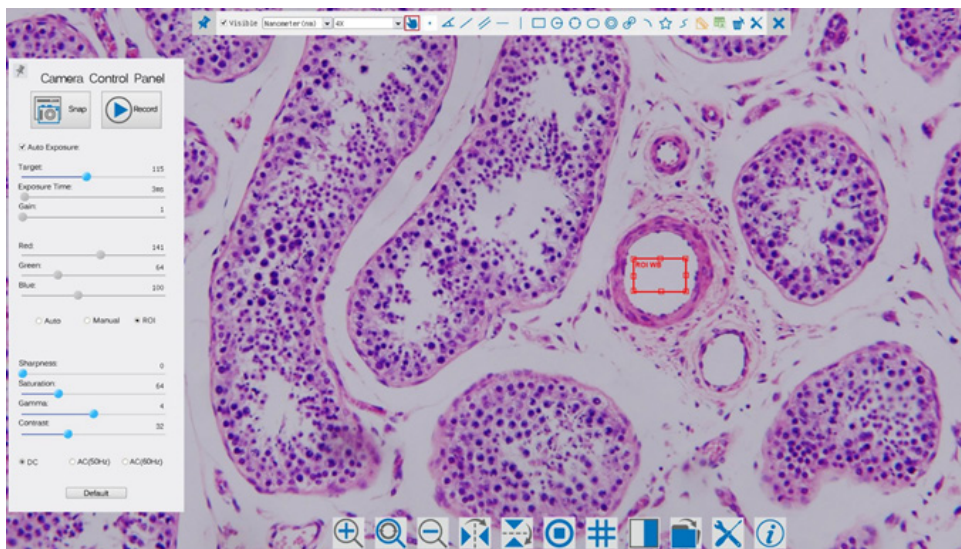
Figure 2 - Camera Control User Interface


Notes:

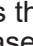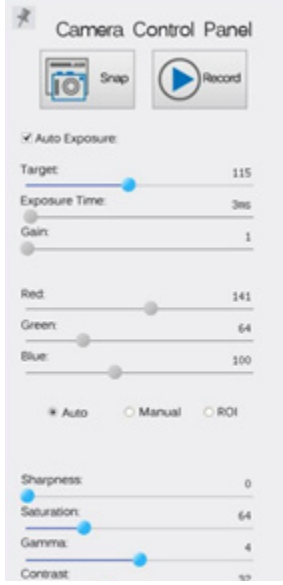1. When users move mouse cursor to the left side of the video window, the *Camera Control Panel* will pop up automatically.
2. When user moves mouse cursor to the bottom of the video window, the *Camera Control Toolbar* will pop up automatically.
3. Move the mouse cursor to the upper side of the video window: a *Measurement Toolbar* will pop up for the calibration and measurement operations. When user left-clicks the *Float/Fixed* button  on the Measurement Toolbar, the Measurement Toolbar will be fixed. In this case the Camera Control Panel will not pop up automatically even if users move mouse cursor to left side of the video windows. Only when user left-clicks the  button on the Measurement Toolbar to exit from measuring procedure will they be able to do other operations on the *Camera Control Panel*, or *Synthesis Camera Control Toolbar*. During the measuring process, when a specific measuring object is selected an *Object Location & Attributes Control Bar*  will appear for changing location and properties of the selected objects.

6.1 The Camera Control Panel

The *Camera Control Panel* controls the camera to achieve the best image quality according to the specific applications. It will pop up automatically when mouse cursor is moved to the left side of the video window (in measurement status, the Camera Control Panel will not pop up. Only when measurement process is terminated will the Camera Control Panel pop up by moving mouse cursor to the left side of the video window).

Left-clicking button  to achieve *Display/ Auto Hide* switch of the Camera Control Panel.

Control Panel	Function	Description	
	Snap	Capture image from the current video window	
	Record	Record video from the current video window	
	Auto Exposure	When Auto Exposure is checked, the system will automatically adjust exposure time according to the value of exposure compensation	
	Target	Available when Auto Exposure is checked. Slide to left or right to adjust Target according to the current video brightness to achieve proper brightness value	
	Exposure Time	Available when Auto Exposure is unchecked. Slide to left or right to reduce or increase exposure time, adjusting brightness of the video	
	Gain	Adjust Gain to reduce or increase brightness of video. The Noise will be reduced or increased accordingly	
	Red	Slide to left or right to decrease or increase the proportion of Red in RGB on video	
	Green	Green is base for reference and cannot be adjusted	
	Blue	Slide to left or right to decrease or increase the proportion of Blue in RGB on the video	
	White Balance	Auto:	White Balance adjustment according to the window video
		Manual:	Slide the Red or Blue to manually set the video White Balance
		ROI:	Set the White Balance according to the ROI. The ROI can be resized and moved
	Sharpness	Adjust Sharpness level of the video window	
	Saturation	Adjust Saturation level of the video window	
	Gamma	Adjust Gamma level of the video. Slide to the right side to increase gamma and to the left to decrease gamma	
	Contrast	Adjust Contrast level of the video. Slide to the right side to increase contrast and to the left to decrease contrast	
	DC	For DC illumination, there will be no fluctuation in light source so no need for compensating light flickering	
	AC(50HZ)	Check AC(50HZ) to eliminate flickering "strap" caused by 50Hz illumination	
AC(60HZ)	Check AC(60HZ) to eliminate flickering "strap" caused by 60Hz illumination		
Default	Set all the settings in the Camera Control Panel to default values		

6.2 Icons and Functions of the Camera Control Toolbar

Icon	Function	Icon	Function
	Zoom In the Video Window		Zoom Out the Video Window
	Reset Zoom to Original 1X		Horizontal Flip
	Vertical Flip		Video Freeze
	Compare Image with the current video		Display Cross Line
	Browse Images and Videos in the SD Card		Settings
	Check the ProView Version		

The  *Setting* function is relatively more complicated than the other functions. Here are more info about it:

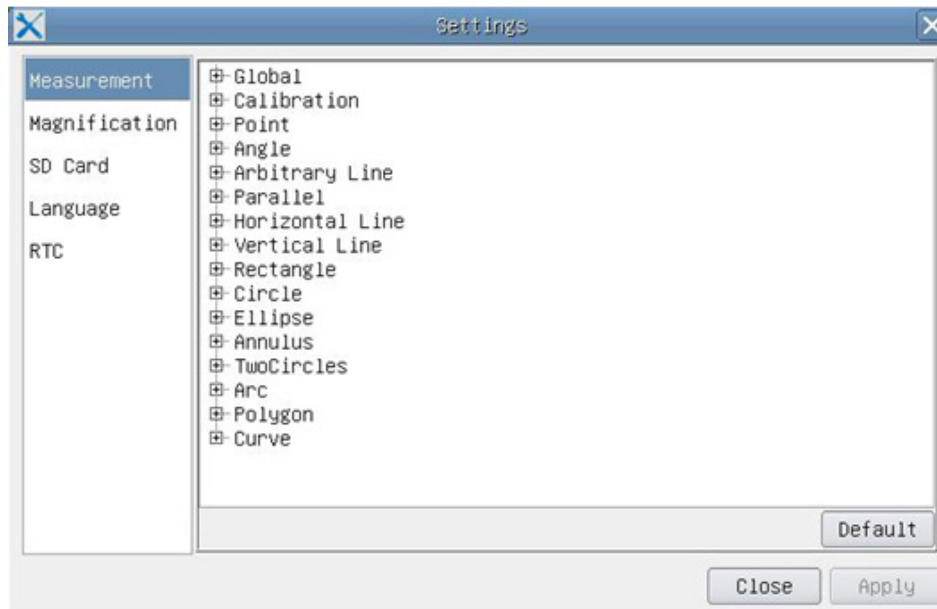
Figure 3 - Measurement Settings Page

- Global:** Used for setting digits behind the decimal point for measurement results;
- Calibration**
 - Line Width** Used for defining width of the lines for calibration;
 - Color** Used for defining color of the lines for calibration;
 - EndPoint**
 - Type:** Used for defining shape of the endpoints of lines for calibration.
 - Null:** Means no endpoints
 - Rectangle:** Means rectangle type of endpoints. It makes alignment more easily

Point, Angle, Line, Horizontal Line, Vertical Line, Rectangle, Circle, Ellipse, Annulus, Two Circles, Polygon, Curve:

Left-click the  beside the measuring patterns mentioned above will unfold the corresponding attribute settings to set the individual property of the measuring objects.

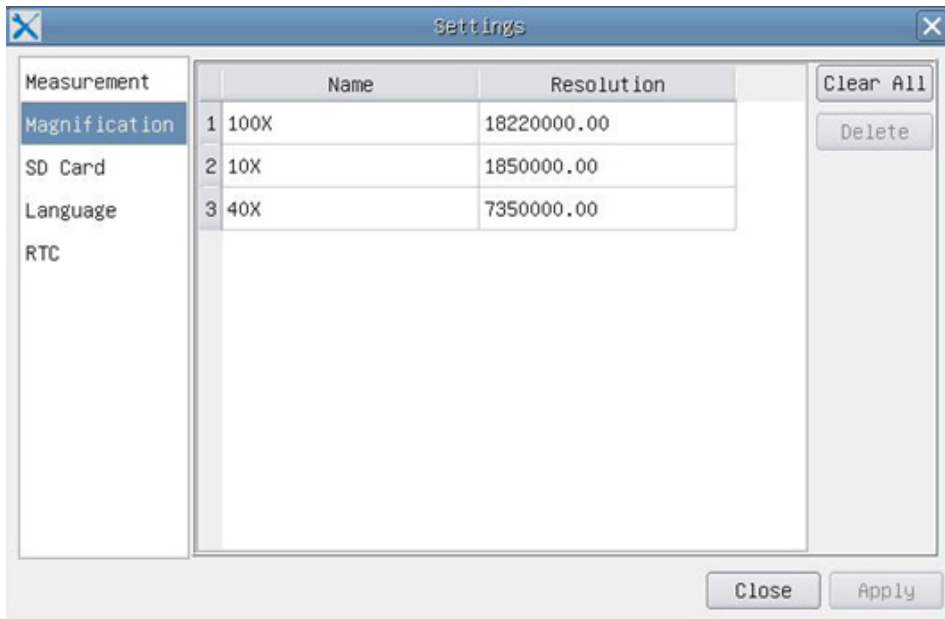
Figure 4 - Magnification Calibration Management Settings Page

Name: Names such as 10X, 40X, 100X are based on magnification of the microscopes. For continuous zoom microscopes, ensure that the selected magnification coincides with the scale alignment line on the microscope zoom knob;

Resolution: Pixels per meter. Devices like microscopes have high resolution value;

Clear All: Click the Clear All button will clear the calibrated magnifications and resolutions;

Delete: Click Delete to delete the selected item for specific resolution.

Figure 5 - SD Card Settings Page

File System FAT32: The file system of SD card is FAT32. The maximum video file size of single file in FAT32 file system is 4GB;

EXFAT: The maximum file EXFAT can store is 2048GB. Use PC to format the SD cards and switch between FAT32 and EXFAT;

Unknown Status: SD card not detected or the file system is not identified.

Figure 6 - Language Selection Settings Page

- English:* Set language of the whole software into English;
- Simplified Chinese:* Set language of the whole software into Simplified Chinese;
- Traditional Chinese:* Set language of the whole software into Traditional Chinese;
- Korean:* Set language of the whole software into Korean;
- Thailand:* Set language of the whole software into Thailand;

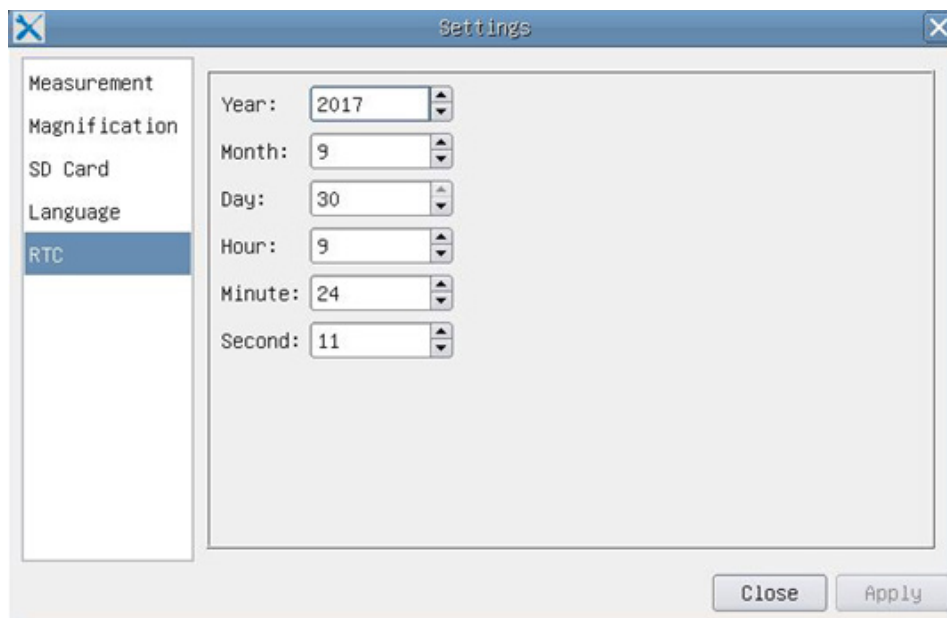
Figure 7 - The Real Time Clock Settings Page

- Year:* Current Year
- Month:* Current Month
- Day:* Current Day
- Hour:* Current Hour
- Minute:* Current Minute
- Second:* Current Second

6.3 The Measurement Toolbar

The *Measurement Toolbar* will pop up when moving mouse cursor to any place near the upper side of the video window.

Figure 8 - The Measurement Toolbar

Icon	Function
	Float/ Fix switch of the Measurement Toolbar
<input checked="" type="checkbox"/> Visible	Define measuring object in Show up/ Hide mode
Nanometer(nm)	Select the desired Measurement Unit
4X	Choose the same Magnification as the microscope to ensure accuracy of measurement result when measurement unit is not in Pixel units
	Object Select
	Point
	Angle
	Arbitrary Line
	Parallel
	Horizontal Line
	Vertical Line
	Rectangle
	Circle
	Ellipse
	Annulus
	Two Circles and Center Distance
	Arc
	Polygon
	Curve
	Make Calibration to determine the corresponding relation between magnification and resolution, this will establish the corresponding relationship between measurement unit and the sensor pixel size. Calibration needs to be done with the help of a micrometer. For detailed steps of carrying out calibration please refer to ProView help manual
	Export the measurement information to CSV file(*.csv)
	Delete All the Measurement Objects

	Settings
	Exit from Current Measurement Mode
	When the measurement ends, left-click on a single measuring object and the Object Location & Properties Control Bar will show up. The icons on the control bar mean Move Left, Move Right, Move Up, Move Down, Color Adjustment and Delete

Notes:

1. When user left-clicks *Display/Hide* button on the *Measurement Toolbar*, the Measurement Toolbar will be fixed. In this case the Camera Control Panel will not pop up automatically even if moving mouse cursor to the left side of the video window. Only when users left click the button on the Measurement Toolbar to exit from the measurement mode they will be able to doing other operations in the Camera Control Panel, the Auto Focus Control Panel or the Synthesis Camera Control Toolbar.
2. When a specific measuring object is selected during the measuring process, the *Object Location & Attributes Control Bar* will appear for changing the object location and properties of the selected objects.

7. C-HE Basic features

C-HE is a HDMI + SD card CMOS camera.
HDMI is used as the data transfer interface to HDMI display.

For HDMI output, the Camera Control Panel+ Measurement Toolbar and Camera Control Toolbar are overlaid on the HDMI screen when the mouse move to the related region.

The USB mouse is used to set the camera.

The C-HE's basic features are as follows:

For HDMI output:

- All in 1(HDMI + SD card) C-mount camera with Aptina high sensitivity CMOS sensor;
- Resolution 2MP (720p);
- Built-in mouse control;
- Built-in image capture to SD card;
- Built-in camera control panel, including exposure (manual/auto) / gain, color adjustment, sharpness control;
- Built-in toolbar including zoom, mirror, freeze, crosshair.

7.1 Package content

- ① Camera
- ② HDMI cable
- ③ Wireless mouse

- ④ SD card 16GB
- ⑤ Power supply

7.2 Quick Instructions for C-HE camera

Before turning on the camera, please connect the standard C-mount camera to the camera adapter and insert it into the microscope's trinocular port.

Figure 8 - C-HE Rear Cover

- | | |
|-----------------------|-----------------|
| ① HDMI connector | ④ SD card slot |
| ② USB mouse connector | ⑤ LED indicator |
| ③ Power supply plug | |

7.2.1 HDMI Mode

1. Plug the HDMI cable into the HDMI port ① to connect the C-HE camera to HDMI display;
2. Plug a USB mouse into USB Mouse port ② to get control of the camera by using built-in software LiteView;
 - **The camera cannot be connected to a PC in any way. The available USB port is used to connect the mouse.**
3. Insert SD card into SD Card Slot ④ for saving captured images;
4. Plug 12V/1A power adapter into DC 12V slot ③ to supply power for the camera. The LED Indicator ⑤ will turn into blue;
5. Move mouse cursor to the left side of the video window, a *Camera Control Panel* will appear. It includes Manual/ Automatic Exposure, White Balance, Sharpness and other functions, please refer to 8.1 for details;
6. Move mouse cursor to the bottom of the video window and a *Camera Control Toolbar* will appear. Operations like Zoom In, Zoom Out, Flip, Freeze, Cross Line, and etc. can be achieved. Please refer to 8.2 for details.

8 C-HE User Interface (UI) and Its Functions

The C-He UI shown in Fig. 9 includes a *Camera Control Panel* on the left side of the video window and a *Camera Control Toolbar* on the bottom of the video window.



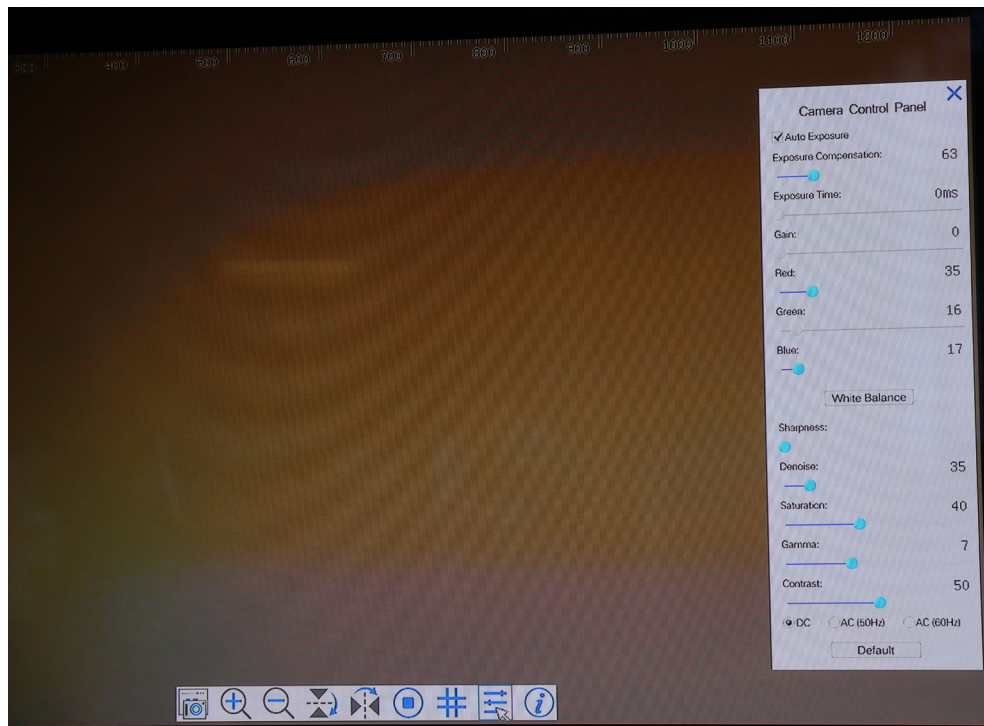
Figure 9 - Camera Control User Interface

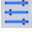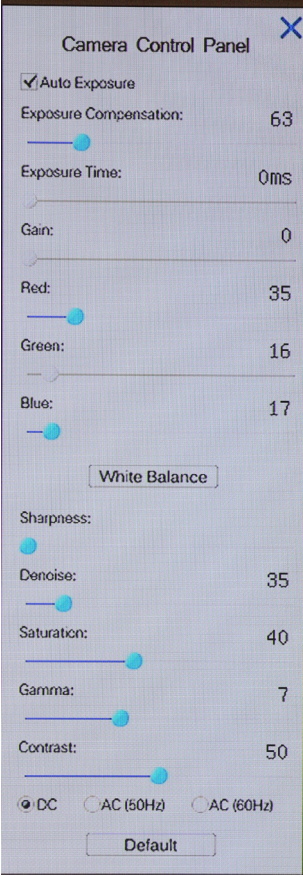Notes:

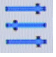
1. When users click on the  button, the *Camera Control Panel* will pop up automatically.
2. When users move mouse cursor to the bottom of the video window, the *Camera Control Toolbar* will pop up automatically.

8.1 The Camera Control Panel

The Camera Control Panel controls the camera to achieve the best image quality according to the specific applications. It will pop up automatically when mouse cursor is moved to the left side of the video window.

Camera Control Panel	Function	Description
	Auto Exposure	When Auto Exposure is checked, the system will automatically adjust exposure time according to the value of exposure compensation
	Exposure Compensation	Available when Auto Exposure is selected. Slide left or right to adjust the exposure compensation according to the current video brightness to get the correct brightness value.
	Exposure Time	Available when Auto Exposure is unchecked. Slide to left or right to reduce or increase exposure time, adjusting brightness of the video
	Gain	Adjust Gain to reduce or increase brightness of video. The Noise will be reduced or increased accordingly
	Red	Slide to left or right to decrease or increase the proportion of Red in RGB on video
	Green	Green is base for reference and cannot be adjusted
	Blue	Slide to left or right to decrease or increase the proportion of Blue in RGB on the video
	White Balance	Adjusting the White Balance according to the monitor display
	Sharpness	Adjust Sharpness level of the video window
	Denoise	Adjust Noise level of the video window
	Saturation	Adjust Saturation level of the video window
	Gamma	Adjust Gamma level of the video. Slide to the right side to increase gamma and to the left to decrease gamma
	Contrast	Adjust Contrast level of the video. Slide to the right side to increase contrast and to the left to decrease contrast
	DC	For DC illumination, there will be no fluctuation in light source so no need for compensating light flickering
	AC(50HZ)	Check AC(50HZ) to eliminate flickering "strap" caused by 50Hz illumination
	AC(60HZ)	Check AC(60HZ) to eliminate flickering "strap" caused by 60Hz illumination
Default	Set all the settings in the Camera Control Panel to default values	

8.2 Icons and Functions of the Camera Control Toolbar

Icon	Function	Icon	Function
	Snap the image and save it to SD card		Horizontal Flip
	Zoom In the Video Window		Video Freeze
	Zoom Out the Video Window		Display Cross Line
	Vertical Flip		Toggle ON/OFF Camera Control Panel

9. C-HPSC / C-HESC

9.1 Overview

9.2 Package content

9.2.1 C-HPSC

- ① C-HP camera
- ② HDMI Monitor
- ③ Mounting bracket
- ④ HDMI cable
- ⑤ USB-Micro USB cable

- ⑥ Wireless mouse
- ⑦ SD card 16GB
- ⑧ Micrometric slide
- ⑨ Power supply (2 pcs)

9.2.2 C-HESC

- ① C-HE camera C-HE
- ② HDMI Monitor
- ③ Mounting bracket
- ④ HDMI cable

- ⑤ Wireless mouse
- ⑥ SD card 16GB
- ⑦ Power supply (2 pcs)

9.3 Assembling

1. Screw the four provided allen screws on the back side of the monitor. (Fig. 10)

Fig. 10

2. Screw the two provided cross screws on the back side of the camera. (Fig. 11)

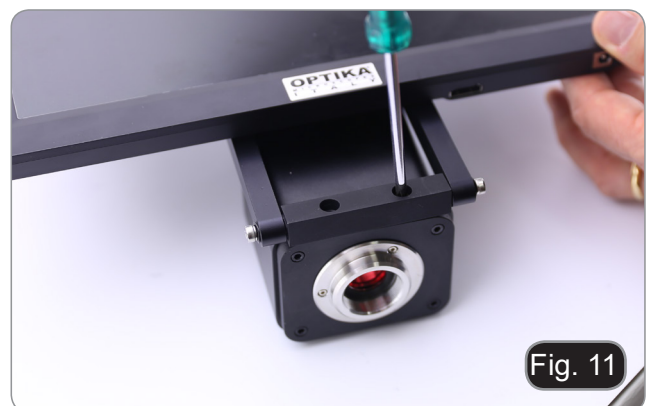
Fig. 11

3. Install the "C" mount and connect the HDMI and power cables to monitor and camera.
4. System is ready.

10. C-HPSC / C-HESC Use of the system

For the operating instructions of the C-HP camera, refer to chapter 6) and for the operating instructions of the C-HE camera, refer to chapter 8).

11. C-WH5 Basic features

C-WH5 is a multiple interfaces (HDMI + WiFi + SD card) CMOS camera. HDMI and WIFI are used as the data transfer interface to HDMI display or computer.

For HDMI output, the *Camera Control Panel + Measurement Toolbar* and *Camera Control Toolbar* are overlaid on the HDMI screen when the mouse move to the related region; in this case, the USB mouse can be used to set the camera, browse and compare the captured image, play the video and perform the measurement.

In the WiFi interface, connect the WiFi adapter to the camera's USB port.

The camera generates a WiFi signal that is recognized by a computer with a WLAN connection or by a mobile device (IOS or Android).

The camera can be managed via a PC using the ProView software. With ProView, you can control the camera, process video and images in the same way as with all other cameras in the OPTIKA series.

Managing the camera via a mobile device requires an App to be downloaded on your device, allowing multiple users to view and capture images from the camera.

The C-WH5's basic features are as follows:

For HDMI output:

- C-mount camera with Sony high sensitivity CMOS sensor;
- Resolution 2MP (1080p);
- HDMI output;
- Built-in mouse control;
- Built-in image capture & video record to SD card;
- Built-in camera control panel, including exposure (manual/auto) / gain, white balance (lockable), color adjustment, sharpness control;
- Built-in video and image measurement;
- Built-in toolbar including zoom, mirror, comparison, freeze, cross, browser function;
- Built-in image & video browsing, display & play;
- Real time clock (RTC).

For WiFi output:

- Ultra-Fine color engine with perfect color reproduction capability;
- Resolution 5MP;
- With advanced video & image processing application ProView, which including professional image processing such as 2D measurement, HDR, image stitching, EDF (Extended Depth of Focus), image segmentation & count, image stacking, color composite and denoising;
- Supports standard UVC protocol for Windows.

11.1 Package content

- ① Camera
- ② WIFI adapter
- ③ HDMI cable
- ④ Wireless mouse
- ⑤ Micrometric slide
- ⑥ SD card 16GB
- ⑦ Power supply

11.2 Quick Instructions for C-WH5 camera

Before turning on the camera, please connect the standard C-mount camera to the camera adapter and insert it into the microscope's trinocular port.

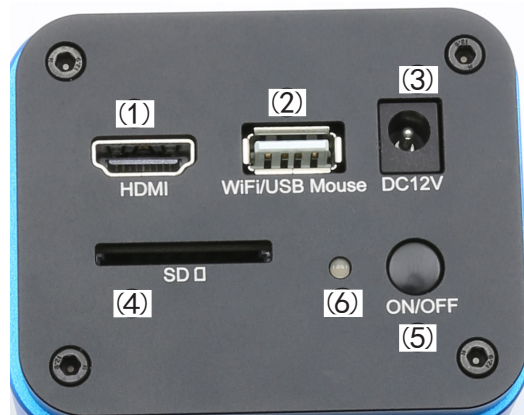
Figure 12 - C-WH5 Rear Cover

- ① HDMI connector
- ② USB mouse/WiFi adapter connector
- ③ Power supply plug
- ④ SD card slot
- ⑤ ON-OFF button
- ⑥ LED indicator

11.2.1 WiFi mode ProView controled

1. Connect the 12V/1A power cable to the DC 12V connector ③ to power the camera. The ⑥ LED indicator will turn red;
2. Connect the provided WiFi adapter to the WIFI/USB Mouse port ②;
3. Press the ON/OFF button ⑤ to turn on the camera; the LED indicator ⑥ will turn blue;
4. Start the computer (equipped with WiFi connection) and in the "WiFi" settings of the computer select **C-WH5-XXXXX**.
 - The password for connecting the camera is "12345678".
5. Start the ProView software, activate C-WH5 by clicking the camera name in the Camera List. For more details, refer to the ProView instruction manual.

11.2.2 WiFi mode App controled

1. Connect the 12V/1A power cable to the DC 12V connector ③ to power the camera. The ⑥ LED indicator will turn red;
2. Connect the provided WiFi adapter to the WIFI/USB Mouse port ②;
3. Press the ON/OFF button ⑤ to turn on the camera; the LED indicator ⑥ will turn blue;
4. In the Settings of your mobile device select the WiFi device **C-WH5-xxxxx**;
 - The password for connecting the camera is “12345678”.
5. Download and launch the TouView App (available on AppStore for IOS and PlayStore for Android);
6. Manage the camera using the App to control all functions (Exposure Control, White Balance, Image or Movies Acquisition).

NOTE: The maximum number of users who can connect simultaneously is limited to 3. This is to ensure that the camera does not lose speed when many users are connected simultaneously. When 3 users are connected, the camera will not allow a fourth user to connect.

11.2.3 HDMI mode

1. Plug the HDMI cable into the HDMI port ① to connect the C-HP camera to HDMI display;
2. Plug a USB mouse into USB Mouse port ② to get control of the camera by using built-in software ProView;
3. Plug 12V/1A power adapter into DC 12V slot ③ to supply power for the camera. The LED Indicator ⑥ will turn into red;
4. Insert SD card into SD Card slot ④ for saving captured images and recorded video;
5. Press ON/OFF Button ⑤ to turn on the camera. The LED indicator ⑥ will turn into blue;
6. Move mouse cursor to the left side of the video window, a *Camera Control Panel* will appear. It includes Manual/ Automatic Exposure, White Balance, Sharpness and other functions, please refer to 12.1 for details;
7. Move mouse cursor to the bottom of the video window and a *Camera Control Toolbar* will appear. Operations like Zoom In, Zoom Out, Flip, Freeze, Cross Line, Comparison and etc. can be achieved. Please refer to 12.2 for details;
8. Move mouse cursor to the upper side of the video window, a *Measurement Toolbar* with calibration and other measurement tools will appear, please refer to 12.3 for details. The measurement data can be output with *.CSV format.

12. C-WH5 User Interface (UI) and Its Functions

The C-WH5 UI shown in Fig. 13 includes a *Camera Control Panel* on the left side of the video window, a *Measurement Toolbar* on the upper side of the video window, a *Camera Control Toolbar* on the bottom of the video window.

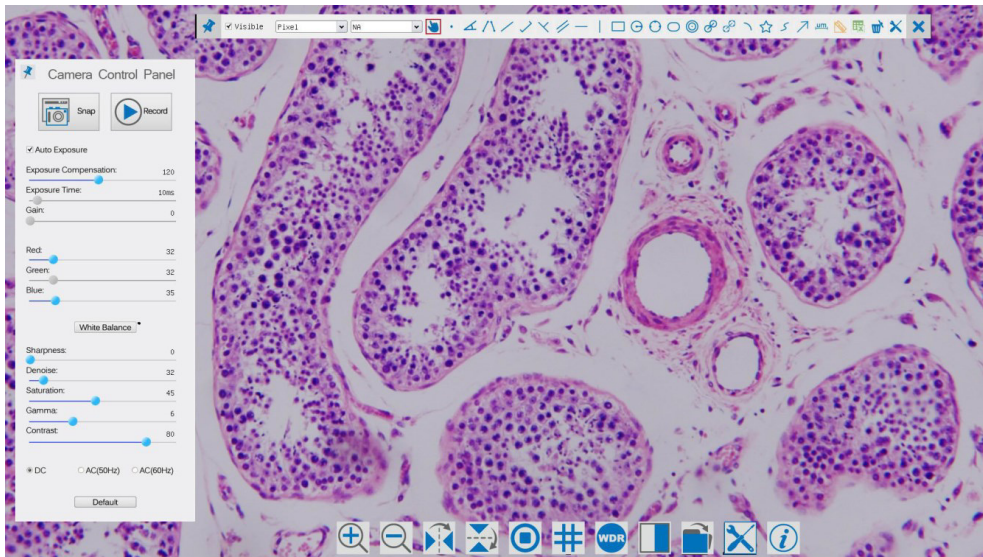
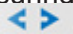
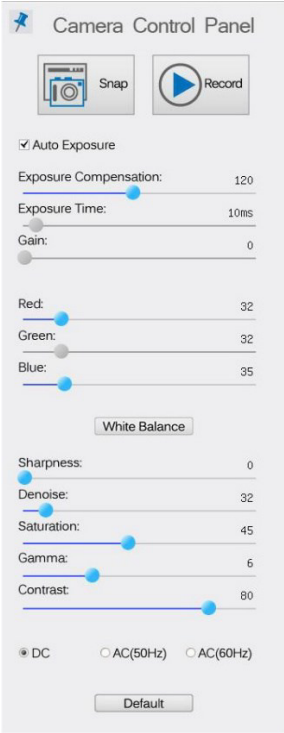
Figure 13 - Camera Control User Interface

Notes:

1. When users move mouse cursor to the left side of the video window, the *Camera Control Panel* will pop up automatically.
2. When user moves mouse cursor to the bottom of the video window, the *Camera Control Toolbar* will pop up automatically.
3. Move the mouse cursor to the upper side of the video window: a *Measurement Toolbar* will pop up for the calibration and measurement operations. When user left-clicks the *Float/Fixed* button  on the Measurement Toolbar, the Measurement Toolbar will be fixed. In this case the Camera Control Panel will not pop up automatically even if users move mouse cursor to left side of the video windows. Only when user left-clicks the  button on the Measurement Toolbar to exit from measuring procedure will they be able to do other operations on the *Camera Control Panel*, or *Synthesis Camera Control Toolbar*. During the measuring process, when a specific measuring object is selected an *Object Location & Attributes Control Bar*  will appear for changing location and properties of the selected objects.

12.1 The Camera Control Panel

The *Camera Control Panel* controls the camera to achieve the best image quality according to the specific applications. It will pop up automatically when mouse cursor is moved to the left side of the video window (in measurement status, the Camera Control Panel will not pop up. Only when measurement process is terminated will the Camera Control Panel pop up by moving mouse cursor to the left side of the video window). Left-clicking button  to achieve *Display/ Auto Hide* switch of the Camera Control Panel.

Control Panel	Function	Description
 <p>The screenshot shows the Camera Control Panel interface with the following settings: Snap and Record buttons; Auto Exposure checked; Exposure Compensation: 120; Exposure Time: 10ms; Gain: 0; Red: 32; Green: 32; Blue: 35; White Balance button; Sharpness: 0; Denoise: 32; Saturation: 45; Gamma: 6; Contrast: 80; DC selected; AC(50Hz) and AC(60Hz) unselected; Default button.</p>	Snap	Capture image from the current video window
	Record	Record video from the current video window
	Auto Exposure	When Auto Exposure is checked, the system will automatically adjust exposure time according to the value of exposure compensation
	Exposure Compensation	Available when Auto Exposure is selected. Slide left or right to adjust the exposure compensation according to the current video brightness to get the correct brightness value.
	Exposure Time	Available when Auto Exposure is unchecked. Slide to left or right to reduce or increase exposure time, adjusting brightness of the video
	Gain	Adjust Gain to reduce or increase brightness of video. The Noise will be reduced or increased accordingly
	Red	Slide to left or right to decrease or increase the proportion of Red in RGB on video
	Green	Green is base for reference and cannot be adjusted
	Blue	Slide to left or right to decrease or increase the proportion of Blue in RGB on the video
	White Balance	Adjusting the White Balance according to the monitor display
	Sharpness	Adjust Sharpness level of the video window
	Denoise	Adjust Noise level of the video window
	Saturation	Adjust Saturation level of the video window
	Gamma	Adjust Gamma level of the video. Slide to the right side to increase gamma and to the left to decrease gamma
	Contrast	Adjust Contrast level of the video. Slide to the right side to increase contrast and to the left to decrease contrast
	DC	For DC illumination, there will be no fluctuation in light source so no need for compensating light flickering
	AC(50HZ)	Check AC(50HZ) to eliminate flickering "strap" caused by 50Hz illumination
	AC(60HZ)	Check AC(60HZ) to eliminate flickering "strap" caused by 60Hz illumination
	Default	Set all the settings in the Camera Control Panel to default values

12.2 Icons and Functions of the Camera Control Toolbar

Icon	Function	Icon	Function
	Zoom In the Video Window		Zoom Out the Video Window
	Vertical Flip		Horizontal Flip
	Video Freeze		Display Cross Line
	Enable Wide Dynamic Range mode		Compare Image with the current video
	Browse Images and Videos in the SD Card		Settings
	Informations on ProView version		

The  *Setting* function is relatively more complicated than the other functions. Here are more info about it:

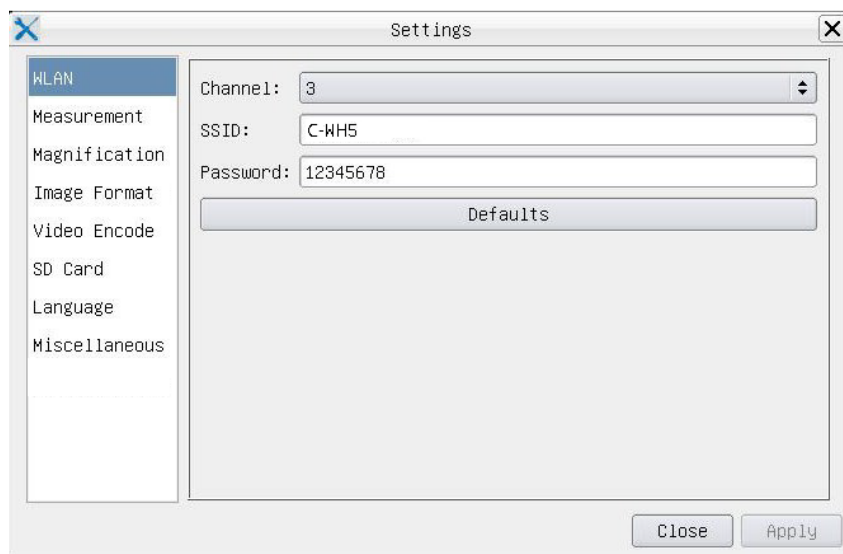
Figure 14 - WIFI Settings Page

Channel: Wireless communication channel, is a data transmission channel with wireless signals as transmission media;

SSID: Camera name sent by WIFI adapter;

Password: WIFI password; default is 12345678.

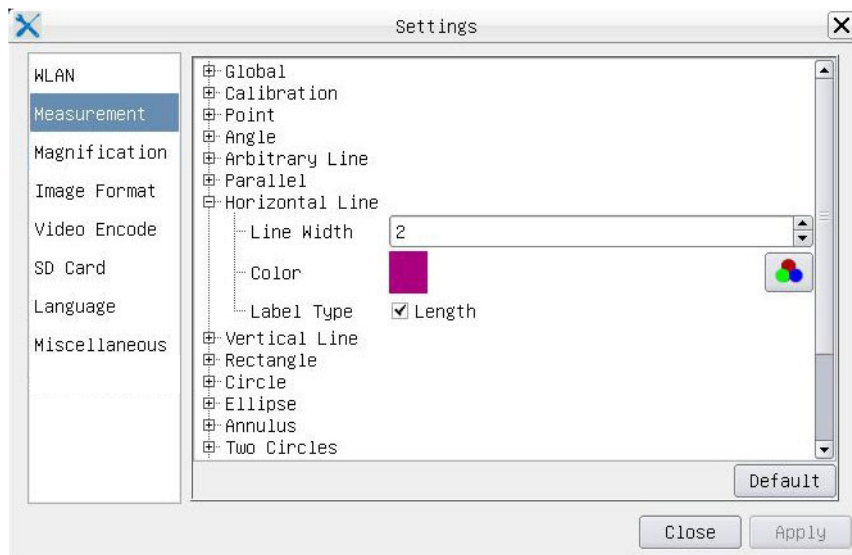
Figure 15 - Measurement Settings Page

Global: Used for setting digits behind the decimal point for measurement results;

Calibration Line Width Used for defining width of the lines for calibration;

Color Used for defining color of the lines for calibration;

EndPoint Type: Used for defining shape of the endpoints of lines for calibration.

Null: Means no endpoints

Rectangle: Means rectangle type of endpoints. It makes alignment more easily

Point, Angle, Line, Horizontal Line, Vertical Line, Rectangle, Circle, Ellipse, Annulus, Two Circles, Polygon, Curve:

Left-click the  beside the measuring patterns mentioned above will unfold the corresponding attribute settings to set the individual property of the measuring objects.

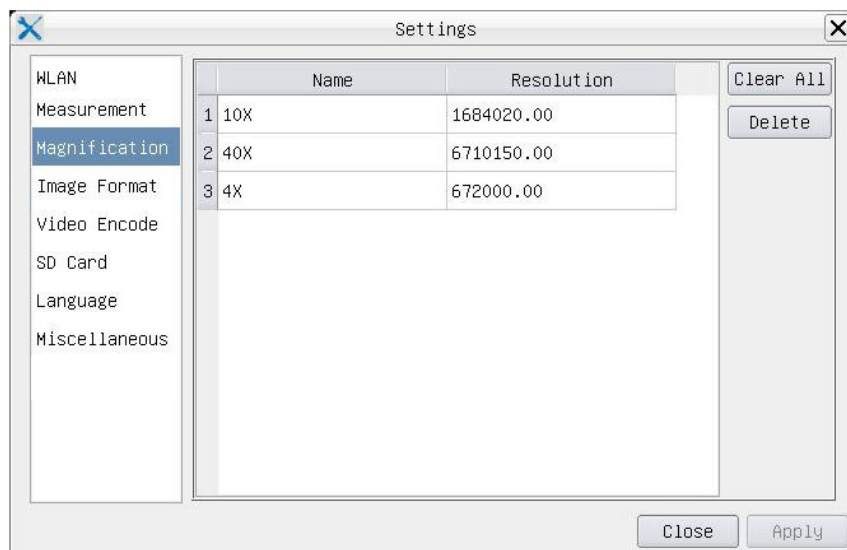
Figure 16 - Magnification Calibration Management Settings Page

Name: Names such as 10X, 40X, 100X are based on magnification of the microscopes. For continuous zoom microscopes, ensure that the selected magnification coincides with the scale alignment line on the microscope zoom knob;

Resolution: Pixels per meter. Devices like microscopes have high resolution value;

Clear All: Click the Clear All button will clear the calibrated magnifications and resolutions;

Delete: Click Delete to delete the selected item for specific resolution.

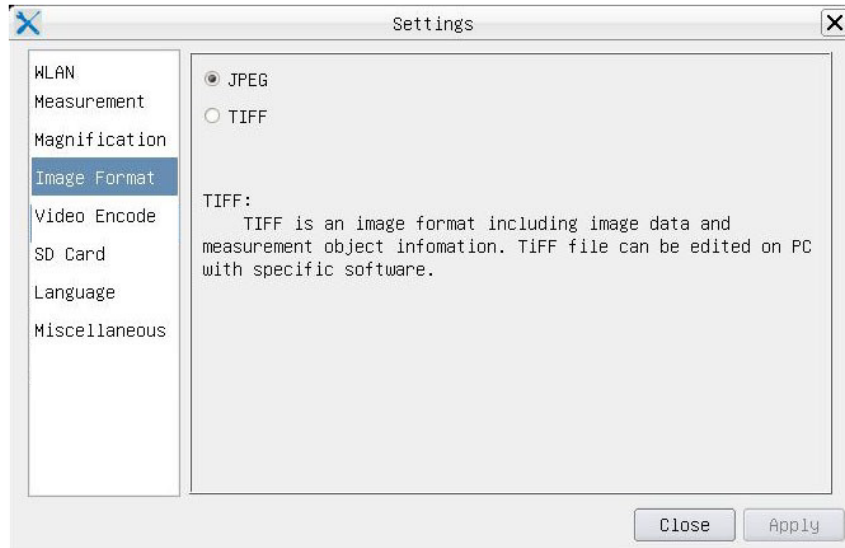
Figure 17 - Image Format Settings Page

JPEG The JPEG file can achieve a very high compression rate and display very rich and vivid images by removing redundant images and color data. In other words, you can get better image quality with minimal disk space. If measuring objects are available, they are burned in the image and the measurement cannot be changed.

TIFF TIFF is an image format that can include both the image data and the measurement object. The measurement objects are saved separately from the image data and can be edited later with specific software.

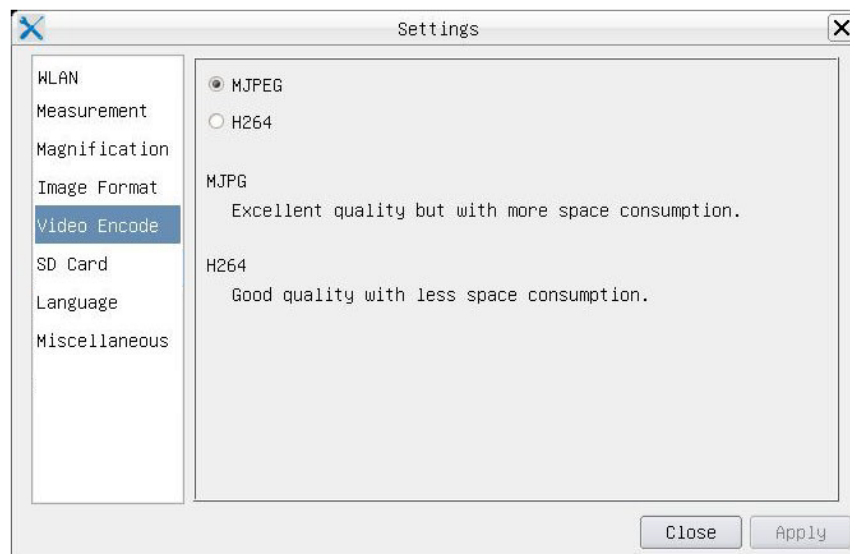
Figure 18 - Video Encode Settings Page

MJPEG Codec format for video is MJPEG. Excellent quality but with more space consumption than H264;

H264 Codec format for video is H264. Good quality with less space consumption.

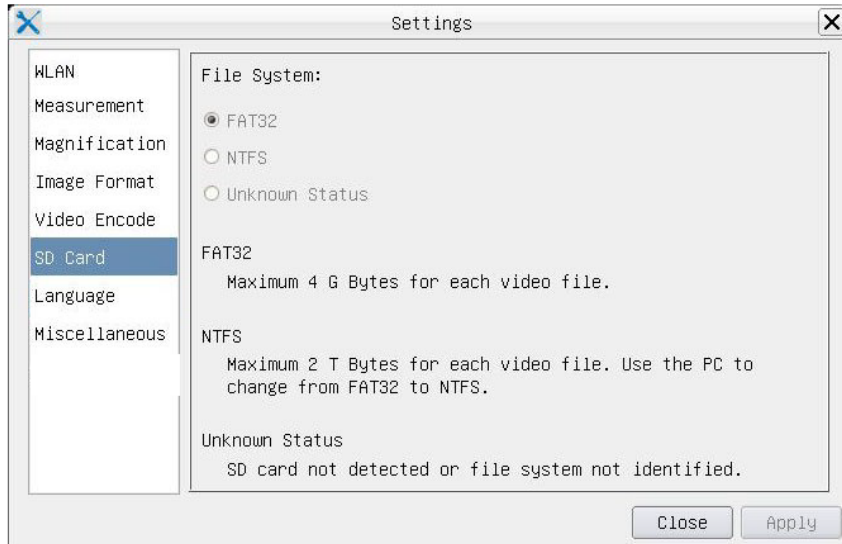
Figure 19 - SD Card Settings Page

- File System FAT32:** The file system of SD card is FAT32. The maximum video file size of single file in FAT32 file system is 4GB;
- NTFS:** The file system of SD card is NTFS. The maximum video file size of single file is 2TB.
Use PC to format the SD cards and switch between FAT32 and NTFS;
- Unknown Status:** SD card not detected or the file system is not identified.

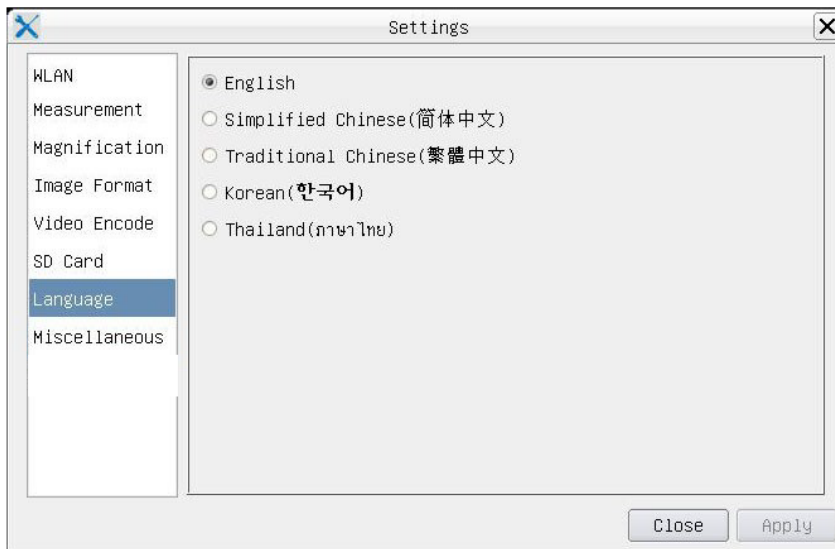
Figure 20 - Language Selection Settings Page

- English:** Set language of the whole software into English;
- Simplified Chinese:** Set language of the whole software into Simplified Chinese;
- Traditional Chinese:** Set language of the whole software into Traditional Chinese;
- Korean:** Set language of the whole software into Korean;
- Thailand:** Set language of the whole software into Thailand.

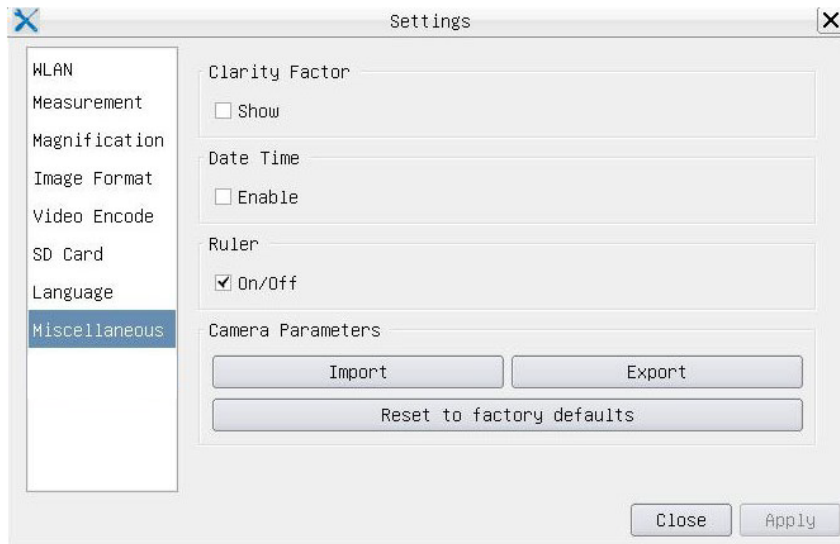
Figure 21 - Miscellaneous Settings Page

- Clarity Factor** The Clarity Factor informs if the sample is in good focus state or not. The larger the Clarity Factor, the better the focus of the sample. Enable *Show* to see if the video is fully focused or not;
- Date Time** For cameras without RTC (Real Time Clock) configuration, users can configure the system date after installation of the camera, which is not saved because there is no built-in RTC;
- Ruler (On/Off)** Enable to show the Ruler on the video window;
- Import** Import parameter settings saved from SD card;
- Export** Export current camera parameters to SD card for later use;
- Reset to factory defaults** Reset camera parameters to factory conditions.

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
	Polygon
	Curve
	Arrow
	Ruler
	Make Calibration to determine the corresponding relation between magnification and resolution, this will establish the corresponding relationship between measurement unit and the sensor pixel size. Calibration needs to be done with the help of a micrometer. For detailed steps of carrying out calibration please refer to ProView help manual
	Export the measurement information to CSV file(*.csv)
	Delete All the Measurement Objects
	Settings
	Exit from Current Measurement Mode
	When the measurement ends, left-click on a single measuring object and the Object Location & Properties Control Bar will show up. The icons on the control bar mean Move Left, Move Right, Move Up, Move Down, Color Adjustment and Delete

- Note:
1. When user left-clicks *Display/Hide* button  on the *Measurement Toolbar*, the Measurement Toolbar will be fixed. In this case the Camera Control Panel will not pop up automatically even if moving mouse cursor to the left side of the video window. Only when users left click the  button on the Measurement Toolbar to exit from the measurement mode they will be able to doing other operations in the Camera Control Panel, the Auto Focus Control Panel or the Synthesis Camera Control Toolbar.
 2. When a specific measuring object is selected during the measuring process, the *Object Location & Attributes Control Bar*  will appear for changing the object location and properties of the selected objects.


13. Micrometric Slide M-005

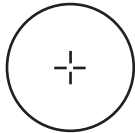Micrometric slide, 26x76mm, with 2 scales
(1mm/100div. for biological microscopes / 10mm/100div. for stereomicroscopes)

1 DIV=0.01mm

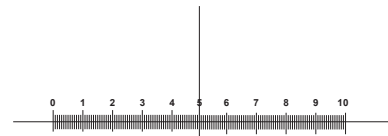For biological microscopes calibration

1 DIV=0.1mm

For stereo microscopes calibration

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.