INSPECTION SYSTEMS Series

INSTRUCTION MANUAL

	Model
	IS-01
	IS-01SMD

Ver. 2.3 2023

Table of contents

1.	Warning	4
2.	Safety Information	4
3.	Package content	5
3.1	IS-01	5
3.2	IS-01SMD	6
4.	Unpacking	7
5.	Intended use	7
6.	Symbols and conventions	7
7.	Instrument description	8
7.1	IS-01	8
7.2	IS-01SMD	9
8.	Assembling	11
9.	Use of the microscope	13
9.1	Light intensity adjustment	13
9.2	Coarse focus tension adjustment	13
9.3	Focus adjustment	13
9.4	Focusing	13
9.5	Magnification change	14
10.	Use of the software	15
10.1	Menu Interface	15
a)	Control	15
b)	Measurement	16
10.1.1	Detailed Control Interface	17
a)	Language Selection	17
b)	Control Panel Lock	17
c)	Exposure	17
d)	White Balance	18
e)	Function	18
f)	Crosshair Tool	18
10.1.2	Image capture	20
a)	USB flash drive insertion	20
b)	Image capture operation	20
c)	Prefix	20
d)	Image Format	20
e)	Image Naming Rules	21
10.1.3	Recording	21
10.1.4	Open Image	22
10.1.5	Image comparison	23
10.1.6	Freeze	23
10.1.7	Save Graphics	23
10.1.8	Flip/ Mirror	24
10.1.9	HDR	24
10.1.10	Anti-screen Flicker	24
10.1.11	Black and White Mode	24
10.1.12	Default	24
10.1.13	Version Number	24
10.2	Detailed Measurement Interface	25
10.2.1	Assist Tool	25
a)	Color Setting	25
b)	Edge Detection	25
10.2.2	Measurement Tool	25
10.2.3	Scan Data	26
10.2.4	Custom Template	26
10.2.5	Measurement Operation	26
a)	Calibration	26
b)	Delete calibration	27
c)	Measuring	27
d)	Delete Measurement Data	28
e)	Select Measurement Data	28
f)	Select Graph	28
g)	Detailed Data	29
h)	Mark Tool	29
11.	Maintenance	30
12.	Troubleshooting	31
	Equipment disposal	32

1. Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.

2. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

3. Package content

3.1 IS-01

- ① Stand
- ② Drop prevention ring
- ③ Focus system
- ④ Microscope body
- ⑤ Ringlight

- ⑥ HDMI cable
- ⑦ Power supply
- ⑧ Wireless mouse
- ⑨ Dust cover

3.2 IS-01SMD

- ① Stand
- ② Drop prevention ring
- ③ Focus system
- ④ Microscope body
- ⑤ Ringlight

- ⑥ 3D attachment
- ⑦ HDMI cable
- ⑧ Power supply
- ⑨ Wireless mouse
- ⑩ Dust cover

4. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

5. Intended use

Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

6. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

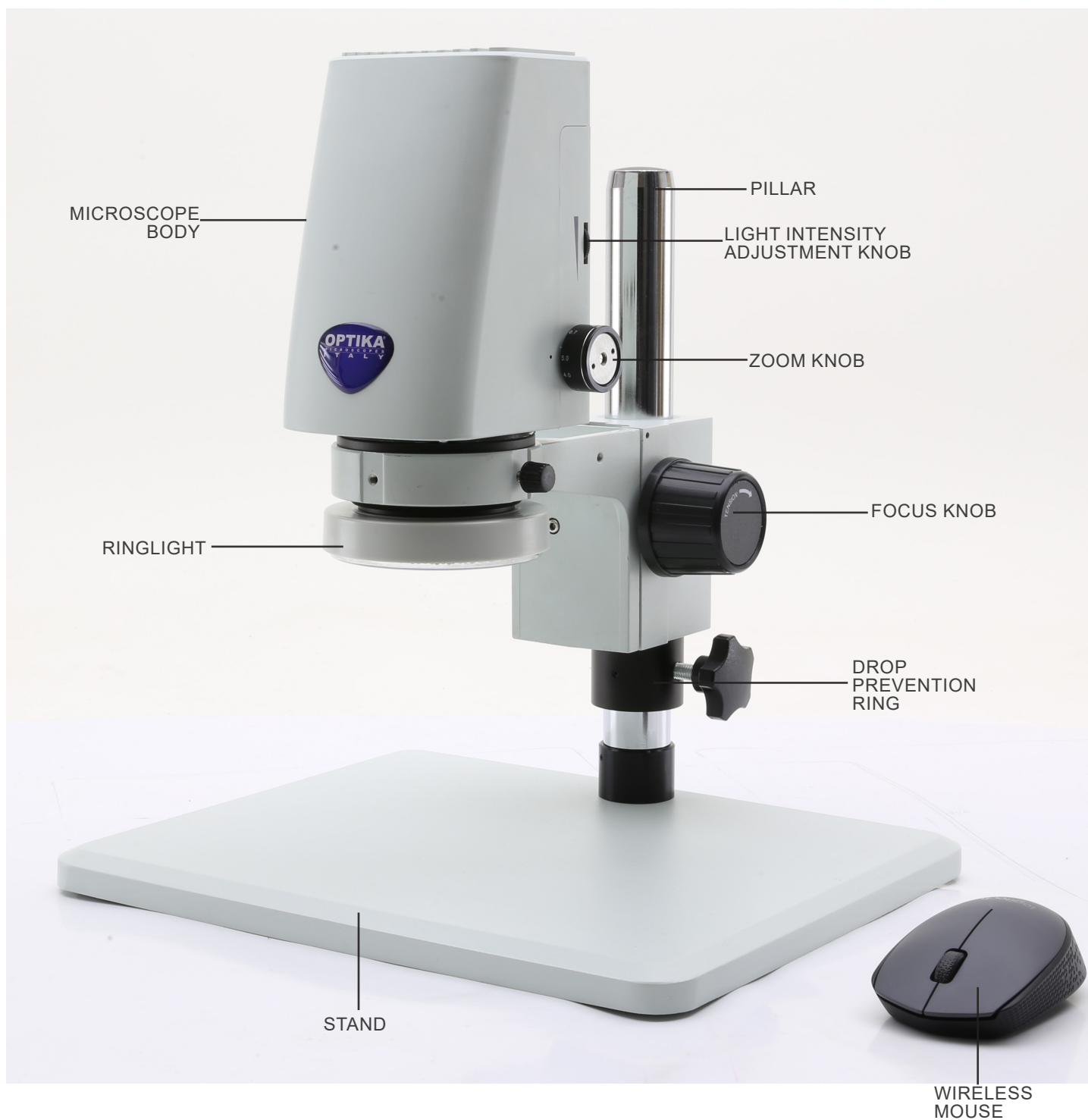
ELECTRICAL SHOCK


This symbol indicates a risk of electrical shock.

7. Instrument description


7.1 IS-01

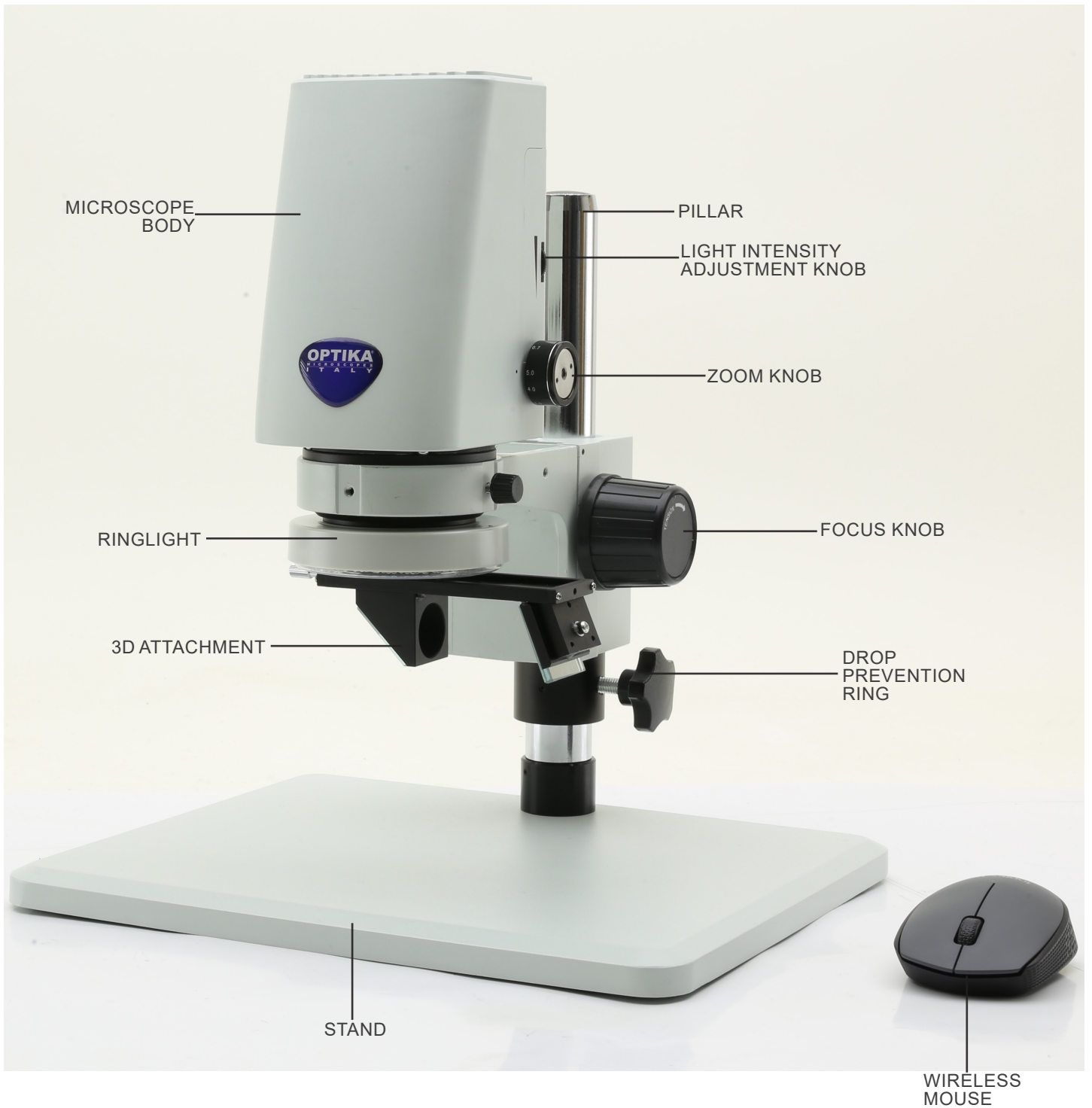7.2 IS-01SMD


Back side

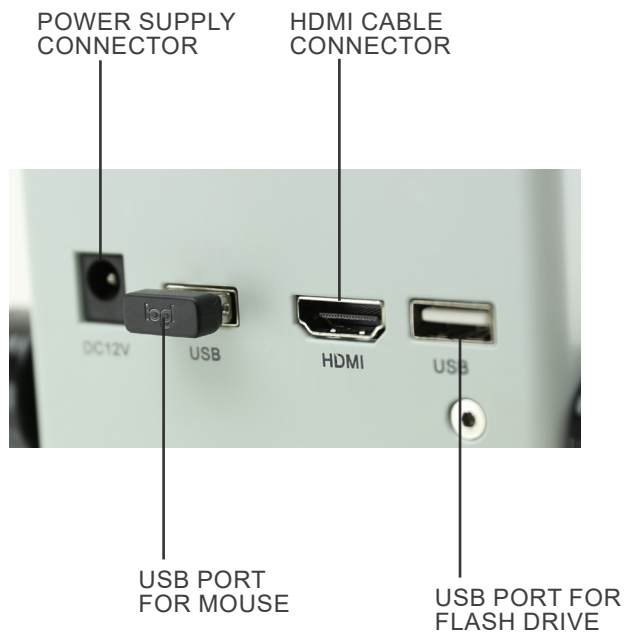8. Assembling

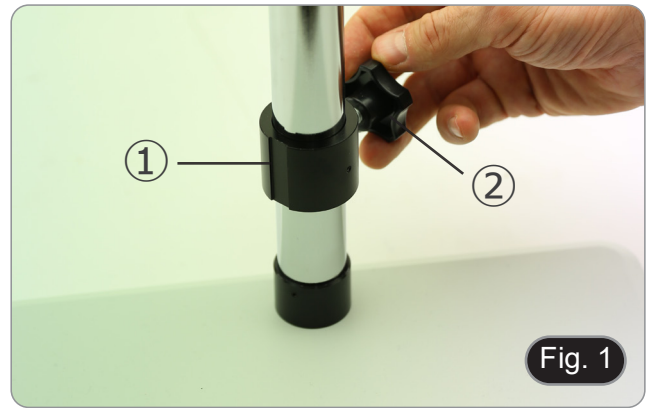1. Insert the drop prevention ring ① in the pillar and lock it with the fixing knob ②. (Fig. 1)


2. Insert the focus system ③ in the pillar and lock it with fixing knob ④. (Fig. 2)

3. Insert the microscope body in the hole of the focus system and lock it with the fixing knob ⑤. (Fig. 3)

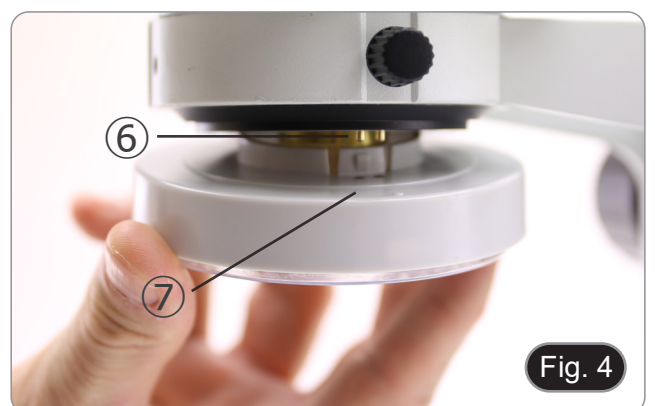4. Align the pins ⑥ of the microscope body with the holes ⑦ on the ringlight illuminator, then gently push the ringlight until it clicks into its position. (Fig. 4)


Only for IS-01SMD

Screw the 3D attachment on the lens thread on the microscope body. (Fig. 5)

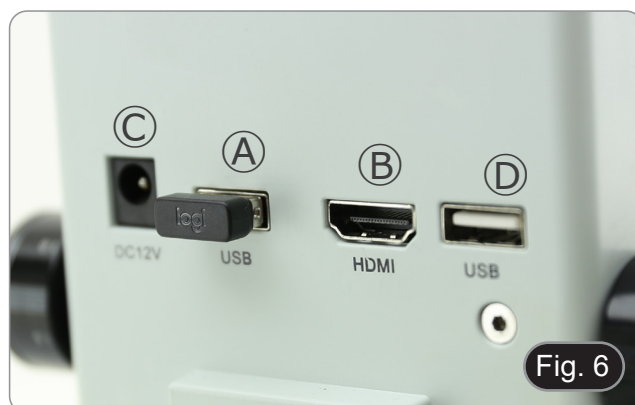5. Connect the wireless mouse receiver in the USB port (A), the HDMI cable in the port (B) and the power supply jack in the connector (C). (Fig. 6)
6. Connect a USB flash disk into the port (D).
 - **There is no ON-OFF switch on the microscope: when the power supply is connected, the system starts up automatically and begins projecting images onto the monitor.**
 - **It takes approximately 10-15 seconds for the system to start up before the image is displayed.**


9. Use of the microscope

9.1 Light intensity adjustment

1. Rotate the light adjustment ring ① on the right side of the microscope body. (Fig. 7)

As an alternative:

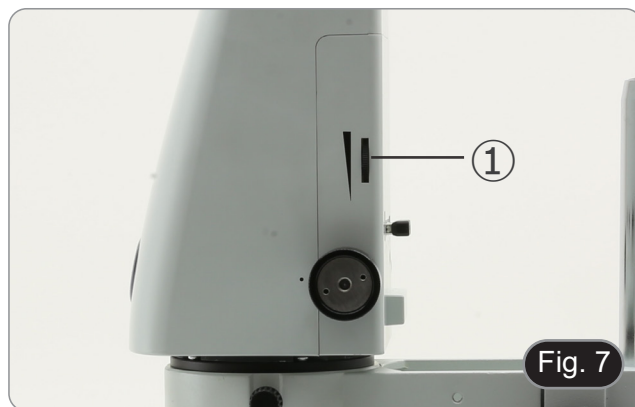1. Move the mouse button on the upper left corner of the screen.
2. The Main Control Panel appears (see chapter 10.1 a).
3. Adjust the illumination using the Camera Control Panel of the built-in software.

9.2 Coarse focus tension adjustment

1. To adjust the focusing tension, hold the left knob ② and rotate the right knob ③. (Fig. 8)
- **Clockwise rotation increases the tension while counter-clockwise rotation loosens it.**
 - Adjust the focusing tension to appropriate position, to prevent the microscope body dropping during observation, and also make focusing stable.

9.3 Parfocality adjustment

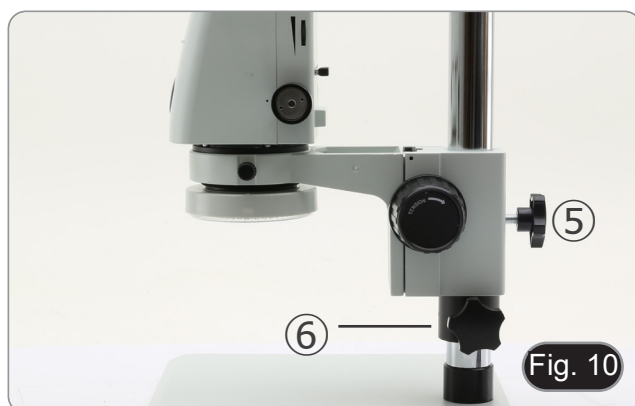
1. Rotate the zoom adjustment ring ④ to the minimum magnification. (Fig. 9)
2. While observing on the monitor, focus the specimen.
3. Rotate the zoom adjustment ring ④ to the maximum magnification.
4. Observe the image on the monitor: if it is not clear, rotate the focusing knob to make it sharp.
5. Return zoom adjustment ring to the minimum magnification. Now parfocality adjustment should be set.

9.4 Focusing

Put the sample to be observed on the stand plate and focus the sample using the focusing knobs.

- If necessary adjust the height of the microscope head along the pillar: unlock the fixing knob ⑤, move the microscope to the desired height and lock the fixing knob ⑤ after setting the height of the microscope. (Fig. 10)
- **In order to prevent accidental lowering of the system due to imperfect locking of the fixing knob, it is advisable to also use the drop prevention ring ⑥.**

9.5 Magnification change

1. Rotate the zoom adjustment ring ④ to change magnification to the desired level. (Fig. 9)
- **While rotating, the zoom ring clicks into fixed positions. This is an indication that the standard magnification positions are obtained.**

Only for IS-01SMD

- **By sliding the 3D attachment, the system allows a quick switching between Straight view and 3D view. (Fig. 11)**
 - **Once inserted into the light path, 3D attachment can be rotated by 360° in order to obtain a oblique observation of the whole specimen.**
1. While looking on the monitor, rotate the 3D attachment until the desired image is achieved.


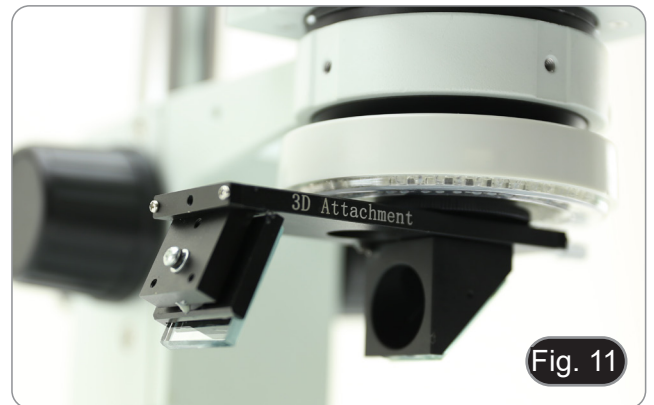
Fig. 11

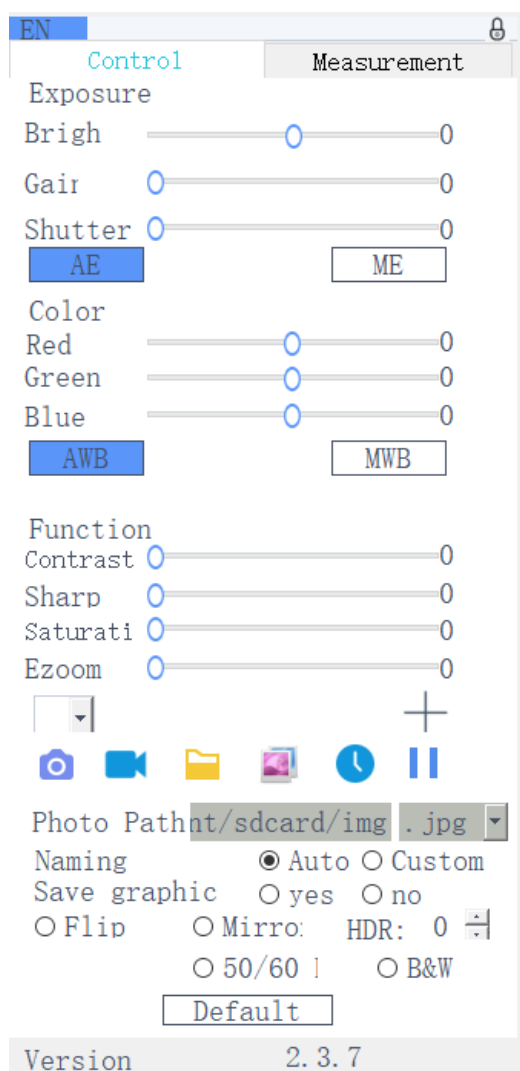10. Use of the software

10.1 Menu Interface

a) Control

b) Measurement

EN Control Measurement

Assist Tools  

Main Tool

Calibration

Calibrate default mm

Scan Data

Custom Template

Name default   

List Show Export Data

ID	Item	Result
----	------	--------

Version 2.3.7

10.1.1 Detailed Control Interface


a) Language Selection


Click the upper left corner of the menu to select the language.
Simplified Chinese, English, Traditional Chinese, Japanese and Russian are available.

b) Control Panel Lock

Click the lock icon in the upper right corner of the menu to switch the menu display status.

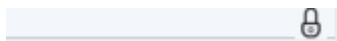
When “*Locked*”, always show menu.

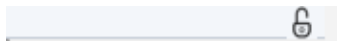When “*Unlocked*”, the menu will be hidden/shown when the mouse is moved to the top left corner of the monitor.

c) Exposure

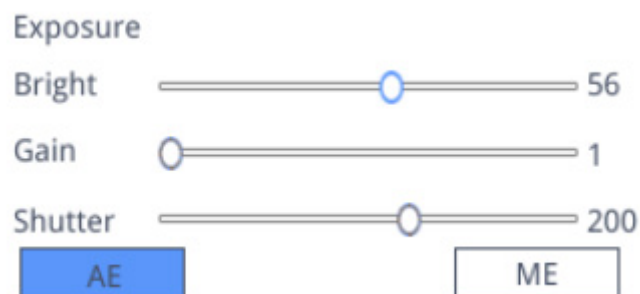
Exposure setting can be *Auto* or *Manual*.

When “*AE*” is selected, user can drag the “*Bright*” slider to adjust the brightness.
When “*ME*” is selected, the “*Gain*” slider can be moved to adjust the gain. Move the “*Shutter*” slider bar to adjust the shutter.
The default value is *AE* (Auto Exposure).

d) White Balance

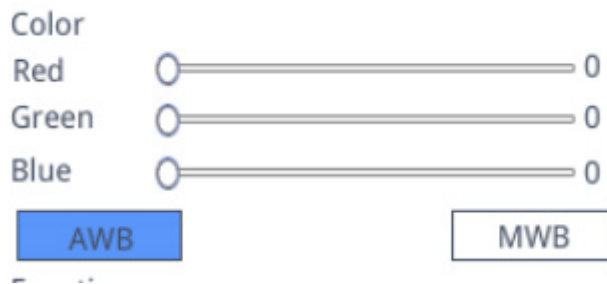White Balance setting can be *Auto* or *Manual*.

AWB: automatic white balance can be achieved under different color temperatures.

MWB: user can move "Red / Green / Blue" sliders to adjust the "Red / Green / Blue" color value.

e) Function

Function setting can be *Contrast* / *Sharp* / *Saturation* / *Ezoom*.

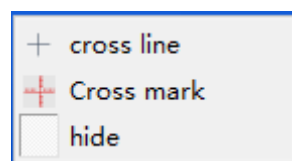
Contrast: adjustable value 0-15. Default value 0.

Sharp: adjustable value 0-15. Default value 4.

Saturation: adjustable value 0-254. Default value 128.

Ezoom: adjustable value 1-6. Default value 1.0.

f) Crosshair Tool


Cross Line

1. Click Crosshair tool .
2. Select "Cross line" to set the number of horizontal and vertical lines.

3. After the setting is complete, select "Display" to show the set crosshair in the image display area, or select "Delete" to not show crosshair.

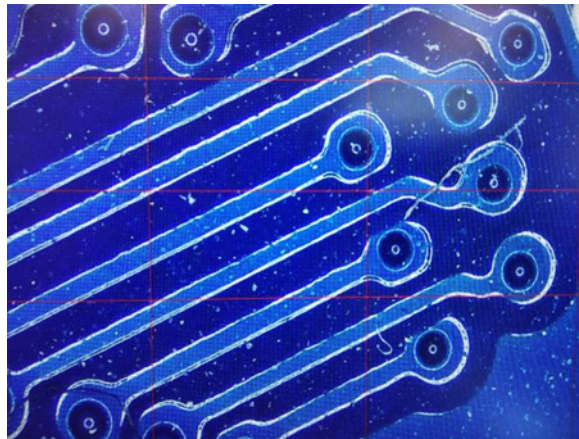Display crosshair:

While selecting one line, hold the left mouse button: user can drag the crosshair.

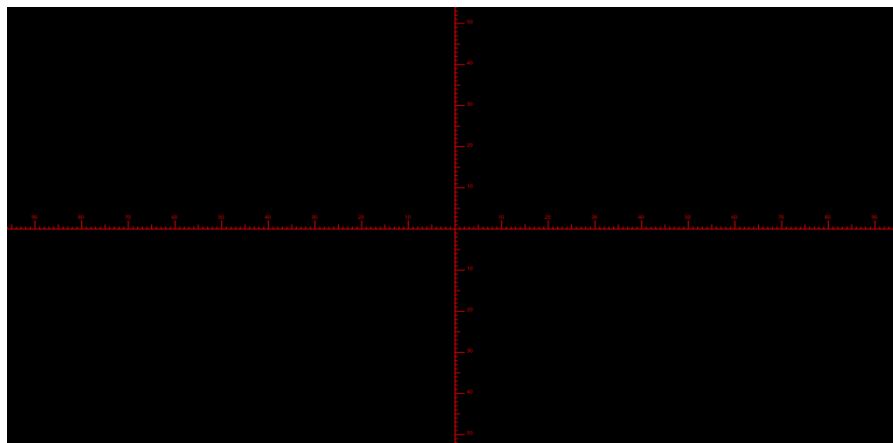
Note: Do not press and hold at the intersection of the horizontal and vertical lines.

Double-click a single crosshair to set its color and position, center and delete.

Cross Mark

1. Click Crosshair tool .
2. Select "Cross mark".
3. The image area will show a crosshair with notches.

Cross Hide

1. Click Crosshair tool .
2. Select "Hide". All cross lines or the crosshair with notches will be hidden in the image area.

Line color

1. Click on the set crosshair color tool .
2. Pull down to set the crosshair color. After the color is set, draw the crosshair again to take effect.
 - **Note: the color of the previously drawn crosshair does not change.**

10.1.2 Image capture

a) USB flash drive insertion

First confirm whether the USB flash drive is successfully inserted.

1. Insert the USB flash drive into the USB slot on the back side of the microscope body.
2. After inserting the USB flash drive, the icon of the successful insertion of the USB flash drive  is displayed for a few seconds in the upper right corner of the display.
3. Now it is possible to take an image.

b) Image capture operation

1. Click the "Capture"  icon.
2. The icon  will be displayed in the center of the screen. After about 3 seconds, the icon disappears and the picture is automatically saved in the `/mnt/sdcard/img` folder of the USB flash drive.
 - **Note: the USB flash drive must be formatted in FAT 32 mode. Any other formatting mode will not allow images to be saved correctly.**


c) Prefix

When the "Naming" is "Auto", the prefix of the file name can be added, and the file is named and saved with "prefix + serial number".

Photo path

d) Image Format

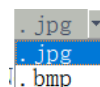
Click on the image format drop-down list to select the image format. There are two formats to choose from: .jpg or .bmp.

e) Image Naming Rules

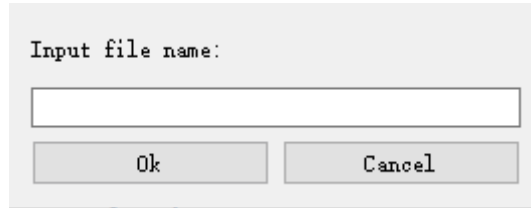
Naming Auto Custom

1. Select "Auto" to name the image according to the time.
2. Select "Custom" to enter a name for saving.

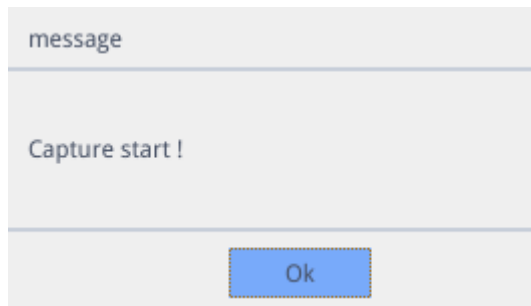10.1.3 Recording

1. Click on "Record"  icon.
2. The dialog box "Capture Start!" appears.

3. Click "OK" to start recording.
4. During recording, the icon  in the lower right corner of the screen flashes.
5. Click on "Record"  icon again.
6. The dialog box "Capture Stop!" appears.

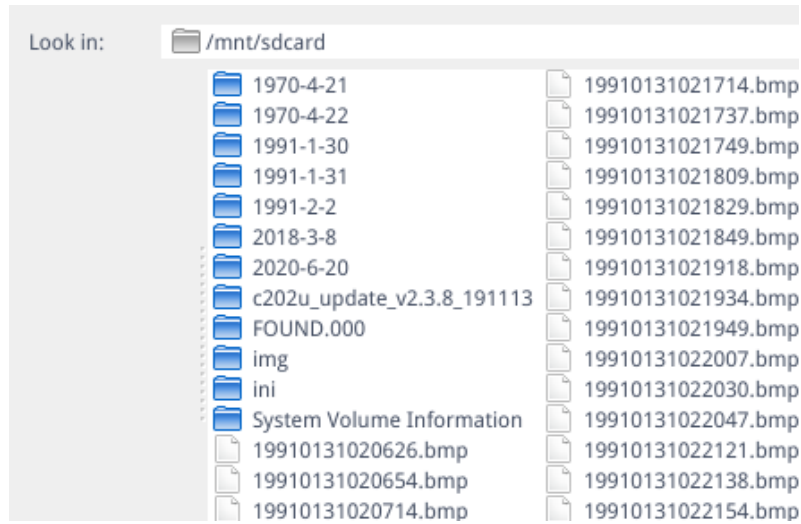
7. Click "OK" to stop recording.

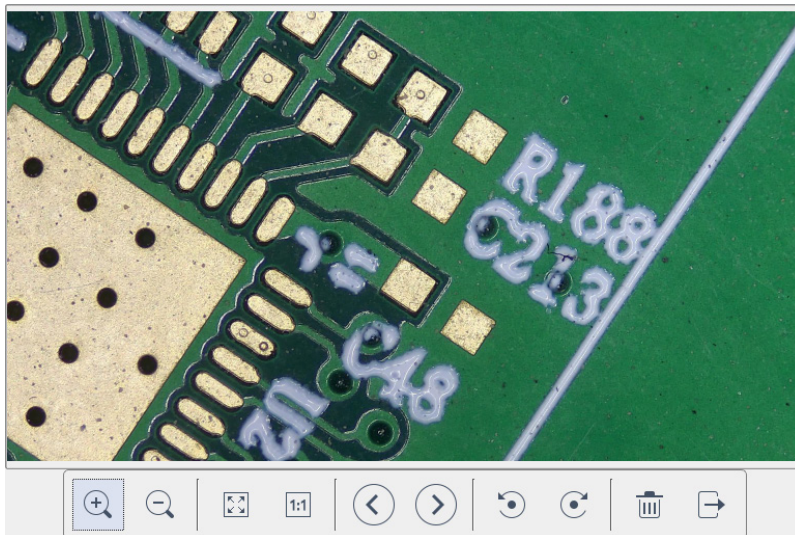
Note: The recorded videos are named according to the date, and are automatically saved in the flash drive with the format of .H264. The maximum length of a single recording is about 2 hours and the size is about 4G. If it exceeds 2 hours, it will automatically stop recording.

10.1.4 Open Image

1. Click on “Open”  icon.
2. The dialog box opens to select the desired image.

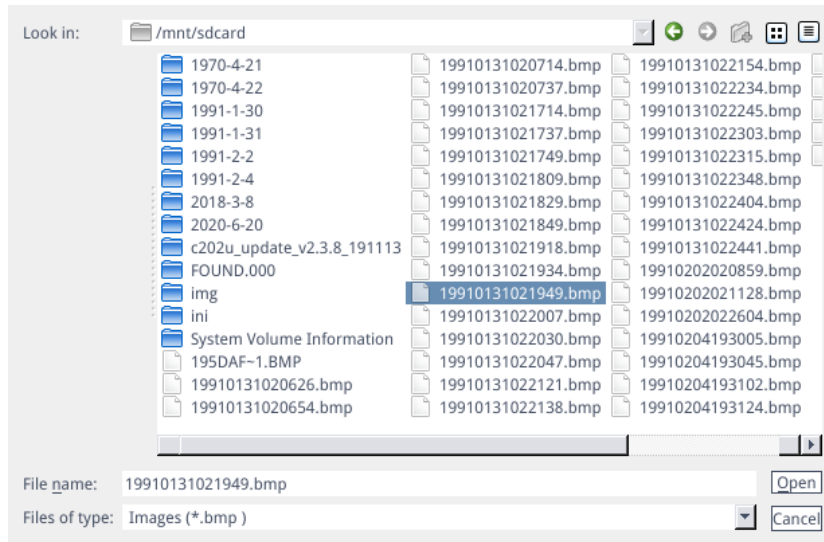
3. Select the image you want to open. The image is opened.

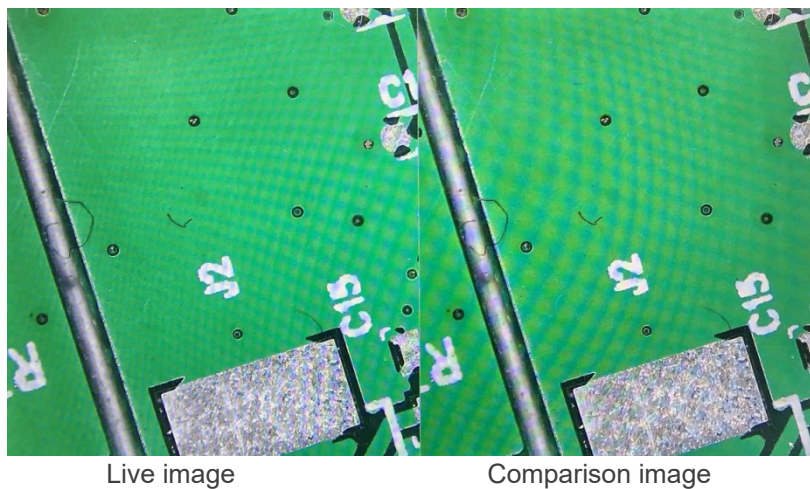
- Click “Scaling”  . User can zoom in and out in the image.
- Click “Full Screen” . The image interface displays the whole image.
- Click “1:1” . The image interface displays part of the picture. Press and hold the left mouse button to move the picture position.
- Click “Browse”   to browse the previous and next images.
- Click “Rotate”  . User can change the image rotation clockwise or counterclockwise.
- Click “Delete”  to delete current image.
- Click “Return”  to close the current image interface.

10.1.5 Image comparison

1. Click the “Compare Picture”  icon.
2. The dialog box opens to select the desired image.

3. Select the picture you want to compare and click “Open”.
4. The image comparison interface appears. The left side of the screen is the real-time video interface, and the right side is the selected image.

5. Click “Compare Picture”  icon. Image comparison function closes.
- Note: Only images in .bmp format are supported for image comparison.**

10.1.6 Freeze

1. Click “Freeze”  icon.
2. The current image is frozen, and the image window remains static, which makes it easy to observe the measured object.
3. After freezing, the icon changes into .
4. Click icon  to restore the real-time display.

10.1.7 Save Graphics

Save graphic yes no

1. When “Yes” is checked, the upper measurement line is saved in the picture.
2. When “No” is checked no measurement is saved.

10.1.8 Flip / Mirror

Flip Mirror

1. Check "Flip": the vertical direction of the image on the screen is the opposite of the current.
2. Check "Mirror": the horizontal direction of the image on the screen is the opposite of the current.

10.1.9 HDR (High Dynamic Range)

HDR: 0 

HDR: adjustable value 0-3. Default value 1.

10.1.10 Anti-screen Flicker

50/60 HZ

Anti-screen flicker effect is achieved after ticking.

10.1.11 Black and White Mode

B&W

Check for black and white mode, default is color mode.

10.1.12 Default

Click "Default" to restore various parameters to factory settings.

10.1.13 Version Number

You can view the current software version number.

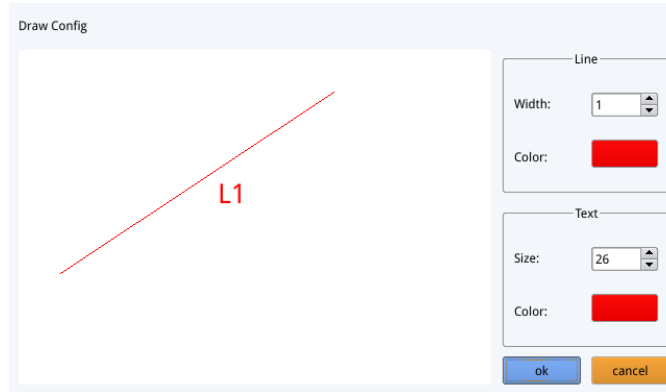
10.2 Detailed Measurement Interface

Click “*Measurement*” to carry out the measurement operation.
 After the measurement operation is completed, if you want to display the menu, click the right mouse button in the image area and move the mouse to the upper left corner of the screen.

10.2.1 Assist Tool

a) Color Setting

1. Click “*Color setting*”  icon.
2. Setting dialog box appears.

- In the “*Line*” field user can set width and color of the line.
- In the “*Text*” field user can set the font size and color of the label.

b) Edge Detection

1. Click “*Edge Detection*”  icon.
2. When selected, the icon turns blue .
3. When measuring an image, the software will automatically find the edge of the object to be measured.

10.2.2 Measurement Tool

Icon	Function	Icon	Function
	Arbitrary line		Arc
	Horizontal line		Circle
	Vertical line		Concentric circles
	Rectangle		Double circle
	Parallel lines		Line circle
	Perpendicular		Perpendicular circle
	Angle		Text
	Polygon		

10.2.3 Scan Data

This function requires a scanner.

After the scanner scans the barcode, the scanned data box displays the scanned data. A folder named after this data is automatically saved in the USB flash drive, and the picture is saved in this folder.

10.2.4 Custom Template

1. After drawing the primitives in the image area, click "New"  icon.
2. The dialog "Whether to convert a primitive to a template?" appears.

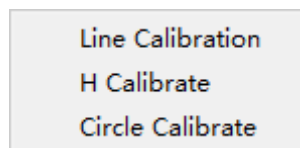
3. Click "OK" to create a new template.
 4. After entering a name in the dialog box, click "Save"  icon to save the template.
- Click "Delete"  icon to delete this template.
 - Click "Edit"  icon to modify the template.
 - Once modification has done, click "Save"  icon again to save the modified template.
 - Click "Back"  icon: this will not save the modified template.

10.2.5 Measurement Operation

a) Calibration

1. Click on the calibration tool drop down.

2. Line Calibration, H calibrate or Circle Calibrate can be chosen.
3. Drag the marking line to align with the engraved line of the scale, enter calibration name (eg. 4x) and length; select the unit, and complete the calibration.

Name Length Unit

 mm ▼

RealPixel 449.755 pixel

RealRuler

Ok Cancel

7	8	9	Del
4	5	6	
1	2	3	Clear
0	.	X	

b) Delete calibration

1. Click the drop-down calibration list, select the calibration group to be deleted, and click the delete  icon to delete this group of calibrations.

c) Measuring

The measurement tool includes a variety of stitch measurements. Depending on the specific application, user must first select the appropriate calibration and unit.

1. Select the appropriate tool for measurement; when measuring, user only needs to click the points of the object to be measured.
2. After you click and release the mouse, the measurement data is displayed in the measurement data list area.


ID	Item	Result
1	Circle	Rad=136.400pixel CLen=857.027pixel Area=58449.156pixel

3. After the measurement is completed, you can export the data in the measurement data list.
4. Click "Export Data", and the data will be automatically saved in the .csv file in the USB flash drive. The user can input the file name.

Input file name:


Ok Cancel

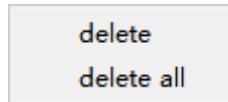d) Delete Measurement Data

1. In the measurement data list, select the data you want to delete.
2. Right-click with the mouse button: user will be prompted to delete a group of data or delete all data.
3. Select according to your needs.

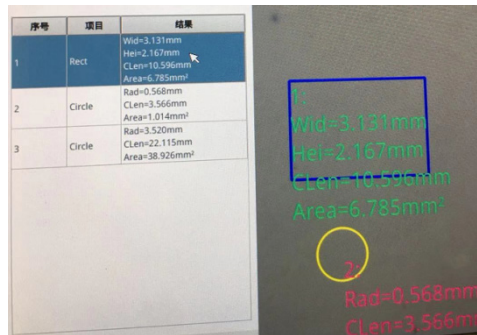e) Select Measurement Data

1. Click the data in the measurement data list with the left mouse button.
2. The selected graph data is displayed in shades of gray.
 - The color of the selected graph in the image display area becomes the reverse of the set color.

f) Select Graph

1. Double-click the data in the measurement data list with the left mouse button.
2. Selected graph data will pop up a dialog box.
3. User can modify the color and line width of the line, and lock the selected graph.
 - The locked graph can no longer be operated.

g) Detailed Data

Double-click the “*List Show*” List Show icon with the left mouse button to pop up the detailed data dialog box to view all the current detailed measurement data.

name	distance	perimeter	area	angle	radian	width	height	radius

h) Mark Tool

1. Click “*Text*” **T** icon.
2. Draw the position to be marked in the image area, enter the comment content in the pop-up window, and click “*OK*” to complete the marking.

Please enter the comment content:

Ok Cancel

11. Maintenance

Microscopy environment

This microscope is recommended to be used in a clean, dry and shock free environment with a temperature of 5°-40°C and a maximum relative humidity of 85 % (non condensing). Use a dehumidifier if needed.

To think about when and after using the microscope

- The microscope should always be kept vertically when moving it and be careful so that no moving parts, such as the eyepieces, fall out.
- Never mishandle or impose unnecessary force on the microscope.
- Never attempt to service the microscope yourself.
- After use, turn off the light immediately, cover the microscope with the provided dust-cover, and keep it in a dry and clean place.

Electrical safety precautions

- Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off-position.
- Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

Cleaning the optics

- If the optical parts need to be cleaned try first to: use compressed air.
- If that is not sufficient: use a soft lint-free piece of cloth with water and a mild detergent.
- And as a final option: use the piece of cloth moistened with a 3:7 mixture of ethanol and ether.
- **Note: ethanol and ether are highly flammable liquids. Do not use them near a heat source, near sparks or near electric equipment. Use these chemicals in a well ventilated room.**
- Remember to never wipe the surface of any optical items with your hands. Fingerprints can damage the optics.
- Do not disassemble objectives or eyepieces in attempt to clean them.

For the best results, use the OPTIKA cleaning kit (see catalogue).

If you need to send the microscope to Optika for maintenance, please use the original packaging.


12. Troubleshooting

Review the information in the table below to solve operating problems.

PROBLEM	CAUSE	SOLUTION
I. Optical Section:		
LED operates, but field of view remains dark.	Power supply is unplugged.	Connect
	Brightness is too low	Set brightness to a proper level
Dirt or dust is visible in the field of view.	Dirt/dust on the specimen	Clean the specimen
	Dirt/dust on the lenses	Clean the lenses
The image is not shown on the screen	Power supply is unplugged.	Connect
II. Mechanical Section:		
Coarse focus knob is hard to turn	Tension adjustment ring is too tight	Loosen tension adjustment ring
Focus is unstable	Tension adjustment ring is too loose	Tighten tension adjustment ring
III. Electrical Section:		
LED doesn't turn on.	Power supply not connected	Check for proper connection
Brightness is not enough	Brightness setting is too low	Adjust brightness
Light blinks	Power supply not well connected	Check for proper connection

Equipment disposal

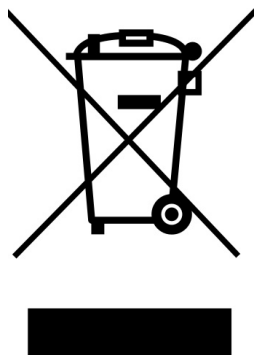Art.13 Dlsg 25 July 2005 N°151. “According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal.”

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.