Solar Battery Pack

# INSTRUCTION MANUAL

<b>Model</b>
M-069

Version: 6  
Issued: 11, 07, 2017

---

## Table of Contents

**Safety information**

**Package contents**

**Symbols and conventions**

**Instructions for use**

*Product Specifications*

*Overview*

*Description of Buttons and Switches*

**Connection guide**

**Operation**

**Troubleshooting**

**Equipment disposal**

---

## Safety Information

### Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position.

Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

## Package Contents

DESCRIPTION	QUANTITY
Solar battery pack unit	1
USB / Micro-USB cable	1
USB / Jack cable	1

## Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

### CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

### ELECTRICAL SHOCK


This symbol indicates a risk of electrical shock.

## Instructions for use

### Product Specifications

<b>Battery</b>	polymer lithium-ion
<b>Capacity</b>	2600 mAh
<b>Input (Micro-USB port)</b>	5 Vdc
<b>Charging time (through USB)</b>	about 5 hours
<b>Output (USB port)</b>	5.5 Vdc / 1 A
<b>Solar panel power</b>	0.7 W max
<b>Solar panel output current</b>	80 mA (at 1000 W/m <sup>2</sup> light)
<b>Working temperature</b>	-10°C ~ +55°C

### Overview

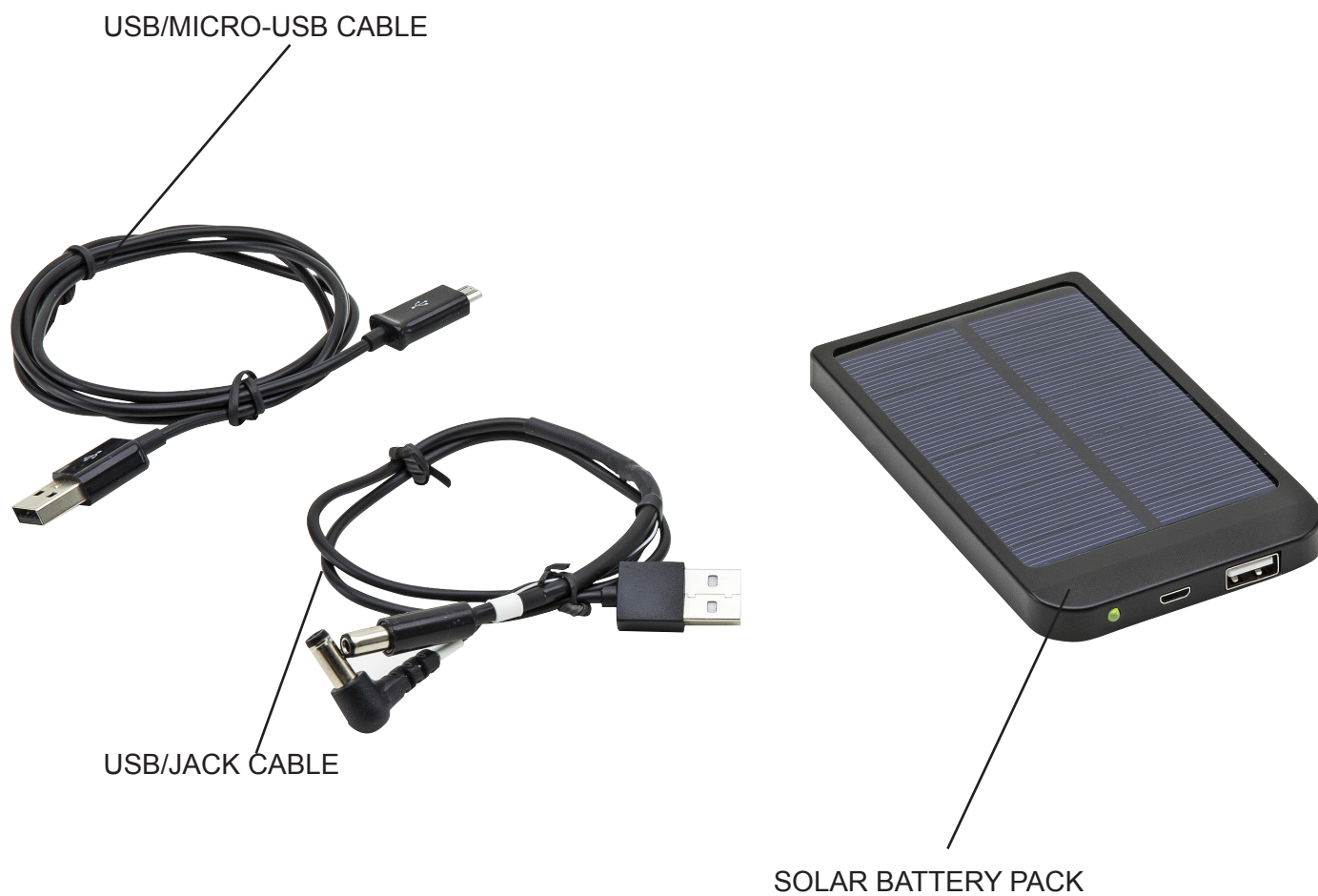---

## Description of Buttons and Switches

## Connection guide

Plug your microscope to the battery pack using the provided USB/Jack cable:

The cable has 2 versions of output jack: 2.1 mm and 2.5 mm internal diameter.  
Check the diameter of the original power supply of your microscope, and use the jack accordingly.

## Operation

**Fully recharge the battery via USB cable and 5 V power supply when using it for the first time.**

Connect the USB/Jack cable to the USB output on the battery pack and to the input Jack of the microscope.

Recharging

- a) Expose the front solar panel to sun, the LED indicator will blink.
- b) Connect via USB/micro-USB cable to an external USB power supply (not included).
- c) Connect via USB/micro-USB cable to a PC USB port.

## Troubleshooting

Review the information in the table below to troubleshoot operating problems.

PROBLEM	CAUSE	SOLUTION
<b>AFTER CHARGING FOR LONG TIME, INDICATORS DON'T SHOW FULL CAPACITY.</b>	Charging time is not enough	Charge it 8 hours continuously
	Power supply or battery pack is faulty	Refer to technical assistance
<b>MICROSCOPE LIGHT IS NOT WORKING.</b>	No power left	Recharge the battery
	Something wrong with cables and connections	Check that USB is well inserted into the output connector of the battery. Check that the jack connector is well inserted into the input of the microscope.
	Wrong settings of microscope switches	Check that the main power switch of the microscope is in ON position. Check that the brightness adjustment knob of the microscope is set to the desired value.

---

## Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection.

The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment.

Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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**OPTIKA® S.r.l.**