ACCESSORIES Series

# INSTRUCTION MANUAL

Model
M-1037

Ver. 1.3 2024

---

## Table of contents

1.	Warning	3
2.	Package content	3
3.	Unpacking	3
4.	Intended use	3
5.	Analysis principle	4
6.	Assembling	5
7.	Use of the system	6
7.1	Light extinction	6
7.2	Gout analysis	6

## 1. Warning

This system is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.

## 2. Package content

- ① Gout attachment
- ② Polarizer

- ③ Analyzer
- ④ Tint plate "1<sup>st</sup> Order Red"

## 3. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

## 4. Intended use

### Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

### IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

---

## 5. Analysis principle

Gout is a metabolic disease, characterized by recurrent attacks of arthritis (pain, redness and swelling in the joints) caused by the deposit of uric acid crystals in the joints.


Gout occurs when uric acid crystals are deposited in the joints, causing inflammation and intense pain.

The crystals are formed due to the high concentration of uric acid in the blood. Usually, uric acid in the blood is eliminated through the kidneys; excessive accumulation (hyperuricemia) can result from the fact that the body produces too much uric acid or that the kidneys cannot dispose of enough uric acid through the urine.

The diagnosis is made by analysis of the joint fluid: a sample of the fluid (arthrocentesis) can be taken by the doctor under local anesthesia using a needle, to be subjected to microscopic analysis.

## 6. Assembling

1. Insert the gout attachment ① on the frame and lock the locking screw ② with the provided Allen wrench. (Fig. 1)

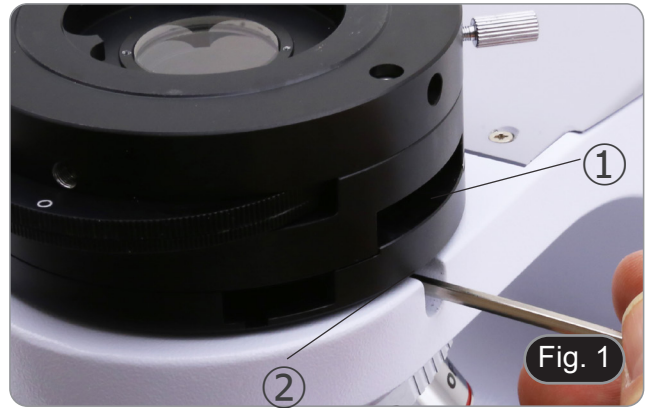2. Insert the optical head on the gout attachment lens and lock the locking screw with the provided Allen wrench. (Fig. 2)

3. Insert the analyzer into the gout attachment. (Fig. 3)

4. Place the polarizer ④ on the light exit at the base of the microscope. (Fig. 4)

## 7. Use of the system

### 7.1 Light extinction

1. Remove the slide from the light path and insert 10X objective.
2. Insert rotatable analyzer in the light path, loosen the locking screw ① and put the vibration scale on 0° ②, then lock the locking screw ①. (Fig. 5)
3. While looking inside the eyepieces, rotate the polarizer to obtain total extinction (total dark in the eyepieces).
4. Once the dark is achieved ("extinction" or "Crossed Nicol" position) it is possible to begin the observation.

### 7.2 Gout analysis

1. Insert 40X objective.
  2. Put a specimen on the stage and focus.
  3. Insert in the light path the "1<sup>st</sup> Order Red" tint plate ③ putting it into the dedicated slot ④. (Fig. 6)
  4. Rotate the specimen until the longitudinal direction of the crystals is parallel to the direction of the axis of the tint plate.
  5. Begin observation.
- If the crystals appear yellow: SODIUM URATE (gout crystals).
  - If the crystals appear blue: POTASSIUM PYROPHOSPHATE (pseudo gout crystals).

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**OPTIKA® S.r.l.**