ACCESSORIES Series

# INSTRUCTION MANUAL

<b>Model</b>
M-870

Ver. 2.1 2022

---

## Table of Contents

1.	Warning	3
2.	Safety Information	3
3.	Package content	4
4.	Unpacking	5
5.	Intended use	5
6.	Symbols and conventions	5
7.	Instrument description	6
8.	Assembling	7
8.1	B-1000MET	7
8.2	IM-5MET	7
9.	DIC observation	8
9.1	DIC observation principle	8
9.2	B-1000MET	9
9.3	IM-5MET	10

---

## 1. Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.

## 2. Safety Information

### Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

---

### 3. Package content

① DIC Prism

---

## 4. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

## 5. Intended use

### Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

### IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

## 6. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

### CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

### ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

## 7. Instrument description

- The DIC prism has a lever that is used to lower or raise the position of the prism within the device.
- This position varies depending on which type of objective is being used.
- In the case of standard objectives the lever must be kept in, while in the case of LWD objectives the lever must be pulled out.
- This is necessary to optimize observation in DIC.
- An incorrect prism position could lead to poor image quality or interference fringes within the image.

DIC PRISM USED WITH  
STANDARD OBJECTIVES

DIC PRISM USED WITH  
LWD OBJECTIVES

## 8. Assembling

- The DIC prism can be used on both the B-1000MET series upright metallurgical microscope and the IM-5MET inverted metallurgical microscope.
- The insertion direction of the prism is different for the B-1000 and the IM-5MET.

### 8.1 B-1000MET

1. Loosen the clamping knob ① in the nosepiece. (Fig. 1)
2. Remove the dummy slider from the nosepiece.


3. Insert the DIC slider, with the prism lever ② facing right. (Fig. 2)
4. Clamp the prism by fixing the clamping knob in the nosepiece ①.

### 8.2 IM-5MET

1. Loosen the clamping knob in the nosepiece ③. (Fig. 3)
2. Remove the dummy slider from the nosepiece.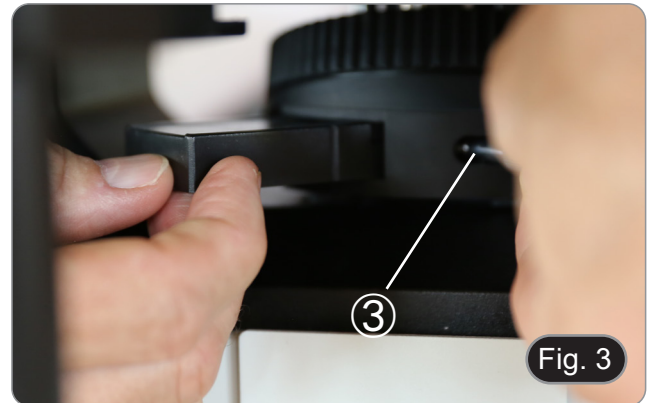

3. Insert the DIC slider, with the prism lever ② facing left. (Fig. 4)
4. Clamp the prism by fixing the clamping knob in the nosepiece ③.

## 9. DIC observation

### 9.1 DIC observation principle

The DIC (Differential Interference Contrast) observation method combines interference and polarization effects and provides more contrasted images with a three-dimensional effect. The use of polarizer and analyzer are needed.

Light (Fig. 5), coming through the polarizer (a), is deflected by the brightfield cube (b) and enters the DIC prism (c). The prism serves as beam splitter and divides the rays (d) into pairs vibrating at 90 degrees to each other with one set of rays traveling at a slower speed than the other.

The light then goes to the specimen (e).

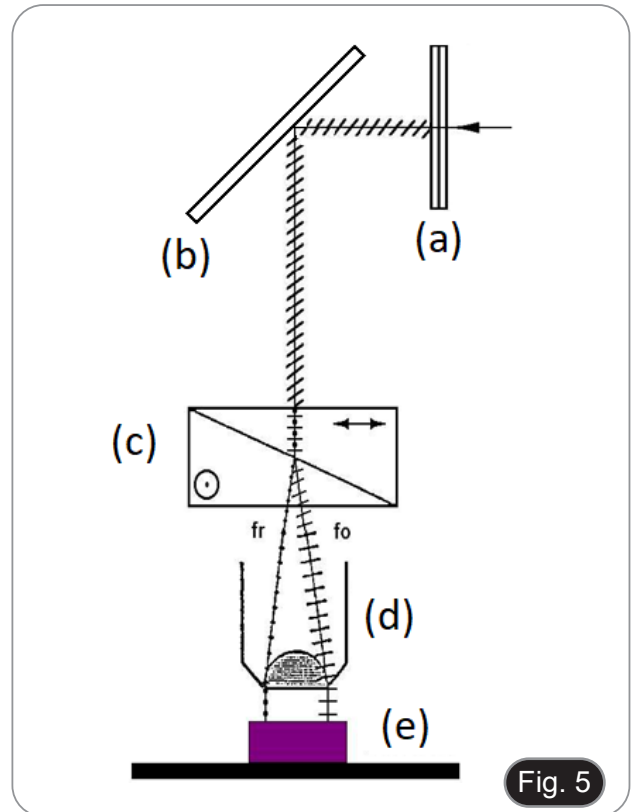
Fig. 5

The light is reflected back into the objective (Fig. 6) and once more passes into the DIC prism (c). The net result is differences in optical path, depending on the original path difference, the specimen's effect and the effect of the prism serving as a beam recombiner in the upward path.

The beams emerging from the prism cannot interfere with one another until they are brought into the same plane of vibration by means of the analyzer (g) which is in place above the brightfield cube (b) as well as above the prism.

The net result of all of this is to produce a so-called "Nomarski-type image" which appears three-dimensional with one side of individual features appearing less bright than the other. This pseudo-3-dimensionality results from differences in light intensity caused by the optical path differences described above.

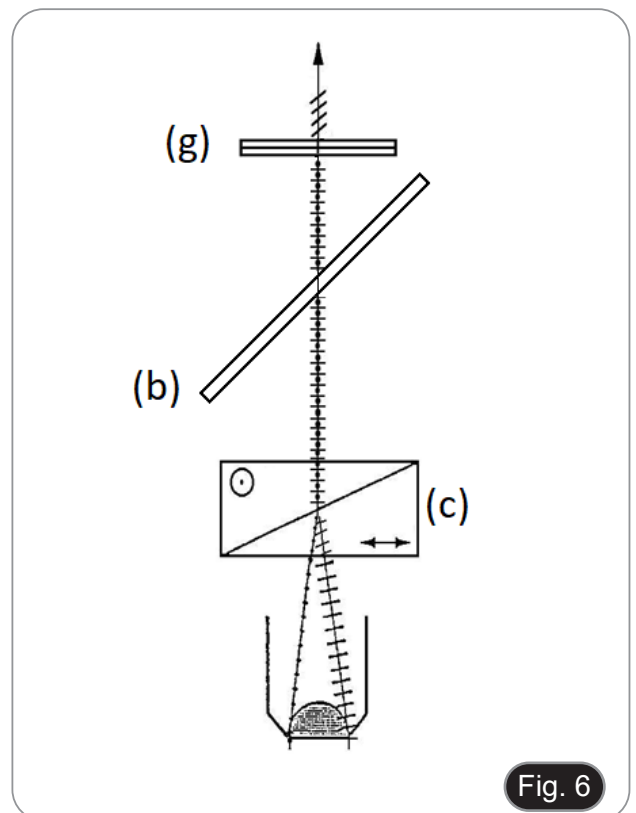
Fig. 6

## 9.2 B-1000MET

- As with observation in polarized light, DIC observation cannot be performed when the BF/DF switch on the microscope is in the 'DF' position.
1. Move the selector on the reflected light illuminator in the position "BF".
  2. Insert the polarizer "PO" with the inscription facing the user (before remove the empty slider). (Fig. 7)
  3. Move the polarizer "PO" into the light path until the slider clicks into its position.
  4. Insert analyzer "AN" with the inscription facing up (first removing the covering plates on the illuminator).
  5. Move the rotatable analyzer "AN" into the light path until the slider clicks into its position. (Fig. 8-9)

6. Put a flat mirror on the stage and get into approximate focus the mirror surface.
  7. While observing into the eyepieces, rotate the scale of the analyzer ① until you achieve the darkest image. (Fig. 10)
- Now the light extinction (so-called "crossed Nicol position") is achieved.

8. Remove the mirror, place the specimen on the stage and adjust the focus.
9. Adjust the background contrast by turning the prism movement knob ② on the DIC slider. (Fig. 11)
10. When the prism movement knob ② on the DIC slider is turned, the interference color in the background varies continuously from the gray to magenta (from -100 to 600 nm). Set the interference color which can provide best contrast with respect to the specimen.

- If the background color is dark, the observation is similar to darkfield.
- If the background color is grey, three-dimensional high contrast images can be achieved rich in grey shades.
- If the background color is close to magenta, even small phase variation can be observed as a change in color.

Fig. 11

### 9.3 IM-5MET

- **As with observation in polarized light, DIC observation cannot be performed when the BF/DF switch on the microscope is in the 'DF' position.**

1. Move the lever in "BF" position for brightfield illumination.
2. Insert the polarizer "PO" with the inscription facing the user. (Fig. 12)
3. Move the polarizer "PO" into the light path until the slider clicks into its position.

Fig. 12

4. Insert analyzer "AN" with the inscription facing up. (Fig. 13)
5. Move the rotatable analyzer "AN" into the light path until the slider clicks into its position.
6. Put a flat mirror on the stage and get into approximate focus the mirror surface.
7. While observing into the eyepieces, rotate the scale of the analyzer ① until you achieve the darkest image. (Fig. 14)
- Now the light extinction (so-called "crossed Nicol position") is achieved.
8. Remove the mirror, put a specimen on the stage and focus the specimen.
9. Adjust the background contrast by turning the prism movement knob on the DIC slider. (Fig. 11)
10. When the prism movement knob ② on the DIC slider is turned, the interference color in the background varies continuously from the gray sensitive color to magenta sensitive color (from -100 to 600 nm). Set the interference color which can provide best contrast with respect to the specimen.

Fig. 13

- If the background color is dark, the observation is similar to darkfield.
- If the background color is grey, three-dimensional high contrast images can be achieved rich in grey shades.
- If the background color is close to magenta, even small phase variation can be observed as a change in color.

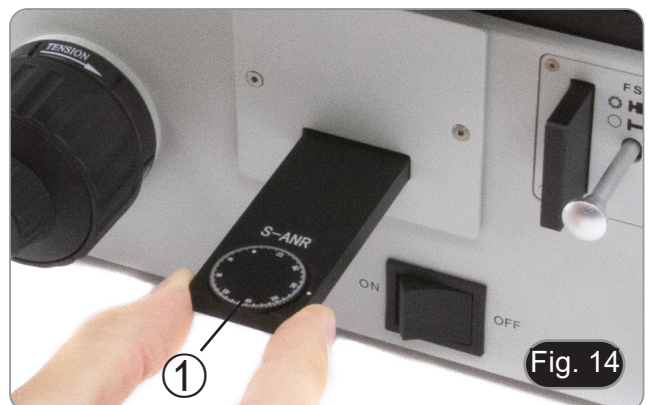
Fig. 14

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**OPTIKA® S.r.l.**