Software Series

# INSTRUCTION MANUAL

Models
Optika Vision Lite

Ver. 1.0 2019

---

## Table of Contents

<b>1.</b>	<b>How to install the software</b>	<b>3</b>
1.1	TB-3W, TB-5W, B-190TB, B-290TB	3
1.2	B-150DM, B-150DB, SFX	3
<b>2.</b>	<b>How to start the software</b>	<b>6</b>
<b>3.</b>	<b>How to select the camera</b>	<b>6</b>
<b>4.</b>	<b>Graphical User Interface (GUI)</b>	<b>7</b>
<b>5.</b>	<b>Tool Bar</b>	<b>7</b>
5.1	Open Report	8
5.2	Save Report as	8
5.3	Print	8
5.4	Activate Live Image	8
5.4.1	Setting the Resolution	9
5.4.2	Setting Camera parameters	10
5.4.3	Snap image	10
5.4.4	Time-Lapse Acquisition	11
5.4.5	Movie Recording	11
5.4.6	Stop recording	12
5.4.7	Full Screen	12
5.5	Import image from file	12
5.6	Export image to file	13
5.7	Flip/Rotate	13
5.8	Measurements	13
5.9	Indicate objects	14
5.10	Measurement line color	14
5.11	Grid	14
5.12	Select calibration	16
<b>6.</b>	<b>Menu bar</b>	<b>17</b>
6.1	File	17
6.1.1	Stop live video	17
6.1.2	Open Report	17
6.1.3	Save Report as	17
6.1.4	Select source and snap image	17
6.1.5	Import image from file	18
6.1.6	Export	19
6.1.7	Print	19
6.1.8	Exit	19
6.2	Tools	20
6.2.1	Flip/Rotate image	20
6.2.2	Indicate objects	20
6.2.3	Linear measurement	21
6.2.4	Calibration	21
6.3	Options	24
6.3.1	Change Language	24
6.3.2	Change print logo	25
6.4	Help	25
6.4.1	Help Index	25
6.4.2	About	25

---

# 1. How to install the software


## 1.1 TB-3W, TB-5W, B-190TB, B-290TB

The software is pre-installed before shipping from the factory. No installation is therefore required.

## 1.2 B-150DM, B-150DB, SFX

1. Operating system requirements: Windows XP, Windows Vista, Windows 7 or Windows 10.
- If your PC is Windows XP: before the installation of your Camera please check if Service Pack2 is installed. If not, please update your XP from the Microsoft web site. If your PC has a USB port less than 2.0, there may be some problems in capturing image from the camera due to the slow transmission.
2. Please insert your CD and run: *\Optika Vision Lite\ setup.exe*
3. *Optika Vision Lite* icon appears on the desktop.
4. Choose from the folder "*drivers*" the model of your microscope, then choose the next options accordingly to the folders displayed on the screen. Now run *driver.exe*. The installer may show warnings about missing Windows certification: ignore such warnings and select "*Continue*".
- **Note : do NOT cancel the installation.**
5. At the end of the installation process restart the PC if requested.
6. Plug the Camera in your PC: the camera will be identified and Windows will recognize the driver which has just been installed from the CD.
7. Ignore any message about Windows compatibility and select "*Continue*".
- **Note : do NOT cancel the installation.**
8. At the end of the identification process, disconnect your Camera and try to plug it again, in order to check if the Driver was installed correctly (no Windows error messages should appear).

- It could happen that the computer reports an error of this kind:

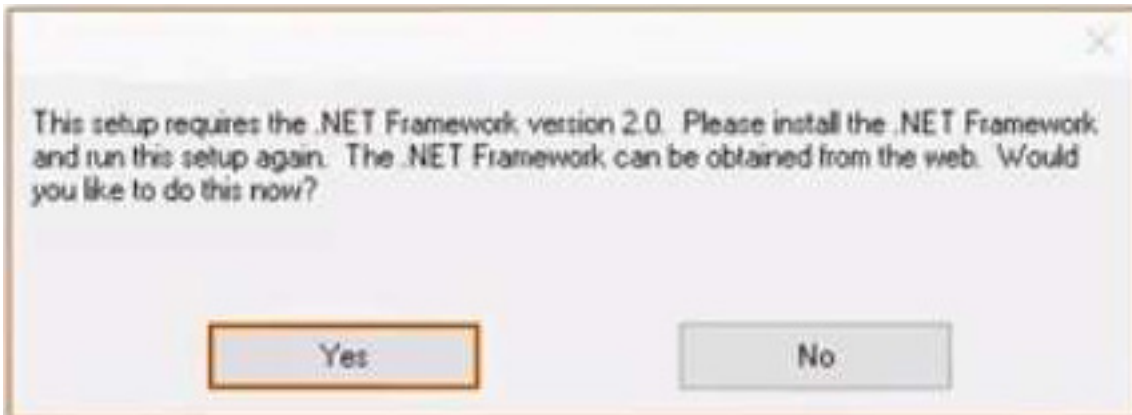
This means that the PC needs to install the .NETframework.

Follow this step-by-step guide to solve the problem:

Click on **“Start”** button


and from there type **“Control Panel”** and select the corresponding item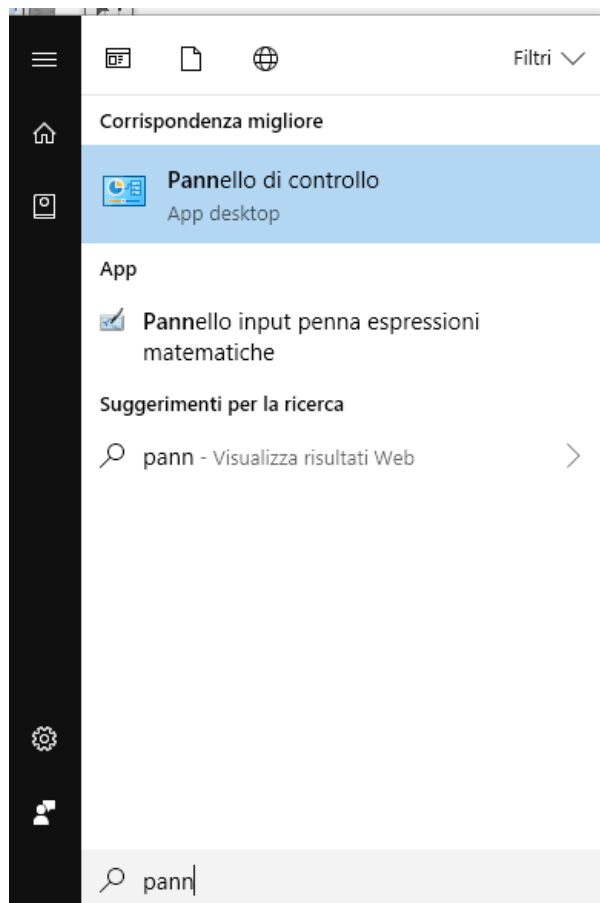


“Control Panel” appears.


Check that the panel display is set to “Category”, then select “Uninstall a program”.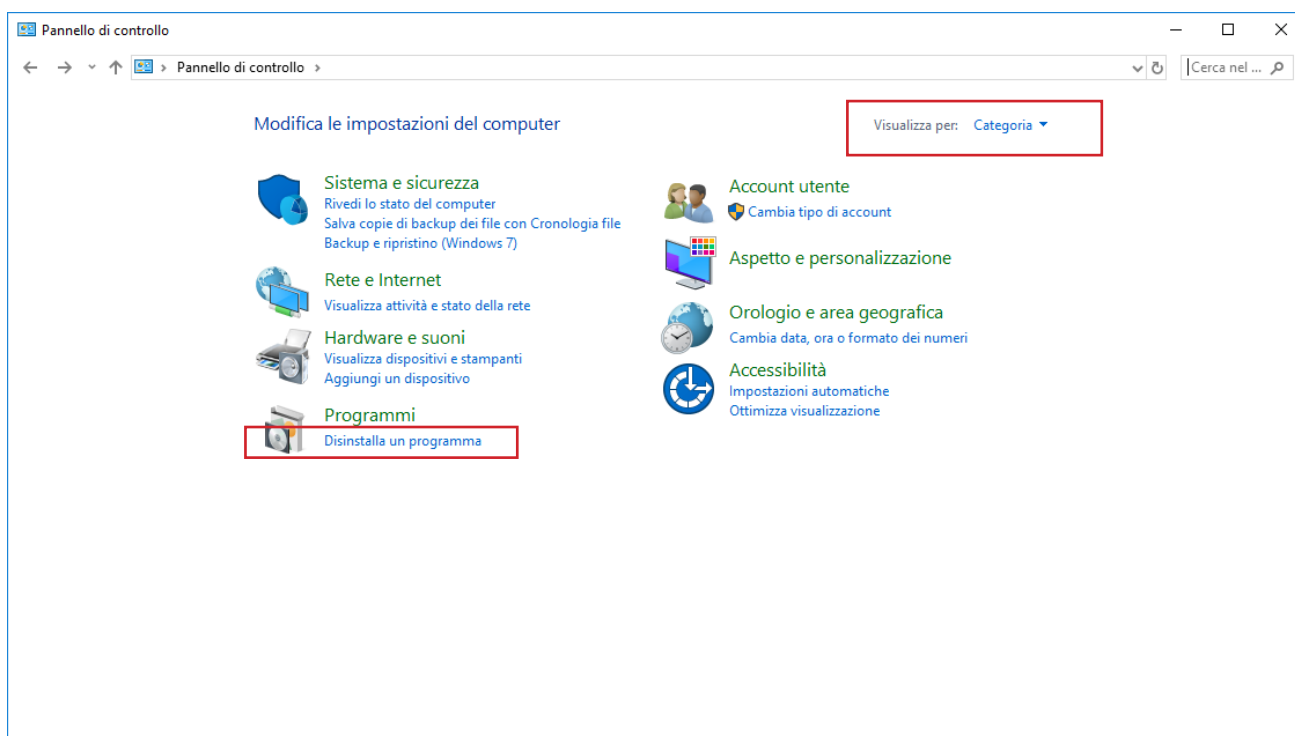

Select “Enabling or disabling Windows functionality”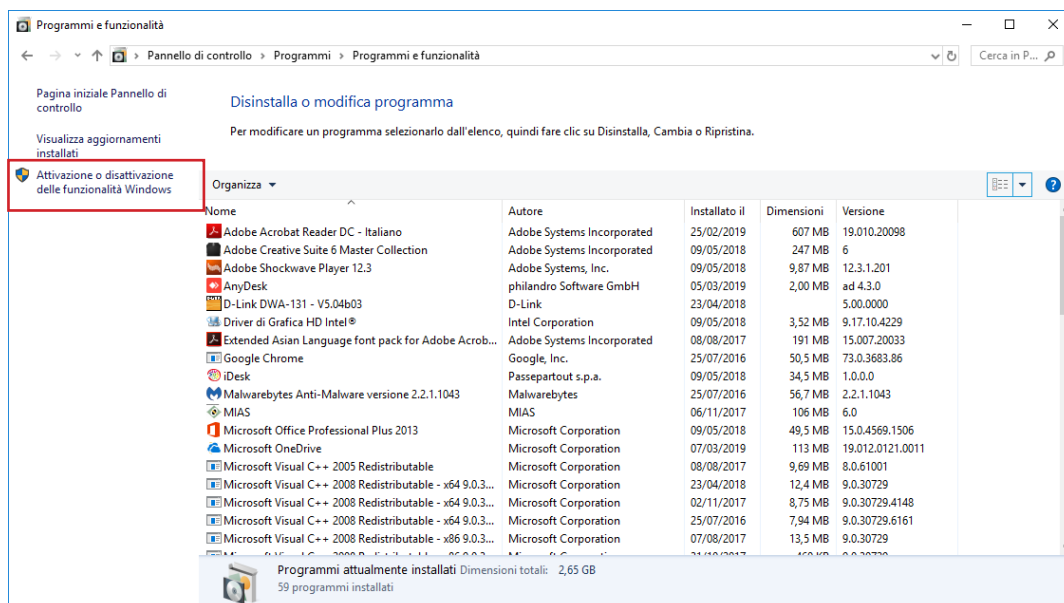

Select the option **“.NET Framework 3.5 (includes .NET 2.0 and 3.0)”** and click on **“OK”**.

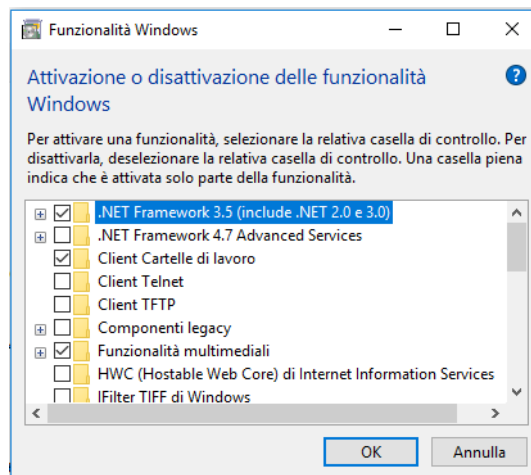Now you start installing the compatibility package. It may happen that the system requires a reboot of the PC.

## 2. How to start the software

From the desktop find the Optika Vision Lite icon, double-click to start the SW.

## 3. How to select the camera


Click on the microscope icon.

A “Device List” dialog will appear listing the various cameras (including any webcam) currently installed on your PC.

Select the one you are interested in and the “live” window will start.

## 4. Graphical User Interface (GUI)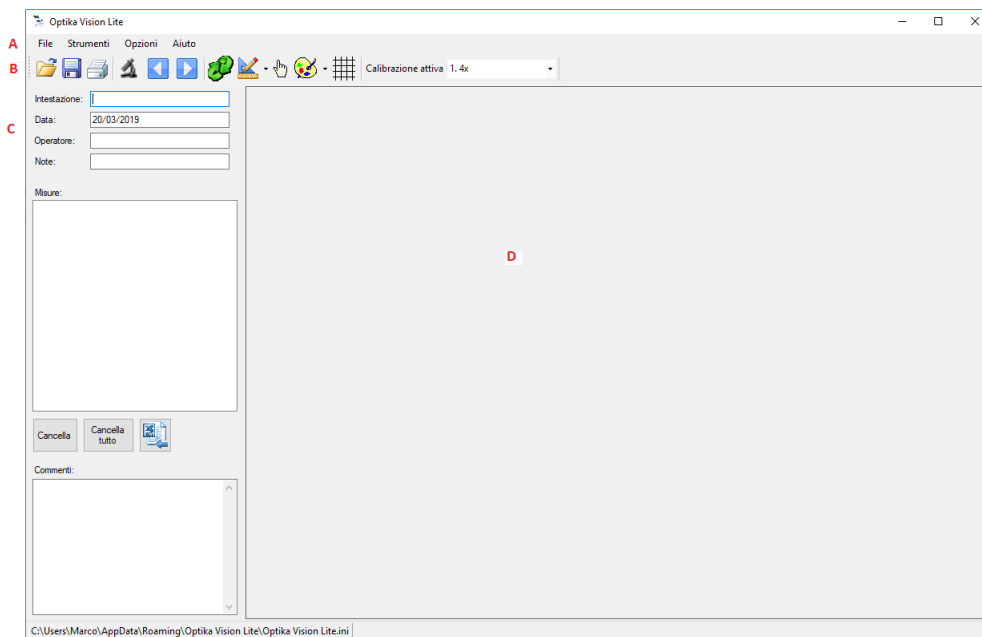

- A: Menu Bar;
- B: Tool Bar
- C: Report Section
- D: Image Window;

## 5. Tool Bar

The Optika Vision Lite toolbar consists of these commands.

- A: Open Report
- B: Save Report;
- C: Print
- D: Open "live" image
- E: Import image from file File
- F: Export image to File
- G: Flip/Rotate Image
- H: Measurement/Calibration
- I: Indicate Object
- J: Color palette for lines
- K: Grid
- L: Calibration list

---

The commands shown in the Toolbar sometimes coincide with those contained in the various menus.

### 5.1 Open Report

A Report is a document in “.asa” format in which a whole series of information has been saved (together with the image).

### 5.2 Save Report as

All the information set in the “Reports” section (Header, Date, Operator, Notes, any Measurements and any Comments) are saved in a report in .asa format.

### 5.3 Print


This command allows you to print a scanned image.

Clicking on this command opens the **Print** dialog box.

### 5.4 Activate Live Image


This command allows you to select the desired image source.

The following dialog opens:

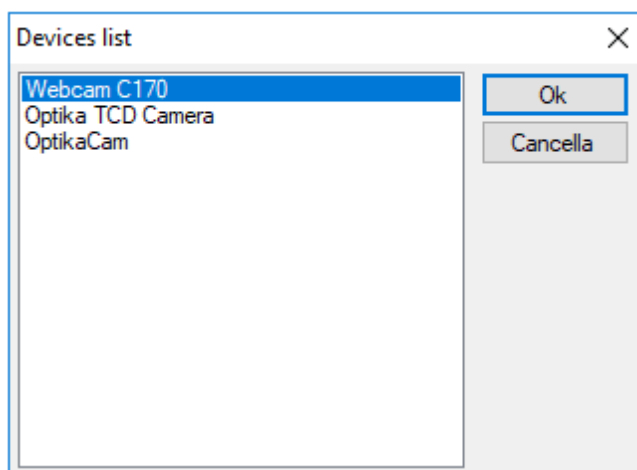Select the desired camera. The live window is activated.

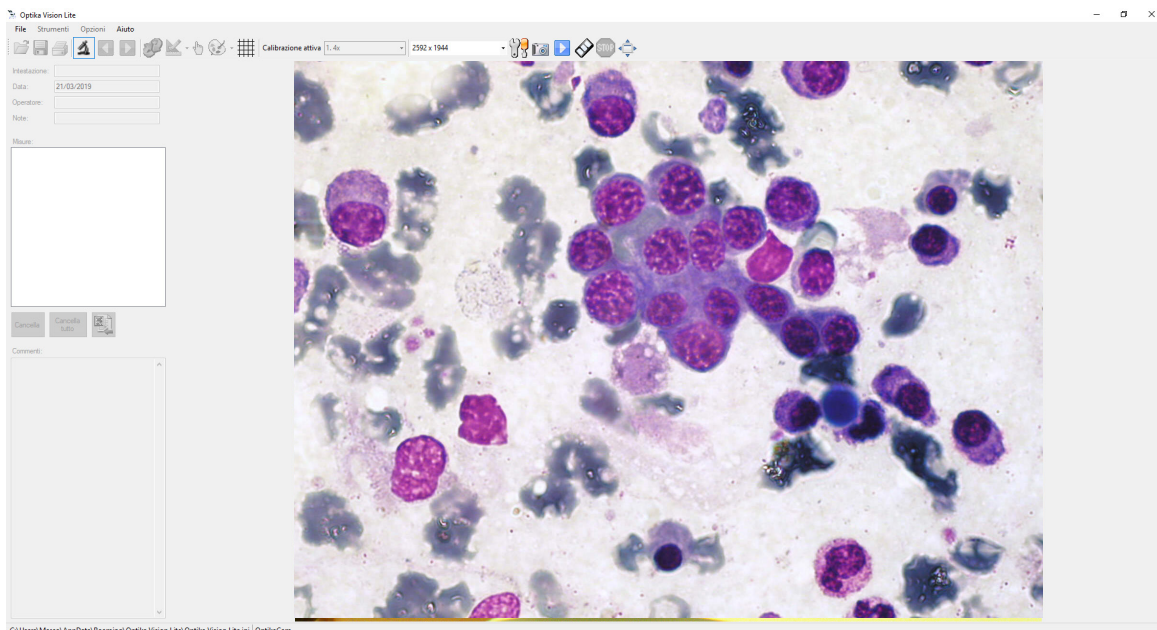
When you activate the “live” image, the toolbar changes by adding a few buttons:

- A: Setting the Resolution;
- B: Setting Camera parameters;
- C: Snap Image;
- D: Recording a movie in “Time-Lapse”;
- E: Recording a movie;
- F: Stop movie recording;
- G: “Full screen” mode


#### 5.4.1 Setting the Resolution

Depending on the type of camera installed, various resolution levels can be selected: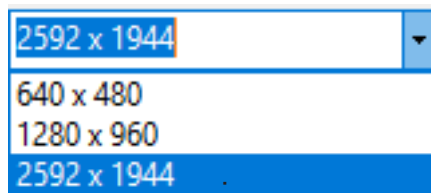

## 5.4.2 Setting Camera parameters

Selecting this command accesses the control panel of all the camera's acquisition functions..

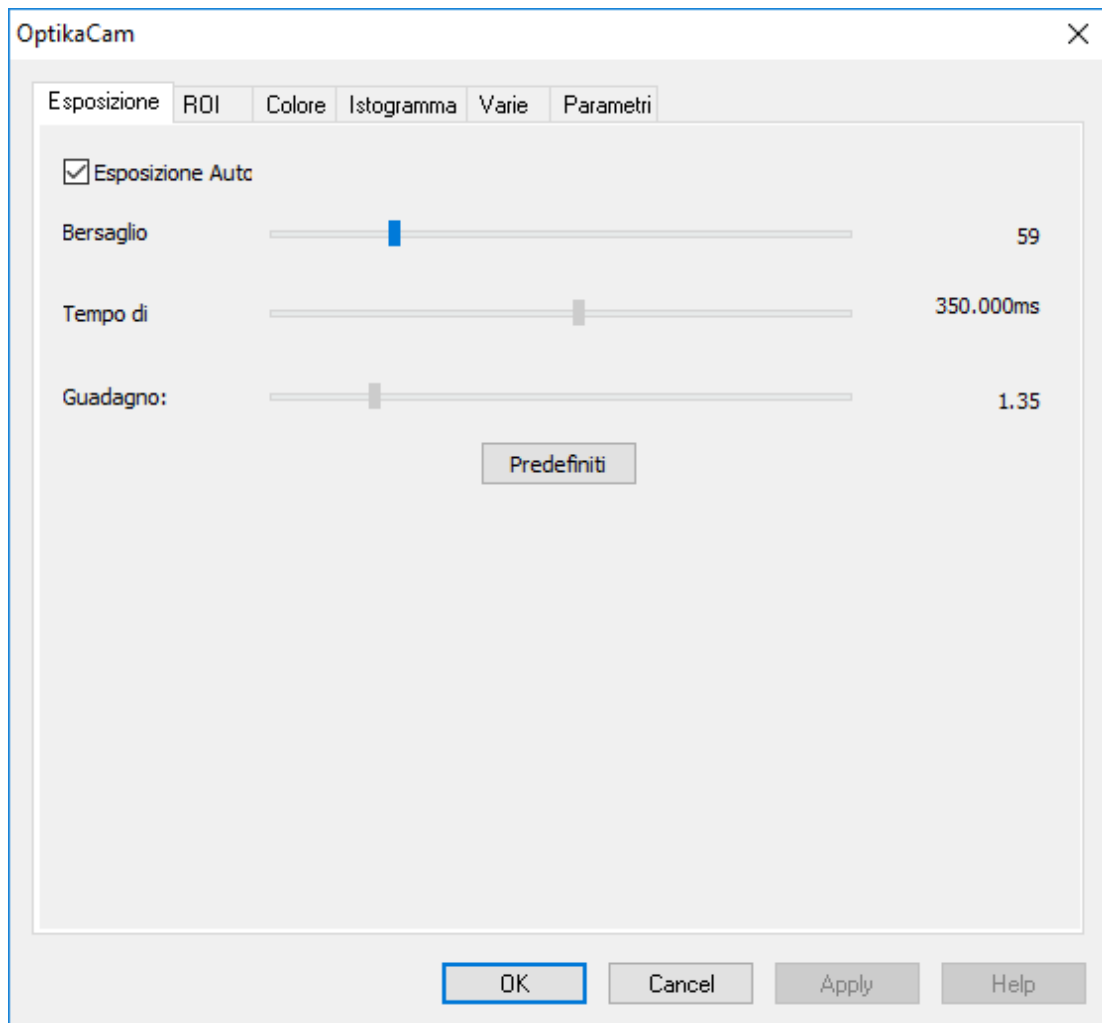Here you can change the parameters of the camera (exposure time, color, etc.) to obtain an image that corresponds to 100% of what you see in the eyepieces in terms of resolution and color fidelity.

## 5.4.3 Snap image

This button is used to acquire the image currently displayed on the live window.


The live preview stops and the newly acquired image is displayed.

To get the live preview again, click on , select the camera and the live view will be reactivated.

#### 5.4.4 Time-Lapse Acquisition


Time-Lapse is a timed acquisition of a number of frames in sequence. Selecting this command accesses the Time-Lapse setting dialog box.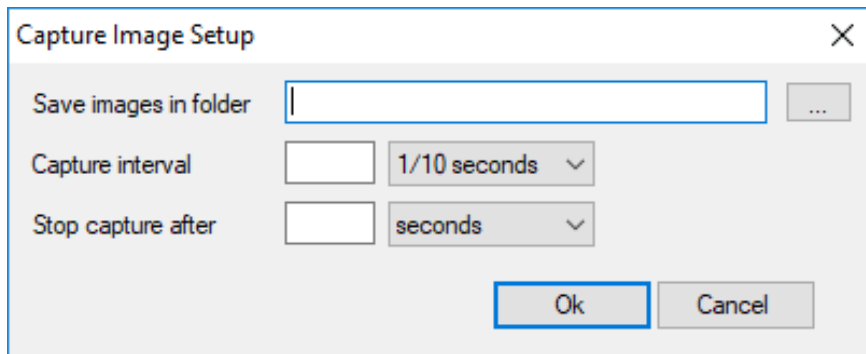


- **Save images in folder:** indicate the destination folder of all images.
- **Capture interval:** indicates the time that elapses between the acquisition of a frame and the next one.
- **Stop capture after:** indicates how long the timed acquisition process must last in total.

Clicking on “OK” starts the process that will automatically stop when the set time expires.

#### 5.4.5 Movie Recording

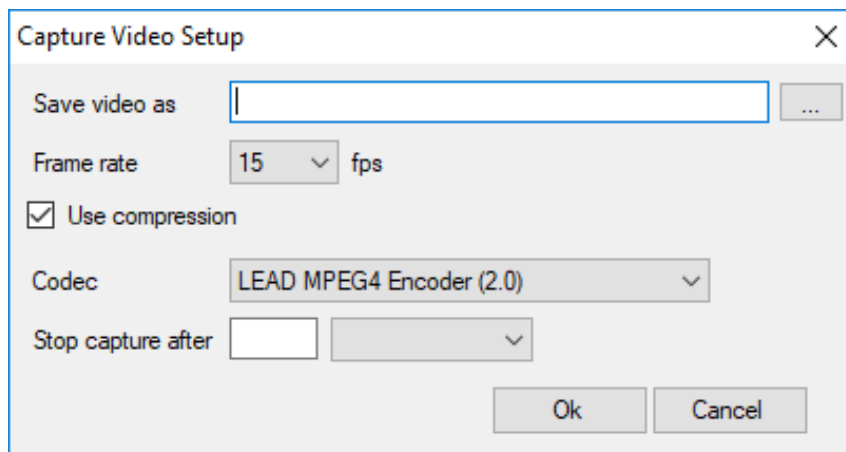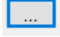- **Save video as:** clicking on the button  next to the editable field, the “Save As” dialog box opens where you can save the file name. The file format will be .avi.
- **Frame Rate:** the number of frames per second (fps) that the camera can capture.
- **Use Compression:** Enabling this option sets which compression protocol you want to use for the movie.
- **Stop capture after:** indicates how long the timed acquisition process must last in total.

Clicking on “OK” starts the process that will automatically stop when the set time expires.

#### 5.4.6 Stop recording

This button is only activated after the acquisition of a movie has been started. To stop the acquisition before the set time has ended, press this key. Movie acquisition stops.

#### 5.4.7 Full Screen


This button allows you to see the live image full screen without displaying the toolbar or menu bar. To return to the standard view, press the ESC key..

#### 5.5 Import image from file

This command allows you to open an image previously saved in any folder on your computer.

The following dialog box opens: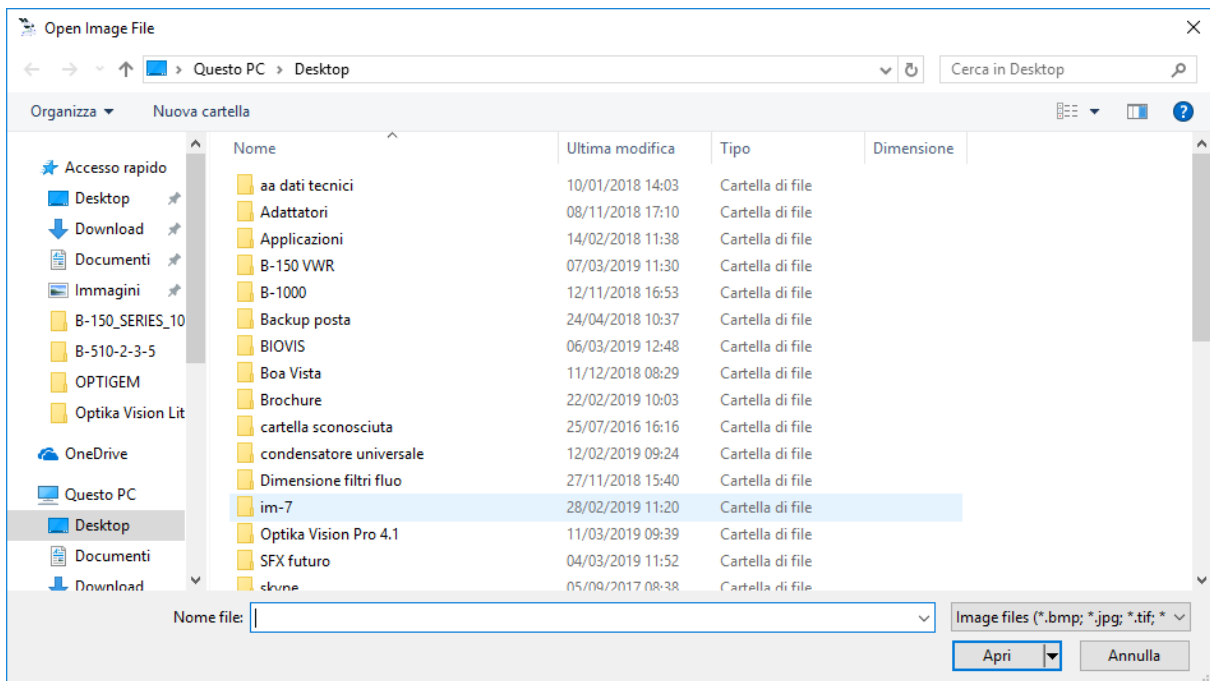


Select the desired file and click on **Open**.

## 5.6 Export image to file

This command allows you to save a captured image to any folder on your computer.

The following dialog box opens: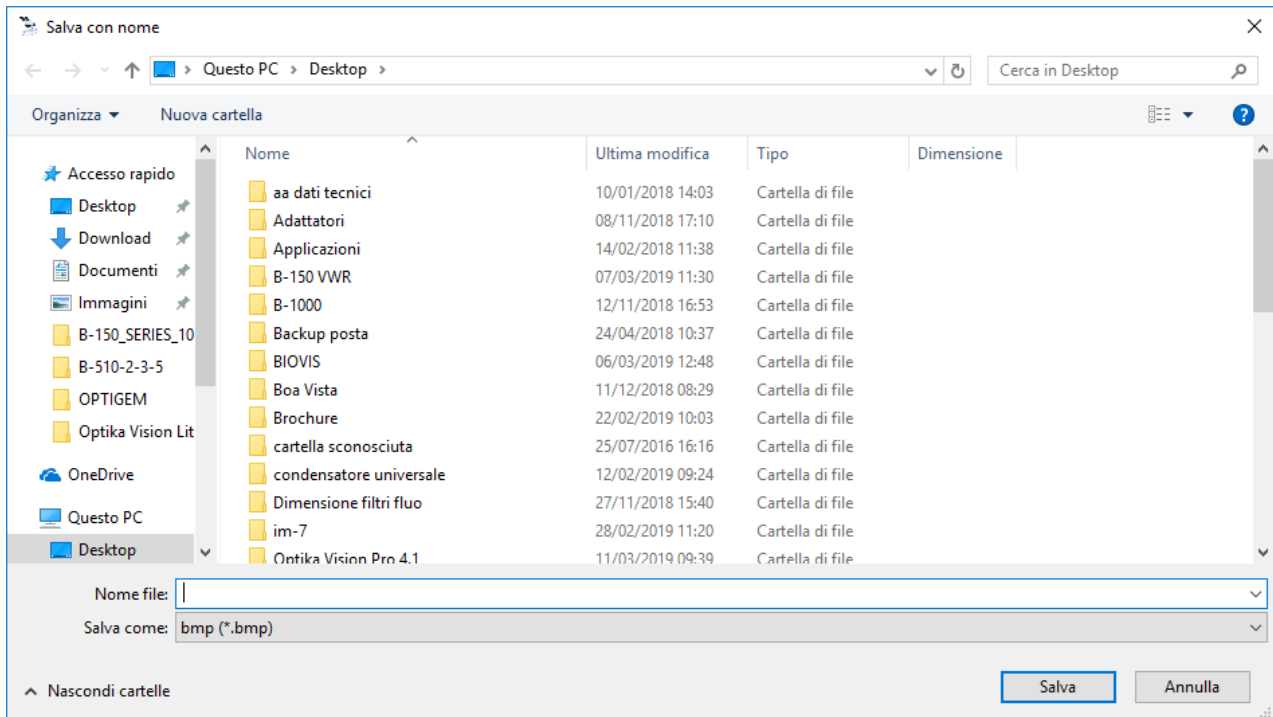


Possible image formats are: .bmp, .jpg, .tif

## 5.7 Flip/Rotate

Select this command to flip the image from the camera horizontally or vertically to make it consistent with the image viewed under the microscope.  
The function acts only on captured images.

## 5.8 Measurements

Select the drop-down menu for this command. Two options are available: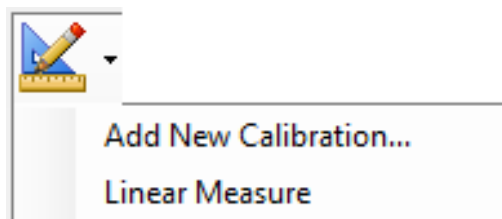

For a detailed explanation of the relevant functions see paragraphs 6.2.3 and 6.2.4.

## 5.9 Indicate objects

For a detailed explanation of the relevant functions see paragraph 6.2.2.

## 5.10 Measurement line color


You can select the color with which to display the lines or grids superimposed on the image. The default color is black. Clicking on the arrow next to the icon opens a drop-down menu with which you can select the color.

The selected color will be immediately applied to all objects present in the overlay on the image.


## 5.11 Grid

You can overlay the image with a grid. This is freely selectable by the user.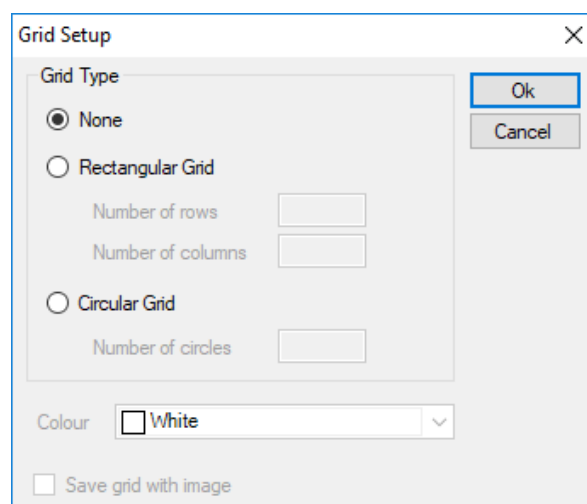


**None:** no overlay grid will be displayed.

**Rectangular grid:** indicate the number of rows and columns you want to draw on the image (e.g. 5x5). The image will be modified as follows: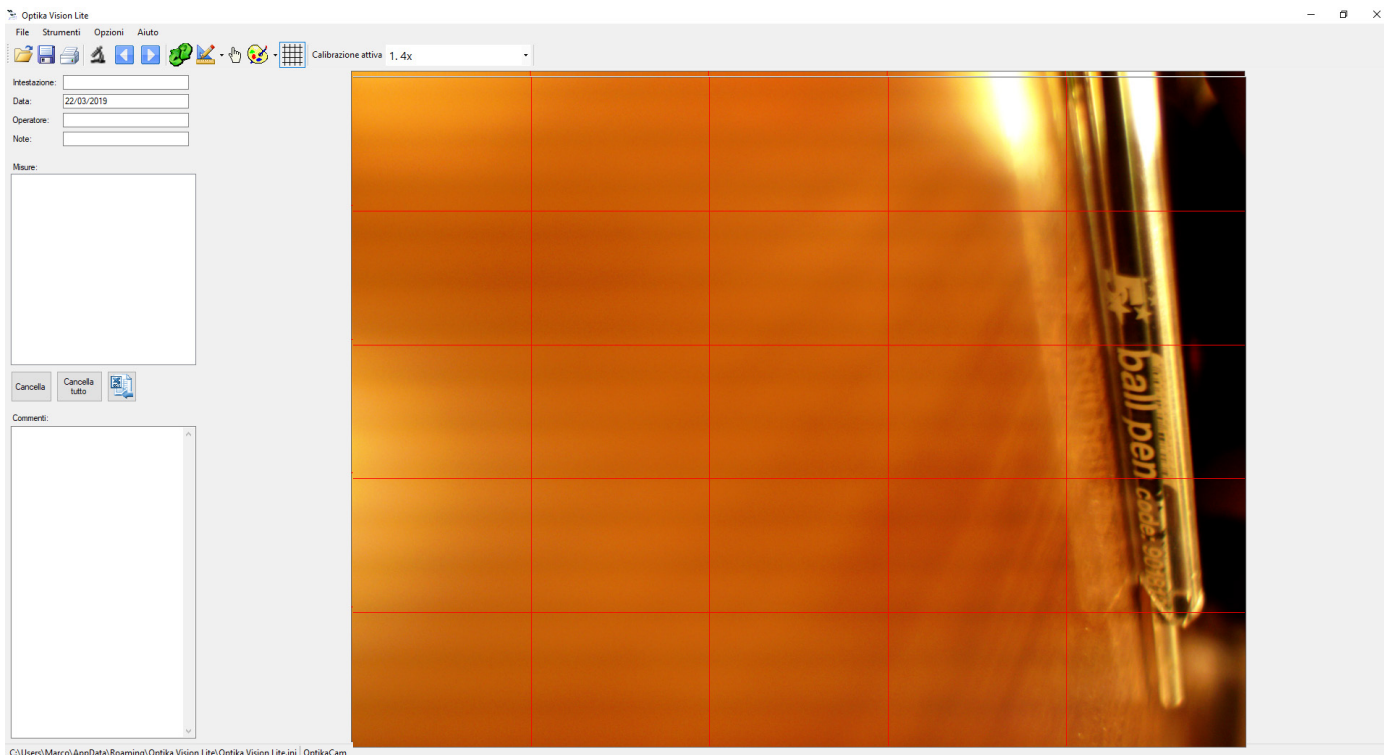

**Circular grid:** indicate the number of circles you want to draw on the image (e.g. 3). The image will be modified as follows.

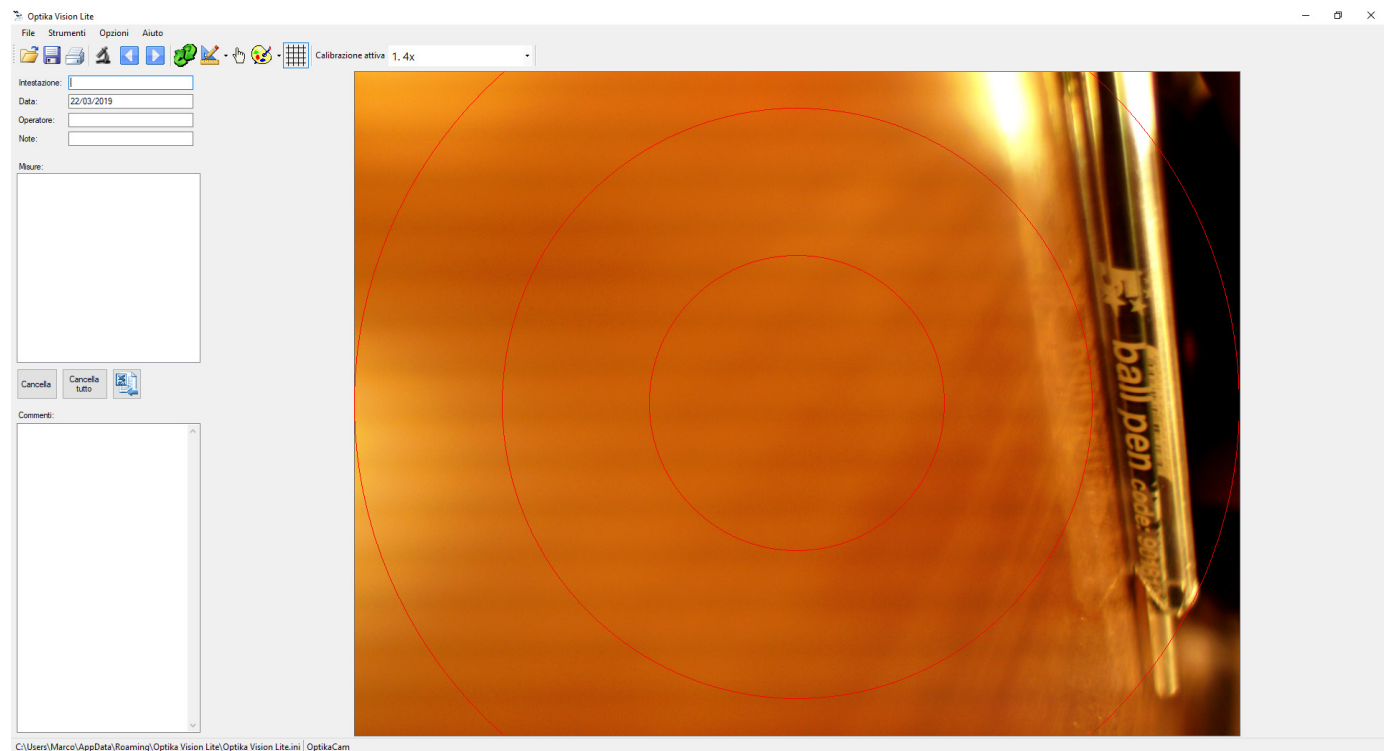
The grid can be drawn both on “live” images and on acquired images.

## 5.12 Select calibration

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In order to make correct measurements (after having calibrated the entire system) it is necessary to select the calibration corresponding to the lens in use.

Click on the arrow next to the field of the active calibration and the list of available calibrations opens.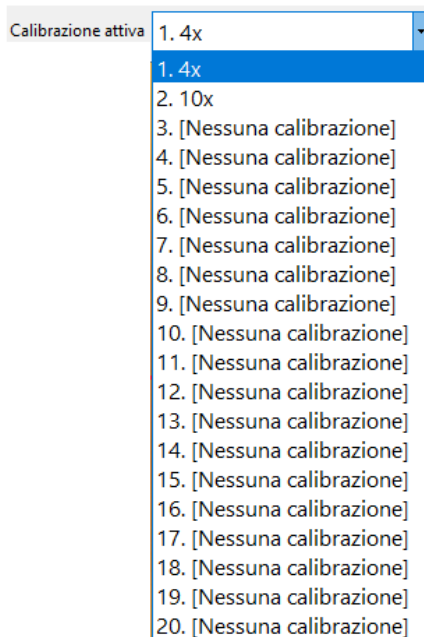

Select the calibration value corresponding to the lens in use.

Correct measurements can now be made.

---

## 6. Menu bar

### 6.1 File

The **File** section contains the following menus:

#### 6.1.1 Stop live video

Select the command **File>Stop live video** to select the camera and activate the live image from the microscope.

This command has the same effect as using the  icon on the Toolbar.

#### 6.1.2 Open Report

A Report is a document in “.asa” format in which a whole series of information has been saved (together with the image).

This command has the same effect as using the  icon on the Toolbar.

#### 6.1.3 Save Report as


All the information set in the “Reports” section (Header, Date, Operator, Notes, any Measurements and any Comments) are saved in a report in .asa format.

This command has the same effect as using the  icon on the Toolbar.


#### 6.1.4 Select source and snap image

This command allows you to manage and acquire images from TWAIN sources.

Clicking on the command **File>Select Source and Capture** this dialog box opens: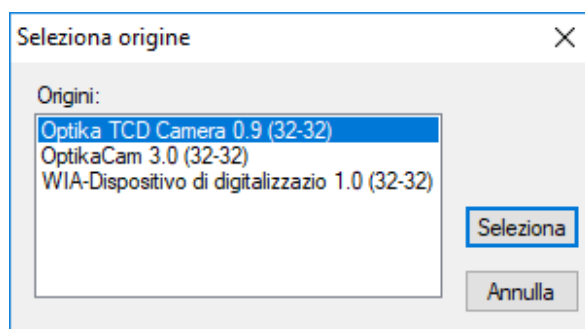


Select the desired camera, click on **Select**: the camera's parameter control window opens..

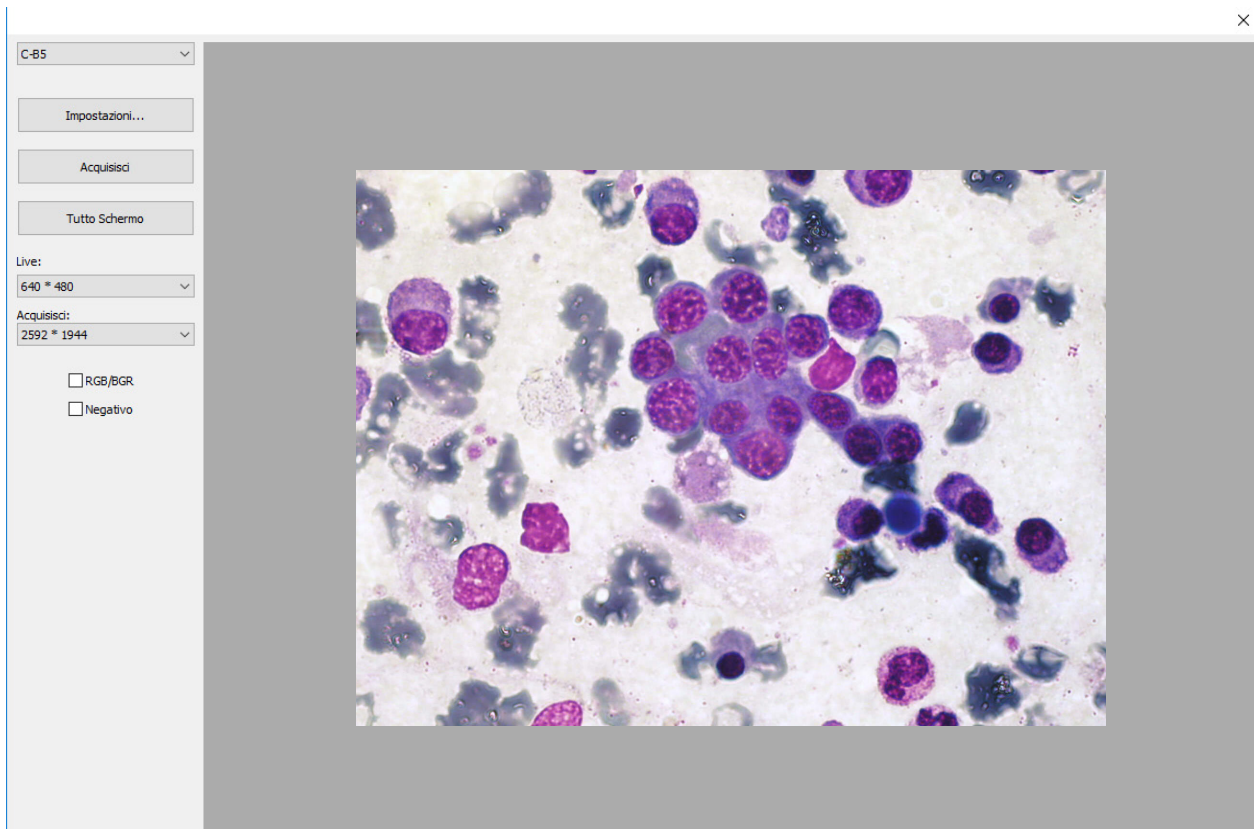Commands vary depending on the selected camera.

### 6.1.5 Import image from file

This command allows you to open a previously saved image.

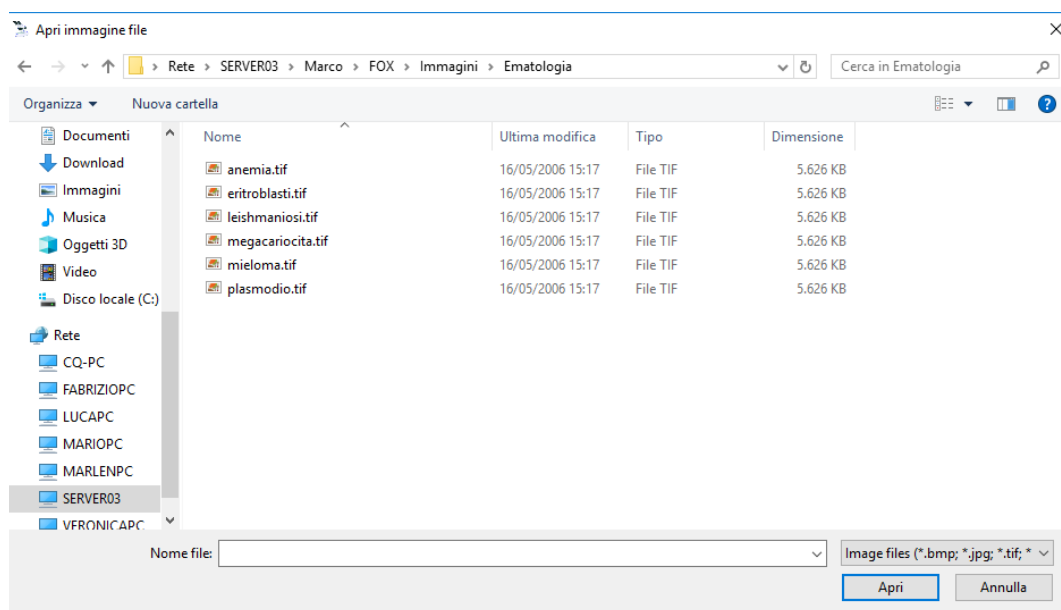
This command has the same effect as using the  icon on the Toolbar.

## 6.1.6 Export

This command allows you to save a captured image or a measurement taken.

Click on **Export>Image to File** to open the Save As dialog box. From here you proceed as a normal Windows save function.

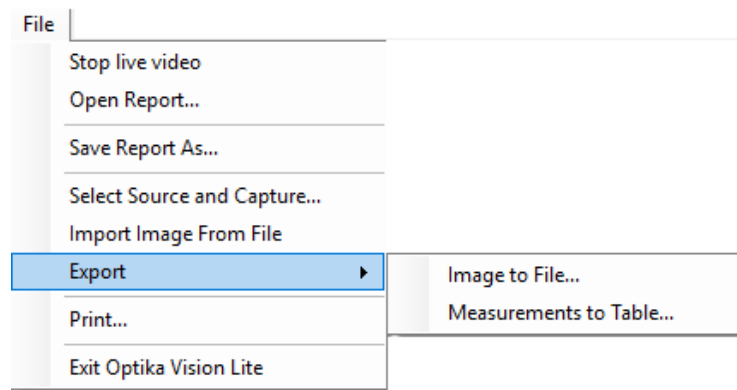Possible image formats are: .bmp, .jpg, .tif

This command has the same effect as using the  icon on the Toolbar.

You can also save measurements made on the image. Click on **Export> Measurements to Table** to save the measurements made on the image.

The file format is .csv.

## 6.1.7 Print

This command allows you to print a captured image.

Clicking on this command opens the **Print** dialog box.

This command has the same effect as using the  icon on the Toolbar.

## 6.1.8 Exit

This command allows you to exit the program.

## 6.2 Tools

The **Tools** section contains the following menus:

### 6.2.1 Flip/Rotate image

Select this command to flip the image from the camera horizontally or vertically to make it consistent with the image viewed under the microscope.

The function acts only on captured images.

This command has the same effect as using the  icon on the Toolbar.

### 6.2.2 Indicate objects

Once the image has been captured, approach the point to be indicated with the mouse and, while holding down the left mouse button, drag the mouse.


You create an arrow whose vertex is the point you want to indicate.

This command has the same effect as using the  icon on the Toolbar.


### 6.2.3 Linear measurement

Draw a line on the object you want to measure. The line you have just drawn will be marked with a progressive number and the measurement you have just made will be added to the left side of the GUI.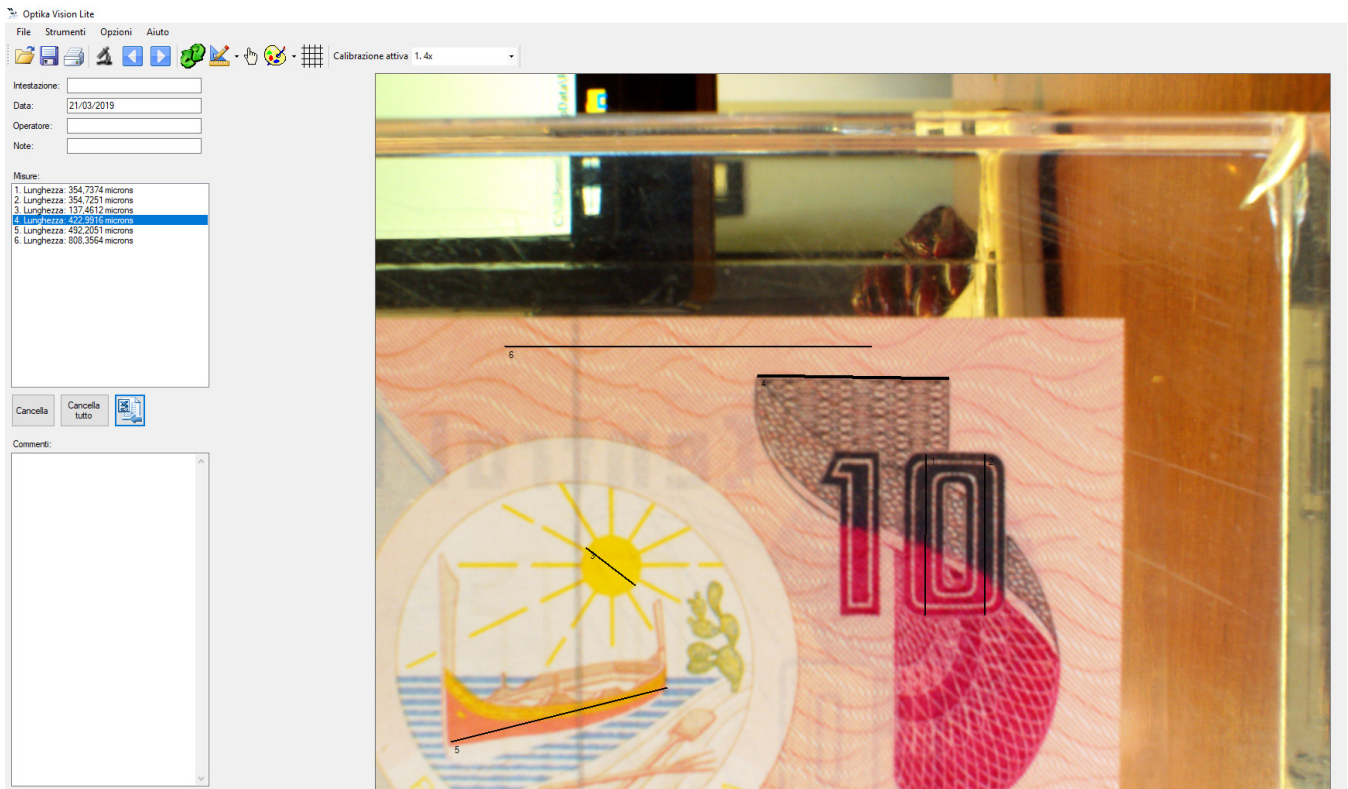

You can delete one or more unwanted measures. Select the measurement and click on **Delete**.

If you want to delete all measurements click on **Delete All**.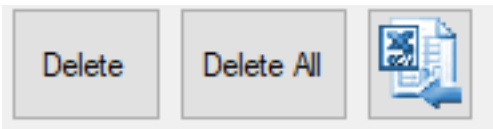


The measurements just made can be exported using the  button or clicking on **File>Export>Measurements to Table**.

### 6.2.4 Calibration


Before making a correct measurement, the system must be calibrated.

The procedure is as follows.


1. Place the object micrometer on the table, insert the lens with the lowest magnification in the optical path and focus on the linear scale.
2. Start the live image of the camera and capture the image of the micrometer scale.
3. Click on **Tools>Calibrate>Add new calibration** or click on the button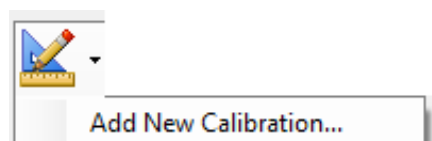


- The dialog box opens

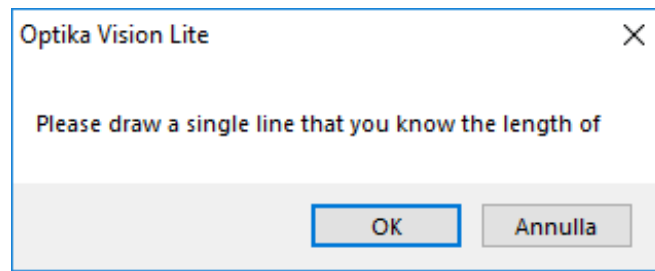
- Draw a line that includes a certain number of notches of the object micrometer and click on "OK".
- The next dialog box opens: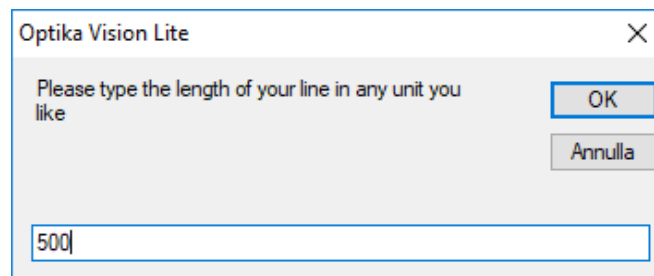


- Enter the length (without any unit of measurement) of the line just drawn: e.g. 500, then click on "OK".
- The dialog box that asks you to indicate the desired unit of measurement (e.g. "micron") opens and then click on "OK".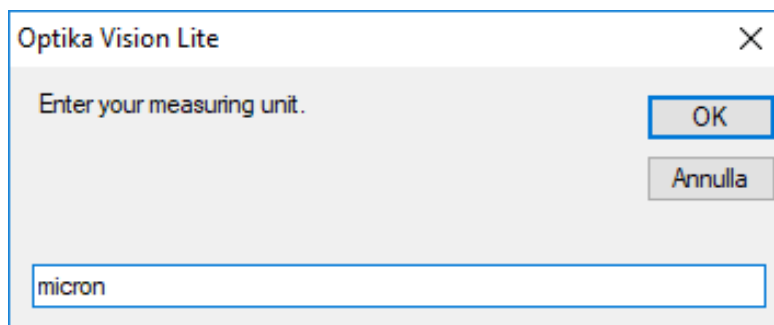


- The dialog box that asks you to name the calibration (e.g. 4x) opens and then click on "OK".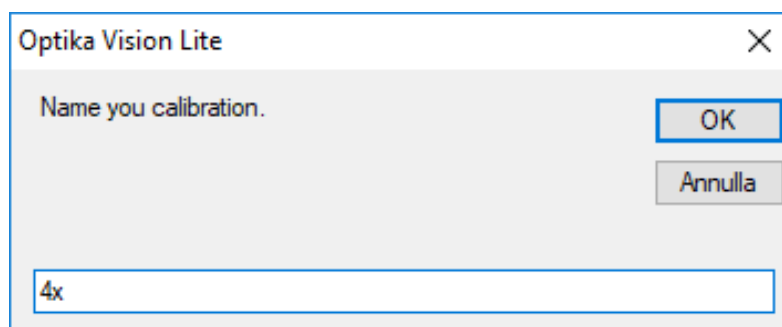

10. The dialog box “**Select position**” opens. This indicates where to save the newly created calibration point. Then click on “**OK**”.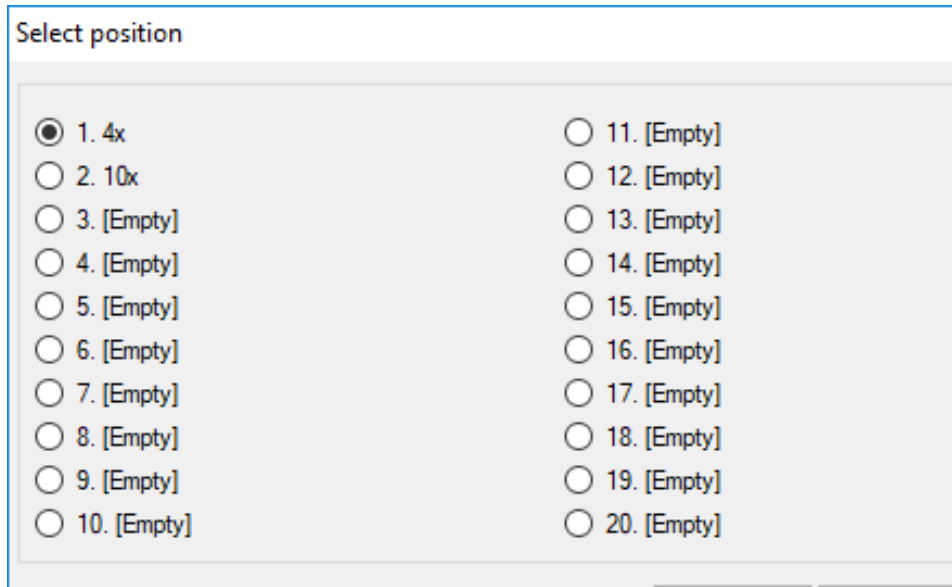


11. Following dialog appears: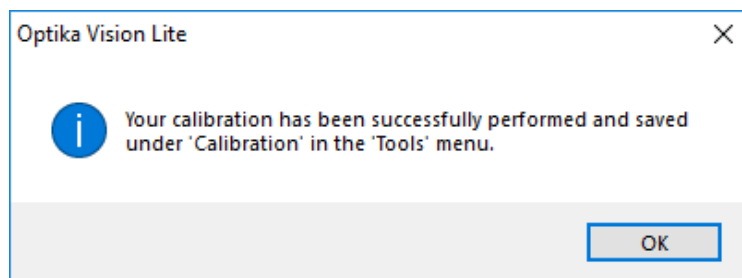

12. Perform the same process for all objectives in use on the microscope.

You can **delete** a calibration point if an error has been made.


Click on **Tools>Calibrate>Remove Calibration**. The dialog box opens: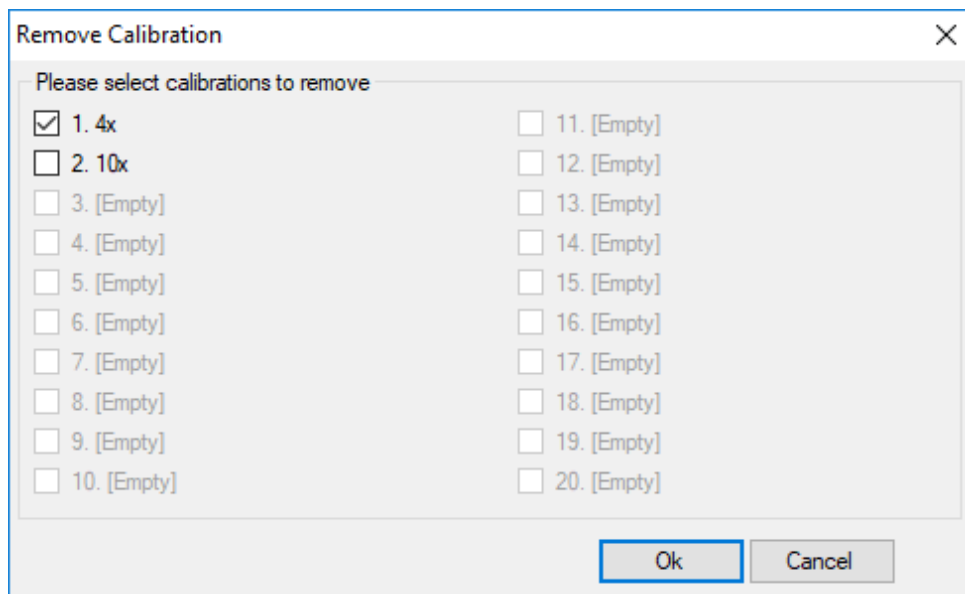

Select the calibration(s) to be deleted and then click on “OK”. The selected values will be deleted from the list.

You can **rename** a calibration point if an error has been made.

Click on **Tools>Calibrate>Rename Calibration**. The dialog box opens: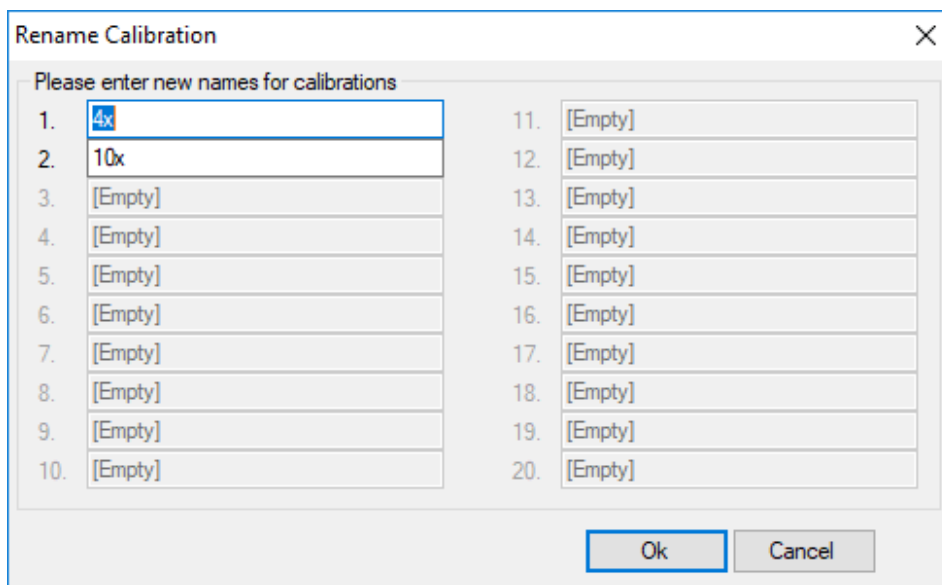

Select the name of the calibration point to be renamed and then press “OK”. The new name will replace the old one.


## 6.3 Options

The **Options** section contains the following menus:

### 6.3.1 Change Language

Select this command to change the language used in the software.

Click on **Options>Change Language**. The dialog box opens:

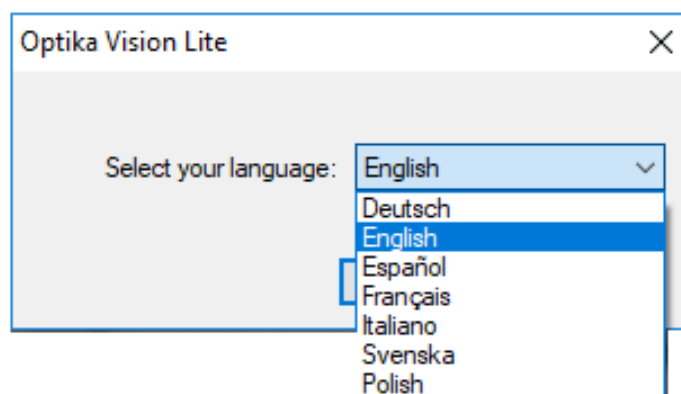
### 6.3.2 Change print logo

You can change the logo set for printing the report.

Click on **Options>Change print logo**. The dialog box opens:

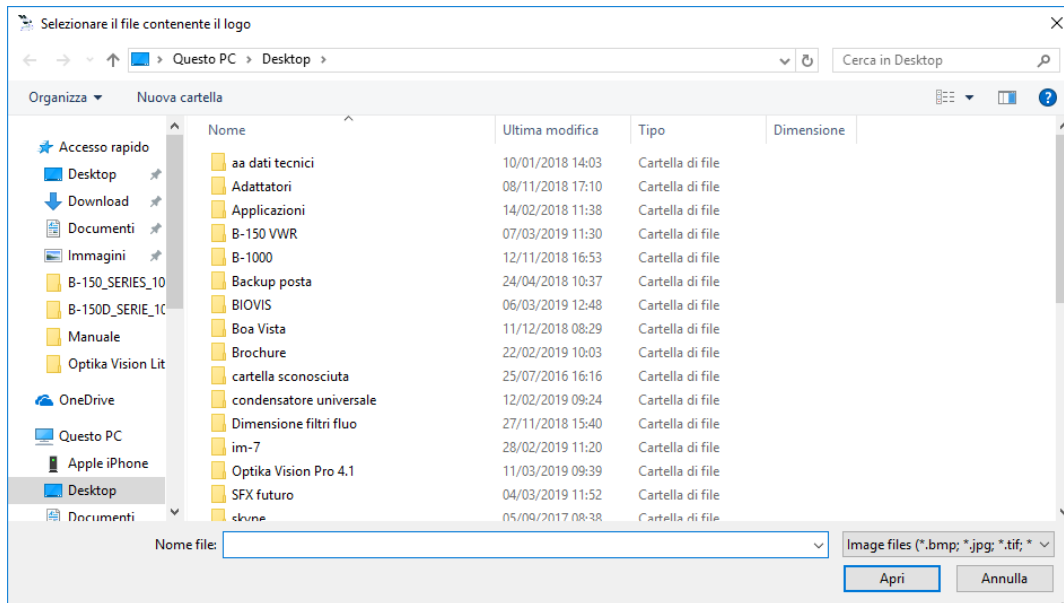Here you select the file that will replace the logo previously set for printing the report, then click on **OK**.

## 6.4 Help

The **Help** section contains the following menus:

### 6.4.1 Help Index

Select this command to open the online summary guide for using the software.

### 6.4.2 About

Select this command to display information about the SW version.

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**OPTIKA® S.r.l.**