ACCESSORIES Series

INSTRUCTION MANUAL

Model
ST-047

Table of Contents

1.	Warning	3
2.	Safety Information	3
3.	Package content	4
4.	Unpacking	5
5.	Intended use	5
6.	Symbols and conventions	5
7.	Instrument description	6
8.	Assembling	7
9.	Use of the stage	8

1. Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.


2. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

3. Package content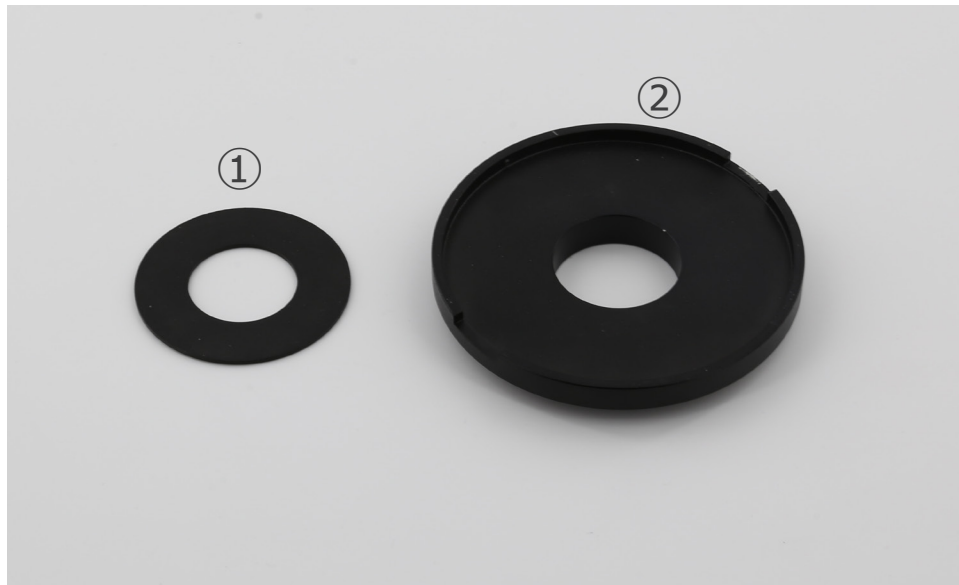

① Fixing plate

② Gliding stage

4. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

5. Intended use

Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

6. Symbols and conventions


The following chart is an illustrated glossary of the symbols that are used in this manual.


CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

7. Instrument description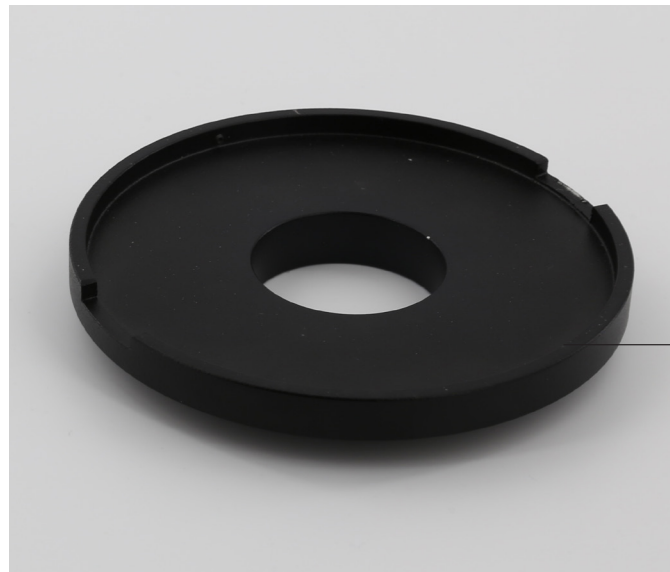


GLIDING STAGE

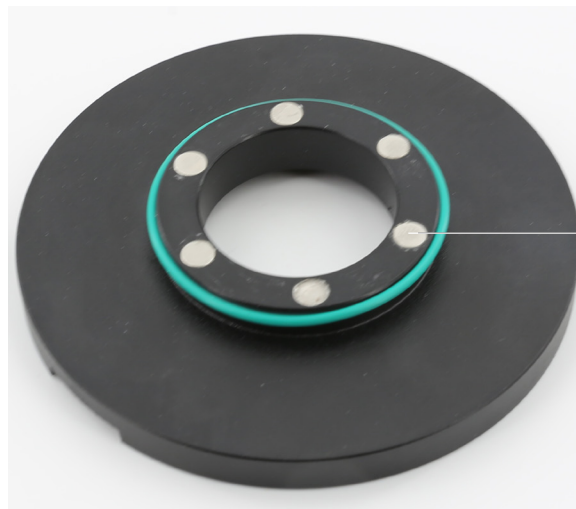
FIXING MAGNETS

8. Assembling

1. Using a Phillips screwdriver (not supplied), remove the two sample clips ①. (Fig. 1)
2. Remove the stage plate ② (glass or plastic) from the base of the microscope.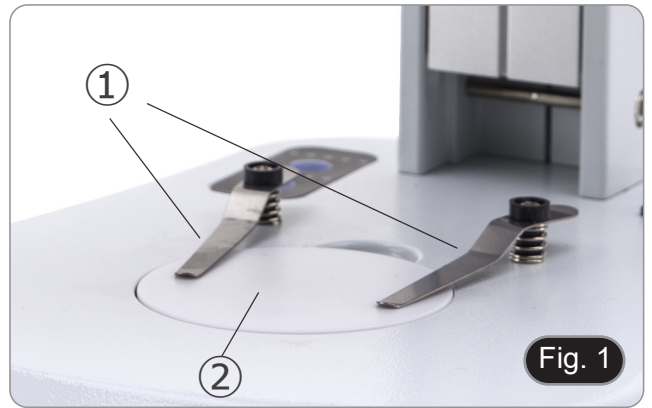

3. Remove the tape protection film from the fixing plate.
4. Insert the fixing plate into the empty hole ③ of the microscope base, making the adhesive part adhere well. (Fig. 2)

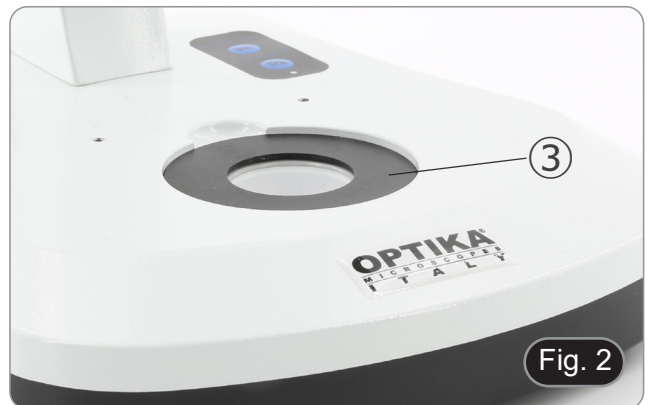5. Attach the gliding stage to the fixing plate using the magnets. (Fig. 3)

9. Use of the stage

1. Place a specimen on the stage and, using both hands, slide the stage to move the specimen to the desired position while observing at the eyepieces. (Fig. 4)

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OPTIKA® S.r.l.