**SZR Series**

# INSTRUCTION MANUAL

Model
SZR-180 + SZR-FLD

Ver. 1.1 2023

---

## Table of Contents

1.	Warning	3
2.	Safety Information	3
3.	Package content	4
4.	Unpacking	5
5.	Intended use	5
6.	Symbols and conventions	5
7.	Instrument description	6
8.	Assembling	9
9.	Using the microscope in Transmitted Light	12
9.1	Adjusting interpupillary distance	12
9.2	Diopter compensation	12
9.3	Magnification	13
9.4	Use of aperture diaphragm	13
9.5	Use of trinocular port	13
9.6	Use of transmitted light illumination	14
9.7	Use of conversion lens	15
10.	Using the microscope in Fluorescence	16
10.1	Use in single fluorescence	16
10.2	Use in multiple fluorescence	18
10.3	Use of the protection shield	19
11.	Simultaneous observation in Transmitted Light + Fluorescence	20
12.	Microphotography	21
12.1	Installing the camera	21
12.2	Parfocality adjustment	21
13.	Maintenance	22
14.	Troubleshooting	23
	Equipment disposal	24

---

## 1. Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use that does not comply with this manual.

## 2. Safety Information

### Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

### 3. Package content

- ① Base
- ② Microscope body
- ③ Objective
- ④ Observation head
- ⑤ Eyepieces
- ⑥ Power supply
- ⑦ Dust cover
- ⑧ Fluorescence epi-illuminator
- ⑨ Fluorescence control box
- ⑩ Fluorescence light source
- ⑪ Light guide
- ⑫ UV protection shield
- ⑬ Power supply + power cord

---

## 4. Unpacking

The microscope is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged. Using both hands (one around the arm and one around the base), lift the microscope from the container and put it on a stable desk.

Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

## 5. Intended use

### Standard models

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

### IVD Models

Also for diagnostic use, aimed at obtaining information on the physiological or pathological situation of the subject.

## 6. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

### CAUTION


This symbol indicates a potential risk and alerts you to proceed with caution.


### ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

## 7. Instrument description
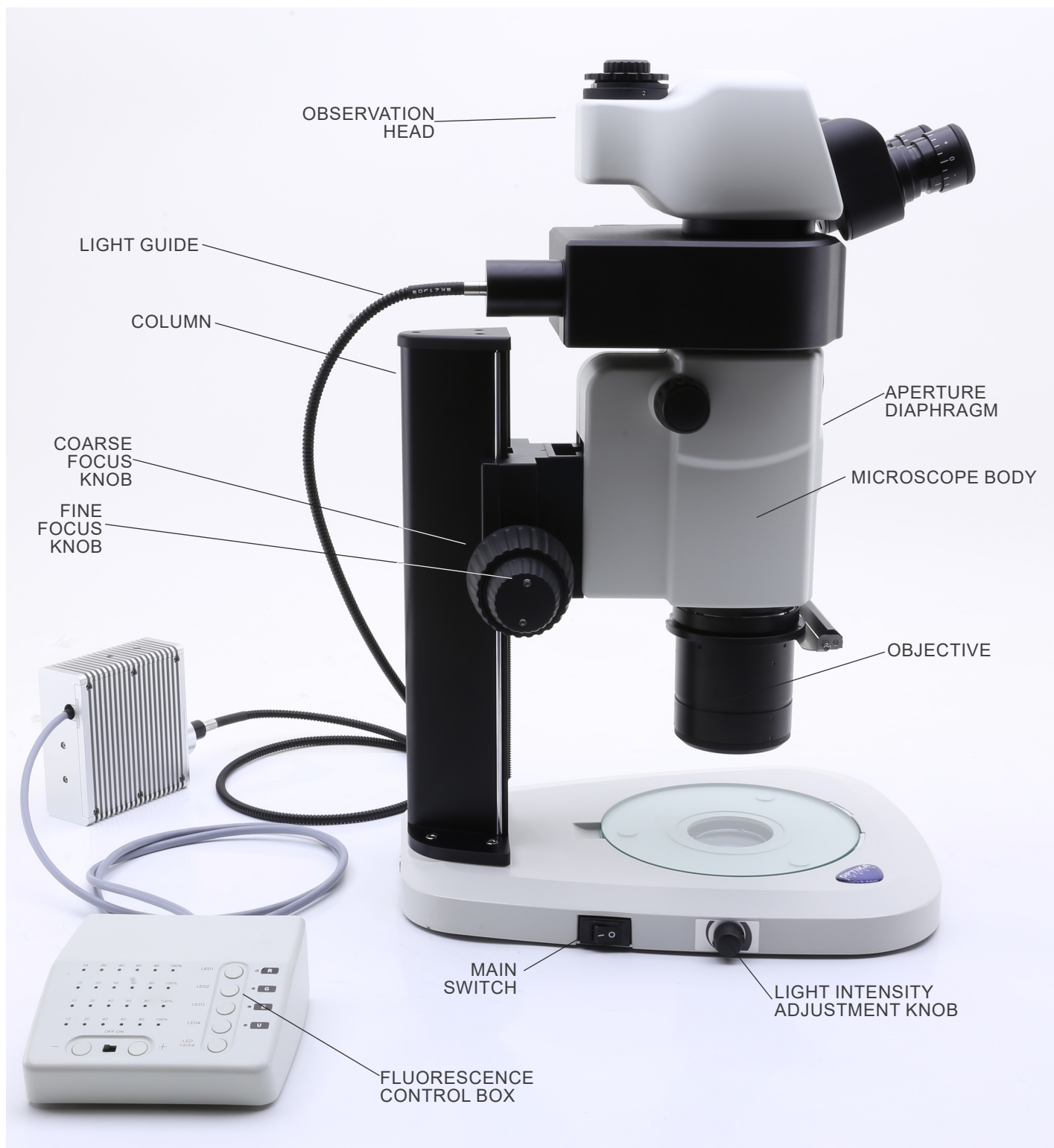


Opposite side

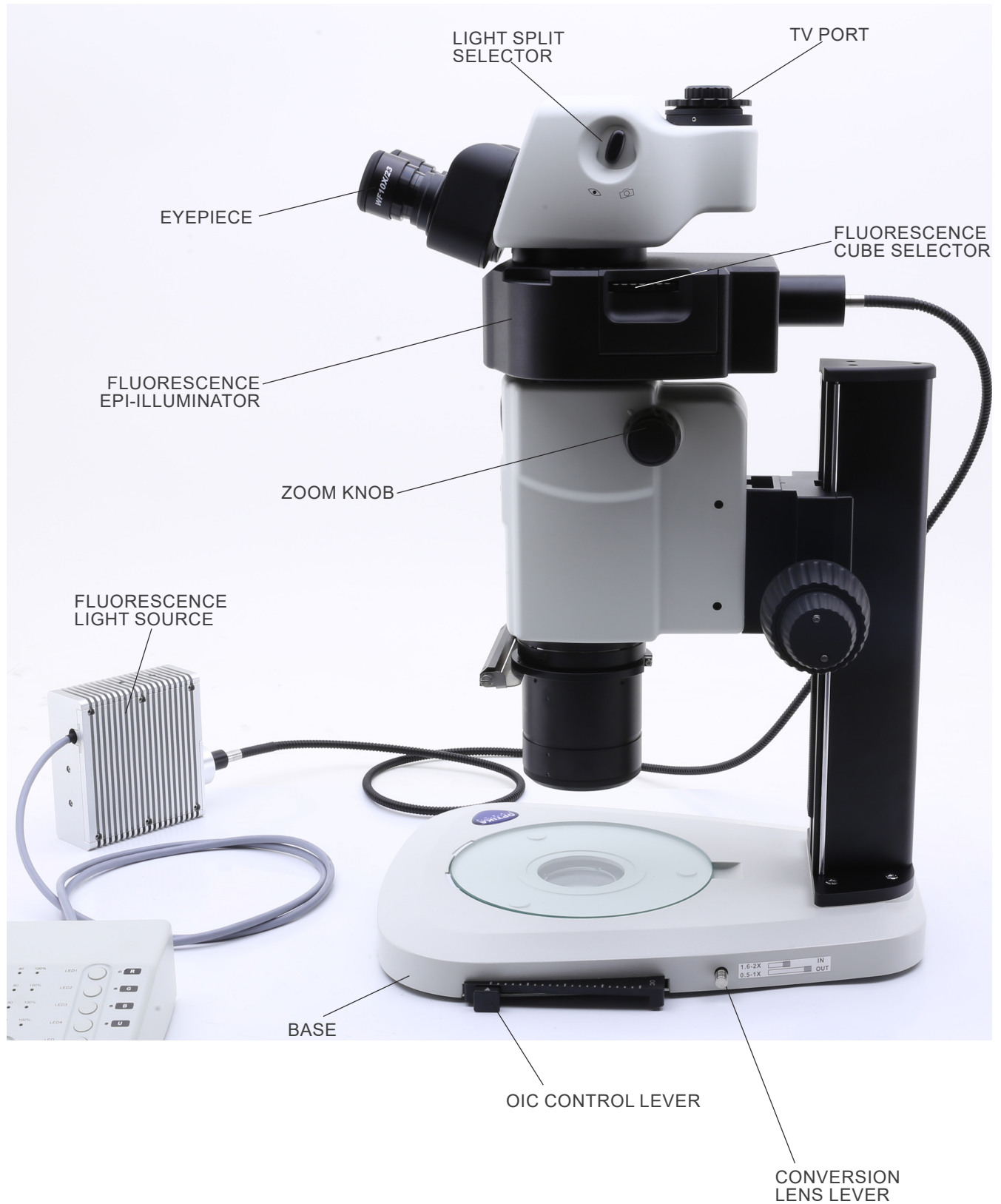
Fluorescence control box

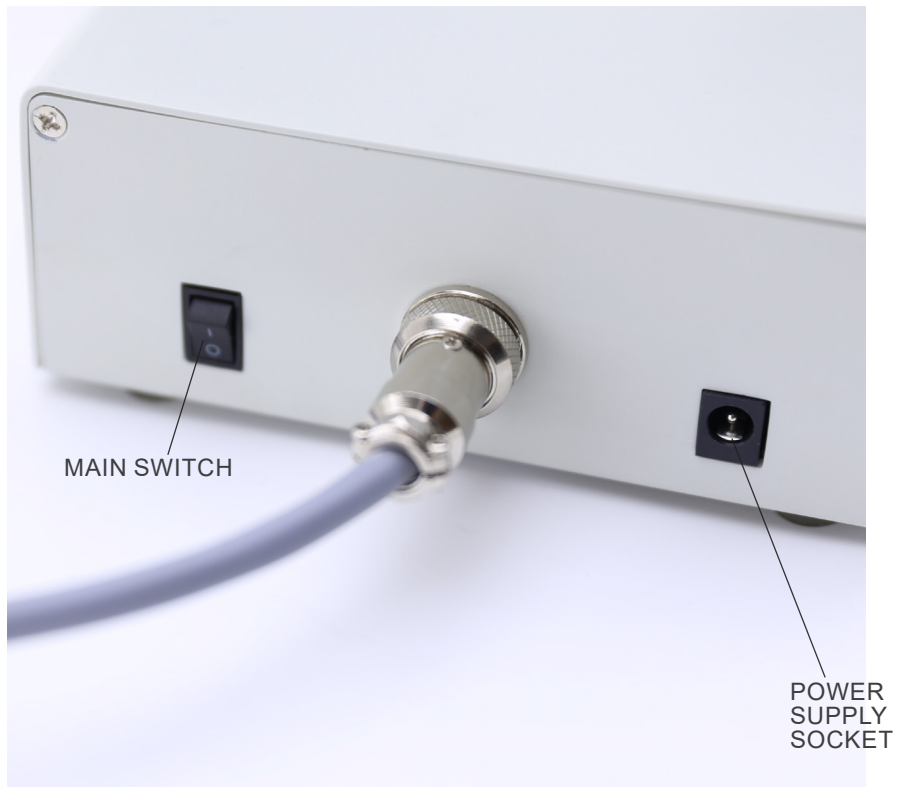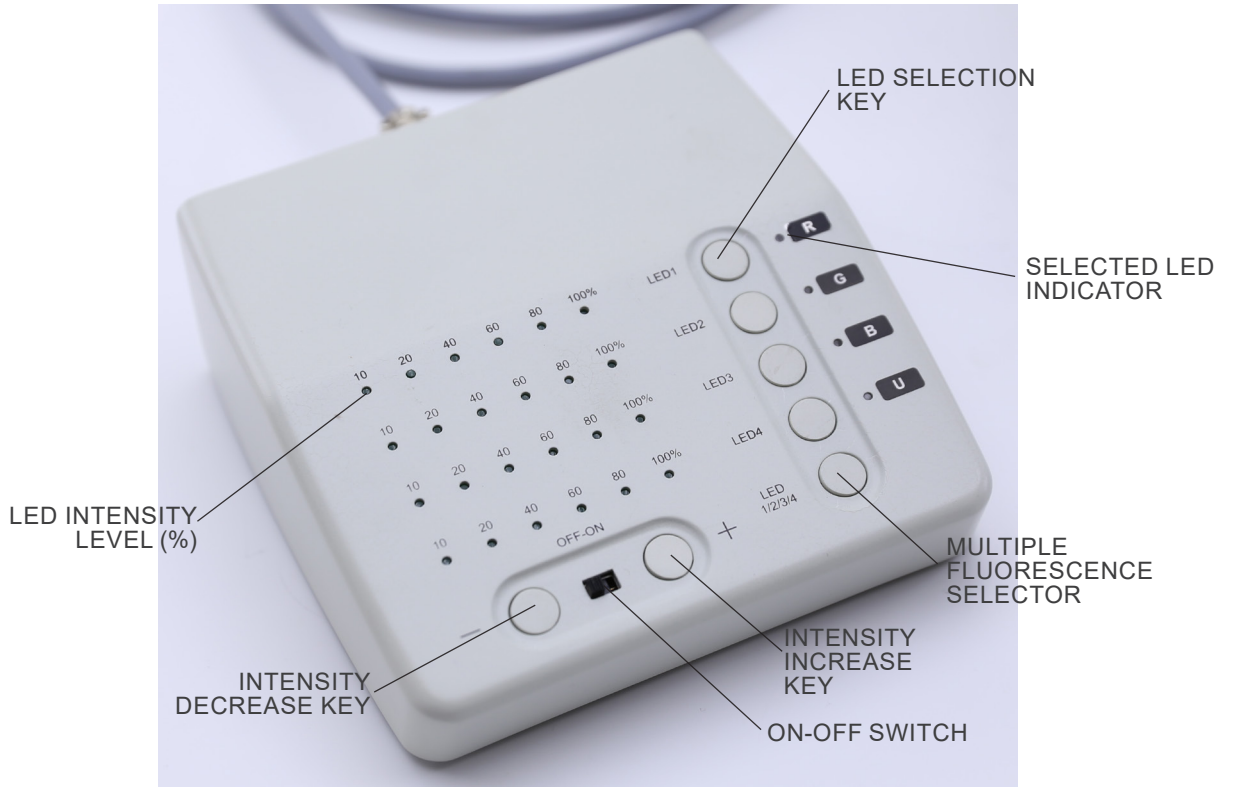
## 8. Assembling

1. Put the microscope base ① flat on the workbench. (Fig. 1)
2. Align the holes of the connecting plate ② of the stand column with the holes of the microscope base.
3. Use the provided fixing screws to connect the base and the column firmly.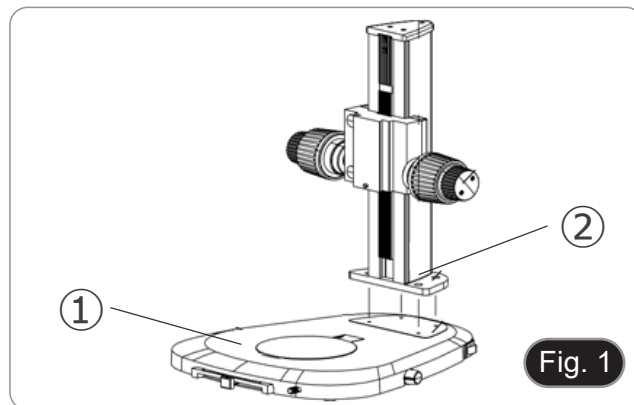

4. Insert the zoom body dovetail ③ into the dovetail of the focus system ④. (Fig. 2).
5. Lock the fixing screws ⑤.

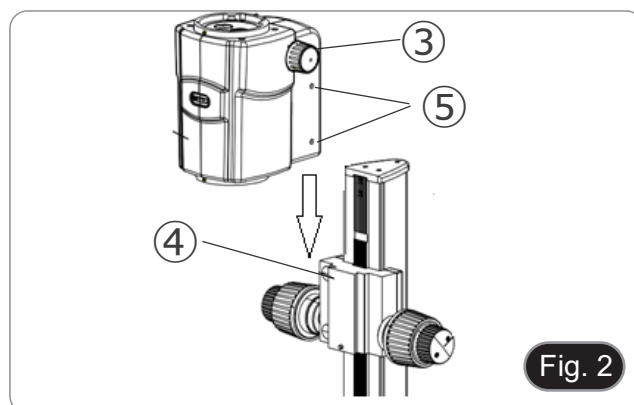
6. Install the fluorescence epi-illuminator by aligning the round dovetail of the illuminator with the hole of the microscope body. (Fig. 3)
7. Lock the fixing screw ⑥.


8. Install the head by aligning the round dovetail of the head with the hole of the fluorescence epi-illuminator. (Fig. 4)
9. Lock the fixing screw ⑦.


10. Insert eyepieces into empty eyepieces sleeves. (Fig. 5)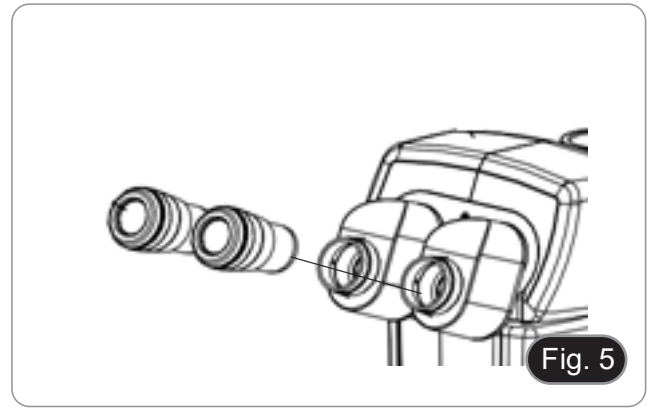

11. Screw the objective in the bottom part of the microscope body. (Fig. 6)  
12. Lock the objective by screwing the fixing screw ⑧.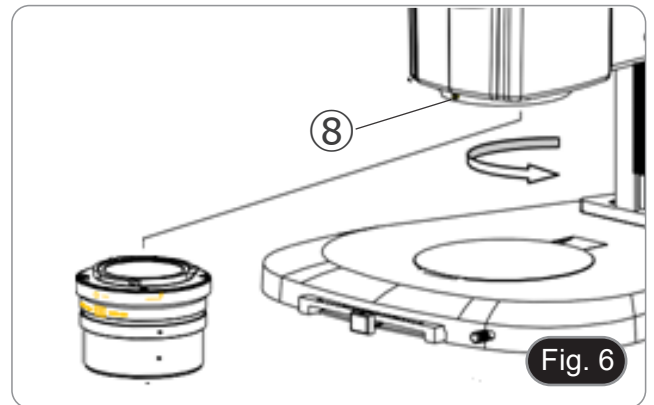

13. Remove the rubber protection caps ⑨ from the light guide. (Fig. 7)

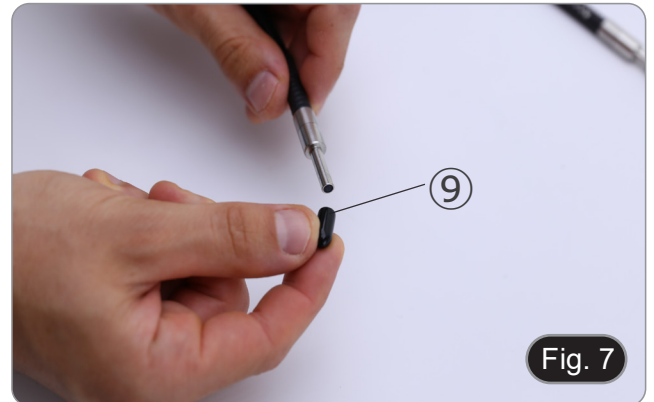14. Insert one end of the light guide into the epi-illuminator. (Fig. 8)  
15. Tighten the fixing screw ⑩.

16. Insert the other one end of the light guide into the light source.  
(Fig. 9)
17. Tighten the fixing screw ⑪.

18. Plug the power cable ⑫ from the light source in the socket ⑬ in the fluorescence control box. (Fig. 10)

19. Plug the jack of the fluorescence power supply ⑭ in the socket of the fluorescence control box. (Fig. 11)
20. Connect the power cord.

21. Plug the jack of the microscope power supply in the socket of the microscope base.
- **Before connecting the power supplies make sure that all switches are in OFF position.**

## 9. Using the microscope in Transmitted Light


### 9.1 Adjusting interpupillary distance

1. Hold the right and left eyepiece tube with both hands and adjust the interpupillary distance by moving the two parts until one circle of light can be seen. (Fig. 12)
- If two circles appear, the interpupillary distance is too big.
- If two overlapped circles appear, the interpupillary distance is too small.

### 9.2 Diopter compensation

- **This compensation makes possible for people wearing glasses to adjust the microscope to their eyes and use the microscope without glasses.**
1. Put the zoom down to the lowest magnification ① and focus the specimen with the focusing knobs ②. (Fig. 13)
  2. Put the zoom to the maximum magnification and repeat the focusing.
  3. Return to the lowest magnification: the specimen will be out of focus.
  4. Adjust the diopter compensation ring of the right eyepiece ③ (Fig. 14) until the image of the right eyepiece is clear and sharp. Repeat the procedure for the left eyepiece ④.
  5. Check the focus of the image for the whole zoom range. It should be perfectly parfocal (focus is maintained during the change of magnification).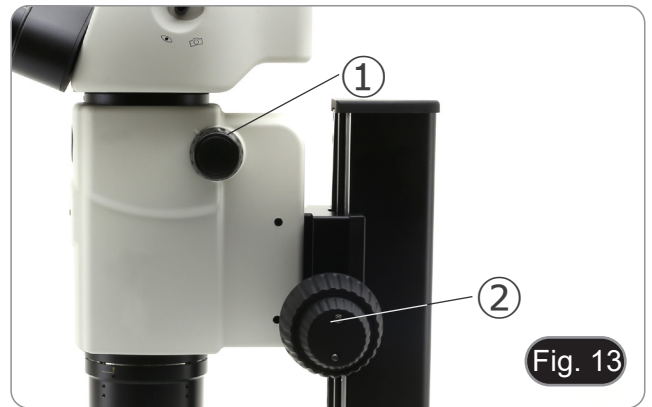

### 9.3 Magnification

1. Select the desired magnification by adjusting the zoom knob ①. (Fig. 15)
- Change the eyepieces if a different magnification range is needed.
  - The microscope body is equipped with a “click-stop” function that allows to obtain a precise setting of the desired magnification.
  - Total magnification used can be calculated as:  
Eyepiece magnification \* Zoom magnification \* Objective lens magnification.

Eyepiece		10x	
Field number (mm)		24	
Objective	Total Mag	F.O.V (mm)	
1X	7.5X - 135X	30.67 - 1.7	

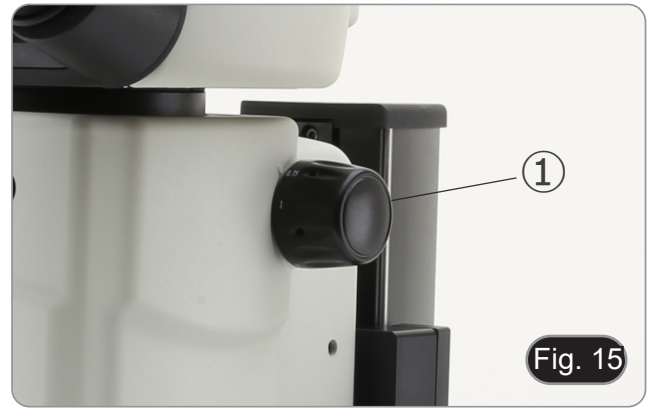
Fig. 15

### 9.4 Use of aperture diaphragm

By varying the diaphragm aperture, the user can adjust the image to achieve better contrast and greater depth of field. (Fig. 16)

1. Move the aperture diaphragm lever ② to the right or left. Rotation towards “○” opens the diaphragm while rotation towards “⊗” closes it.
  2. Adjust the diaphragm during observation to optimize the effect on contrast and depth of field.
- **Do not close the diaphragm too tightly so as not to have a poor image quality and not to have an excessively dark image.**

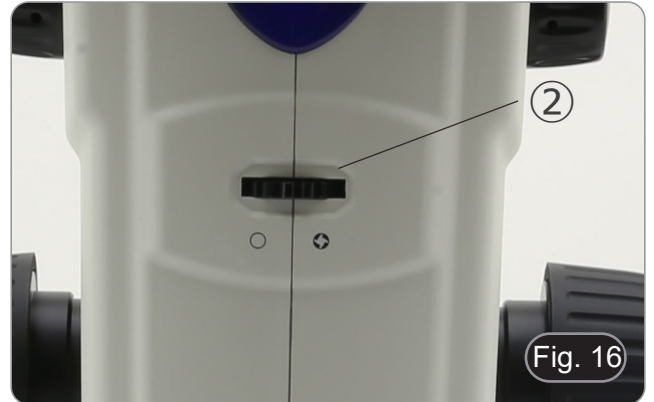
Fig. 16

### 9.5 Use of trinocular port

Observation head is equipped with a light split selector ③. (Fig. 17)

1. Moving the selector to the “eye” position ④ allows the binocular observation.
  2. Moving the selector to the “camera” position ⑤ allows the trinocular observation.
- **When the selector is in the “camera” position the right eyepiece will be completely dark. This is not a defect.**

Fig. 17


## 9.6 Use of transmitted light illumination

- When working in transmitted light make sure that the fluorescence filter holder turret is in an empty position. However, if the turret is full, make sure that the UV (DAPI) position is selected to prevent colors from the fluorescence filters from damaging the image quality in transmitted light.


1. Set the main switch ① to "I" to turn on the base. (Fig. 18)
2. Turn the light intensity adjustment knob ② to adjust the brightness to the sample under observation.

3. Move the lever ③ located on the right side of the base. (Fig. 19)
- This lever controls a black foil which slides horizontally along the light exit.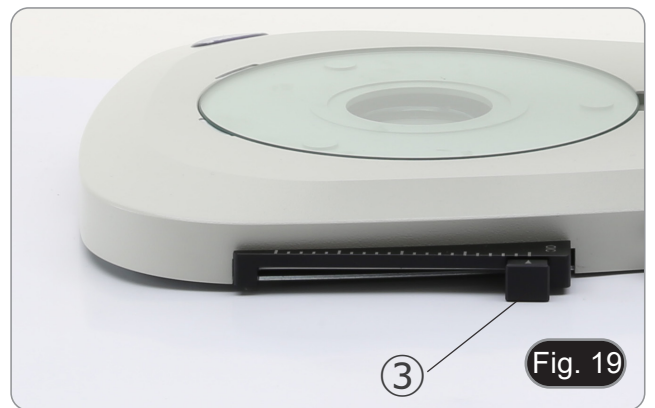

- Using the OIC (Oblique Illumination Contrast) system allows you to modulate the light to obtain three different types of observation.
4. When the foil is outside the optical path (Fig. 20), brightfield (BF) observation is obtained.

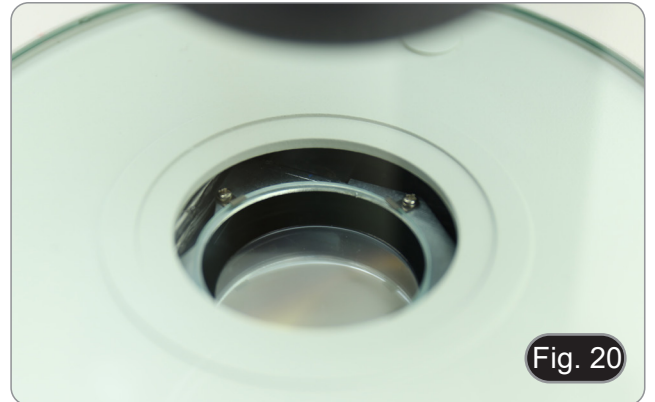5. When the foil is completely inside the optical path (Fig. 21), dark field (DF) observation is obtained.


6. Moving the foil to an intermediate position (Fig. 22) results in the modulation of light called Oblique Illumination (OIC).

### 9.7 Use of conversion lens

- The conversion lens is used to optimize the light intensity according to the objective in use. (Fig. 23)
- 1. Fully insert the lever ① when using 1.6x or 2x objectives.
- 2. Pull lever ① all the way out when using 0.5x or 1x objectives.
- **There are currently no additional objectives available other than the standard 1X objective.**

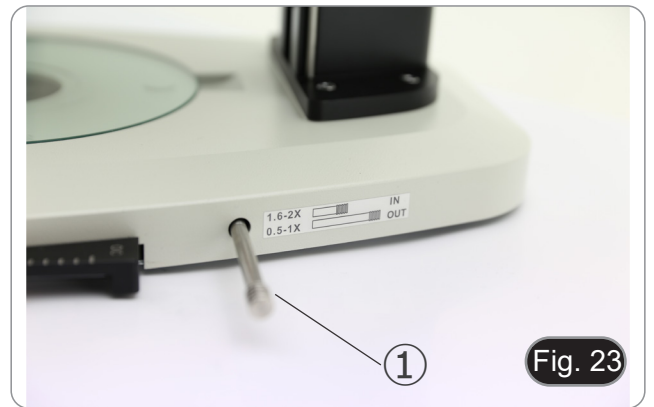
## 10. Using the microscope in Fluorescence

### 10.1 Use in single fluorescence


1. Turn on the main switch ① on the back side of the fluorescence control box. (Fig. 24)

2. Move the ON-OFF switch ② on the front panel of the fluorescence control box in the "ON" position. (Fig. 25)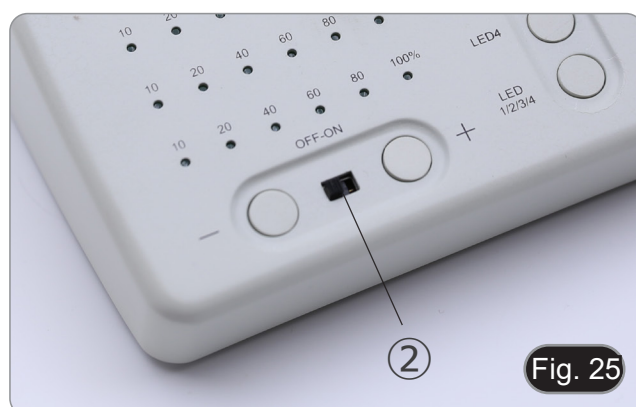

3. Press one of the LED selection keys ③ to activate the desired light. (Fig. 26)
4. Once activated, the indicator LED ④ lights up.
5. The LED light within the fluorescence light source, related to the selected LED (U, B, G or R) lights up.

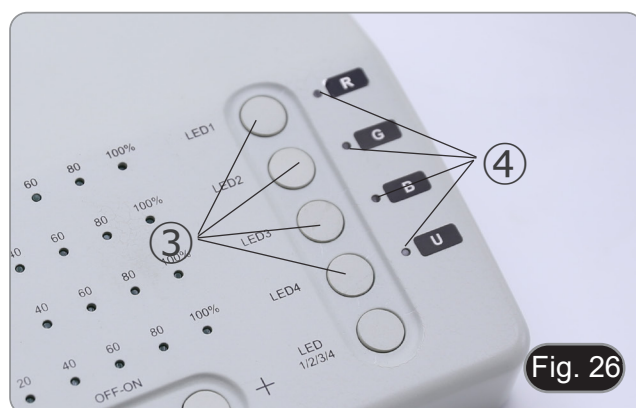
6. The LED light intensity level ⑤ of the selected LED lights up. (Fig. 27)

7. Rotate the fluorescence filter turret ⑥ to insert into the light path the fluorescence filter corresponding to the activated LED. (Fig. 28)

8. Focus the specimen using the focus knobs on the microscope.
9. While observing in the eyepieces increase or decrease the fluorescence light intensity using the "+" or "-" keys ⑦. (Fig. 29).
10. The numbers "10", "20", "40", "60", "80" or "100" ⑧ on will turn on or off according to the selected light intensity. (Fig. 30)
11. Begin the fluorescence observation.

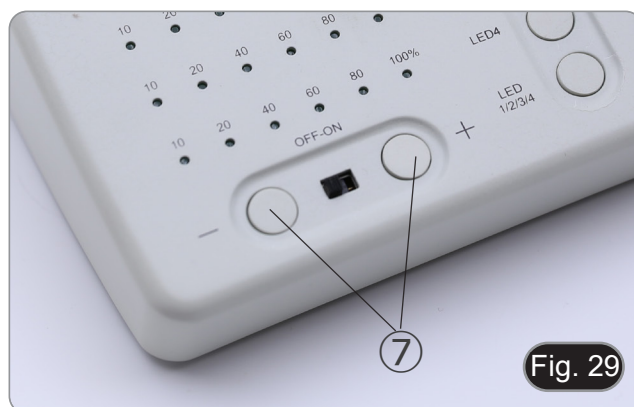## 10.2 Use in multiple fluorescence

- It is possible to work in multiple fluorescence, activating two or more LEDs simultaneously. Of course, fluorescence filter selection will have to be done manually by rotating the filter turret according to the currently active LEDs.
- Press the selection keys ③ of the LEDs you want to activate. (Fig. 26)
  - All selected LED sources within the fluorescence light source light up.
  - Each selected indicator LED ⑨ lights up in green. (Fig. 31)
  - Only one of the illuminated indicator LEDs will be lit red (Fig. 31). This indicates which LED source can be adjusted in intensity by means of the “+” e “-” keys ⑦. (Fig. 29)
  - Each selected LED light intensity level ⑩ of the selected LED lights up. (Fig. 31)
  - Press the “LED 1/2/3/4” key ⑪. (Fig. 32)
  - A single press of the key ⑪ cyclically selects one of the indicator LEDs ⑨, by turning the relevant LED red. (Fig. 31)
  - Rotate the fluorescence filter turret ⑥ to insert into the light path the fluorescence filter corresponding to the activated LED. (Fig. 28)
  - While observing in the eyepieces increase or decrease the fluorescence light intensity using the “+” or “-” keys ⑦. (Fig. 29).
  - The numbers “10”, “20”, “40”, “60”, “80” or “100” ⑩ will turn on or off according to the selected light intensity. (Fig. 31)
  - Begin the fluorescence observation.
  - To change fluorescence: rotate the fluorescence filter turret ⑥.
  - If it is necessary to vary the light intensity of the LED source according to the fluorescence filter inserted, press the “LED 1/2/3/4” key ⑪, (Fig. 32), in order to use the “+” and “-” keys for that specific LED.

Fig. 31


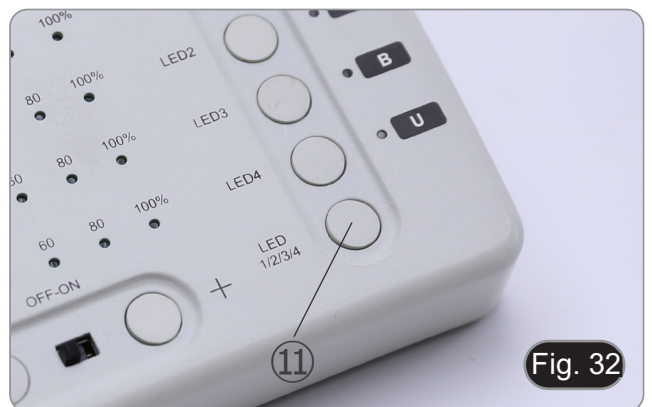
Fig. 32

FILTER CUBE	EXCITATION (nm)	DICHROIC (nm)	EMISSION (nm)	LED (nm)
U	330-385	410	420LP	365
B	460-490	500	510 LP	485
G	510-550	570	575LP	535
R	620-650	660	670-750	660
DAPI	360-390	415	435-485	365
FITC	460-495	505	510-550	485
TRITC	528-553	565	578-633	535

### 10.3 Use of the protection shield

- In order to prevent possible damages from UV radiation, use the orange protection screen.

1. Use the provided screwdriver to loosen the fixing screw ①. (Fig. 33)

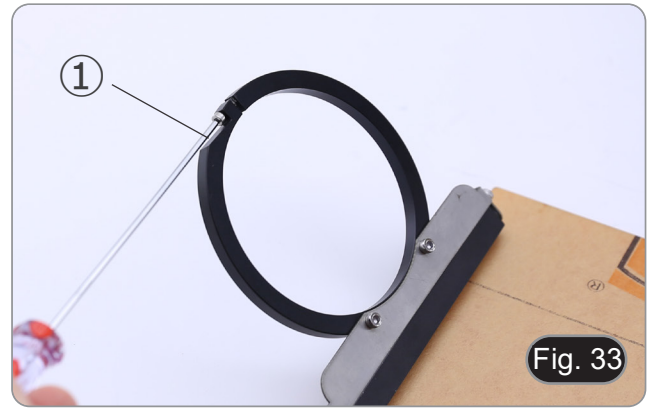2. Place the round ring of the protective screen on the microscope objective. (Fig. 34)

3. Tighten the fixing screw. (Fig. 35)

---

## 11. Simultaneous observation in Transmitted Light + Fluorescence

- **This microscope allows observation in Transmitted Light in combination with Fluorescence. Samples with rapid decay must first be observed in Fluorescence and then in Transmitted Light. The combined observation allows you to easily identify some areas of the sample that emit fluorescence.**
1. Turn on the microscope main switch.
  2. Move the fluorescence filter selector to an empty position or, if the filter holder is full, to the position containing the UV filter.
  3. Turn the transmitted light intensity adjustment knob to adjust the brightness to the sample under observation.
  4. Move the lever located on the right side of the base, to select Brightfield, OIC or Darkfield.
  5. Focus the sample.
  6. Move the fluorescence filter selector to the desired position.
  7. Adjust the light intensity of the reflected light.
  8. To obtain the proper observation of the sample, adjust the light intensity of the transmitted light to modulate the intensity of the fluorescence with the one of the transmitted light (Brightfield, OIC or Darkfield).

## 12. Microphotography

### 12.1 Installing the camera

1. Screw the camera on the provided C-mount adapter.
2. To obtain a perfect alignment of the camera with respect to the eyepieces, loosen the fixing screw ① on the C-mount adapter, rotate the camera until correct alignment is achieved, then tighten the screw. (Fig. 36)

### 12.2 Parfocality adjustment

In order to have the same focus when observing the specimen through the eyepieces and on the screen/monitor, please check that the microscope is properly installed and follow the instructions below.

1. Use a low power magnification and focus the specimen.
2. Reach the highest magnification available using the zoom knob and then focus the specimen again.
3. Turn on the live-view on the camera, without changing the focus on the microscope.
4. While observing the image on the screen/monitor, adjust the focus by turning the knurled knob ② on the C-mount adapter. (Fig. 37)

---

## 13. Maintenance

### Operative environment

This microscope is recommended to be used in a clean, dry and shock free environment with a temperature of 5°-40°C and a maximum relative humidity of 75 % (non condensing). Use a dehumidifier if needed.

### To think about when and after using the microscope

- The microscope should always be kept vertically when moving it and be careful so that no moving parts, such as the eyepieces, fall out.
- Never mishandle or impose unnecessary force on the microscope.
- Never attempt to service the microscope yourself.
- After use, turn off the light immediately, cover the microscope with the provided dust-cover, and keep it in a dry and clean place.

### Electrical safety precautions

- Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off- position.
- Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users do have full responsibility to use this equipment safely.

### Cleaning the optics

- If the optical parts need to be cleaned try first to: use compressed air.
- If that is not sufficient: use a soft lint-free piece of cloth with water and a mild detergent.
- And as a final option: use the piece of cloth moistened with a 3:7 mixture of ethanol and ether.
- **Note: ethanol and ether are highly flammable liquids. Do not use them near a heat source, near sparks or near electric equipment. Use these chemicals in a well ventilated room.**
- Remember to never wipe the surface of any optical items with your hands. Fingerprints can damage the optics.
- Do not disassemble objectives or eyepieces in attempt to clean them.

**For the best results, use the OPTIKA cleaning kit (see catalogue).**

If you need to send the microscope to Optika for maintenance, please use the original packaging.

## 14. Troubleshooting


Review the information in the table below to troubleshoot operating problems.

PROBLEM	CAUSE	SOLUTION
<b>I. Optical Section:</b>		
The illumination is ON, but the field of view is dark	The plug is not connected to the illumination	Connect the cable
	The brightness is too low	Adjust to a proper setting
	Fluorescence filter selector is not in a click stop	Move the selector to a click stop
	Fluorescence filter is not suitable for the specimen	Use a suitable filter
	Fluorescence filter is not matching the LED source	Match LED light and fluorescence filter.
Dust and stains can be seen in the field of view.	There are stains and dust on the specimen	Clean the specimen
	There are stains and dust on the eyepiece	Clean the eyepiece
Poor image quality: <ul style="list-style-type: none"> <li>The image is not sharp</li> <li>The contrast is not high</li> <li>The details are not clear</li> <li>Image glares</li> </ul>	The lenses (additional lens, objective, eyepieces) are dirt	Thoroughly clean all the optical system
One side of the image is out of focus.	The specimen is out of place (tilted)	Place the specimen flat on the stage.
<b>II. Mechanical Section:</b>		
The focus knob is hard to turn.	The tension adjustment collar is too tight	Loosen the tension adjustment collar
The focus is unstable.	The tension adjustment collar is too loose	Tighten the tension adjustment collar
<b>III. Electric Section:</b>		
The LED doesn't turn on.	No power supply	Check the power cord connection
The brightness is not enough	The brightness adjustment is low	Adjust the brightness
The light blinks	The power cord is poorly connected	Check the power cord
<b>IV. Viewing tube assembly:</b>		
The field of view of the two eyes is different	The interpupillary distance is not correct	Adjust the interpupillary distance
	The diopter correction is not right	Adjust the diopter correction
	The viewing technique is not correct, and the operator is straining the eyesight	When look into the eyepieces, do not stare at the specimen but look at the whole field of view. Periodically, move the eyes away to look at a distant object, then back into the eyepieces
<b>V. Microphotography and video:</b>		
The image is unfocused	Incorrect focusing	Adjusting the focus system as in the present manual
Bright patches appear on the image	Stray light is entering the microscope through the eyepieces and through the camera viewfinder	Cover the eyepieces and the viewfinder with a dark cloth

---

## Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

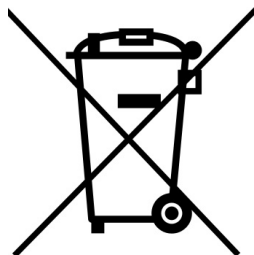The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste. The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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**OPTIKA® S.r.l.**