VC Series

INSTRUCTION MANUAL

Model
VC-05

Version: 2
Issued: 03, 12, 2014

Safety Warning

- 1.0 **Do not install the camera near electric or magnetic fields.**
Install the camera away from radio transmitter, electric motors, audio speakers because the magnetic fields generate from above devices will distort the video image.
- 2.0 **Do not install the camera in the high temperature environment.**
Install the camera away from stoves, or other heat generating devices as the high temperature could cause deformation, discoloration or other damages of the camera.
- 3.0 **Never face the camera toward the sun.**
Never face the camera at the sun or other extremely bright objects whether it is in use or not.
- 4.0 **Cleaning**
Do not touch the surface of CCD sensor with hands directly. Use a soft cloth to remove the dirt from the camera body. Use lens tissue or a cotton tipped applicator to clean the CCD sensor and the camera lens.

Specification

Main Features :

- Colour CDD 1/3" with 420 TV lines
- 12Volts power suppli
- PAL video signal

Micro Camera Specification-Analog Signals Series

VC-05 Analog Camera:

SPECIFICATION	
CONNECT MODE	DIRECTLY PUT INTO EYEPIECE TUBE
OUTPUT SIGNALS	VIDEO SIGNAL
PIXELS	370K PIXELS
TV LINES	420 TV LINES
IMAGE SENSOR	1/3"SENSOR
IMAGE DEFINITON	628(H)×586(V)-PAL 510(H)×496(V)-NTSC
WHITE BALANCE	AUTO
SNR	≥48DB
DYNAMIC RANGE	≥56DB
WORKING VOLTAGE	DC 12V
CONSUMPTION	≤85 MA
WORKING TEMPERATURE	-10 – 50°C


Installation Instructions

Put the specimen on focus , then put off the eyepiece from your microscope and put on the camera, use the 30mm black ring if the hole is bigger than the nosepiece of your camera.

Connect the power supply cable and the TV cable, then turn on the TV and the camera.

Equipment disposal

Art.13 Dlsg 25 july 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

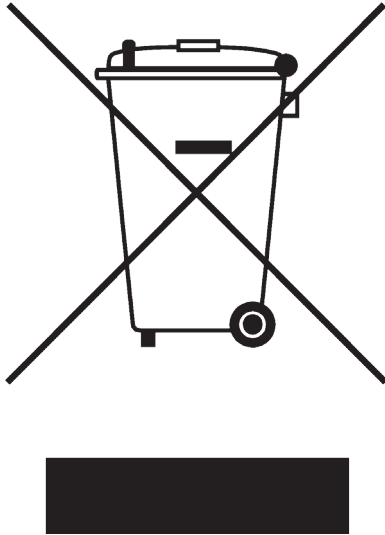The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection.

The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment.

Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.