OPTIKAM B1/B3/B5/B9

INSTRUCTION MANUAL

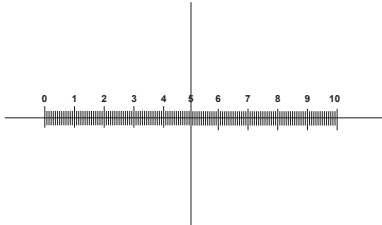
Code	Model
4083.B1	OPTIKAM B1
4083.B3	OPTIKAM B3
4083.B5	OPTIKAM B5
4083.B9	OPTIKAM B9

Version: 10
Issued: 24, 01, 2017

M-005

Micrometer slide for software calibration, 1mm/10um, 10mm/100um.

1 DIV=0.1mm

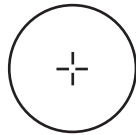
For stereo microscopes calibration

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1 DIV=0.01mm

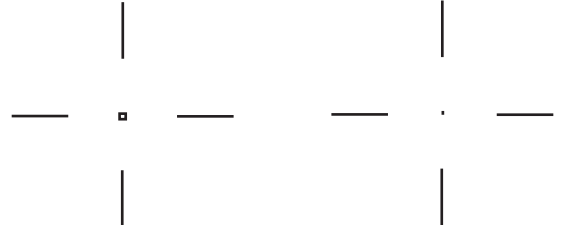
For biological microscopes calibration

For biological microscopes calibration

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Ø 0.15mm

For checking the calibration done by drawing a line or a circle

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Installation of camera drivers and software

Operating system requirements: Windows XP, Windows Vista, Windows 7, Windows 8 or Windows 10. If your PC is Windows XP: before the installation of your Camera please check if Service Pack2 is installed. If not, please update your XP from the Microsoft web site. If your PC has a USB port less than 2.0, there may be some problems in capturing image from the camera due to the slow transmission.

1. Please insert your CD and run:
 \optika vision lite \ setup.exe
2. Now Optika Vision Lite icon appears on your desktop.
3. Please run \ optika IsView \ setup.exe
4. Now the Optika IsView icon appear on your desktop.
5. Choose from the folder “drivers” the model of camera you want to use and run all the “.exe” files.
6. The installer may show warnings about missing Windows certification: ignore such warnings and select “Continue”. Note : do NOT cancel the installation.
7. At the end of the installation process restart the PC if requested.
8. Plug the Camera in your PC: the camera will be identified and Windows will recognize the driver which has just been installed from the CD.
 Then ignore any message about Windows compatibility and select “Continue” .
 Note : do NOT cancel the installation.
9. At the end of the identification process, disconnect your Camera and try to plug it again, in order to check if the Driver was installed correctly (no Windows error messages should appear).
 Should any error message appear on your PC, you will find a “\Troubleshooting” folder on your CD which shows how to solve any problem which Windows could give.

Some notes

You should know that your Windows PC needs to install a Driver before capturing images from the camera: after installing the driver from the CD, Windows needs to recognize the camera before using it.

Often a PC has some ports on the front and some on the back: the ports on the back are usually faster. We suggest to use these ones.

All our cameras, even though they may have an external connection, do not need any external power supply because it comes from the USB cable of your PC.

If you connect the camera into another USB port, Windows needs to recognize its Driver again even if it has

already done this operation for the first USB port.

Generally, every time you use a peripheral device and after that you connect your Camera, then Windows needs to recognize the Driver again.

It may happen sometimes that Windows doesn't recognize a Camera using one of your PC USB ports, while gives no problem using another port.

To prevent any installation problem we have put on the CD all the software and drivers in specific folders with an appropriate names.

How to test your Camera

In order to check if your Camera works correctly, it's not necessary to put it into the microscope: it's enough to connect it into the USB port, run Vision Lite using the icon on your desktop, select the right driver and check whether the shown image is not totally black.

Select the right drivers

It's very easy:

For Optikam B1 Camera you have to select the driver:
OPTIKA TCD 1.21

For Optikam B3 Camera you have to select the driver:
OPTIKA TCB 3.0

For Optikam B5 Camera you have to select the driver:
OPTIKA TCB 5.0

For Optikam B9 Camera you have to select the driver:
OPTIKA TCB 10

With Optika Vision Lite you have to tell the Software which is the Driver by selecting the microscope-shaped icon.

With Optika IsView the live image coming from your camera will automatically appear.

Getting started with Optika Vision Lite, a very easy measuring and documentation software.

Run Optika Vision Lite by clicking on its icon on your desktop.

Select the right Driver for capturing the image. Note that after choosing the driver and before capturing the image you can set the brightness, saturation, gamma ecc..
Then you can capture the image.

The image is now ready for being manipulated using Optika Vision Lite.

Getting started with Optika IsView, a complete software for measuring, adjusting, capturing images and video clip.

Run Optika IsView by clicking on your desktop.

The live image with easy buttons for set your camera live view will automatically appear.

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection.

The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment.

Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA S.r.l.