OPTIKAM HDMI EASY

INSTRUCTION MANUAL

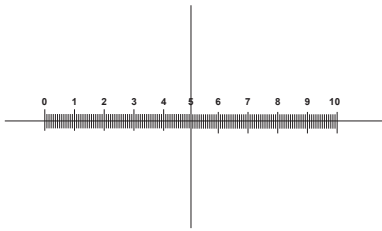
Code	Model
4083.13E	OPTIKAM HDMI EASY

v 1.0 2017

M-005

Micrometer slide for software calibration, 1mm/10um, 10mm/100um.

1 DIV=0.1mm

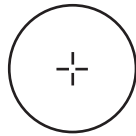
For stereo microscopes calibration

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1 DIV=0.01mm

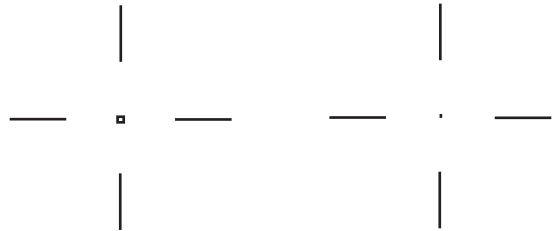
For biological microscopes calibration

For biological microscopes calibration

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Ø 0.15mm

Ø 0.07mm

For checking the calibration done by drawing a line or a circle

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Installation of camera drivers and software

1. Operating system requirements: Windows Vista, Win 7, Win 8, Win 10
2. Please insert your CD and run:
\\optika Vision Lite \ setup.exe
3. Now the Optika Vision Lite icon appears on your desktop.
4. You don't need to install any driver.
5. Plug the Camera in your PC: the camera will be identified automatically.

How to test your Camera

In order to check if your Camera works correctly, it's not necessary to put it into the microscope: it's enough to connect it into the USB port, run Vision Lite using the icon on your desktop, select the right driver and check whether the shown image is not totally black.

Select the right drivers

It's very easy, you have to select the driver: YW 2310 CAMERA

With Optika Vision Lite you have to tell the Software which is the Driver by selecting the microscope-shaped icon.

Getting started with Optika Vision Lite, a very easy measuring and documentation software.

Run Optika Vision Lite by clicking on its icon on your desktop.

Select the right Driver for capturing the image.

The image is now ready for being manipulated using Optika Vision Lite.

Camera interface

This HDMI Easy camera has this interface on the back :

1. USB port
2. Micro-SD port
3. HDMI port
4. Power Indicator : Red Led
5. Recording Indicator : Green Led
6. Power Adapter port
7. Keyboard

The keyboard has six buttons as below :

1. Power ON/OFF, Freeze/Unfreeze
2. UP, Zoom In
3. OK, Snapshot , Record
4. Menu
5. DOWN, Zoom Out
6. Model, AE

Operations

1.1 Power on your device

1. Plug the power adapter in.
2. Press the ON/OFF button, the Red power indicator lights on.

1.2 Live image preview

Connect the camera to a TV monitor with HDMI input, The TV Monitor will show the live preview image 1280 x 720 resolution.

The camera starts by default in the snapshot mode.

Common operations

1.1 Snapshot

In preview mode Press the "OK" button, a 5 M Pixel picture will be saved in the Micro-SD Card.

1.2 Recording

Press the "MODEL" button for video recording , a recorder ICON appears.

Now you can press the "OK" button for saving your video and press it again to stop the recording.

When you are recording a flickering red circle appears too.

1.3 Playback

Press the "MODE" button to playback mode, either a triangle for the images or a film ICON appears.

Press the "UP" or "DOWN" buttons to navigate among them and press "OK" to play.

Press the "MODE" button again for returning to the preview mode.

1.4 Digital Zoom

Press the "UP" or "DOWN" buttons in the preview image , the magnification number appears on the screen.

Press and hold one of the two buttons for a quick activation of the zoom.

1.5 Freeze/Unfreeze

In preview mode press the button “ON/OFF” for freezing the preview image.

1.6 Auto/Manual Exposure

In preview mode press and hold the “MODEL” button to switch the exposure from Auto to Lock mode.

1.7 Exposure

Press the “Menu” button and then choose “Exposure”, now press “UP” or “DOWN” buttons to change the EV offset for increasing or decreasing the exposure.

1.8 Image Color

Press the “Menu” button and choose “Color”, now press “UP” “DOWN” buttons to change the image effect from Color to Gray or Sepia.

1.9 Line set

your HDMI Easy camera provides eight lines, all of them can be shown horizontally or vertically and the position of them can be changed, into details you can set as follows:

1.9.1 Line set into details

Press the “Menu” button and choose “Line Set”, then choose “Line”, now you can press “MODE” or “OK” button to the number of line you want to show.

- You can choose the line direction by changing the “Line DIR” parameter.
- You can change the position of the line by changing the “Line Position” parameter.
- You can change the color of the line by changing the “Line Color” parameter.
- You can change the size of the line by changing the “Line Weight” parameter.

1.10 Crosshair

Press the “Menu” button and choose “Line Set”, now select “Cross Hair” and press “MODE” or “OK” button for activating the cross.

1.11 Language

In preview mode select the “Menu” button and choose “Language”, the languages interfaces appears with ten available languages.

1.12 Light Frequency

In preview mode press the “Menu” button and choose “Frequency”, now you can select the right current frequency in your country, this reduces the image flicker behaviour if appears.

1.13 Format Micro-SD Memory Card

If the Micro-SD card is not working properly or it needs to be formatted please press the “Menu” button and choose “Format”, then choose “OK”, now the Micro-SD card will be formatted.

PAY ATTENTION : All the data on your Card will be erased.

1.14 Default Setting

Press the “Menu” button and choose “Default Setting”, now choose “OK” to get the default parameters.

How to use your HDMI with the PC

If you connect the camera to a computer with the USB Cable the TV preview will be disconnected, the USB connection has the priority.

Please also note :

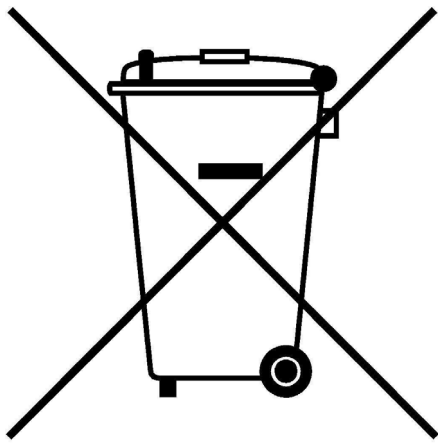
1. The USB camera needs the power adapter plugged in and the camera itself turned up
2. If there is a plugged in Micro-SD card the USB cable will be used to control the Micro-SD, so for using the USB camera in live view on the PC please remove the Micro-SD card before.

Power off your HDMI camera

Press and hold the “ON/OFF” button for more than 5 seconds, then the camera saves all the parameters and turns off.

Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection.

The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment.

Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA S.r.l.