WiFi Camera

INSTRUCTION MANUAL

Model
OPTIKAM WiFi

Version: 4 2019

Table of Contents

1. Warning	3
2. Symbols and conventions	3
3. Safety Information	3
4. Intended use	3
5. Overview	4
6. Unpacking	4
7. Assembling	4
8. Connecting the Camera	5
8.1 Use with PC	5
8.2 Use with WiFi	5
9. Installation of camera drivers and software on your PC	5
10. Starting the Software	5
11. Selecting the right driver	6
13.1 Camera connection to a Browser	9
13.2 Camera connection to IOS	10
13.3 Camera connection to Android	14
14. Advanced settings	14
Equipment disposal	15

1. Warning

This device is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use. We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users. We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.

2. Symbols and conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

CAUTION

This symbol indicates a potential risk and alerts you to proceed with caution.

ELECTRICAL SHOCK

This symbol indicates a risk of electrical shock.

3. Safety Information

Avoiding Electrical Shock

Before plugging in the power supply, make sure that the supplying voltage of your region matches with the operation voltage of the equipment and that the lamp switch is in off position. Users should observe all safety regulations of the region. The equipment has acquired the CE safety label. However, users have full responsibility to use this equipment safely. Please follow the guidelines below, and read this manual in its entirety to ensure safe operation of the unit.

4. Intended use

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

5. Overview

6. Unpacking

The device is housed in a moulded Styrofoam container. Remove the tape from the edge of the container and lift the top half of the container. Take some care to avoid that the optical items (objectives and eyepieces) fall out and get damaged.

 Do not touch with bare hands optical surfaces such as lenses, filters or glasses. Traces of grease or other residuals may deteriorate the final image quality and corrode the optics surface in a short time.

7. Assembling

Once opened the camera box, the parts are the following:

- ① Camera body + antenna
- ② Power supply
- ③ USB cable
- ④ Stage micrometer

- ⑤ Projection lens
- ⑥ 30 mm reduction ring
- ⑦ 30,5 mm reduction ring
- ⑧ "C" mount additional ring

8. Connecting the Camera

The camera can be used as a normal PC camera or as a WiFi camera. Please refer to specific connection guide (below) in order to ensure a proper operation of the device.

8.1 Use with PC

Plug the USB cable into the USB socket. Camera power will be ensured by the USB cable and therefore there is no need to plug the provided power supply.

8.2 Use with WiFi

Plug the jack of the power supply into the socket of the camera. Using the camera as a WiFi unit doesn't require the connection of the USB cable.

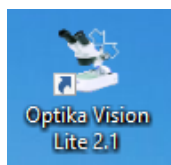
NOTE: the connection of the power supply automatically disables the USB connection. If the user wants to use the camera under PC, he must not plug the power supply.

9. Installation of camera drivers and software on your PC

1. Operating system requirements: all the Windows versions.
2. Insert the provided CD and run: `\\optika vision lite\ setup.exe`
3. Now the Optika Vision Lite icon appears on your desktop.
4. You don't need to install any driver.
5. Plug the Camera in your PC: the camera will be automatically identified.


10. Starting the Software

Run **Optika Vision Lite** by clicking on its icon on your desktop:


11. Selecting the right driver

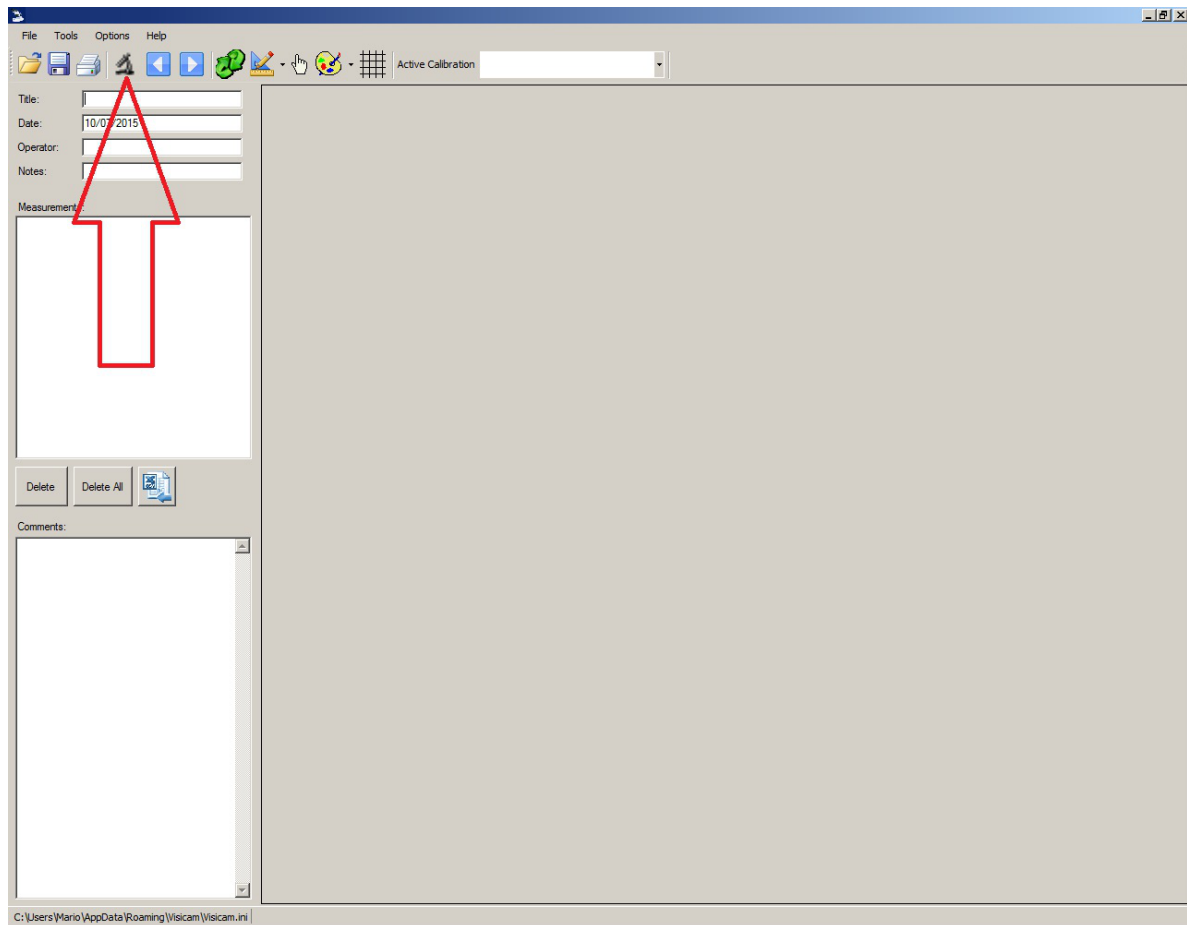Click on microscope icon

A drop-down list appears:

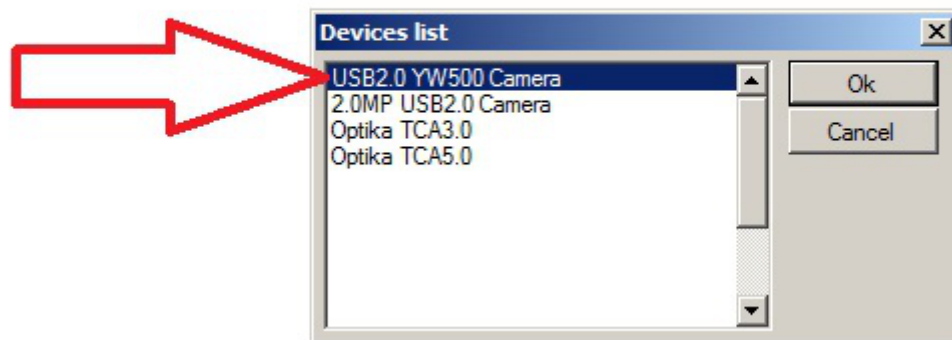Select **USB2.0 YW500 Camera**

Now the live image appears.

Choose the best resolution you need accordingly to a Video capture or Image capture.

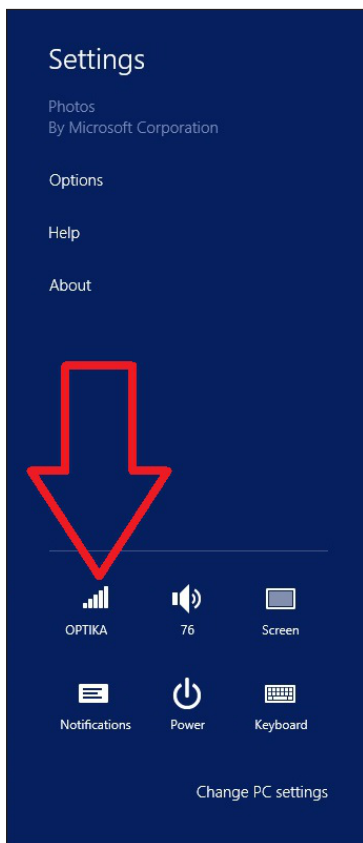
For a complete explanation of the Optika Vision Lite functions, please refer to the relevant Instruction Manual.

13. WiFi Camera connection

Please connect the WiFi camera to the power supply and wait a minute, allowing the PC to recognise the new WiFi device.

The connection can be done on any device having a Wifi connection.

Example for Windows connection: from **Start** menu select **Settings** and choose the **WiFi** signal.

You should find **Netcam012** or **OpenWrt_C12C5F** (see below). Connect to the hotspot:

It doesn't matter if the WiFi signal is found as **Limited**.

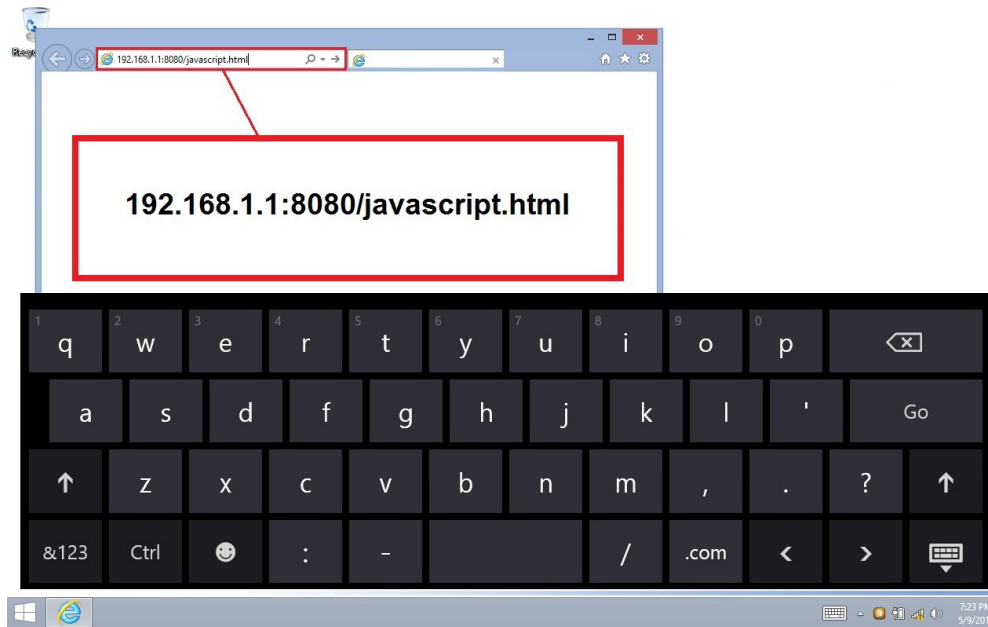
Now the WiFi connection has been established.

13.1 Camera connection to a Browser

These settings can be used for PC, Laptop or Tablet (under Windows O.S.) or any phone (Windows, IOS or Android).

Run Google Chrome and type in the IP address: **192.168.1.1:8080/javascript.html**

- **NOTE: The application has been specially designed for Google Chrome; with other browsers it may not be working properly.**

Immediately, the Wifi camera connects to Windows and sends all the informations to get a live preview:

From this page it is very easy to save videos or images; it can be done by selecting **Snapshot** or by keeping your finger pressed on the touch screen and then choosing **Save video** or **Save image**.

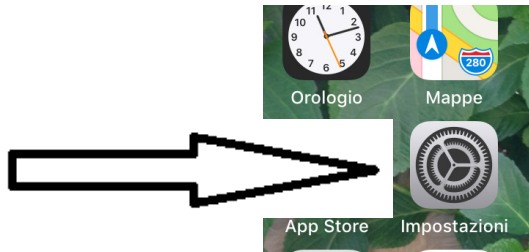
13.2 Camera connection to IOS

To connect IOS device please download from the APP store the following app:

IPCam Viewer lite

<https://itunes.apple.com/us/app/ip-cam-viewer-lite/id397322405?mt=8>

1. From your IOS device tap on **Setting** icon:

2. Select **WiFi**:

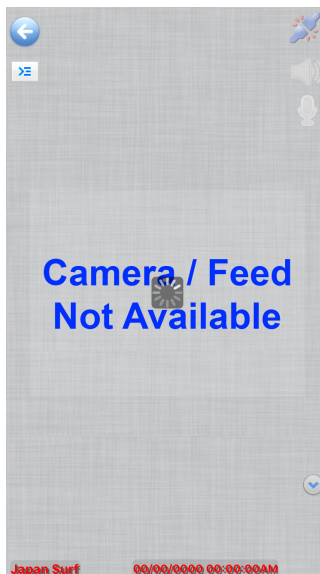
3. Now select **Netcam012**. WiFi connection is established.
4. Exit from **Settings** menu.


5. Launch **IPCam Viewer** app. The following screen appears:


6. Tap on any of the windows **Camera Feed not Available**. The dialog box changes as follows:

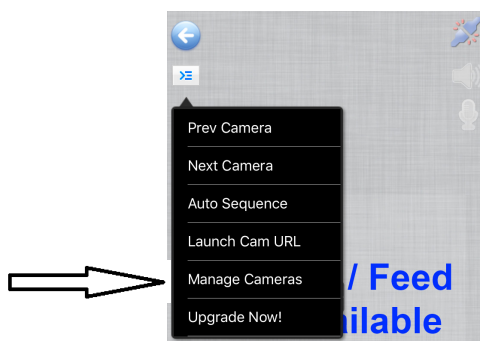
7. Tap on the icon  on the upper left side of the screen. A new dialog appears:

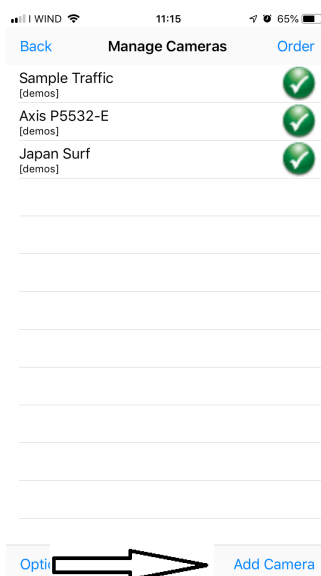8. Tap on **Manage Cameras** to select the camera you want to connect. The dialog changes as follows:

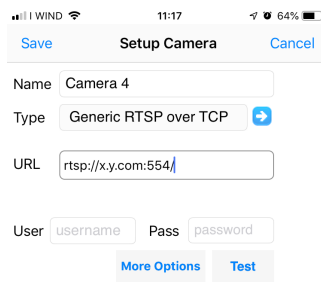9. Tap on **Add Camera** on the lower left corner of the screen. A new dialog appears:


Enter the full RTSP over TCP URL (eg. `rtsp://x.com:554/video.mp4`). If output is gray, enable Prefer Quality over Speed in More Options.
Click 'More Options' to invert pan, tilt, image, etc.

10. Here you can rename the camera using the name you prefer.

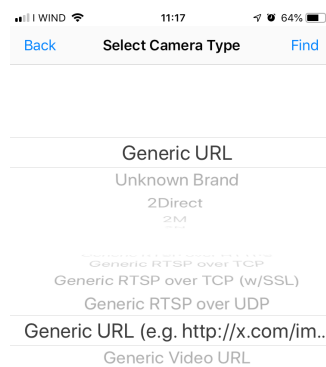
11. Select the camera type: tap on **Generic URL**.


12. In the drop-down list below the camera type select **Generic URL (e.g. http://x.com/....)**.

13. Then tap on **Back**.

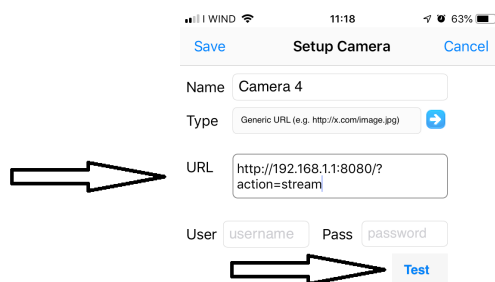
14. You are switched back to the previous dialog. Here you must type in the URL address for the proper connection to the camera: **`http://192.168.1.1:8080/?action=stream`**.

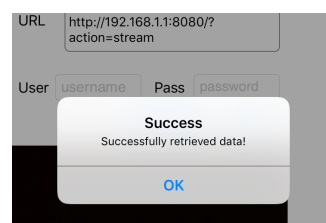
15. Tap on **Test**. If the connection succeeded, you will get this message:

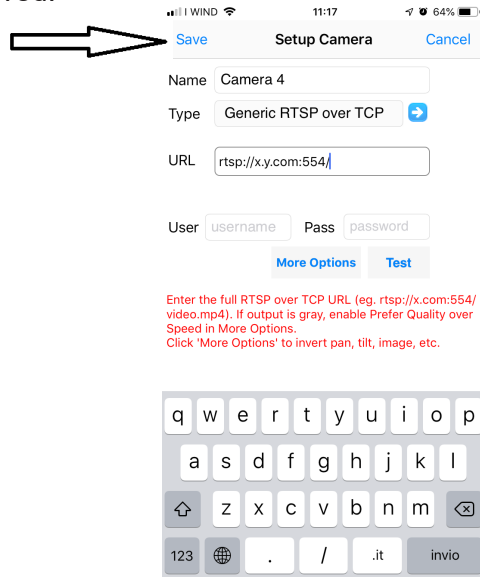
16. Tap on **OK**. Then tap on **Save**.
17. Connection will be saved.

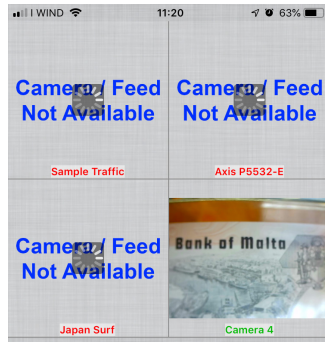18. The initial dialog box appears but now with a new camera option available.

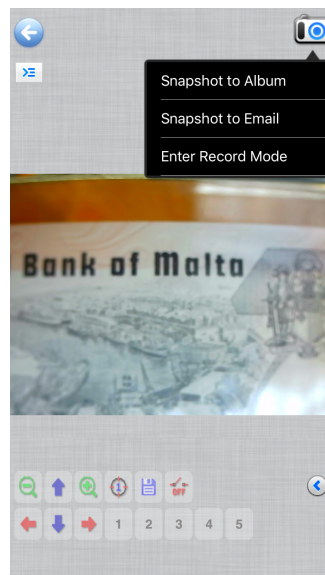19. Tap on **Camera 4**. The live image coming from the WiFi camera is now available.


13.3 Camera connection to Android

To connect Android device please download from the Google Play the following app:

SimpleMjpegView

<https://play.google.com/store/apps/details?id=com.tk.android.simplemjpeg>

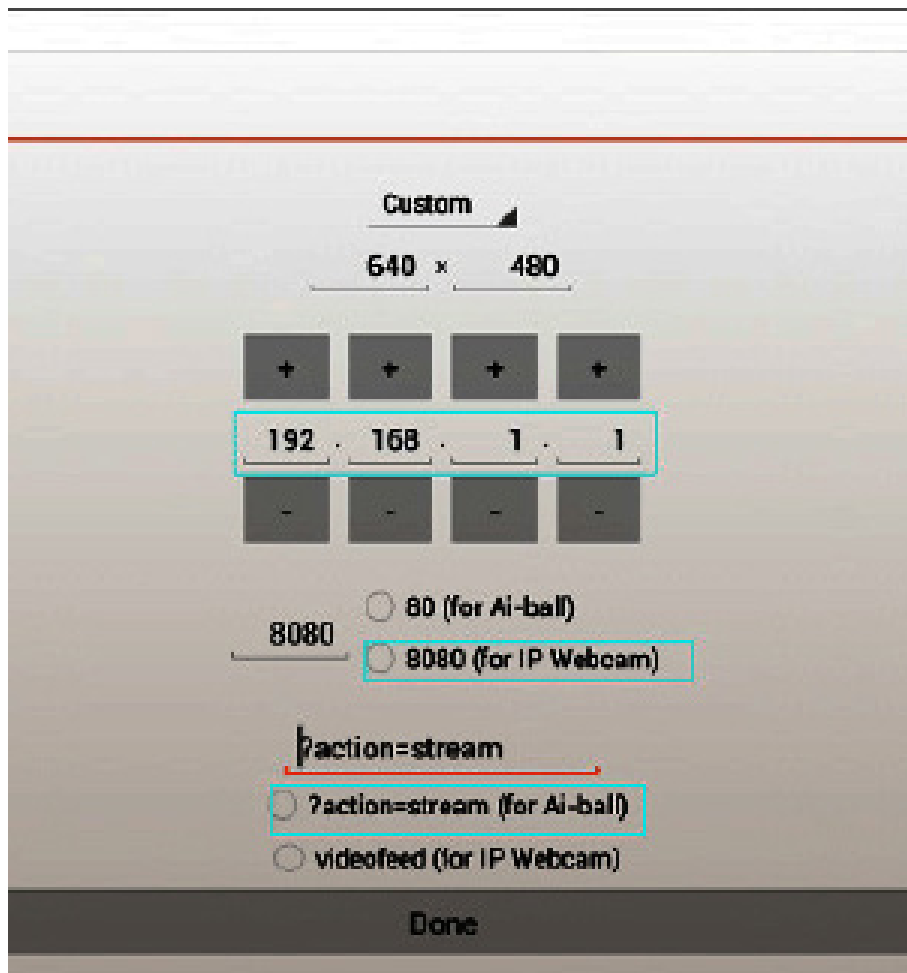Please note that the camera can be detected by using this parameters only:

14. Advanced settings

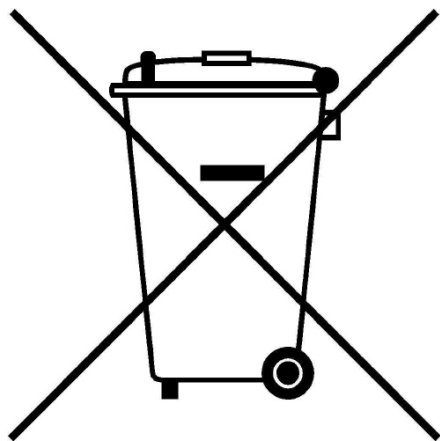
The camera is shipped from the factory with default settings for the WiFi connection. In order to ensure a fast frame rate, when used under WiFi condition, the resolution is set at 640x480.

However it is possible to change these default settings, although it is not advisable to do this.

Should the user needs to change settings accordingly to personal needs, please contact our Technical Dept. for the operative instructions.

Equipment disposal

Art.13 Dlsg 25 july 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection.

The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment.

Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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OPTIKA® S.r.l.