OPTISCAN

INSTRUCTION MANUAL

Model	Code
OPTISCAN10	4083.SC10

Version: 2
Issued: 06, 02, 2017

Table of Contents

Warning
Package Contents
Overview
Set up
How to scan microscope slides
There are four major parts of the user interface

Warning

Check the package content before getting started.

Important! Save the original box, receipt and packing material for future shipping needs.

- Only use the supplied USB-Cable. Connect the computer to the scanner directly using the USB 2.0 cable.
- Do not use a Hub or extension cable. Only use the USB ports on the rear of desktop computers.
- If there are any other devices (multifunction device, printers, scanners, webcam) connected to the computer through USB disconnect them.
- A minimum of 1GB RAM (Windows/Mac) is required.
- **Warning! Using any other power adapter may severely damage the scanner and will void the product warranty.**

- Operating Ambient Temperature Range 50° to 104°F (10° to 40°C)
- Operating Ambient Humidity Range 15% to 76% relative humidity (non-condensing)
- **Note:** If the scanner is stored or to be transported in cold temperature and then brought into a warm / humid environment condensation may occur.
- **Note:** Do not turn on the scanner unless it is within this range. If the scanner has been stored or transported outside of this range, allow it to return to within this range before turning it on.

Package Contents

DESCRIPTION	QUANTITY
Power Supply	1
Slide Scanner	1
USB 2.0 Cable	1
SciView Software Cd	1

Overview

Set up

Please read this section completely before using your scanner with the SciView software. Always set up your scanner on a smooth surface, such as a desk or table.

DURING INSTALLATION, A WARNING DIALOGUE BOX MAY APPEAR, IF THIS IS THE CASE PLEASE SELECT "INSTALL THIS DRIVER ANYWAY".

1. In "Found New Hardware Wizard" window, select "Search automatically or Browse computer for driver software."
2. In "Completing the Found New Hardware Wizard" window, click "Finish".

NOTE: Not all PCs will display the "Found New Hardware" window, some will install in the background.

INSTALLING SOFTWARE AND DRIVER

1. Double click the "SciView" icon to begin the installation process. Follow the onscreen prompts to complete the installation process.

INSTALLING HARDWARE

1. Locate the power switch on the back side of the scanner. Make sure the switch is in OFF position.
2. Plug the AC adapter into an AC power socket, before plugging it into the scanner.
3. Attach the USB cable to the computer using the standard USB adapter end. Attach the USB adapter end of the cable to back panel of the scanner.

Please Note: When installing your new scanner onto a Windows system that already has another USB Flatbed scanner or All-in-one Scan/Copier/Printer device installed, it may be necessary for you to disconnect the other scanner from the USB port for the scanner to work properly. This does not apply to Wi-Fi Printing/scanning devices.

How to scan microscope slides

Note: Before using the scanner, make sure you have properly prepared your slides according to the slide making instructions.


STEP 1 – Open SciView Application

Open SciView Application from the Desktop icon or Start (Windows icon) → All Programs → SciView

STEP 2 – Select Scanning Mode

Microscope Slides

Tissue Sections
Common Things
Aquatic Organisms

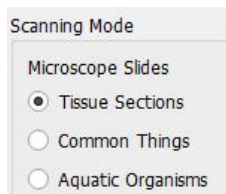
STEP 3 – Place Slide into the scanning area (see picture)

Once the slide is secure in the slide holder of the scanner, close the lid gently.


STEP 4 – Set Basic Scan Preferences

1. In the SciView software click on Preferences...
2. Configure the settings for this session (for Custom 1-5 settings see Step 4a)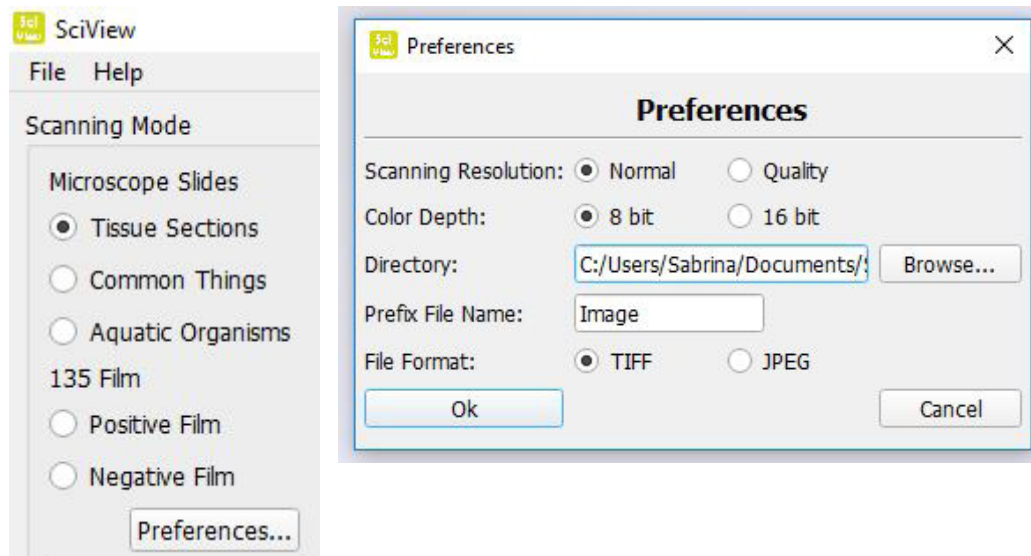

STEP 4a – Set Basic Scan Preferences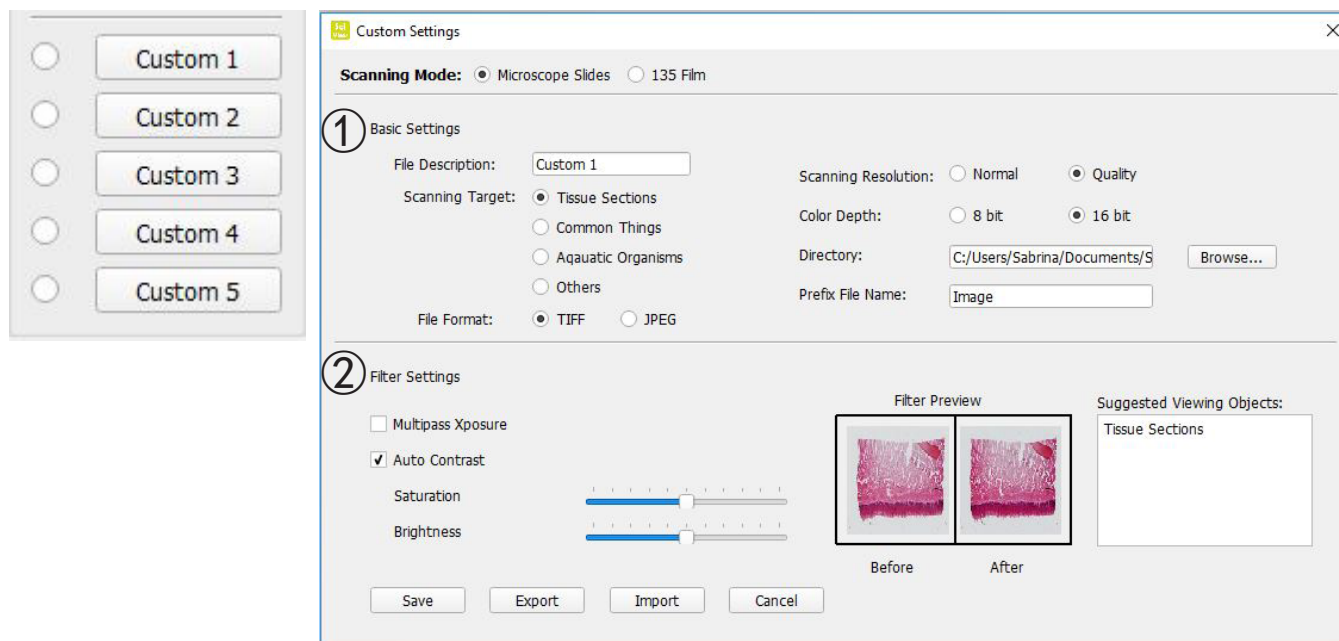

Once in Custom Settings you can select your scanning mode
Microscope Slides

① Basic Settings:

1. **File Description:** change the name on the Custom button
2. **Scanning Target:** tissue Sections; Common Things; Aquatic Organisms; Other
3. **File Format:** TIF (uncompressed) or JPEG
4. **Scanning resolution:** Quality or Normal
5. **Color Depth:** 8 or 16 bit color
6. **Directory:** change where the files are saved to
7. **Prefix File Name:** the name of the scanned images when saved

② Filter Settings:

1. Automatic adjustments applied to the scans
2. Multipass Exposure
3. Auto Contrast
4. Saturation
5. Brightness

Save / Export / Import


Save: apply these settings to the current scan session

Export: saves these settings to a directory of your choice for future recall

Import: recalls previously exported custom settings

STEP 5 – Prescan

Select the “Prescan “ to prescan the slide currently loaded into the scanner.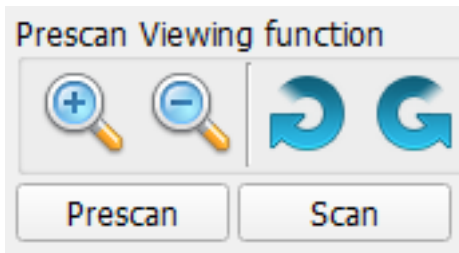

Pre-Scanning should be used to manually adjust color/edit settings for each slide before doing the final scan. It's possible to set different settings for each slide to be scanned.

1. Once the whole image shows up in the Prescan window, you may use the Rotate and Flip functions to adjust how your images will be presented after scanning.
2. After settings have been applied, press and hold left mouse button at one corner of the area where you want to scan, and then slide diagonally to the opposite corner of the area to highlight the area before the final Scan. See Fig 1. You may also move the highlight box to a new area you want to scan by pressing and holding left mouse button at the center of the highlight box after an arrow appears. Fig. 2 & 3.

Note: You may delete the highlight by using the delete button on the keyboard.

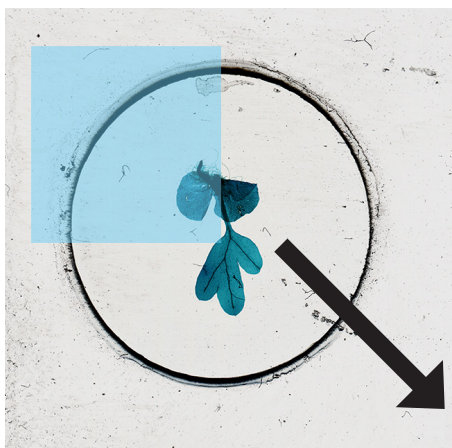
Fig. 1


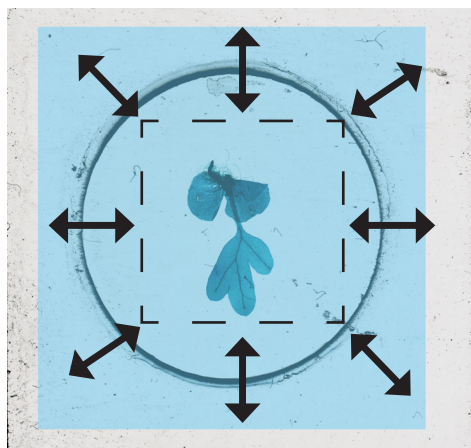
Fig. 2


Fig. 3

STEP 6 – Scan / Exit

Select Scan to scan and save the file.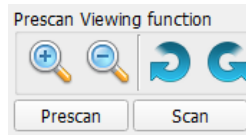

Once you have finished scanning Exit the program by clicking on File on the menu then Close.

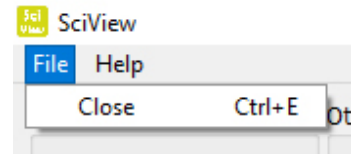One Button Scan

1. Open the SciView (see above to acquire the SciView driver program)
2. Configure Scan Settings (see above)
3. Press the one touch scan button to begin the scan to file process.

Suggestions for Making Your Own Prepared Microscope Slides

Always make sure clean slides and cover glass are used.
Handle slides carefully by the edges to avoid fingerprints.
If you have forceps or tweezers, use it to carefully handle cover glass.
The thin cover glass will break if too much pressure is applied!

It is highly recommended to use a blade for cutting specimens. Keep the blade covered and store in a safe place when not in use.

Making a Temporary “Wet” Mounted Slide

• Tissue sections:

1. Use the blade to carefully cut the specimen as thin as possible.
2. Place the specimen on the slide.
3. Use the plastic dropper to place one drop of water on the specimen. Note: Steps 2 & 3 can be reversed. Find the best protocol for each specimen.
4. Carefully place a cover glass over the specimen, allowing the water to spread under the glass.
5. The cover slip should be flattened against the specimen and slide as much as possible, removing air bubbles if any.
6. Blot any excess water carefully with paper tissues.
7. Carefully place slide on the scanner holder (cover glass side up) and scan.
8. Remember to properly name your scanned images in the SciView software.

• Tip:

1. Slides may be carefully washed and re-used for future specimens. Cover glass is extremely fragile and not recommended for re-using. Carefully discard after use.
2. With scanner, you now have the ability to observe a wide range of different objects, such as tissue sections, common things, and aquatic organisms. Below are some more ideas for observation.
3. Many specimens are transparent. You may try staining your tissue specimens in food color for a couple of minutes before mounting it on the slide. Use the tweezers to pick up the “stained” specimen.
4. Scanning fresh specimens that are several-cells thick with pigments/colors, such as plant leaves and black hair, will likely return with dark black images without showing any detail of the specimens.

Some More Ideas for Observation

- **Crystals:**

1. Mix salt/sugar with some warm water in a small cup.
2. Use the plastic dropper to place a couple drops on a slide. Let it dry and observe the re-crystallized salt/sugar without the cover slip.
3. Choose Tissue Sections Mode for your first scan.

- **Flower pollens:**

1. Shake some grains of pollen from a flower onto a blank slide.
2. Follow the Wet Mount protocol for the rest steps.
3. You may also try observing the pollens directly without making a wet mounted slide.
4. Choose Tissue Sections Mode for your first Scan.

- **Hairs in white or light colors:**

1. Use human or pet hair, or pieces of feathers collected from outside.
2. Use a small piece of tape at each end of the hair to hold it on the slide or follow the wet mounted procedure to prepare your slides.
3. Try comparing hairs from various animals.
4. Choose Common Things Mode for your first scan.

- **Fibers :**

1. Find different kinds of paper, a newspaper or printing paper, or fibers from different types of fabric.
2. Tear paper/fabric to obtain few threads of fibers from paper or fabric.
3. Follow steps for hairs to hold fibers on the slides.
4. Compare natural fibers such as cotton and wool to synthetics like polyester.
5. Choose Common Things Mode for your first scan.

- **Water from a pond that contains live organisms:**

1. Place a drop onto a clean slide.
2. Cover with a cover slip then observe.
3. You may need to mix the sample with couple drop of 95% ethanol to paralyze those moving organisms for easy observation.
4. Choose Aquatic organisms Mode for your first scan


- **Very small insects or parts such as legs or wings of larger bugs.**

1. Always handle insects with tweezers and wash hands after working with your specimens.
2. Use a small piece of tape at each end of the specimen to hold it on the slide.
3. Choose Common Things Mode for your first scan.


- **Mold on old bread or other food**

1. You may make your own mold by leaving a small piece of bread in a dish for several days add a small amount of moisture to encourage the mold growth.
2. Add couple drops of water on the center of slide.
3. Use tweezers to pick a tiny amount of mold from bread then spread it in the water.
4. Cover with a cover slip then observe.
5. Choose Tissue Sections Mode for your first scan.

There are four major parts of the user interface: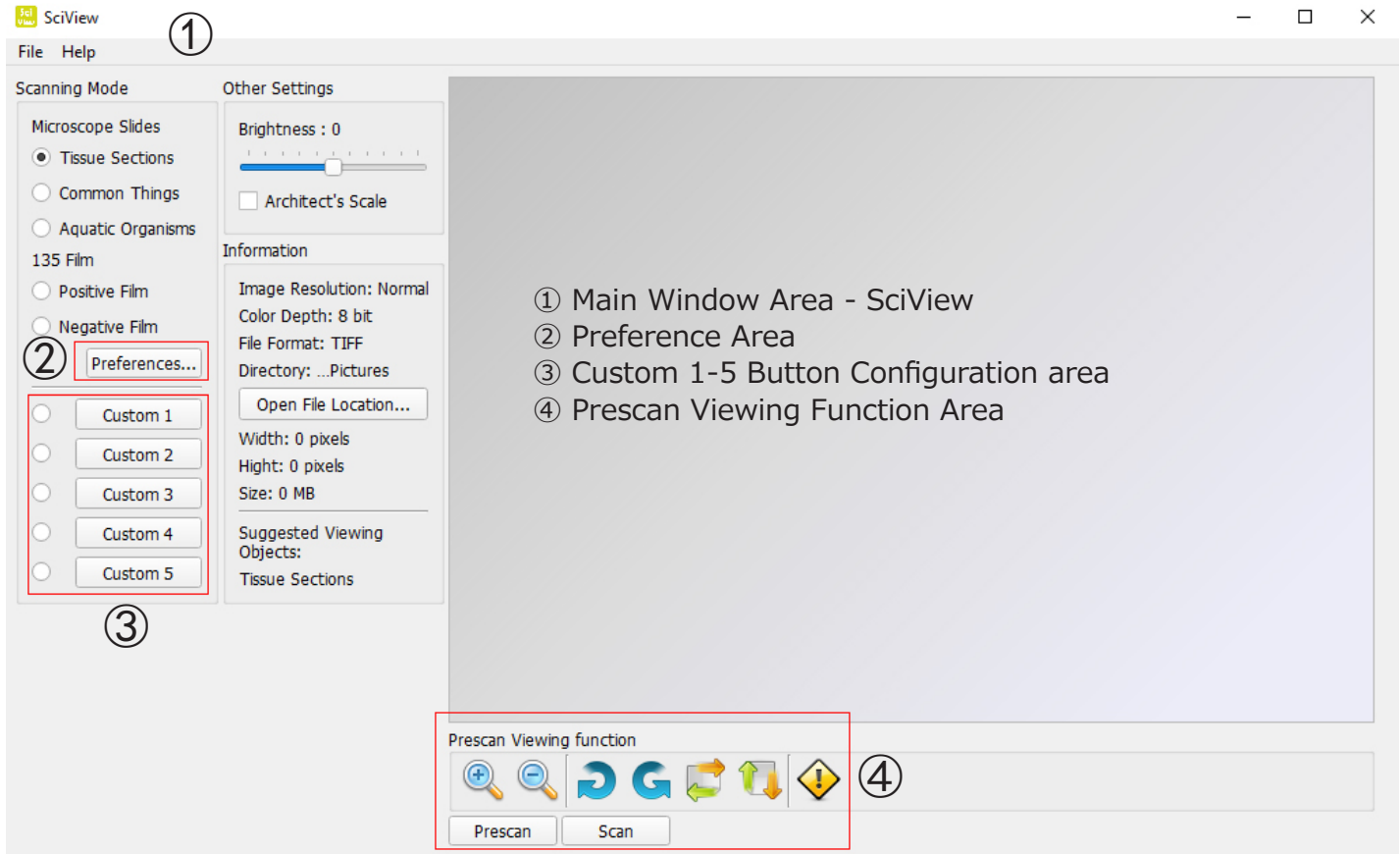

① Main Window Area:

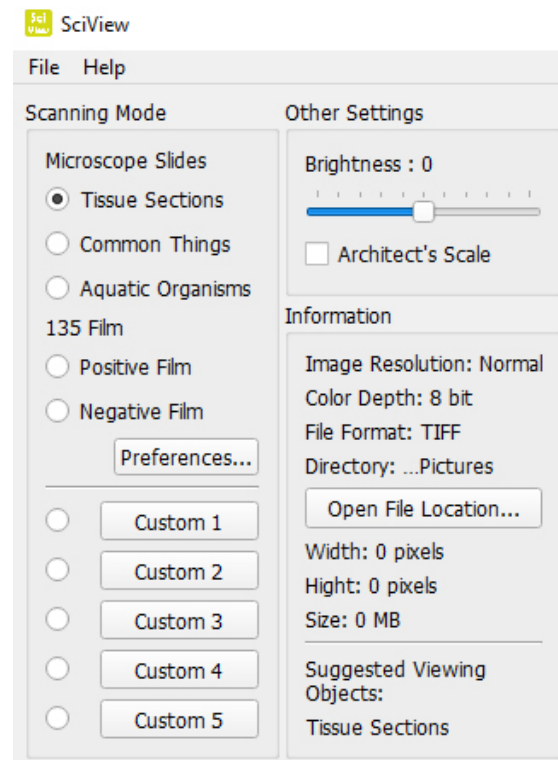
Scanning Mode

Microscope Slides:



- Tissue Sections
- Common Things
- Aquatic Organisms
- Other

File Menu:
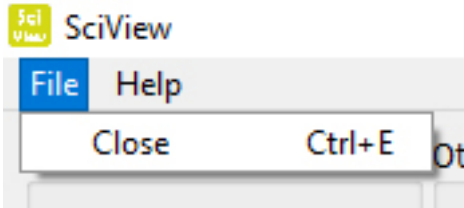


User Manual - Opens this SciView manual (PDF)
Firmware Update - Update the scanner hardware
About – Software and hardware information

Help Menu: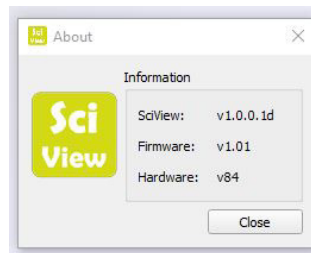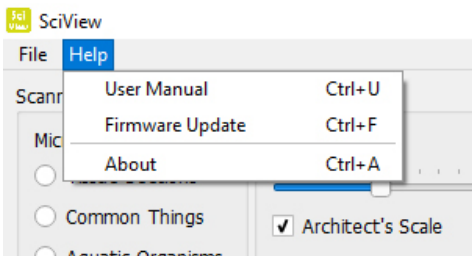

② Preference Section:

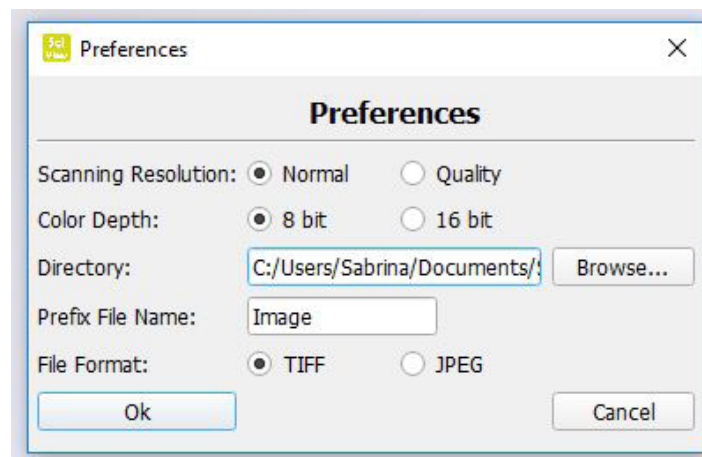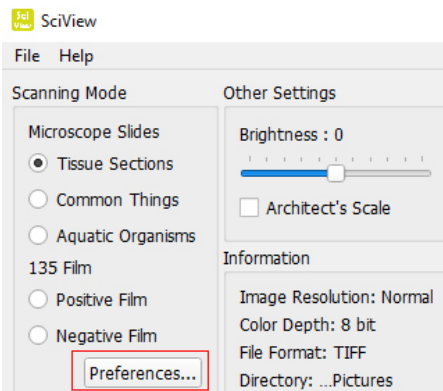Main Preference setting area

Scanning Resolution: Quality and Normal

Color Depth: 8 or 16 bit color options – creating 24 or 48 bit color final images

Directory: Choose the location by browsing to a folder of your choice


Prefix File Name: Select the default name of the files

File Format: TIFF or JPEG – TIFF is an uncompressed larger file format, JPEG is for use in slide shows and presentations.



③ Custom Settings for buttons 1-5:

These buttons are used to save custom scans settings for future use. When you prescan an image configure the settings the way you like then Export these settings to a file, you can then use those same settings by selecting the Import command during future sessions. There are separate sets of functions for both Microscopy and 135mm film formats.

Microscopy Format (Set up explained in STEP 4a)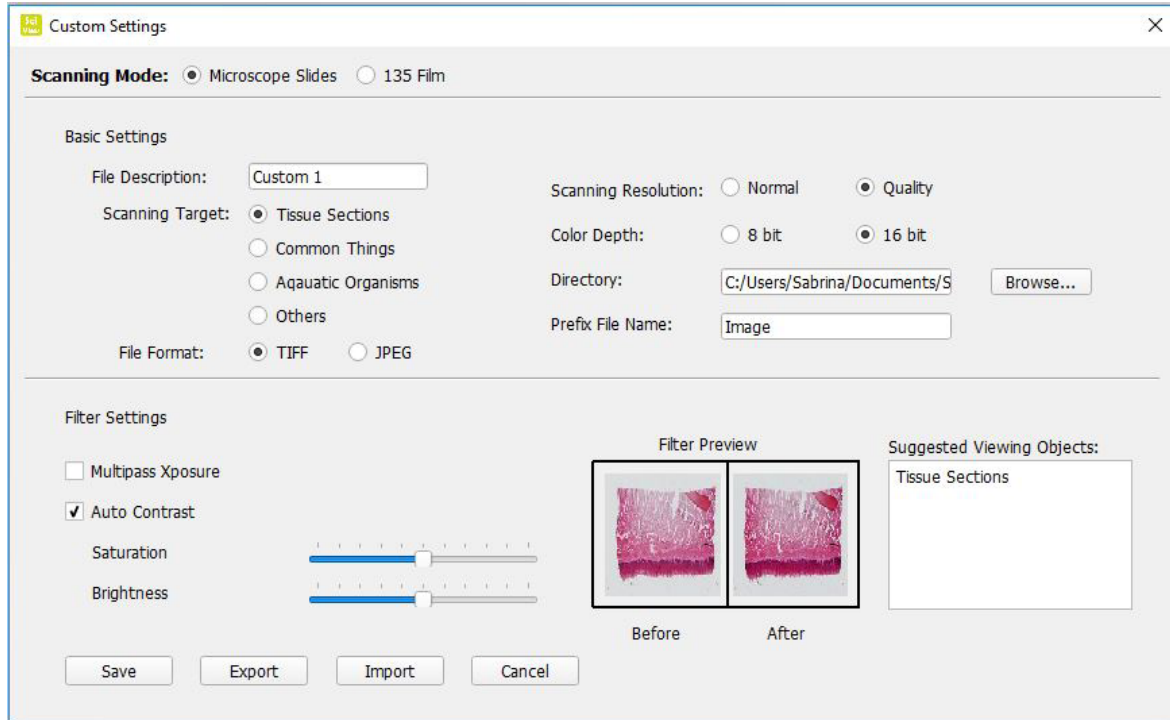


④ Prescan Viewing Function:

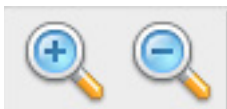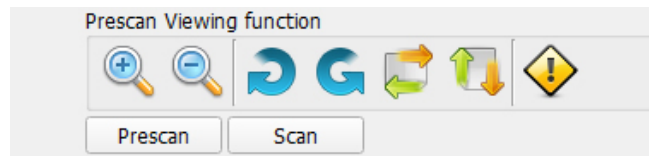Zooms in and out on Prescanned image

Rotates Image Left or Right

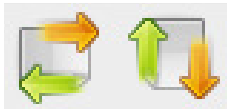Flips Image Horizontally/Vertically

Clears the Prescanned image

Prescan - Previews the slide or film in the scanner

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OPTIKA® S.r.l.